

Quantitative Determination of Urinary COOH-THC for Forensic Toxicology: Comparison to a Validated Reference Method

Xin Zhang,¹ Jonathan Danaceau,¹ Lawrence Andrade,² Kelli Demers,² and Erin Chambers¹

¹Waters Corporation, Milford, MA; ²Dominion Diagnostics, North Kingstown, RI

APPLICATION BENEFITS

- Rapid extraction of COOH-THC from urine
- Elimination of evaporation and reconstitution steps, minimizing the risk of analyte loss from non-specific binding
- Direct sample concentration on the SPE device
- Excellent recovery and minimal matrix effects
- Excellent correlation with a fully validated quantification method

WATERS SOLUTIONS

[Oasis® PRiME HLB μElution™ Plate](#)

[ACQUITY UPLC® BEH C₁₈ Column](#)

[ACQUITY UPLC I-Class System](#)

[Xevo® TQ-S Mass Spectrometer](#)

KEYWORDS

COOH-THC, SPE, Oasis PRiME HLB, cannabis, tetrahydrocannabinol, forensic toxicology

INTRODUCTION

Cannabis continues to be the most widely abused recreational drug in the United States. In addition, the growing number of states legalizing cannabis for medical and/or recreational purposes means that there is a growing need for analytical methods for the quantification of THC and its metabolites. We have recently developed a simple, rapid, and clean SPE extraction method for these analytes using Oasis PRiME HLB.¹ While this method demonstrated excellent accuracy over a wide calibration range, a side by side comparison with a fully validated method from an external laboratory is a key component of method validation. This application note details an inter laboratory comparison of COOH-THC quantification using Oasis PRiME HLB.

MATERIALS AND METHODS

Authentic urine samples were obtained from an outside lab that conducts toxicology tests. Samples were prepared as follows: 25 μL of internal standard (100 ng/mL) was added to 350 μL of urine. 320 μL of 50% KOH was added and the samples were incubated for 2 hours at 65 °C to hydrolyze the COOH-THC glucuronide. 75 μL of the pretreated sample was added directly to the wells of an Oasis PRiME HLB μElution Plate ([p/n 186008052](#)). All samples were washed with 2 x 300 μL of 25:75 MeOH:H₂O and eluted with 2 x 25 μL of 60:40 ACN:IPA. The samples were then diluted with 100 μL of 40% ACN in water. 5 μL was injected onto the UPLC-MS/MS system. The SPE procedure is shown in Figure 1.

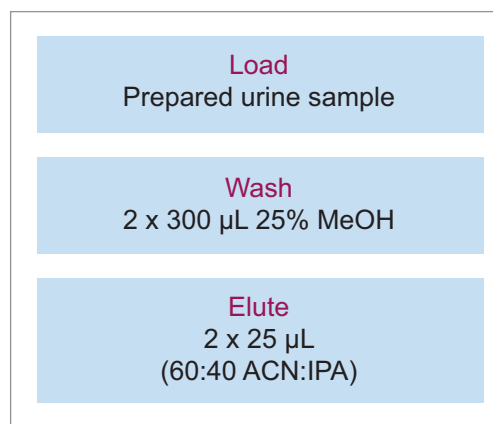


Figure 1. SPE procedure for the extraction of COOH-THC from urine samples using Oasis PRiME HLB μElution Plates.

Mobile phases A and B consisted of 0.1% formic acid in water (MPA) and ACN (MPB). The LC gradient is shown in Table 1. Separation was achieved on an ACQUITY UPLC BEH C₁₈, 1.7 µm, 2.1 x 50 mm Column ([p/n 186002350](#)) at a temperature of 40 °C. Detection was achieved with a Xevo TQ-S. MRM parameters are listed in Table 2.

RESULTS AND DISCUSSION

25 authentic urine samples were analyzed and compared to the validated comparison method. The samples ranged in concentration from 6.70–458 ng/mL, covering nearly the entire linear range of the reference method (5.00–500 ng/mL). A Deming regression (Figure 2) had a slope of 0.995 demonstrating parallelism between the two methods. The correlation (R) of 0.998 indicated an excellent correlation between the results obtained by the two laboratories. The Bland-Altman plot for the Oasis PRiME HLB method for the analysis of COOH-THC in authentic urine samples vs. reference method indicated that the two sets of data are similar at 95% limits of agreement (Figure 3). Table 3 details the results obtained by the two methods. 78% of the sample results are within 20% of each other, exceeding the FDA-GLP specification of 67% for incurred sample reanalysis.² Most results showed a slight negative bias not seen in the standards or QCs. Since the standards and QC samples were prepared in surrogate matrix (Surine), it is possible that the combination of different SPE methods and different chromatographic conditions differentially remove or chromatographically resolve an endogenous substance from the urine samples causing slight signal suppression during ionization. Despite the fact that the samples were subject to different extraction procedures as well as different LC-MS/MS conditions, the results show excellent agreement and indicate that the simplified SPE methodology, which eliminates conditioning and equilibration, gives equivalent results for authentic urine samples.

Time (min)	Flow (mL/min)	%A	%B
0	0.6	50	50
1.0	0.6	50	50
3.0	0.6	5	95
3.5	0.6	5	95
3.6	0.6	50	50
4.0	0.6	50	50

Table 1. Mobile phase gradient. The compositions of MPA and MPB are listed in the Materials and Methods section.

Analyte	RT (min)	MRM transitions (m/z)	Cone voltage (V)	Collision energy (eV)
THC-COOH	1.92	345.3>327.3	50	20
		345.3>299.3	50	25
THC-COOH-d3	1.92	348.3>330.3	50	20

Table 2. Mass spectral parameters for all analytes and internal standards.

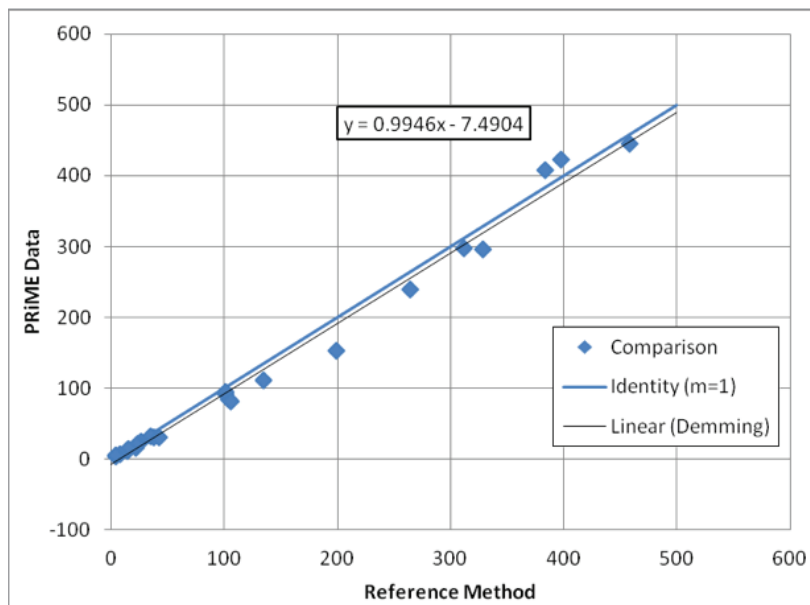


Figure 2. Correlation between the reference method and the Oasis PRiME HLB method for the analysis of COOH-THC in authentic urine samples. The blue line indicates the theoretical identity line of a perfect correlation. The black line is the plotted Deming regression with the equation listed on the chart.

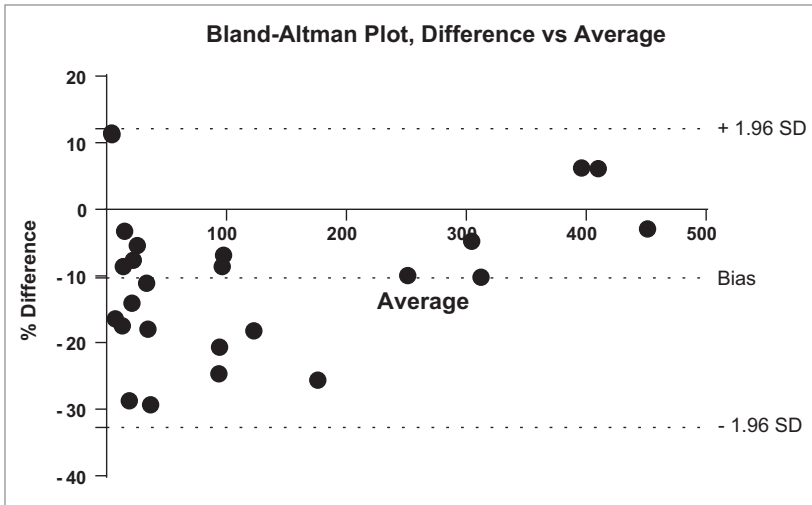


Figure 3. Bland-Altman plot at 95% limits of agreement for the Oasis PRiME HLB method for the analysis of COOH-THC in authentic urine samples vs. reference method.

Sample number	Urine THC concentration (ng/mL) original Oasis MAX reference method	Urine THC concentration (ng/mL) PRiME HLB method	Mean concentration (ng/mL)	%BIAS
1	BQL*	BQL	N/A	N/A
2	BQL	BQL	N/A	N/A
3	7.90	6.70	7.30	-16%
4	14.3	12.0	13.2	-17%
5	14.6	13.4	14.0	-9%
6	15.5	15.0	15.3	-3%
7	21.9	16.4	19.2	-29%
8	22.8	19.8	21.3	-14%
9	23.1	21.4	22.3	-8%
10	26.5	25.1	25.8	-5%
11	35.2	31.5	33.4	-11%
12	37.6	31.4	34.5	-18%
13	42.2	31.4	36.8	-29%
14	101	92.4	96.6	-9%
15	101	94.4	97.8	-7%
16	104	84.5	94.3	-21%
17	105	82.1	93.7	-25%
18	134	112	123	-18%
19	199	154	176	-26%
20	264	239	251	-10%
21	312	297	304	-5%
22	328	297	312	-10%
23	384	409	396	6%
24	398	423	410	6%
25	458	445	451	-3%
			AGREEMENT	78%

Table 3. COOH-THC concentrations from the reference method and the Oasis PRiME HLB method detailed in this application note. *BQL: Below Quantifiable Limit.

CONCLUSIONS

This work demonstrates the applicability of a simplified SPE method using Oasis PRiME HLB μ Elution Plates for the analysis of COOH-THC in authentic urine samples. Correlation was excellent with a validated method from Dominion Diagnostics that used a different SPE protocol and LC-MS/MS analysis procedure. Previous work had shown that Oasis PRiME HLB had excellent recovery with very low variability and could nearly eliminate matrix effects in urine samples. These data demonstrate that the differences between different samples do not affect the quantification of COOH-THC. The use of Oasis PRiME HLB enables a clean extraction without the need for conditioning or equilibration and gives quantitative results equivalent to a fully validated method.

References

1. Zhang, X., J.P. Danaceau, E. Chambers: Quantitative Analysis of THC and Metabolites in Urine with a Simple, Fast, and Clean Oasis PRiME HLB μ Elution Plate. Waters Application Note: [720005556EN](#), 1–5 (2015).
2. Booth, B., M.E. Arnold, B. DeSilva et al.: Workshop Report: Crystal City V – Quantitative Bioanalytical Method Validation and Implementation: The 2013 Revised FDA Guidance. *The AAPS Journal* 17(2), 277–288 (2015).

For forensic toxicology use only.

Waters

THE SCIENCE OF WHAT'S POSSIBLE.®

Waters, The Science of What's Possible, ACQUITY UPLC, Oasis, and Xevo are registered trademarks of Waters Corporation. μ Elution is a trademark of Waters Corporation. All other trademarks are property of their respective owners.

©2016 Waters Corporation. Produced in the U.S.A. December 2016 720005869EN AG-PDF

Waters Corporation

34 Maple Street

Milford, MA 01757 U.S.A.

T: 1 508 478 2000

F: 1 508 872 1990

www.waters.com