

917 Coulometer



Manual

8.917.8004EN / 2022-04-30



Metrohm AG
CH-9100 Herisau
Switzerland
+41 71 353 85 85
info@metrohm.com
www.metrohm.com

917 Coulometer

Program version 5.917.0025

Manual

Technical Communication
Metrohm AG
CH-9100 Herisau

This documentation is protected by copyright. All rights reserved.

This documentation has been prepared with great care. However, errors can never be entirely ruled out. Please send comments regarding possible errors to the address above.

Disclaimer

Deficiencies arising from circumstances that are not the responsibility of Metrohm, such as improper storage or improper use, etc., are expressly excluded from the warranty. Unauthorized modifications to the product (e.g. conversions or attachments) exclude any liability on the part of the manufacturer for resulting damage and its consequences. Instructions and notes in the Metrohm product documentation must be strictly followed. Otherwise, Metrohm's liability is excluded.

Melody for the BEEP command: excerpt from "En Altfrentsche", with kind permission of the Laseyer Quartett, Appenzell.

Table of contents

1	Introduction	1
1.1	Instrument description	1
1.1.1	Titration and measuring modes	1
1.1.2	Connectors	2
1.1.3	Intended use	2
1.2	About the documentation	2
1.2.1	Symbols and conventions	3
2	Safety instructions	4
2.1	General notes on safety	4
2.2	Electrical safety	4
2.3	Tubing and capillary connections	5
2.4	Flammable solvents and chemicals	5
2.5	Recycling and disposal	6
3	Overview of the instrument	7
3.1	Front of the instrument	7
3.2	Rear of the instrument	8
4	Installation	9
4.1	Setting up the instrument	9
4.1.1	Packaging	9
4.1.2	Checks	9
4.1.3	Location	9
4.2	Setting up the titration cell for coulometry	10
4.2.1	Inserting the titration cell	10
4.2.2	Preparing the titration cell	10
4.2.3	Mounting the addition and aspiration tube	14
4.3	Connecting the coulometer to the power supply	15
4.4	Connecting sensors	16
4.4.1	Connecting generator electrodes	16
4.4.2	Connecting indicator electrodes	18
4.4.3	Connecting a temperature sensor	20
4.5	Connecting MSB devices	21
4.5.1	Connecting a dosing device	22
4.5.2	Connecting an additional stirrer or titration stand	23
4.5.3	Connecting a Remote Box	23
4.6	Connecting USB devices	24
4.6.1	General	24

4.6.2	Connecting a USB hub	24
4.6.3	Connecting a printer	24
4.6.4	Connecting a balance	25
4.6.5	Connecting a PC keyboard	26
4.6.6	Connecting a barcode reader	26
4.6.7	Connecting the 885 Compact Oven SC	27
4.7	Connecting the Coulometer to a network	28
5	Titration	29
5.1	Coulometric water content determination according to Karl Fischer (KFC)	29
5.2	Bromine index determination (BRC)	30
6	Operation	31
6.1	Switching the instrument on and off	31
6.2	Fundamentals of operation	32
6.2.1	Touch-sensitive screen	32
6.2.2	Display elements and controls	33
6.2.3	Status display	34
6.2.4	Entering text and numbers	35
7	System settings	37
7.1	General system settings	37
7.1.1	Selecting the dialog language	37
7.1.2	Setting the date, time and local time	38
7.2	System-specific dialog options	39
7.3	User administration	42
7.3.1	Editing the user configuration	44
7.3.2	Creating an identification profile	47
7.3.3	Defining login options	48
7.3.4	Password options	50
7.3.5	Modification options	52
7.3.6	Reasons	52
7.3.7	Audit Trail	53
7.4	Measured value display	55
7.5	Acoustic signals	55
8	Reagents	56
8.1	Editing reagent data	56
8.2	Reagent monitoring	57
9	Sensors	60
9.1	Adding a new sensor	61
9.2	Editing the sensor data	61

9.3	Monitoring the working life	62
10	Device manager	64
10.1	Adding a new device	65
10.2	Configuring the instrument	65
10.3	Coulometer	66
10.3.1	E-mail	67
10.3.2	PC/LIMS report	68
10.3.3	Shared memory	69
10.3.4	TCP/IP settings	71
10.4	Metrohm control instruments	72
10.4.1	Properties – Measuring input	73
10.4.2	Properties – MSB connector	74
10.4.3	Properties – Peripheral devices	75
10.5	885 Compact Oven SC	75
10.5.1	Properties - 885 Compact Oven SC	76
10.5.2	Properties - Tower	77
10.5.3	Properties - Oven module	78
10.6	Printer	79
10.6.1	PDF settings	80
10.6.2	Network printer	81
10.6.3	More options	83
10.7	Balance	83
10.8	USB/RS-232 adapter	85
10.9	PC keyboard	87
10.10	Barcode reader	88
11	File manager	91
11.1	Managing files	91
11.1.1	Copying a file	93
11.1.2	Renaming a file	93
11.1.3	File properties	94
11.2	External storage medium	95
11.3	Creating backups / Restoring data	97
11.3.1	Restoring data	97
12	GLP manager	99
12.1	Automatic system test	100
12.2	Test tools	100
12.3	GLP tests for measurement and titration	101
12.3.1	Parameter description	102



12.4 System validation	105
12.4.1 Parameter description	106
12.5 System monitoring	109
12.5.1 Service interval	109
12.5.2 Backup interval	109
13 Common variables	111
13.1 Editing common variables	112
13.2 Properties of common variables	113
13.3 Monitoring validity	114
13.4 Assigning a result automatically to a common variable	115
14 Templates	117
14.1 Sample data	117
14.1.1 Sample identification list	118
14.1.2 Sample assignment table	119
14.2 Custom result templates	121
14.2.1 Editing result templates	122
14.3 Input lines	125
14.3.1 Editing the input signal	126
14.4 Output lines	127
14.4.1 Editing the output signal	129
14.5 Report header	130
15 Methods	132
15.1 Creating a new method	132
15.2 Saving a method	133
15.3 Loading a method	134
15.4 Editing a method	135
15.4.1 Inserting a command	136
15.5 Method options	137
15.5.1 Start options	138
15.5.2 Stop options	140
15.5.3 Sample data	141
15.5.4 Method properties	144
15.5.5 Note	146
15.5.6 Saving a determination automatically	146
16 Electronic signatures	149
16.1 Signing methods/determinations electronically	149
16.2 Deleting electronic signatures	150

17 Control	151
18 Favorites	155
18.1 Creating favorites	156
18.1.1 Editing favorites	156
19 Sample data	159
19.1 Entering sample data in the main dialog	159
19.2 Requesting sample data at the start of the determination	160
20 Sample table	162
20.1 General	162
20.2 Editing sample data	165
20.3 Properties	167
21 Determination run	170
21.1 Carrying out a single KFC determination	170
21.2 Performing a sample series	171
21.3 Canceling determinations manually	172
22 Live modifications	173
22.1 Editing the sample data of the running determination	173
22.2 Editing the sample table while a determination is running	174
22.3 Live display	175
22.4 Live parameters	176
23 Results and more determination data	178
23.1 More determination data	179
23.1.1 Details	180
23.2 Messages	182
23.3 Local common variables	182
23.4 Determination properties	183
23.5 Loading a determination	185
23.5.1 Determination list	186
23.6 Saving a determination	190
23.7 Curves	191
23.8 Recalculation and reevaluation	193

28.2.2	Titration parameters	241
28.2.3	Stop conditions	242
28.2.4	Conditioning	242
28.2.5	Cell	244
28.2.6	Sensor	245
28.2.7	Stirrer	246
28.3	Measurements (MEAS)	247
28.3.1	Measuring parameters	247
28.3.2	Sensor	249
28.3.3	Stirrer	251
28.4	Evaluations (EVAL)	252
28.4.1	Fixed endpoint evaluation (EVAL FIX-EP)	253
28.4.2	Minimum and maximum evaluation (EVAL MIN/MAX)	254
28.4.3	Rate evaluation (EVAL RATE)	255
28.5	Calculations	256
28.5.1	Calculations (CALC)	256
28.5.2	Calculations (CALC LIVE)	262
28.5.3	Formula editor	264
28.6	Reports (REPORT)	271
28.6.1	General report options	271
28.6.2	Settings of the individual reports	272
28.6.3	List of reports	273
28.7	Dosing and Liquid Handling	276
28.7.1	Preparing an exchange or dosing unit (PREP)	276
28.7.2	Emptying a dosing unit (EMPTY)	277
28.7.3	Dosing a specified volume (ADD)	278
28.7.4	Liquid Handling (LQH)	282
28.8	Communication	285
28.8.1	Scanning remote lines (SCAN)	285
28.8.2	Setting remote lines (CTRL)	287
28.8.3	Scanning the RS-232 interface (SCAN RS)	288
28.8.4	Defining RS-232 commands (CONTROL RS)	289
28.9	Automation	289
28.9.1	Rotating sample rack (MOVE)	289
28.9.2	Moving the lift (LIFT)	291
28.9.3	Resetting the rack (RACK)	292
28.9.4	Defining the sample variable (SAMPLE)	292
28.9.5	Creating a subsequence (SUBSEQ)	293
28.9.6	Controlling the heating (HEATER)	296
28.9.7	Controlling the gas flow (FLOW)	296
28.9.8	Starting the measurement for temperature and gas flow (MEAS T/F ON)	297
28.9.9	Exiting the measurement for temperature and gas flow (MEAS T/F OFF)	298
28.10	Miscellaneous commands	298
28.10.1	Controlling a stirrer (STIR)	298

28.10.2	Pausing the method run (WAIT)	299
28.10.3	Scan data (REQUEST)	299
28.10.4	Defining an acoustic signal (BEEP)	300
28.10.5	Signing a determination (SIGN)	301
28.10.6	Canceling the method run (END)	301
29	Operation and maintenance	302
29.1	System initialization	302
29.2	Maintenance	303
30	Troubleshooting	304
30.1	Editing methods	304
30.2	Sample table	304
30.3	Results/Statistics	305
30.4	Printing	305
30.5	Manual control	306
30.6	Titration stand with pump	306
30.7	Coulometric Karl Fischer titration	306
31	Appendix	310
31.1	Dosing unit	310
31.1.1	Maximum dosing and filling rate	310
31.1.2	Default parameters for preparing (PREP) and emptying (EMPTY)	310
31.2	Exchange unit	311
31.2.1	Maximum dosing and filling rate	311
31.2.2	Default parameters for preparing (PREP)	311
31.3	Stirring rate	311
31.4	Balance	312
31.5	Result variables as parameter setting	313
31.6	Using AuditTrailViewer	314
31.6.1	Installing AuditTrailViewer	314
31.6.2	Opening the Audit Trail	315
31.6.3	Contents of the Audit Trail	316
31.6.4	Filtering the Audit Trail	317
31.6.5	Exporting the Audit Trail	318
31.7	Diagnosis	318
31.7.1	LCD test	319
31.7.2	Formatting an external storage medium	319
31.7.3	Removing an external storage medium	320
31.7.4	Adjusting the touch screen	320
31.7.5	Testing the touch screen	321

31.7.6	Software update (loading program versions and language files)	322
31.7.7	Service	325
32	Technical specifications	326
32.1	Touch screen	326
32.2	Measuring inputs	327
32.2.1	Generator electrode	327
32.2.2	Indicator electrode	327
32.2.3	Temperature sensor	327
32.2.4	Determination range	328
32.2.5	Resolution	328
32.2.6	Reproducibility	328
32.3	Internal stirrer	328
32.4	Pump	328
32.5	Interfaces and connectors	329
32.6	Power supply	329
32.7	Declarations of Conformity	329
32.8	Ambient temperature	330
32.9	Reference conditions	330
32.10	Dimensions	330
32.11	Storage capacity	331
33	Accessories	332
	Index	333



Table of figures

Figure 1	Front	7
Figure 2	Rear	8
Figure 3	Inserting the titration cell	10
Figure 4	Filling the adsorber tube	11
Figure 5	Equipping the titration cell	12
Figure 6	Mounting the addition and aspiration tube	14
Figure 7	Connecting the power supply unit	15
Figure 8	Unscrewing the protective cap from the generator electrode	17
Figure 9	Screwing the electrode cable to the generator electrode	17
Figure 10	Connecting a generator electrode	18
Figure 11	Unscrewing the protective cap from the indicator electrode	18
Figure 12	Screwing the electrode cable to the indicator electrode	19
Figure 13	Connecting an indicator electrode	19
Figure 14	MSB connections	21
Figure 15	MSB connector	22
Figure 16	Directory structure on the external storage medium	96
Figure 17	Evaluation of minimum and maximum	254
Figure 18	Rotational speed depending on the stirring rate	312

1 Introduction

1.1 Instrument description

The 917 Coulometer is a compact titration system for coulometric Karl Fischer titration. This titrator combines in a single device the touch-sensitive color monitor for convenient and efficient operation, the magnetic stirrer and the membrane pump for adding and aspirating solvents. The upper side of the housing offers space for the titration cell. Thanks to its compact construction, you can use the 917 Coulometer in a small space as a stand-alone titrator.

You manage reagents, sensors, methods, etc. conveniently in the internal memory of the coulometer. You can also save your files externally through the USB connector, e.g. on a USB flash drive. On this storage medium you can not only store your methods and determinations, but also create a backup together with all of the data and settings of your system.

The integrated Ethernet connector is available to you should you wish to connect your 917 Coulometer to a network. The network connection offers you the following advantages:

- Saving data to a computer within the network
- Printing reports on a network printer
- Sending displayed messages as e-mails

1.1.1 Titration and measuring modes

The 917 Coulometer supports the following titration and measuring modes.

- **KFC**
Coulometric water content determination according to Karl Fischer with voltametric measurement.
- **BRC**
Coulometric bromine index determination. Determining the amount of double bonds in e.g. mineral oils.
- **MEAS**
Measuring modes:
 - **Ipol** (voltametric measurement with selectable polarization current)
 - **Upol** (amperometric measurement with selectable polarization voltage)
 - **T** (temperature measurement)



1.1.2 Connectors

The 917 Coulometer is equipped with the following connectors:

- **Power connection**
For connecting to the power grid using the power supply unit provided.
- **Two MSB connectors (Metrohm Serial Bus)**
For connecting dosing devices, one additional stirrer or a Remote Box.
- **USB connector**
For connecting peripheral devices (printer, PC keyboard, etc.), a USB flash drive, a Sample Processor or a USB hub.
- **Sensor connectors**
One connector each for:
 - Generator electrode
 - Indicator electrode
 - Temperature sensor (Pt1000 or NTC)
- **Ethernet connector**
For connecting the Coulometer to a network.
- **Two connection nipples**
For connecting tubing for replacing reagent using the built-in pump.

1.1.3 Intended use

The 917 Coulometer is designed for usage as a titrator in analytical laboratories. Its main application field is coulometric Karl Fischer titration.

This instrument is suitable for processing chemicals and flammable samples. Therefore, the use of the instrument requires the user to have basic knowledge and experience in handling toxic and caustic substances. Knowledge regarding the application of fire prevention measures prescribed for laboratories is also mandatory.

1.2 About the documentation










CAUTION

Please read through this documentation carefully before putting the instrument into operation. The documentation contains information and warnings which the user must follow in order to ensure safe operation of the instrument.

1.2.1 Symbols and conventions

The following symbols and formatting may appear in this documentation:

(5-12)	Cross-reference to figure legend The first number refers to the figure number, the second to the instrument part in the figure.
1	Instruction step Perform the steps one after the other.
Method	Dialog text, parameter in the software
File ► New	Menu or menu item
[Continue]	Button or key
	WARNING This symbol draws attention to a possible life-threatening hazard or risk of injury.
	WARNING This symbol draws attention to a possible hazard due to electrical current.
	WARNING This symbol draws attention to a possible hazard due to heat or hot instrument parts.
	WARNING This symbol draws attention to a possible biological hazard.
	WARNING Warning of optical radiation
	CAUTION This symbol draws attention to possible damage to instruments or instrument parts.
	NOTICE This symbol highlights additional information and tips.



2 Safety instructions

2.1 General notes on safety



WARNING

Operate this instrument only according to the information contained in this documentation.

This instrument left the factory in a flawless state in terms of technical safety. To maintain this state and ensure non-hazardous operation of the instrument, the following instructions must be observed carefully.

2.2 Electrical safety

The electrical safety when working with the instrument is ensured as part of the international standard IEC 61010.



WARNING

Only personnel qualified by Metrohm are authorized to carry out service work on electronic components.



WARNING

Never open the housing of the instrument. The instrument could be damaged by this. There is also a risk of serious injury if live components are touched.

There are no parts inside the housing which can be serviced or replaced by the user.

Supply voltage



WARNING

An incorrect supply voltage can damage the instrument.

Only operate this instrument with a supply voltage specified for it (see rear panel of the instrument).

Protection against electrostatic charges



WARNING

Electronic components are sensitive to electrostatic charges and can be destroyed by discharges.

Do not fail to pull the power cord out of the power socket before you set up or disconnect electrical plug connections at the rear of the instrument.

2.3 Tubing and capillary connections



CAUTION

Leaks in tubing and capillary connections are a safety risk. Tighten all connections well by hand. Avoid applying excessive force to tubing connections. Damaged tubing ends lead to leakage. Appropriate tools can be used to loosen connections.

Check the connections regularly for leakage. If the instrument is used mainly in unattended operation, then weekly inspections are mandatory.

2.4 Flammable solvents and chemicals




WARNING

All relevant safety measures are to be observed when working with flammable solvents and chemicals.

- Set up the instrument in a well-ventilated location (e.g. fume cupboard).
- Keep all sources of flame far from the workplace.
- Clean up spilled liquids and solids immediately.
- Follow the safety instructions of the chemical manufacturer.



2.5 Recycling and disposal

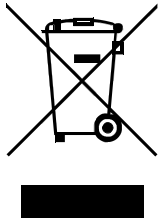

This product is covered by European Directive 2012/19/EU, WEEE – Waste Electrical and Electronic Equipment.

The correct disposal of your old instrument will help to prevent negative effects on the environment and public health.

More details about the disposal of your old instrument can be obtained from your local authorities, from waste disposal companies or from your local dealer.

3 Overview of the instrument

3.1 Front of the instrument

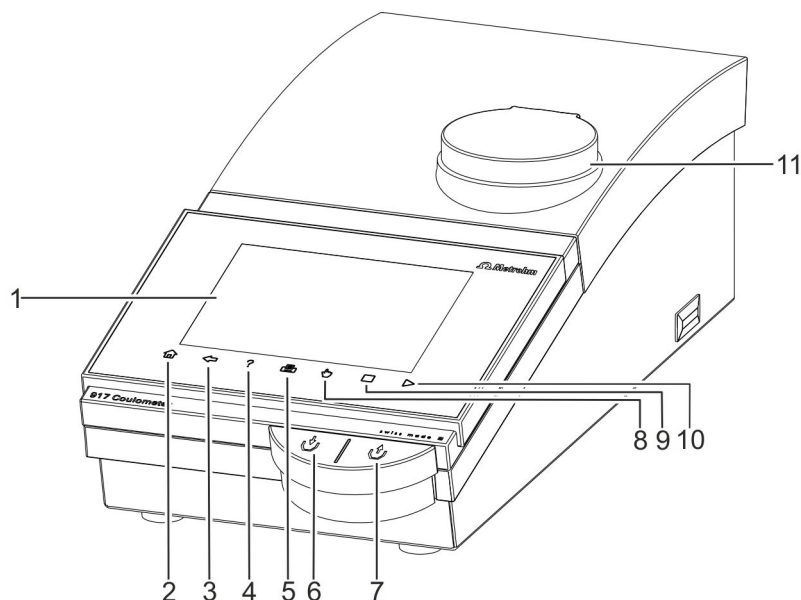


Figure 1 Front

<p>1 Display Touch screen.</p>	<p>2 Fixed key [Home] Opens the main dialog.</p>
<p>3 Fixed key [Back] Saves the entry and opens the next-higher dialog page.</p>	<p>4 Fixed key [Help] Opens the online help for the dialog displayed.</p>
<p>5 Fixed key [Print] Opens the print dialog.</p>	<p>6 Key Pressing the key pumps air into the solvent bottle. The overpressure in the solvent bottle pushes solvent into the titration cell.</p>
<p>7 Key Pressing the key aspirates air out of the aspiration bottle. The vacuum in the aspiration bottle aspirates the liquid out of the titration cell and into the aspiration bottle.</p>	<p>8 Fixed key [Manual] Opens the manual control.</p>
<p>9 Fixed key [STOP] Cancels the running determination.</p>	<p>10 Fixed key [START] Starts a determination.</p>
<p>11 Magnetic stirrer For placing the titration vessel holder.</p>	

3.2 Rear of the instrument

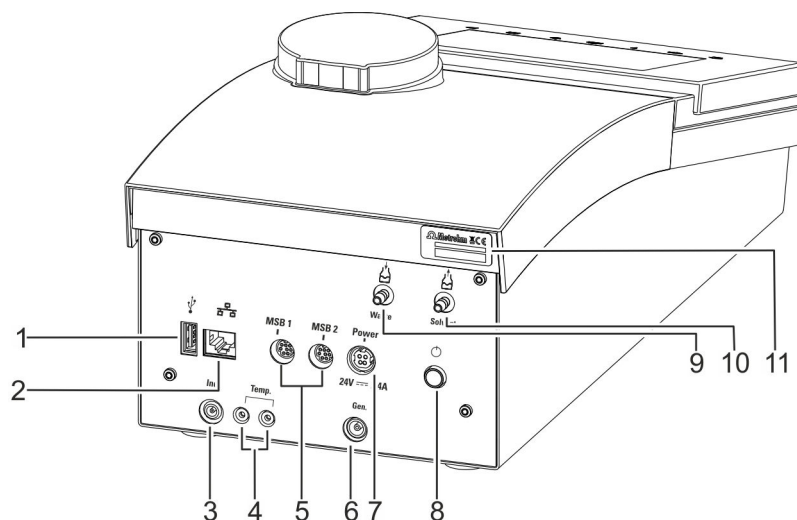


Figure 2 Rear

1 USB connector (type A)

For connecting printers, USB flash drives, USB hubs, 885 Compact Oven SC, etc.

3 Indicator electrode connector (Ind.)

For connecting the indicator electrode.

5 MSB connector (MSB 1 and MSB 2)

Metrohm Serial Bus. For connecting external dosing devices, stirrers or a Remote Box. Mini DIN, 8-pin.

7 Power socket (Power)

For connecting the external power supply unit.

9 Connection nipple for PVC tubing

For aspirating the contents of the titration cell.

11 Type plate

Contains the serial number.

2 Ethernet connector (RJ-45)

For connecting to a network.

4 Temperature sensor connector (Temp.)

For connecting temperature sensors (Pt1000 or NTC). Two B sockets, 2 mm.

6 Generator electrode connector (Gen.)

For connecting the generator electrode.

8 Power switch

Switches the instrument on/off.

10 Connection nipple for PVC tubing

For aspirating solvent.

4 Installation

4.1 Setting up the instrument

4.1.1 Packaging

The instrument is supplied in protective packaging together with the separately packed accessories. Keep this packaging, as only this ensures safe transportation of the instrument.

4.1.2 Checks

Immediately after receipt, check whether the shipment has arrived complete and without damage by comparing it with the delivery note.

4.1.3 Location

The instrument has been developed for operation indoors and may not be used in explosive environments.

Place the instrument in a location of the laboratory which is suitable for operation and free of vibrations and which provides protection against corrosive atmosphere and contamination by chemicals.

The instrument should be protected against excessive temperature fluctuations and direct sunlight.

4.2 Setting up the titration cell for coulometry

4.2.1 Inserting the titration cell

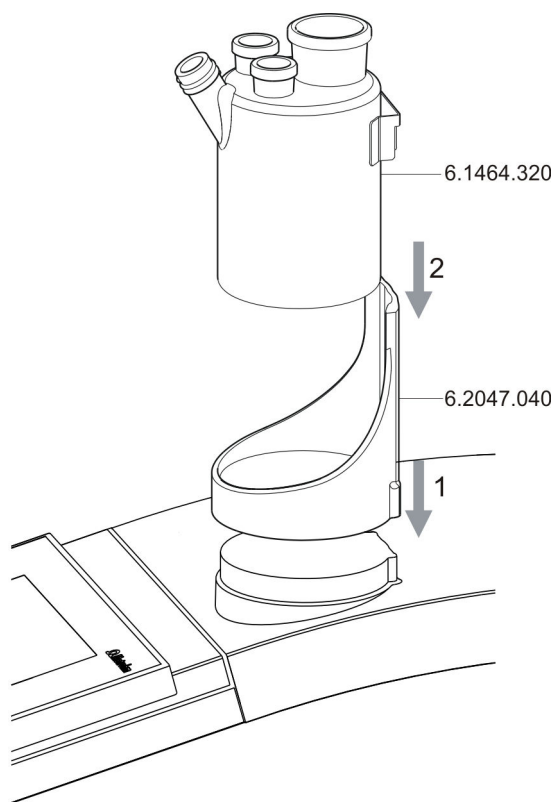


Figure 3 Inserting the titration cell

Set up the titration cell on the magnetic stirrer as follows:

- 1 Attach the 6.2047.040 titration vessel holder to the magnetic stirrer.
- 2 Insert the 6.1464.320 titration cell into the titration vessel holder.

4.2.2 Preparing the titration cell

Filling the adsorber tube

Before setting up the titration cell, the 6.1403.030 adsorber tube has to be filled with 6.2811.000 molecular sieve. Proceed as follows:

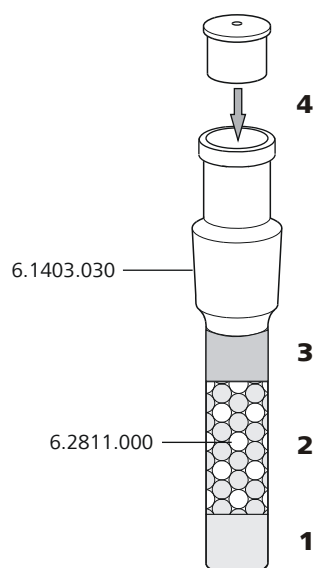


Figure 4 Filling the adsorber tube

- 1** Insert a small cotton plug into the bottom of the adsorber tube. Do not pack the cotton too tightly.
- 2** Fill the molecular sieve up to $\frac{3}{4}$ of the height.
- 3** Place a small cotton plug on the molecular sieve. Do not pack the cotton too tightly.
- 4** Seal the adsorber tube with the appropriate cover.



NOTICE

Note that the molecular sieve must be replaced at regular intervals. Each time you refill the adsorber tube with molecular sieve, you can, for example, write the date directly on the adsorber tube.

Equipping the titration cell

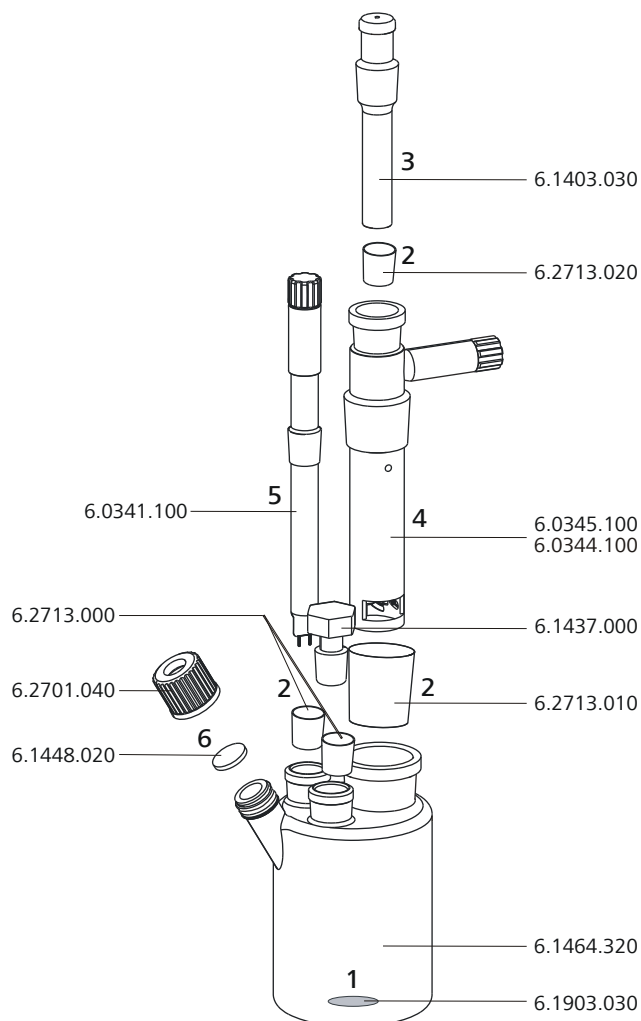


Figure 5 Equipping the titration cell

Equip the titration cell as follows:

- 1** Place the 6.1903.030 stirring bar in the titration cell.
- 2** Cut the 6.2713.0x0 ground-joint sleeves to the correct length and attach them to the ground joints of the inserts (electrodes, adsorber tube, etc.).
Take care to ensure that the edges of the ground-joint sleeves are cut to size cleanly and that there are no fringes. The ground-joint sleeves must not protrude at the lower edge of the ground-joint opening.
- 3** Insert the 6.1403.030 adsorber tube into the generator electrode.

- 4 Insert the 6.0345.100 generator electrode without diaphragm or the 6.0344.100 generator electrode with diaphragm together with the adsorber tube into the large ground-joint opening at the rear.
- 5 Insert the 6.0341.100 indicator electrode into the left ground-joint opening.
- 6 Place the 6.1448.020 septum on the front opening of the titration cell and screw it shut with the 6.2701.040 screw cap.
Tighten the screw cap only enough so that it seals. The septum must not bend.

Filling the titration cell (generator electrode without diaphragm)

Proceed as follows when using a generator electrode without a diaphragm:

- 1 Fill approximately 100 mL of reagent into the titration cell using the 6.2738.000 funnel.
- 2 Close the remaining ground-joint opening on the right with the 6.1437.000 ground-joint stopper (with ground-joint sleeve attached).

Filling the titration cell (generator electrode with diaphragm)

Proceed as follows when using a generator electrode with a diaphragm:

- 1 Fill approximately 5 mL of catholyte into the generator electrode.
- 2 Fill approximately 100 mL of anolyte into the titration cell using the 6.2738.000 funnel. The level of the anolyte should be roughly 1 - 2 mm above the level of the catholyte.
- 3 Close the remaining ground-joint opening on the right with the 6.1437.000 ground-joint stopper (with ground-joint sleeve attached).



4.2.3 Mounting the addition and aspiration tube

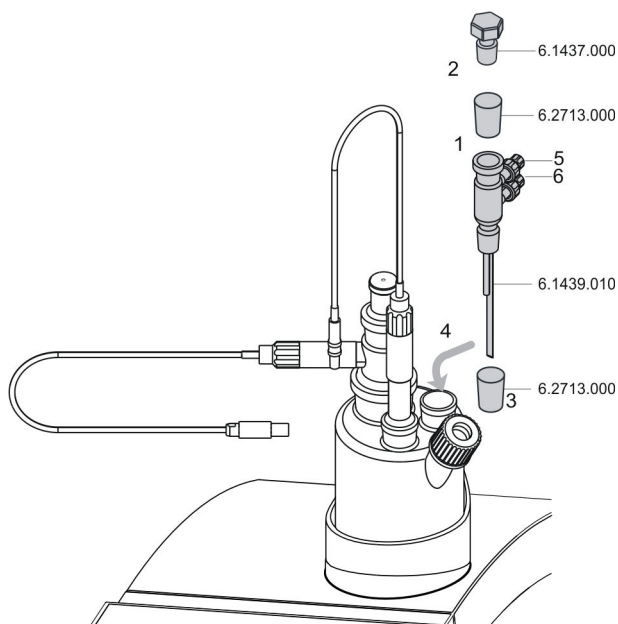


Figure 6 Mounting the addition and aspiration tube

Insert the addition and aspiration tube as follows into the titration cell:

- 1** Attach the 6.2713.000 ground-joint sleeve that has been cut to size to the ground joint of the 6.1437.000 stopper.
- 2** Insert the stopper into the 6.1439.010 addition and aspiration tube.
- 3** Attach the 6.2713.000 ground-joint sleeve that has been cut to size to the ground joint of the addition and aspiration tube.
- 4** Insert this assembly into the ground-joint opening.
- 5** Connect the tubing for the reagent addition at the upper connector of the addition and aspiration tube (5).
- 6** Connect the tubing for the aspiration of the titration cell at the lower connector of the addition and aspiration tube (6).

Details regarding how to connect the addition tubing and the aspiration tubing can be found in the tutorial for the 917 Coulometer.

4.3 Connecting the coulometer to the power supply

Connecting the power supply unit

The 917 Coulometer has an external power supply unit for a 24 V power supply (DC). This is connected to the power socket of the coulometer.



WARNING

An incorrect supply voltage can damage the instrument.

Operate the instrument only with the supply voltage specified for it. Use the supplied power supply unit exclusively, see chapter *Technical specifications*.

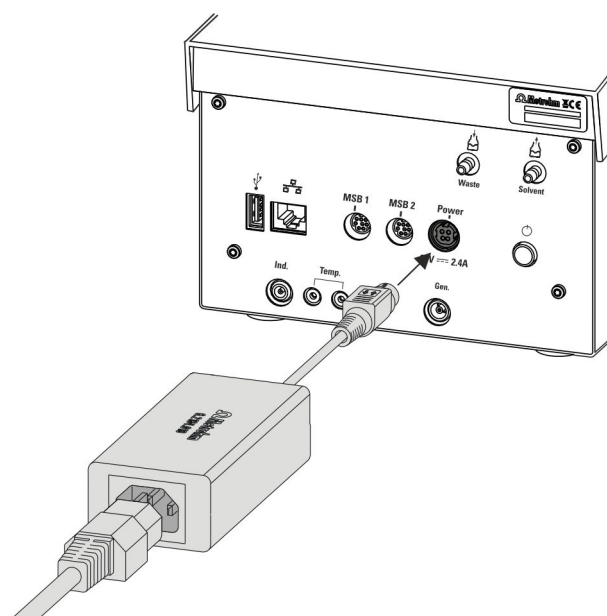


Figure 7 Connecting the power supply unit

Proceed as follows:

- 1 Connect the plug of the external power supply unit with the power socket of the coulometer.

**NOTICE**

The plug of the power supply unit is protected against accidental disconnection of the cable by means of a pull-out protection feature. If you wish to pull out the plug, you first need to pull back the outer plug sleeve marked with arrows.

- 2 Connect the power cord with the external power supply unit of the coulometer and with the power grid.

**CAUTION**

Switch off the Coulometer correctly by pressing the power switch before you disconnect the electricity supply. If this is not done, then there is a danger of data loss.

4.4 Connecting sensors

The measuring interface contains the following measuring inputs:

- **Gen.** for a generator electrode
- **Ind.** for a double Pt electrode
- **Temp.** for a temperature sensor of the Pt1000 or NTC type

**CAUTION**

Under all circumstances, avoid mixing up the electrode cable from the indicator electrode with the one from the generator electrode. Mark the screw heads of the cables accordingly.

4.4.1 Connecting generator electrodes

Screwing the electrode cable to the generator electrode

- 1 Unscrew the protective cap from the generator electrode.

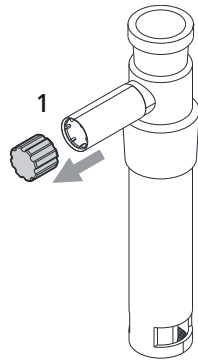


Figure 8 Unscrewing the protective cap from the generator electrode

- 2 Tighten the 6.2104.120 electrode cable to the generator electrode.

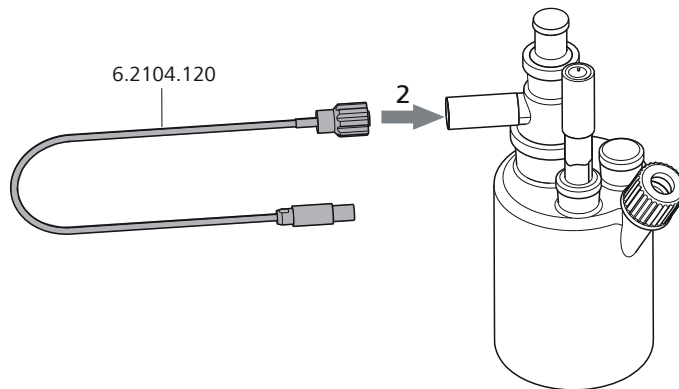


Figure 9 Screwing the electrode cable to the generator electrode

Connecting the electrode cable to the coulometer

- 1 Plug the electrode plug into the **Gen.** socket of the coulometer.

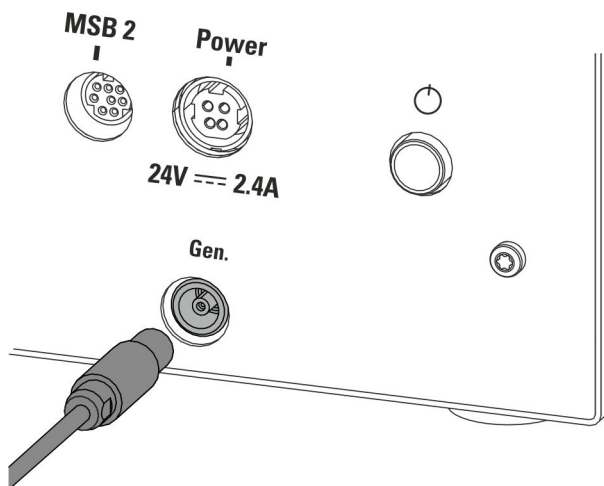


Figure 10 Connecting a generator electrode



NOTICE

The electrode cable is protected against accidental disconnection of the cable by means of a pull-out protection. If you wish to pull out the plug again, you first need to pull back the outer plug sleeve.

4.4.2 Connecting indicator electrodes

Screwing the electrode cable to the indicator electrode

- 1 Unscrew the protective cap from the indicator electrode.

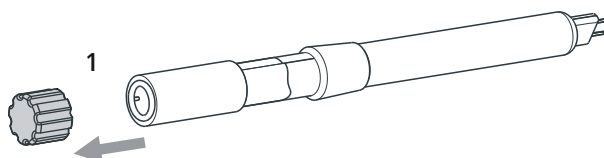


Figure 11 Unscrewing the protective cap from the indicator electrode

- 2 Tighten the 6.2104.020 electrode cable to the indicator electrode.

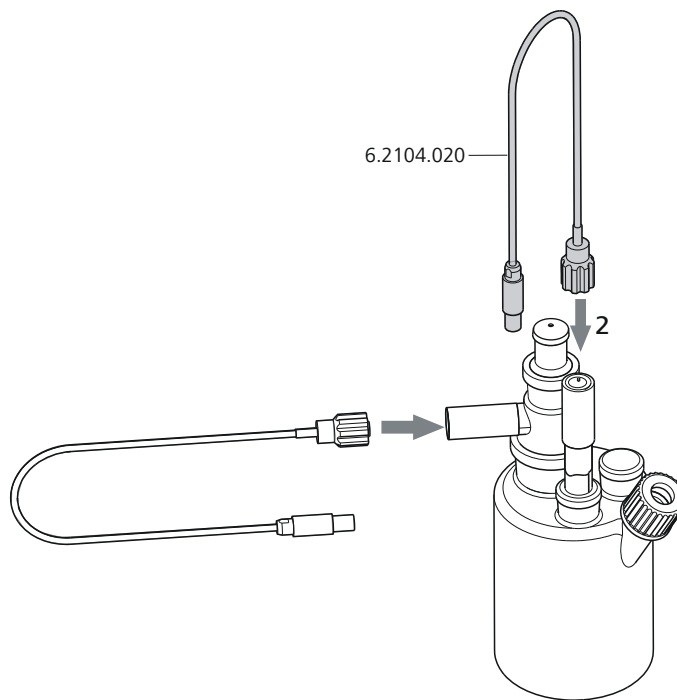


Figure 12 Screwing the electrode cable to the indicator electrode

Connecting the electrode cable to the coulometer

- 1 Plug the electrode plug into the **Ind.** socket of the coulometer.

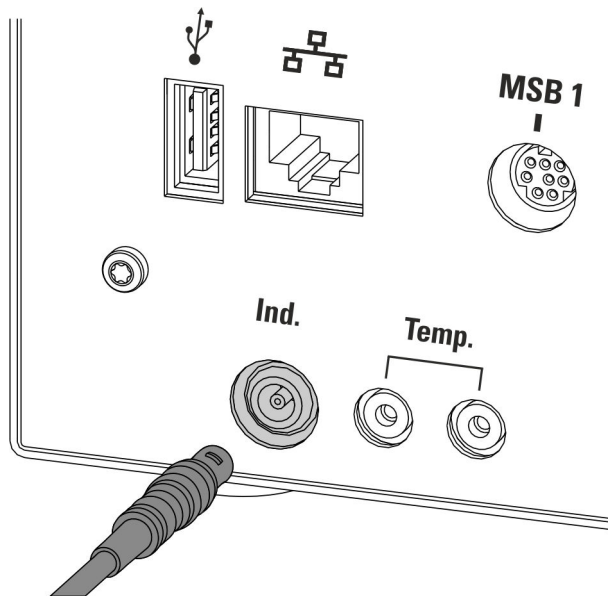


Figure 13 Connecting an indicator electrode

**NOTICE**


The electrode cable is protected against accidental disconnection of the cable by means of a pull-out protection. If you wish to pull out the plug again, you first need to pull back the outer plug sleeve.

4.4.3 Connecting a temperature sensor

A temperature sensor of the Pt1000 or NTC type can be connected to the **Temp.** connector.

Connect the temperature sensor as follows:

- 1 Insert the plugs of the temperature sensor into the **Temp.** sockets of the coulometer.

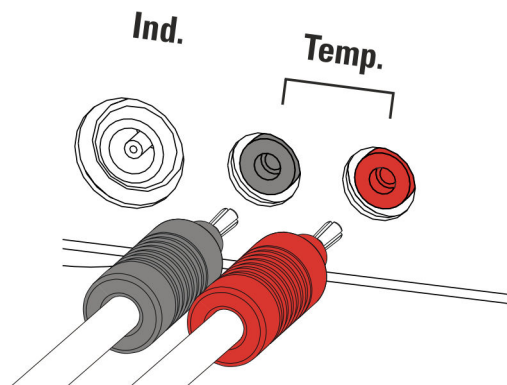**NOTICE**

Always insert the red plug into the red socket. This is the only way that shielding against electrical interference can be ensured.

4.5 Connecting MSB devices

In order to connect MSB devices, e.g. dosing device or Remote Box, the Coulometer has two connectors at what is referred to as the *Metrohm Serial Bus* (MSB). Various peripheral devices can be connected in sequence (daisy chain) at a single MSB connector (8-pin Mini DIN socket) and be controlled simultaneously by the Coulometer. In addition to the connection cable, stirrers and the Remote Box are each equipped with their own MSB socket for this purpose.



NOTICE

The additional stirrer must not be connected to the MSB 1!

The following figure provides an overview of the instruments that can be connected to an MSB socket, along with a number of different cabling variations.

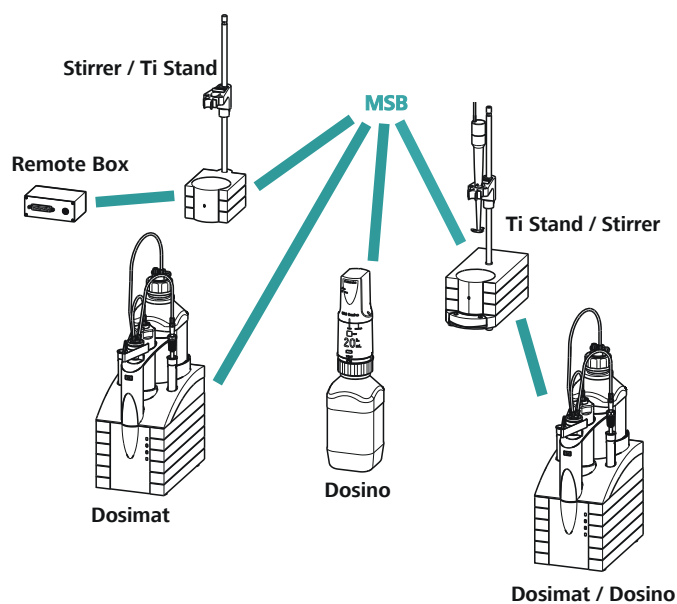


Figure 14 MSB connections



NOTICE

When connecting MSB devices together, the following must be observed:

- No additional stirrer may be connected to the MSB 1!
- Only one device of the same type may be used at a single MSB connector at one time.
- When making the connection, take care to ensure that the flat part of the MSB plug marked with arrows is pointing in the direction of the marking on the MSB connector (see figure 15, page 22).

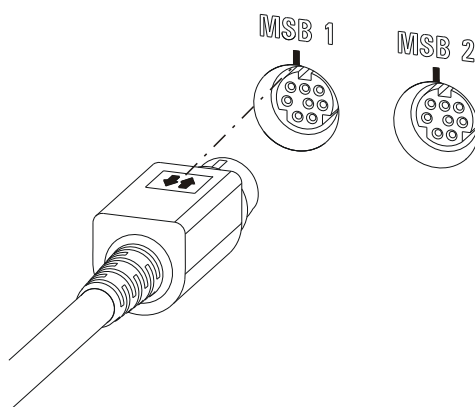


Figure 15 MSB connector



CAUTION

Switch off the Coulometer before you plug in MSB devices. When it is switched on, the Coulometer automatically recognizes which device is connected to which MSB connector. The connected MSB devices are entered automatically in the device manager.

MSB connections can be extended with the 6.2151.010 cable. The maximum connection length permitted is 6 m.

4.5.1 Connecting a dosing device

You can connect two dosing devices to the Coulometer.

The types of dosing devices that are supported are:

- 800 Dosino
- 805 Dosimat

Proceed as follows:

- 1 Switch off the Coulometer.
- 2 Connect the dosing device connection cable to an MSB connector on the rear side of the Coulometer.
- 3 Switch on the Coulometer.

4.5.2 Connecting an additional stirrer or titration stand

Proceed as follows:

- 1 Switch off the Coulometer.
- 2 Connect the connection cable of the magnetic stirrer or of the titration stand to MSB 2 on the rear of the Coulometer.
- 3 Switch on the Coulometer.

4.5.3 Connecting a Remote Box

Instruments that are controlled via remote lines and/or that send control signals via remote lines can be connected via the 6.2148.010 Remote Box. In addition to Metrohm, other instrument manufacturers also use similar connectors that make it possible to connect different instruments together. These interfaces are also frequently given the designations "TTL Logic", "I/O Control" or "Relay Control" and they generally have a signal level of 5 volts.

Control signals are understood to be electrical line statuses or brief (> 200 ms) electrical pulses which display the operational state of an instrument or which trigger or report an event. Sequences on a variety of instruments can thus be coordinated in a single complex automation system. However, no exchange of data is possible.

Proceed as follows:

- 1 Switch off the Coulometer.
- 2 Connect the Remote Box connection cable to an MSB connector on the rear side of the Coulometer.
- 3 Switch on the Coulometer.

You will find precise information concerning the pin assignment of the interface on the Remote Box in the *appendix*.

4.6 Connecting USB devices

4.6.1 General

The 917 Coulometer has a USB connector (type A socket) for peripheral devices with a USB interface and for the 885 Compact Oven SC. If you wish to connect more than one device to the USB, you can use a commercially available USB hub.



NOTICE

We recommend that the Coulometer be switched off while you set up or disconnect connections between the devices.

4.6.2 Connecting a USB hub

Use a USB hub with its own power supply.

Connect the USB hub as follows:

- 1 With the help of the 6.2151.030 cable (length 0.6 m) or the 6.2151.020 cable (length 1.8 m), connect the USB connector of the Coulometer (Type A) with the USB connector of the hub (Type B, see manual for the USB hub).

The USB hub is recognized automatically.

4.6.3 Connecting a printer

Printers connected to the 917 Coulometer must meet the following requirements:

- Printer languages: HP-PCL, HP-PCL-GUI, Canon BJI Commands or Epson ESC P/2
- Paper size: A4 or Letter.

Connect the printer as follows:

- 1 With the aid of the 6.2151.020 cable, connect the USB connector of the Coulometer (type A) with the USB connector of the printer (type B, see manual for the printer).
- 2 Configure the printer in the device manager of the Coulometer (*see chapter 10.6, page 79*).

4.6.4 Connecting a balance

If you want to connect a balance to the Coulometer you need a USB/RS-232 adapter (6.2148.050).

The following table offers an overview of the balances that you can use together with the Coulometer and of which cable you will need for connection to the RS-232 interface:

Balance	Cables
AND ER, FR, FX with RS-232 interface (OP-03)	6.2125.020 + 6.2125.010
Mettler AB, AG, PR (LC-RS9)	In the scope of delivery for the balance
Mettler AM, PM, PE with interface option 016 or Mettler AJ, PJ with interface option 018	6.2146.020 + 6.2125.010 Also from Mettler: ME 47473 adapter and either ME 42500 hand switch or ME 46278 foot switch
Mettler AT	6.2146.020 + 6.2125.010 Also from Mettler: ME 42500 hand switch or ME 46278 foot switch
Mettler AX, MX, UMX, PG, AB-S, PB-S, XP, XS	6.2134.120
Mettler AE with interface option 011 or 012	6.2125.020 + 6.2125.010 Also from Mettler: ME 42500 hand switch or ME 46278 foot switch
Ohaus Voyager, Explorer, Analytical Plus	Cable AS017-09 from Ohaus
Precisa balances with RS-232-C interface	6.2125.080 + 6.2125.010
Sartorius MP8, MC, LA, Genius, Cubis	6.2134.060
Shimadzu BX, BW	6.2125.080 + 6.2125.010



Connect the balance as follows:

- 1** Connect the USB plug of the USB/RS-232 adapter with the USB connector of the Coulometer (Type A).
The USB/RS-232 adapter will be recognized automatically and entered in the device manager of the Coulometer.
- 2** Connect the RS-232 interface of the USB/RS-232 adapter with the RS-232 interface of the balance (see table for cable).
- 3** Switch on the balance.
- 4** If necessary, switch on the RS-232 interface of the balance.
- 5** Configure the RS-232 interface of the USB/RS-232 adapter in the device manager of the Coulometer (*see chapter 10.8, page 85*).
- 6** Enter and configure the balance in the device manager of the Coulometer (*see chapter 10.7, page 83*).
- 7** Make sure that the parameters of the USB/RS-232 adapter configured in the device manager match those of the balance.

4.6.5 Connecting a PC keyboard

The PC keyboard is used as an aid for text and numerical input.

Connect the PC keyboard as follows:

- 1** Connect the USB plug of the keyboard with the USB connector of the Coulometer (Type A).
- 2** Enter and configure the keyboard in the device manager of the Coulometer (*see chapter 10.9, page 87*).

4.6.6 Connecting a barcode reader

The barcode reader is used as an aid for text and numerical input. You can connect a barcode reader with USB interface.

Connect the barcode reader as follows:

- 1** Connect the USB plug of the barcode reader with the USB connector of the Coulometer (Type A).

- 2 Enter and configure the barcode reader in the device manager of the (see chapter 10.10, page 88).

Settings on the barcode reader:

Program the barcode reader as follows (also see manual for the barcode reader):

- 1 Switch the barcode reader to programming mode.
- 2 Specify the desired layout for the keyboard (USA, Germany, France, Spain, German-speaking Switzerland).
This setting must match the setting in the device manager.
- 3 Make sure that the barcode reader is set in such a way that Ctrl characters (ASCII 00 to 31) can be sent.
- 4 Program the barcode reader in such a way that the ASCII character 02 (STX or Ctrl B) is sent as the first character. This first character is normally referred to as the "Preamble" or "Prefix Code".
- 5 Program the barcode reader in such a way that the ASCII character 04 (EOT or Ctrl D) is sent as the last character. This last character is normally referred to as the "Postamble", "Record Suffix" or "Postfix Code".
- 6 Exit the programming mode.

4.6.7 Connecting the 885 Compact Oven SC



NOTICE

In order for the 885 Compact Oven SC to be controlled with the 917 Coulometer, the following firmware versions must be installed on the instruments as a minimum requirement:

- 917 Coulometer 5.917.0025
- 885 Compact Oven SC 5.885.0013

The 885 Compact Oven SC is connected via the USB interface. The 6.2151.110 USB cable is required for this.

If the USB interface on the 917 Coulometer is also needed for other USB devices (USB flash drive, USB keyboard, USB/RS-232 adapter for balance), use a USB hub with its own power supply.



Connect the 885 Compact Oven SC as follows:

- 1** Switch off the 917 Coulometer.
- 2** Connect the 885 Compact Oven SC to the Coulometer using a USB cable.
- 3** Switch on the 885 Compact Oven SC.
- 4** Switch on the 917 Coulometer.



NOTICE

When the 917 Coulometer is started, the 885 Compact Oven SC is added to the device manager and the settings from the device manager are transferred to the instrument.

The switch-on sequence must be strictly observed with the 885 Compact Oven SC.

- 5** Configure the 885 Compact Oven SC in the device manager (*see chapter 10.5, page 75*).

4.7 Connecting the Coulometer to a network

The 917 Coulometer has a network connection (Ethernet). This can be used to integrate your Coulometer in your network. You can, for example, store data on a PC within the network or print reports on a network printer. In *Chapter 10, page 64*, you will find information as to which settings are necessary for the network connection.

Connecting the Coulometer to a network

5 Titrations

5.1 Coulometric water content determination according to Karl Fischer (KFC)

The coulometric Karl Fischer titration is a variation of the classic water content determination method according to Karl Fischer.

The conventional method works with a methanolic solution of iodine, sulfur dioxide and a base as buffer substance. When titrating a sample containing water several reactions take place, which can be summarized by the following equation:



According to the equation above, the I_2 reacts quantitatively with H_2O . This chemical relation is the base for the water content determination.

With the coulometric Karl Fischer titration, the necessary iodine is directly and electrochemically generated in the submitted electrolyte ("electronic buret"). Between the amount of electric charge and the amount of generated iodine, there is a strictly quantitative relationship, which is used for high-precision dosing of the iodine. No titer needs to be determined as the coulometric Karl Fischer method is an absolute determination. It must only be ensured that the reaction generating the iodine runs with a 100% current efficiency. This is the case with all the reagents available at present.

The endpoint indication is effected voltametrically by modulating an alternating current of constant strength to a double Pt electrode. This produces a voltage difference between the Pt wires, which significantly declines when lowest amounts of free iodine are present. This circumstance is used for detecting the endpoint of the titration.

5.2 Bromine index determination (BRC)

The bromine index is a measure for the content of olefins in a sample.

The basis for the bromine index determination is the addition of bromine to the double bond of olefins. A bromine molecule is used for every carbon-carbon double bond.

With the bromine index determination, the necessary bromine is directly and electrochemically generated in the submitted electrolyte ("electronic buret"). Between the amount of electric charge and the amount of generated bromine, there is a strictly quantitative relationship, which is used for high-precision dosing of the bromine. No titer needs to be determined as the bromine index determination is an absolute determination.

The bromine index is usually indicated in mg bromine per 100 g sample.

6 Operation

6.1 Switching the instrument on and off

Switching on the instrument



CAUTION

Peripheral devices (e.g., printers, 885 Compact Oven SC, etc) must be connected and switched on before you switch on the 917 Coulometer.



NOTICE

English is set as the default dialog language when the instrument is switched on for the first time.


Proceed as follows:

- 1 ▪ Press the power switch on the left-hand side of the back panel of the 917 Coulometer.
The 917 Coulometer is initialized. A system test is performed. This process takes some time.



NOTICE

If a buret unit is connected, then a request appears to carry out the **Prepare** function:

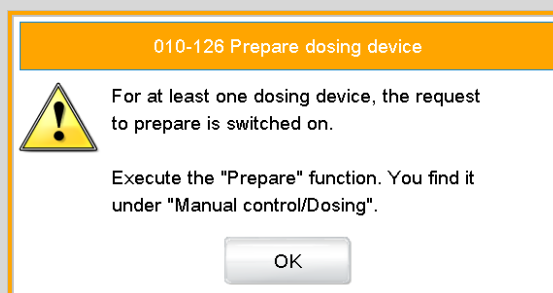

All tubings and the cylinder are rinsed with the **Prepare** function. The preparing of the buret unit is described in *chapter 27.3.3, page 221*.

- Confirm the message with **[OK]**.

The main dialog is displayed:

Switching off the instrument



CAUTION

The 917 Coulometer must be switched off by pressing the power switch on the rear of the instrument before the electricity supply is interrupted. If this is not done, then there is a danger of data loss.


Proceed as follows:

- 1 Press the power switch on the left-hand side of the back panel of the 917 Coulometer.

The current data is saved and the system is shut down. This process takes just a short time. At the same time, all other instruments connected to the 917 Coulometer via a USB cable are also being switched off (except the 885 Compact Oven SC).

6.2 Fundamentals of operation

6.2.1 Touch-sensitive screen

The entire 917 Coulometer user interface is touch-sensitive. Simply touch a few of the buttons on the interface to learn how a touch-sensitive screen reacts. You can always return to the main dialog by touching [].

In order to enable an element on the 917 Coulometer user interface, just touch the screen with your fingertip, the eraser of a pencil or a stylus (special tool for operating instruments with touch-sensitive screens).




CAUTION

Never touch the touch screen with a pointed or sharp object such as a ballpoint pen.





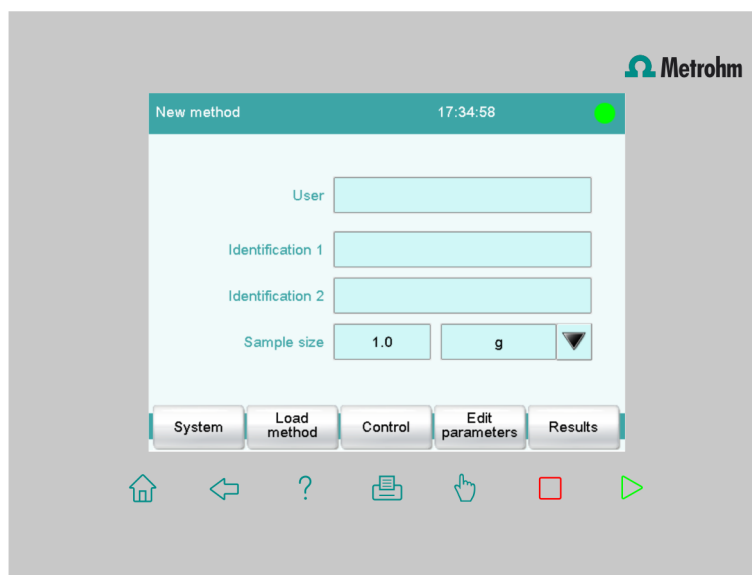


In the default setting, the software is configured in such a way that an acoustic signal will be generated every time an active control is touched. This setting can be deactivated in the system settings (*see chapter 7.5, page 55*).

6.2.2 Display elements and controls



The following display elements and controls are available:



Table 1 Fixed keys which are always available


	[Home] always opens the main dialog.
	[Back] saves the entry and opens the next-higher dialog page.
	[Help] opens the online help for the dialog displayed.
	[Print] opens the printing dialog.
	[Manual] opens the manual control.
	[Stop] cancels the running determination.
	[Start] starts a determination.

The file name of the loaded method, the time and the system status are displayed in the main dialog in the **Title bar**.


In the other dialogs, the title bar shows the headings of the next upper level and of the displayed dialog. This is an aid for orientation during navigation through the user dialog.

Table 2 Screen elements

	Buttons open a new dialog when they are tapped.
	

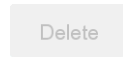

Inactive buttons with gray lettering indicate that the respective function is not available at the moment.

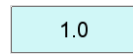

Input fields open an input dialog when tapped.



Tapping on the **selection symbol** opens a selection list.



A **check box** can also be activated or deactivated by tapping on it.

6.2.3 Status display

The current status of the system is displayed in the upper right-hand corner of the title bar.

Table 3 Status displays



The instrument is in normal status.



The working medium is being conditioned.



Conditioning has been paused.



The working medium is conditioned.



A method has been started.



A method has been paused.



An action has been started in manual control.

6.2.4 Entering text and numbers

In the editing dialog for text or numerical input, enter the individual characters by tapping in the input field. The following functions are available:

Text editor

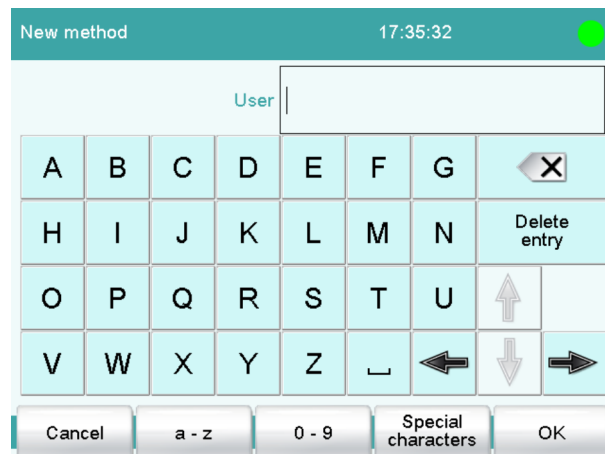


Table 4 Editing functions

Editing function	Description
[OK]	The modification is applied and the editing dialog is exited.
[Cancel]	The editing dialog is exited without applying the modification.
[Delete entry]	The content of the input field is deleted completely.
[□]	The character in front of the cursor is deleted.
[←]	The cursor within the input field is shifted to the left by one character.
[⇒]	The cursor within the input field is shifted to the right by one character.
[a...z]	The lower-case letters are displayed. The label changes to [A...Z]. The upper-case letters are displayed again by tapping.
[0-9]	Numbers and mathematical characters are displayed.
[Special characters]	Special characters are displayed. You can use the button [More] to navigate through all available characters.

Number editor

Edit command / Stop conditions

Stop time s

Input:
1 ... 999999

Default value:
off

7	8	9	off
4	5	6	R1 ▼
1	2	3	
0	+/-	.	

Cancel Delete entry OK

Table 5 Editing functions

Editing function	Description
[OK]	The modification is applied and the editing dialog is exited.
[Cancel]	The editing dialog is exited without applying the modification.
[Delete entry]	The content of the input field is deleted completely.
[off]	If not only numbers but also special values (e.g., off) can be entered, then the corresponding buttons will be shown to the right of the numerical keypad.
[R1]	For many parameters, a result previously defined in the method can also be entered in place of a number (see chapter 31.5, page 313). You can select the result variable by touching [R1] .




NOTICE

A commercially available USB keyboard can be connected to make text and numerical input easier. The key assignment is described in *chapter 10.9, page 87*.

7 System settings

Main dialog: **System ► System settings**

This chapter describes the various system settings and configurations.

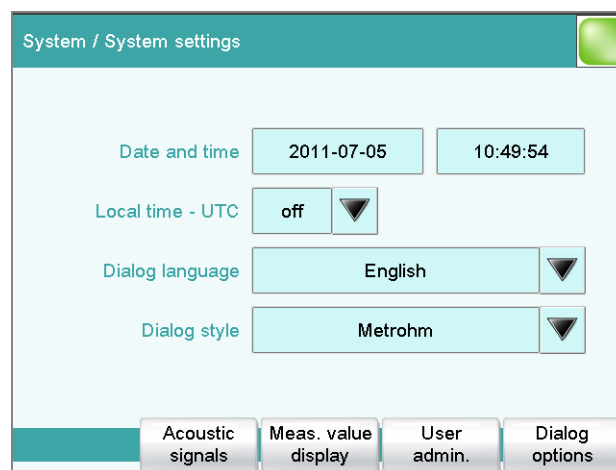

- Selecting the dialog language (see chapter 7.1.1, page 37).
- Setting the date and time (see chapter 7.1.2, page 38).
- Defining system-specific dialog options (see chapter 7.2, page 39).
- User administration (see chapter 7.3, page 42).
- Defining settings for the measured value display (see chapter 7.4, page 55).
- Configuring settings for acoustic signals (see chapter 7.5, page 55).

7.1 General system settings

7.1.1 Selecting the dialog language

The user interface is available in several languages. In addition to the two default dialog languages *English* and *German*, additional languages can be selected.

Proceed as follows to select the dialog language:

1 Opening the system settings

- In the main dialog, tap on **[System]**.
- Tap on **[System settings]**.

The **System / System settings** dialog is displayed.



2 Selecting the dialog language

- Tap on the **Dialog language** list box and select the desired language.

3 Saving the settings

Tap on the fixed keys [] or [].

The main dialog is displayed in the respective dialog language.

7.1.2 Setting the date, time and local time

The Coulometer displays the date and time in accordance with ISO standard 8601.

Proceed as follows to set the date and time:

1 Opening the system settings

- In the main dialog, tap on **[System]**.
- Tap on **[System settings]**.

The **System / System settings** dialog is displayed.

2 Entering the date

- Tap on the input field for the date.
The editor opens.
- Enter the current date in the format **YYYY-MM-DD** and confirm with **[OK]**.
The arrow keys [] and [] are used to move the cursor to the left or to the right by one character.

The input is saved and the editor is closed.

3 Entering the time

- Tap on the input field for the time.
The editor opens.
- Enter the current time in the format **hh:mm:ss** (24-hour format) and confirm with **[OK]**.
The arrow keys [] and [] are used to move the cursor to the left or to the right by one character.

The input is saved and the editor is closed.

4 Enter the local time

- Tap on the list box **Local time - UTC** and select the difference from the UTC (Coordinated Universal Time).
The selection **off** means that the time is saved with no difference from the UTC.


5 Saving the settings

Tap on the fixed keys [] or [].

The time settings are saved.

7.2 System-specific dialog options

Main dialog: **System ► System settings ► Dialog options**

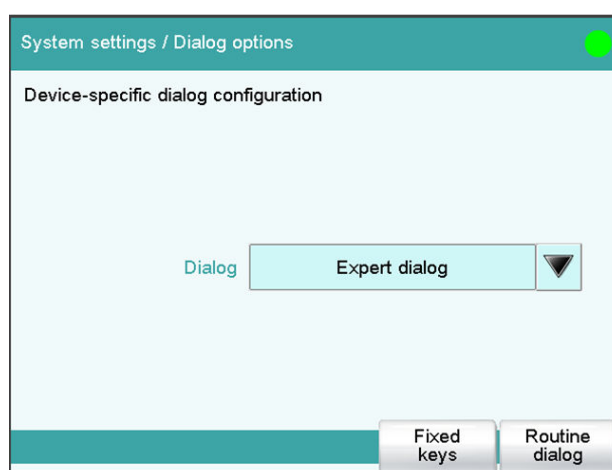

If you work without a login function, in this dialog you can define whether the system should generally be operated in expert mode or in routine mode. If you work with the login function activated, you must define this setting separately for each user (*see chapter 7.3.1, page 44*).

Dialog

Dialog mode in which the user may operate the system.

Selection	Expert dialog Routine dialog
Expert dialog	All functions that are supported by the system are available.
Routine dialog	The user dialog can be limited for routine operations. Only selected functions are available (<i>see Configuring the routine dialog, page 41</i>).

[Fixed keys]

Block unneeded fixed keys (see "*Blocking unneeded fixed keys*", page 40).

[Routine dialog]

Configure functions for the routine dialog (see "*Configuring the routine dialog*", page 41).

Selecting the dialog mode

Proceed as follows to change the dialog mode:

1 Select the dialog mode

Open the selection list **Dialog** and select either **Expert dialog** or **Routine dialog**.

2 Saving the settings

Tap on the fixed keys [] or [].

The setting will apply to all dialogs.



NOTICE

If you have selected **Routine dialog** and if the routine dialog was configured in such a way that the dialog **System settings / Dialog options** is blocked, then you can switch back over to the expert dialog as follows:

- Operation without login function:
In the main dialog, enter **User = Metrohm**.
- Operation with login function:
A user who works with expert dialog must log in.

Blocking unneeded fixed keys

This following configurations apply for **both** dialog modes: routine dialog and expert dialog.

Blocking fixed keys

Proceed as follows to block unneeded fixed keys:

1 Display fixed keys which can be blocked


Tap on the **[Fixed keys]** button.

All fixed keys which can be blocked are displayed.

2 Deactivate fixed keys

Deactivate those fixed keys which are not permitted to be used.

3 Saving the settings

Tap on the fixed keys [] or [].

Deactivated fixed keys cannot be used.

Configuring the routine dialog

A suitable **Standard configuration** has already been saved for routine operations:

- Methods can only be loaded, but not modified or created.
- Determinations cannot be recalculated.

You can readjust this standard configuration by disabling additional functions or re-enabling disabled functions.



NOTICE


The configuration of the routine dialog applies for all routine users. You also have the option of defining user-specific routine settings. To do this, you must work with the login function activated and create an identification profile for each user (*see chapter 7.3.2, page 47*).

Proceed as follows to modify the configuration for the routine dialog:

1 Opening the dialog

Tap on the [**Routine dialog**] button.

The list of all buttons in the main dialog, in the manual control, etc. is displayed:

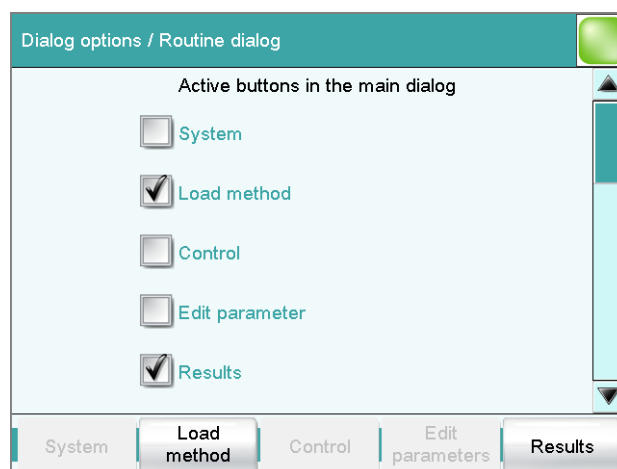

2 Deactivate buttons

Deactivate those buttons which are not permitted to be used.

All deactivated buttons will appear grayed-out, i.e. they are inactive.

3 Deactivate other functions

Many additional buttons and parameters can be disabled in the dialogs of **[System]**, **[Load method]**, **[Control]**, **[Edit parameters]** and **[Results]**. The corresponding option must be activated in order to enable these buttons.

4 Saving the settings

Tap on the fixed keys [] or [].

All deactivated functions will appear grayed-out, i.e. they are inactive.

7.3 User administration

Main dialog: **System ► System settings ► User admin.**

All of the functions of the user administration are described in this chapter. No user administration is mandatory for the simple operation of a titration system. If, however, you wish to be in compliance with *FDA Guidance 21 CFR Part 11*, then you must use the functions of the user administration. Additional information regarding the requirements of the FDA Guidance can be found in the document *Compliance Guide Touch Control*.



NOTICE

If you work with the login function activated, then the user administration is accessible only for users with administrator rights. This means that you must ensure that at least two users have administrator rights so at least one of them will be available. Keep the access rights for a user with administrator rights in a safe place so that they are accessible in an emergency.

System settings / User administration		
User	Dialog	Status
Administrator	Expert dialog	active
ak	Routine dialog	active
ga	Routine dialog	active
jb	Expert dialog	active
wl	Expert dialog	active

The following data is displayed in the user list for each user:

- Name
- Dialog mode in which the user may operate the system.
- Status

You can use the user list two different ways:

- Operation with login function:
If you work with the login function activated, i.e. if each user must log into the system before starting work, then only those users entered in the list can log in. The user who is currently logged in is shown in the main dialog.
- Operation without login function:
If you work without the login function, the users whose names are entered in the list can be selected in the main dialog or a user name can be entered. This makes it possible to document who has operated the titration system.

[Login options]

Define the settings for the login, Audit Trail, etc. (see chapter 7.3.3, page 48).

[Create ID profile]

Create an identification profile for the selected user on a storage medium (see chapter 7.3.2, page 47).

[New]

Adding a new user to the list (see chapter 7.3.1, page 44).

[Delete]

Delete the selected user from the list.

**NOTICE**

Once users have worked with the login function and password protection, they can no longer be deleted, even if the password protection is disabled again. The status of these users must be set to **inactive** (requirement of FDA Guidance 21 CFR Part 11).

The last user with administrator rights cannot be deleted.

[Edit]

Editing the data of the selected user (*see chapter 7.3.1, page 44*).

7.3.1 Editing the user configuration

User list: **User** ► **New** / **Edit**

User

The designation of the user is used for unambiguous identification, e.g. the company internal shorthand symbol or the personal number. The user name is printed out in all reports containing determination data and stored in the determination file. Each file always contains the name of the user who created it and the name of the last user to edit it.

Entry	24 characters maximum
Default value	empty

Full name

Complete name of the user.

Entry	24 characters maximum
Default value	empty

Dialog

Dialog mode in which the user may operate the system.

Selection	Expert dialog Routine dialog
-----------	---------------------------------------

Expert dialog

All functions that are supported by the system are available.

Routine dialog

The user dialog can be limited for routine operations. Only selected functions are available (see *Configuring the routine dialog*, page 41).

Status

Status of the user. Users can be deactivated. This function is useful, for instance, if the user is no longer authorized to operate the system or no longer works for the company.

Selection	active inactive
Default value	active

active

The user is authorized to operate the system.

inactive

The user is not authorized to operate the system and can no longer log in.

Admin. rights

on | off (Default value: **off**)

If this option is activated, then the user has administrator rights.



NOTICE

The last user with administrator rights cannot be deleted anymore.

[Favorites]

Saving methods and sample tables as user-specific favorites (see *chapter 18, page 155*).

[Signature method]

Assign the permission to use and sign methods (see *"Edit user / Signature method" dialog*, page 45).

[Signature determ.]

Assign the permission to use and sign determinations (see *"Edit user / Signature determination" dialog*, page 46).

"Edit user / Signature method" dialog

User: **Edit ► Signature method**

Authorizations for using and signing methods can be defined in this dialog. These settings are only effective if you are working with the login function enabled and password protection.

Use only released methods

on | off (Default value: **off**)

If this option is activated, then the user may only start methods that have been released.

Review methods (signature level 1)

on | off (Default value: **off**)

If this option is activated, then the user may only sign methods which have the status **saved**. The method is assigned the status **reviewed**.

Release methods (signature level 2)

on | off (Default value: **off**)

If this option is activated, then the user may only sign methods which have the status **reviewed**. The method is assigned the status **released**.



NOTICE

If the two options **Review methods** and **Release methods** are selected for a user, then this user can sign different methods at either level 1 or level 2, but not the same method at both level 1 and level 2.

Delete signatures

on | off (Default value: **off**)

If this option is activated, then the user may delete the signatures of a method which has been released. The method is assigned the status **saved**. The signatures of a method can only be deleted if it has the status **released**.

"Edit user / Signature determination" dialog

User: **Edit** ► **Signature determ.**

Authorizations for signing determinations can be defined in this dialog. These settings are only effective if you are working with the login function enabled and password protection.

Review determinations (signature level 1)**on | off** (Default value: **off**)

If this option is activated, then the user may only sign determinations at the first level. The determination is assigned the status **reviewed**.

Release determinations (signature level 2)**on | off** (Default value: **off**)

If this option is activated, then the user may only sign determinations at the second level. The determination is assigned the status **released**.

Delete signatures**on | off** (Default value: **off**)

If this option is activated, then the user may delete the signatures of a determination which has been released. The signatures of a determination can only be deleted if it has the status **released**.

7.3.2 Creating an identification profileUser list: **User ► Create ID profile**

If you plan to carry out the login with an identification profile (*see chapter 7.3.3, page 48*), then you must first create an identification profile on a storage medium for each user. A check can then be made at the time of login as to whether or not the user does exist and whether or not he or she is working in the expert dialog or in the routine dialog. After a successful login the routine dialog settings stored on the card are loaded.

**NOTICE**

In addition to the user name, the current routine dialog settings are also stored in this identification profile. This means you can define user-specific routine dialog settings for each user. However, you must configure them (*see "Configuring the routine dialog", page 41*), **before** you create the identification profile.

Before you create the identification profile, check whether the desired routine dialog settings are active.

1 Plug in a storage medium

Plug in the USB storage medium.

2 Select a user

In the user list, select the user for whom the profile is to be created.

3 Create the identification profile

Tap on **[Create ID profile]**.

The user configuration and the current routine dialog settings are saved.

7.3.3 Defining login options

Main dialog: **System ► System settings ► User admin. ► Login options**

There are a number of different ways to log onto the system:

- Without login
- Login via user name
- Login via user name and password
- Login via identification profile
- Login via identification profile and password

**NOTICE**

If you work with the login function activated, then the user administration is accessible only for users with administrator rights. This means that you must ensure that at least two users have administrator rights so at least one of them will be available. Keep the access rights for a user with administrator rights in a safe place so that they are accessible in an emergency.

**NOTICE**

If you exit this dialog with [] or [], and if you have selected one of the login variants **Login via user name** or **Login via identification profile**, then the login dialog will open automatically and you must also log in to the system.

This means that you must make sure that you have first defined all the users and created the identification profiles before you activate the login function.

User administration / Login options

Login via user name

Login via identification profile

Password required

Logout automatically

after min

Login only for the same user

Audit Trail Reasons Modific. options Password options

Login via user name

on | off (Default value: **off**)

If this option is activated, then the user must log in with his or her unambiguous identification.

Login via identification profile

on | off (Default value: **off**)

If this option is activated, then the login will take place via USB storage medium with the identification profile stored on it.

Password required

on | off (Default value: **off**)

If this option is activated, then the user must enter a password in addition to his or her user name or identification profile.

Logout automatically

on | off (Default value: **off**)

If this option is activated, then the user will be logged out automatically after the specified time.

Input range	1 to 60 min
-------------	--------------------

Login only for the same user

on | off (Default value: **off**)

If this option is activated, then only the same user may log in again after he or she has logged out. Users with administrator rights can, however, log in at any time.

[Audit Trail]

Define the settings for recording an Audit Trail (*see chapter 7.3.7, page 53*).



[Reasons]

Predefining a list of reasons which can be selected for the modification/signature of a method or determination (see chapter 7.3.6, page 52).

[Modific. options]

Defining the modifications for which a reason is required (see chapter 7.3.5, page 52).


[Password options]

Define the settings for the password, see following chapter.

7.3.4 Password options

Main dialog: **System ► System settings ► User admin. ► Login options ► Password options**

You can make various settings for password entry in the password options.

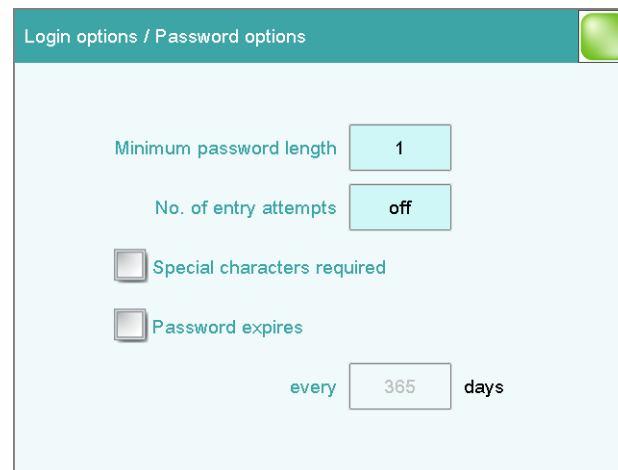

Minimum password length

Minimum number of characters of the passwords.

Input range	1 to 10
Default value	1

No. of entry attempts

If the user has logged in incorrectly this many times, then it will automatically be deactivated. It can only be reactivated by a user with administrator rights.

Input range	2 to 5
Selection	off
Default value	off

Special characters required

on | off (Default value: **off**)

If this option is activated, then the password must contain one of the following special characters: ° § + | @ * # ç % & ¬ () = ' ^ ` ~] [] { - _ : . ; , > < £ !

Password expires

on | off (Default value: **off**)

If this option is activated, then the user must define a new password after the time specified. A password that has already been used cannot be used again.

Input range	1 to 999 days
Default value	365 days

Forgotten password



CAUTION

If a user has forgotten his password, a new user name must be defined. The same user name can only be used again after a re-installation and re-creating the user list.

Proceed as follows:

1 Create a backup

Create a backup (*see chapter 11.3, page 97*).

2 Executing a system initialization

Switch off the Coulometer and execute a system initialization (*see chapter 29.1, page 302*).

3 Restore backed-up data

Use the function **Restore** to reload the data from the backup into your system (*see chapter 11.3.1, page 97*).

Deactivate the options **User list** and **System settings / User admin..**

4 Switch on the Coulometer

Switch the Coulometer back on again after a few seconds.

5 Restore the user list

Create the user list again and redefine the login options.

7.3.5 Modification options

Main dialog: **System ► System settings ► User admin. ► Login options ► Modific. options**

In the dialog **Login options / Modification options**, you can define for which actions a reason must be entered. These reasons are documented in the Audit Trail (*see chapter 7.3.7, page 53*) together with the modification. The reason for the last modification is shown in the properties of the method or determination.



NOTICE

The reasons are only requested when working with activated login function and password.

Saving modified method

on | off (Default value: **off**)

If this option is activated, then a reason must be given when saving a method modification.

Recalculating determination

on | off (Default value: **off**)

If this option is activated, then a reason must be given when determinations are being recalculated.

7.3.6 Reasons

Main dialog: **System ► System settings ► User admin. ► Login options ► Reasons**

In the dialog **Login options / Reasons**, you can create a selection list containing reasons from which a selection can be made when signing and modifying methods and determinations. Some reasons have already been provided.

**[New]**

Add a new reason to the list.

[Delete]

Delete the selected reason from the list.

[Edit]

Change the designation of the selected reason.

7.3.7 Audit Trail

Main dialog: **System ► System settings ► User admin. ► Login options ► Audit Trail**

An Audit Trail is an automatically generated log of all user activities. An Audit Trail contains precise logs of user actions (date, time, user, action, etc.). Recording an Audit Trail is important for compliance with *FDA Guidance 21 CFR Part 11* when using PC programs. A step-by-step description of how you must proceed to remain in compliance with the FDA Guidance is included in the document *Compliance Guide Touch Control*. The Audit Trail is saved to the internal memory.

You can also use the Audit Trail function specifically to record the data which are of interest to you.

You can view, filter and export the Audit Trail on a PC with the supplied software program *AuditTrailViewer*. Details concerning utilization of the *AuditTrailViewer* can be found in the Appendix (*chapter 31.6, page 314*).

If necessary, you can delete all of the entries in the Audit Trail with **[Delete Audit Trail]**. Do not fail to create a backup before doing so, however.

7.4 Measured value display

Main dialog: **System ► System settings ► Meas. value display**

This setting refers exclusively to the display of the measured values in the live display and in the manual control. The values will, however, always be stored with their full accuracy.

7.5 Acoustic signals

Main dialog: **System ► System settings ► Acoustic signals**


You can define acoustic signals in order to direct attention to particular events. You can define signals for the following events:

- **Wrong manipulation**
An acoustic signal will sound each time an invalid action is attempted (e.g. pressing [?] again when Help is open).
- **Display a message**
A short beep will sound each time a message appears on the display. This informs the user that the message must be confirmed.
- **Button contact**
Each time a button is touched on the touch screen, this will be confirmed by an acoustic signal.
- **External data input**
An acoustic signal will provide confirmation each time data is received from external devices (e.g. balance, barcode reader).

8 Reagents

Main dialog: **System ► Reagents**

This chapter describes how you can create a list of the reagents used in the system.

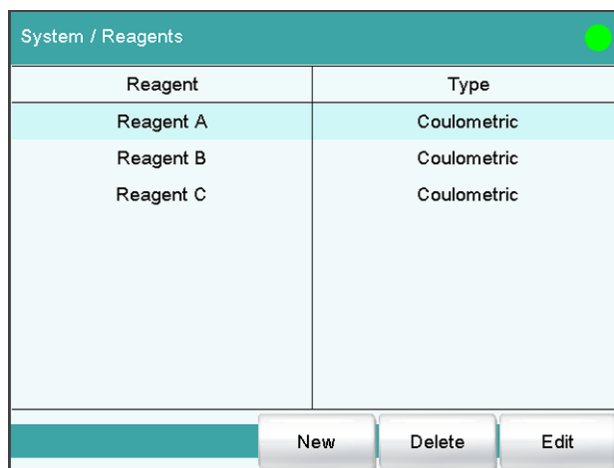

Reagent	Type
Reagent A	Coulometric
Reagent B	Coulometric
Reagent C	Coulometric

The list of reagents gives the designation and type of each reagent.

[New]

Add a new reagent to the list, see following chapter.

[Delete]

Delete the selected reagent from the list.

[Edit]

Edit the data of the selected reagent, see following chapter.

8.1 Editing reagent data

Main dialog: **System ► Reagents ► New / Edit**

Reagent

The designation of the reagent is used for unambiguous identification.

Entry **24 characters maximum**

Comment

Entry **24 characters maximum**

[Reagent monitoring]

Set the parameters for the reagent monitoring, see following chapter.

8.2 Reagent monitoring

The conditions for the monitoring of the reagent are defined in the dialog **Edit reagent / Reagent monitoring**.

If one of the following values is reached, then the reagent must be replaced. The values are checked in the following cases:

- at the start of the determination.



- at the end of the determination.

Number of determ.

The number of determinations to be carried out with a certain amount of reagent depends on the type of sample and its amount.

Input range	1 to 999
Selection	off
Default value	off

Working life

Working life of the reagent.

Input range	1 to 999 days
Selection	off
Default value	off

Reagent capacity

This parameter is only visible with coulometric reagents.

Water capacity of the reagent.

Input range	1 to 9999 mg
Selection	off
Default value	off

Drift

This parameter is only visible with coulometric reagents.

During conditioning the measured drift has to be in the following range for 2 min: '**specified drift value + 50 µg/min**'.

Input range	1 to 999 µg/min
Selection	off
Default value	off

[Status]

Display the status overview of the current values of the reagent monitoring.

[Reagent replacement]

Edit the parameters for the reagent exchange.

"Reagent monitoring / Status" dialog

The current reagent monitoring values are displayed in this dialog.

[Reset]

Reset the values to zero.

Dialog "Reagent monitoring / Reagent replacement"

The parameters for the reagent replacement are defined in this dialog.

Reagent replacement

The reagent can either be exchanged manually or automatically.

Selection	manual auto
Default value	manual

manual

If a monitored parameter has reached the limit set, a message is being displayed. Then the reagent has to be exchanged manually.

auto

If a monitored parameter has reached the limit set, the method defined below is started automatically.

Memory

This parameter can only be edited with **Reagent replacement = auto**.

Memory location the method is loaded from. All memory locations are selectable, even if they are currently not accessible.

Selection	Internal memory External memory 1 External memory 2 Shared memory
Default value	Internal memory

Shared memory

Shared directory in the network.

Method

This parameter can only be edited with **Reagent replacement = auto**.

Method used for emptying the titration cell.



NOTICE


Make sure that the memory is accessible.

Entry	32 characters maximum
Selection	Selection of stored methods

9 Sensors

Main dialog: **System ► Sensors**

This chapter describes how you can create a list of the sensors used in the system.

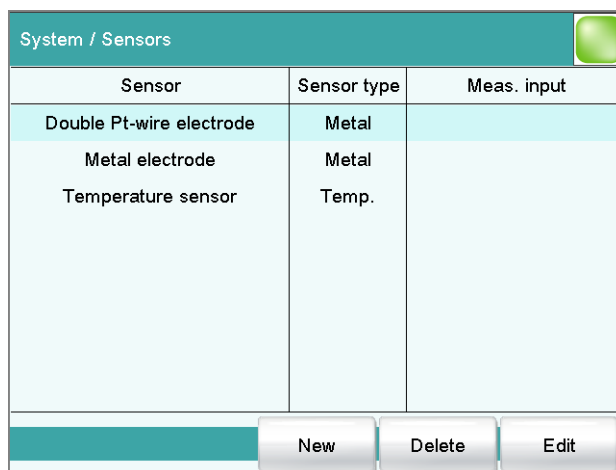

Sensor	Sensor type	Meas. input
Double Pt-wire electrode	Metal	
Metal electrode	Metal	
Temperature sensor	Temp.	

Two standard sensors are defined in the sensor list: **Metal electrode** and **Temperature sensor**. These sensors cannot be deleted or renamed. A maximum of 25 additional sensors can be added to these sensors.

For each sensor, the following data is displayed in the sensor list:

- Designation
- Type
 - **Metal**: Metal electrode
 - **Temp.**: Temperature sensor
 - **Other**: Other sensor

The following sensor data is stored in the list of sensors:

- Name
 - Each sensor in the system is identified by its unambiguous name.
- Working life
- etc.

[New]

Adding a new sensor to the list (*see chapter 9.1, page 61*).

[Delete]

Delete the selected sensor from the list.

[Edit]

Editing the data of the selected sensor (*see chapter 9.2, page 61*).

9.1 Adding a new sensor

Before you can use a sensor, you must add it to the sensor list. To do this, use the button **[New]**.

- The properties dialog is opened after the sensor type has been selected, see following chapter. The following sensor types can be selected:
 - Metal electrode (Pt electrode)
 - Other sensor
 - Temperature sensor

9.2 Editing the sensor data

Sensor list: **Sensor ▶ New / Edit**

All of the data for the selected sensor is displayed in the dialog **Sensors / Edit**.

Sensor

The designation of the sensor is used for unambiguous identification.

Entry	24 characters maximum
-------	------------------------------

Order number

Order number of the sensor.

Entry	24 characters maximum
Default value	empty

Serial number

Serial number of the sensor.



Entry	8 digits maximum
-------	-------------------------

Comment

Entry	24 characters maximum
-------	------------------------------

[Working life]

Defining the working life of the sensor (*see chapter 9.3, page 62*).

9.3 Monitoring the working life

Sensor: **Edit ► Working life**

In the dialog **Edit sensor / Working life**, you can define the time interval after which the sensor must be replaced. If you do not wish to monitor the working life, then you can enter only the date of manufacture for documentation purposes.

Start-up

Date on which the sensor was used for the first time.

Format: YYYY:MM:DD

Monitoring

on | off (Default value: **off**)

If this parameter is activated, then the working life will be monitored.

Working life

If you define a time interval for the working life, then the **Expiry date** will be tracked automatically.

Input range	1 to 999 days
Default value	999 days

Expiry date

If you define an expiry date, then the **Working life** will be tracked automatically.

Format: YYYY:MM:DD

Action

Selection of the action which is carried out when the time interval has expired.

Selection	Display message Document message Cancel determination
Default value	Display message For all three options it is documented in the determination data (see dialog More determination data / Messages), that the time interval has been expired.

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message

In the determination data it will be documented, that the time interval has been expired.


Cancel determination

The determination is stopped.

10 Device manager

Main dialog: **System ► Device manager**

This chapter describes how you can configure the Coulometer and the connected peripheral devices. A detailed description of the hardware installation can be found in the corresponding installation manual.

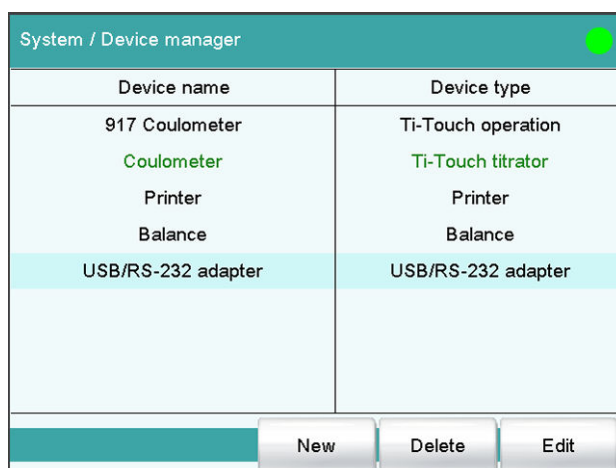

Device name	Device type
917 Coulometer	Ti-Touch operation
Coulometer	Ti-Touch titrator
Printer	Printer
Balance	Balance
USB/RS-232 adapter	USB/RS-232 adapter

The instrument list gives the name and type of every instrument. The **Ti-Touch titrator** is a component of the 917 Coulometer and therefore permanently entered as control instrument in the device manager. If an 885 Compact Oven SC is connected and switched on, it is automatically entered in the list. Peripheral devices (dosing devices, stirrers, etc.) connected to the MSB connector are entered in the list automatically. A USB/RS-232 adapter will also be automatically recognized and entered in the instrument list with default settings. A printer is entered in the instrument list in the default settings. You must enter a PC keyboard, a barcode reader or a balance in the instrument list yourself.

[New]

Add a new device to the list (*see chapter 10.1, page 65*).

[Delete]

Delete the selected device from the list.



NOTICE

Devices which are recognized automatically cannot be deleted from the list while they are still connected.

[Edit]

Configure the selected device (*see chapter 10.2, page 65*).

10.1 Adding a new device

Of the following device types, you can enter one device each in the device list, even if it is not yet connected:

- Sample Processor
- Balance
- Barcode reader
- USB/RS-232 adapter
- PC keyboard

Proceed as follows:

1 Displaying the device selection

Tap on **[New]**.

2 Selecting the instrument

Tap on the button for the desired device.

The new device is entered in the list.

10.2 Configuring the instrument

Instrument list: **Instrument ► Edit**


The data stored for an instrument depends on the type of instrument. You can define an instrument name and a comment for each instrument. The instrument name of the Coulometer is printed out in the standard report header.

The description of the individual instruments can be found in the following specific chapters:

- Coulometer (*see chapter 10.3, page 66*)
- Metrohm control instruments (*see chapter 10.4, page 72*)
- 885 Compact Oven SC (*see chapter 10.5, page 75*)
- Printer (*see chapter 10.6, page 79*)
- Balance (*see chapter 10.7, page 83*)
- PC keyboard (*see chapter 10.9, page 87*)
- USB/RS-232 adapter (*see chapter 10.8, page 85*)
- Barcode reader (*see chapter 10.10, page 88*)

10.3 Coulometer

Instrument list: **917 Coulometer** ▶ Edit

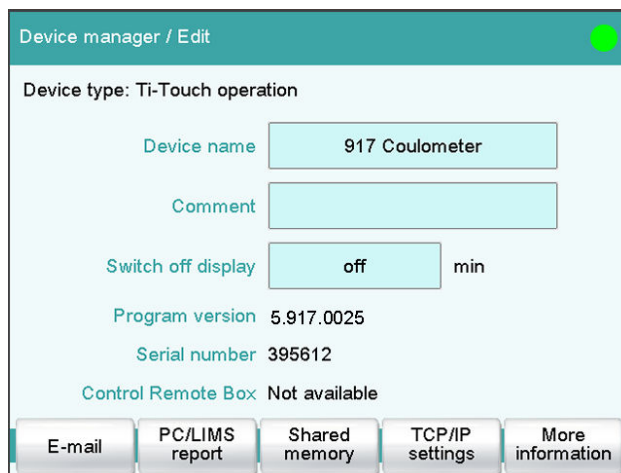

Device name

This designation is used for identification purposes when selecting control devices (command, manual control).

Entry	24 characters maximum
-------	------------------------------

Comment

Entry	24 characters maximum
-------	------------------------------

Switch off display

If this time interval has expired without the Coulometer having been operated, then the display will be switched off. Touching the display will switch it back on again at any time.

Input range	1 to 999 min
Selection	off
Default value	off

Program version

Program version of the software.

Serial number

Shows the serial number of the device.

Control Remote Box

Shows to which control device and MSB connector the Remote Box is connected.



The "Control Remote Box" is the interface via which the system can be started and stopped externally. If multiple Remote Boxes are connected, then the one that is recognized first when the program starts will be used as the "Control Remote Box."

Selection	Name of the control device / Number of the MSB connector Not available
-----------	--

10.3.1 E-mail

917 Coulometer: **Edit ► E-mail**

The system allows you to send displayed messages as e-mails. The Coulometer must be connected to a network for this to function. The following types of messages can be sent:

- : general warning messages
- : error messages

Configuring e-mail dispatch

Proceed as follows so that messages can be sent as e-mails:

1 Activating the option

- In the instrument properties of the 917 Coulometer, tap on the **[E-mail]** button.
- In the **Edit device / E-mail** dialog, activate the **Send the following messages as e-mail:** option.

2 Configuring e-mail addresses



- Tap on the **[E-mail settings]** button. The **E-mail / Settings** dialog is displayed.
- Enter the addresses of the mail server, the sender and the desired recipient.

Parameter description

Send the following messages as e-mail:

on | off (Default value: **off**)

If this parameter is activated, then messages with the following symbols will be sent as e-mails:

- : General warning messages
- : Error messages

Send only during running determination**on | off** (Default value: **on**)

If this parameter is activated, then messages will be sent as e-mails only if a determination is running. Deactivate this parameter if messages are also to be sent in normal status.

Mail server

Address of the mail server for outgoing mail, e.g. mail.metrohm.ch. You can find the address of the mail server either in your e-mail program settings or obtain it from your IT department.

Entry	50 characters maximum
Default value	empty

Sender

E-mail address of the sender. This address must be formatted as an e-mail address, but need not necessarily correspond to an existing e-mail account, e.g. coulometer@metrohm.com.

Entry	50 characters maximum
Default value	empty

Recipient


The messages will be sent to this e-mail address.

Entry	50 characters maximum
Default value	empty

10.3.2 PC/LIMS report

917 Coulometer: **Edit ► PC/LIMS report**

You can generate a machine-readable report with all of the important data concerning a determination, which is referred to as a PC/LIMS report. This report can be saved as a TXT file (according to ISO/IEC 8859-1) or as a UTF-8 file or sent to a terminal program or a LIMS via an RS-232 interface:

- manually with the [] fixed key (*see chapter 26, page 205*).
- automatically at the end of a determination (*see chapter 15.5.6, page 146*).

The file name of the report file is constructed as follows: *PC_LIMS_Report-ID1-YYYYMMDD-hhmmss.txt*. A detailed description of the contents of the PC/LIMS report can be found in the *PC/LIMS Report Guide*.

Memory

Memory location where the PC/LIMS report is stored as a file. The report will be saved in the directory *pc_lims_report*. This directory will be created

the first time a PC/LIMS report is generated. All three memory locations are listed as possible selections, even if they cannot be accessed at the moment.

Selection	off External memory 1 External memory 2 Shared memory
Default value	off

off

The report will not be saved as a file.

Shared memory

The report will be saved in a shared directory on the network. The shared directory is selected in the **Edit device / Shared memory** dialog (see chapter 10.3.3, page 69).

RS-232

The RS-232 interface via which the PC/LIMS report is sent. The interface parameters are adjusted in the **Edit device / Port parameters** dialog (see chapter 10.8, page 85).

Selection	off COM 1 COM 2
Default value	off

off

The report will not be sent via an RS-232 interface.

COM 2

This interface is inactive.

Coding

Format in which the PC/LIMS report is coded and stored.

Selection	ISO 8859-1 UTF-8
Default value	ISO 8859-1

ISO 8859-1

This format is recommended for all languages that use the extended ASCII code (e.g. German, English, Spanish, etc.).

UTF-8

This format is required for all languages that do not use the extended ASCII code (e.g. Russian, Chinese, Korean, etc.).

10.3.3 Shared memory

917 Coulometer: **Edit ► Shared memory**

If you have your Coulometer connected to your network, then you can define in this dialog a shared memory location on a computer within your network for the purpose of saving data (methods, determinations, etc.).



CAUTION

If the computer on which you share a memory location does not have the same subnet as the Coulometer, then a WINS server must be present. The computer must be entered in this server. In order for the communication to work, the SMB protocol version 1 (SMBv1) must be activated. Since this function is not standard anymore in Windows 10, it has to be activated manually. From firmware version 5.917.0024, the SMB protocol version 2 and 3 is also supported.

Computer

Host name of the computer on which a memory location is to be shared. If you are working with Windows, then you will find the host name of the computer as follows: In the Windows **command prompt** window, enter the command **ipconfig -all**. The host name is listed together with other parameters of the computer.

Do not under any circumstances enter an IP address in this input field.

Entry	max. 32 characters
Default value	empty

Share name

Share name of the shared memory location (file directory) on the above-specified computer. Please note that the Share name of a file directory often does not match the name of the file directory. The share name can be found on the Release tab in the Properties dialog of the shared file directory.

Entry	max. 32 characters
Default value	empty

Domain

Network domain in which the above-specified computer is located. If you are working with Windows, then you will find the domain name as follows: In the Windows **command window**, enter the command **ipconfig -all**. The domain is listed together with other parameters of the computer. Leave the field empty if the computer is not located within a domain.

Entry	max. 32 characters
Default value	empty

User

User name of the user authorized to access the shared memory location.

Entry	max. 32 characters
Default value	empty

Password

Password of the user configured on the computer.

Entry	max. 32 characters
Default value	empty

[Connect]

Establish the network connection. If the connection has been set up correctly, then all of the input fields will become inactive and the label switches to **[Disconnect]**. The network connection can be disconnected with this.

10.3.4 TCP/IP settings

917 Coulometer: **Edit ► TCP/IP settings**

If you have connected your Coulometer to your network, then you will have to define the network-relevant settings in this dialog. The Coulometer requires an IP address so that it is identifiable on the network. The Coulometer can acquire the IP address either dynamically from a DHCP server or you can enter the address directly.

Get IP address automatically (DHCP)

on | off (Default value: **on**)

If this parameter is activated, then the Coulometer will obtain its IP address directly from a DHCP server. In this case, the remaining parameters can no longer be edited.

IP address

IP address for the Coulometer. IP addresses are 32-bit numbers and are written as sequences of four decimals, each separated by a period, e.g. "10.157.212.8".

Entry	x.x.x.x
Default value	192.168.10.11
	"x" is a decimal between 0 and 255.

Subnet mask

The net mask or subnet mask, together with the IP address, indicates to which network the device to be connected belongs. Subnet masks are 32-bit numbers and are written as sequences of four decimals, each separated by a period.

Entry	x.x.x.x
Default value	255.255.255.0

- Properties of the measuring inputs (*see chapter 10.4.1, page 73*)
- Properties of the MSB connectors (*see chapter 10.4.2, page 74*)
- Properties of the peripheral devices on the MSB connector (*see chapter 10.4.3, page 75*)

10.4.1 Properties – Measuring input

Instrument list: **Control instrument** ► **Edit** ► **Measuring input 1** ► **Properties**

Device manager / Edit

Coulometer / Measuring input 1

ADC type 3.851.1200

Serial number 27251

Thick film type 3.680.0780

Temperature sensor

Type **Pt1000** ▼ R (25 °C)

B value K

In this dialog, you define the type of temperature sensor you have connected to the selected measuring input.

ADC type

Type of analog-digital converter.

Serial number

Serial number of the measuring interface.

Thick film type

Type of the thick film.

Temperature sensor

Type

The instrument supports the use of two different temperature measurement techniques:

- NTC (Negative Temperature Coefficient)
- Pt1000 (Platinum resistance)

Select here the type that has been connected to the instrument. If an NTC sensor is used, then it is also necessary to enter two sensor characteristics. These characteristics are listed in the specifications of the sensor.

Selection	Pt1000 NTC
Default value	Pt1000

R (25 °C)

This parameter can only be edited with **Type = NTC**.

Nominal resistance of the NTC sensor at 25 °C.

Input range	1000 to 99999 ohm
Default value	30000 ohm The default value applies to Metrohm sensors with an NTC sensor.

B value

This parameter can only be edited with **Type = NTC**.

Material constant of the NTC sensor. B values of NTC sensors are frequently based on different reference temperatures (usually 25 °C and 50...100 °C). When entering the B value the influence of the second reference temperature is negligible in comparison with the measuring accuracy of an NTC sensor.

Input range	1000 to 9999 K
Default value	4100 K The default value applies to Metrohm sensors with an NTC sensor. If no B value is given for your sensor then you can retain the default value.

10.4.2 Properties – MSB connector

Instrument list: **Control instrument ▶ Edit ▶ MSB connector 1/2 ▶ Properties**

Device manager / Edit

Coulometer / MSB 1

When would you like to be asked to prepare the dosing device on MSB 1?

Switch on

Attach an exchange or dosing unit

Time interval h

In this dialog, you can define when the request to carry out the **Prepare** function for connected dosing devices is to be displayed. This setting applies for all dosing devices of the selected MSB connector. The manual

control contains a description of how to prepare the exchange unit and/or the dosing unit (*see chapter 27.3.3, page 221*).

Switch on

on | off (Default value: **on**)

If this parameter is activated, then you will be requested to prepare the dosing device when the Coulometer is switched on.

Attach an exchange or dosing unit

on | off (Default value: **on**)

If this parameter is activated, then you will be requested to prepare the dosing device when the exchange/dosing unit is attached.

Time interval

on | off (Default value: **off**)

Activate this parameter if you wish to receive a regular request to prepare the dosing unit.

Input range	0.1 to 999.9 h
Default value	12.0 h

10.4.3 Properties – Peripheral devices

Instrument list: **Control instrument ► Edit ► Peripheral device ► Properties**

The following data is displayed for the connected dosing devices and stirrers:

- Type
- Serial number


10.5 885 Compact Oven SC



NOTICE

In order for the 885 Compact Oven SC to be controlled with the 917 Coulometer, at least the following firmware versions must be installed on the instruments:

- 917 Coulometer 5.917.0025
- 885 Compact Oven SC 5.885.0013

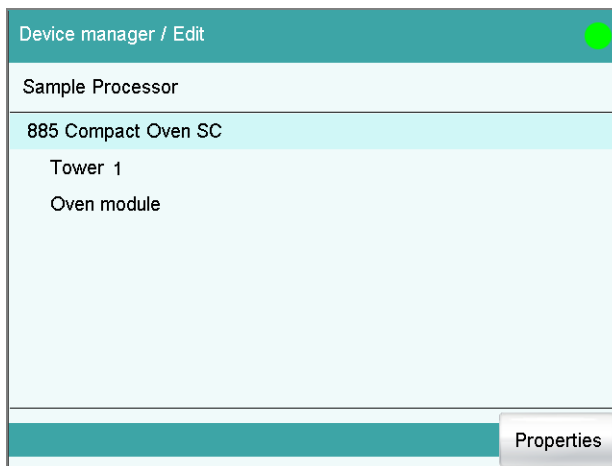

The dialog displays the 885 Compact Oven SC with its tower and the oven module. If the control instrument is not connected, then the properties dialog of the control instrument will be displayed directly (see chapter 10.5.1, page 76).

Properties

Open the Properties dialog of the highlighted entry.

10.5.1 Properties - 885 Compact Oven SC

Instrument list: **885 Compact Oven SC** ▶ **Edit** ▶ **Properties**

Device name

This designation is used for identification purposes when selecting control instruments (command, manual control).

Entry **24 characters maximum**

Comment

Entry **24 characters maximum**

Program version


Program version of the device software.

Serial number

Display of the serial number of the device.

10.5.2 Properties - Tower

Instrument list: **885 Compact Oven SC** ▶ **Edit** ▶ **Tower** ▶ **Properties**

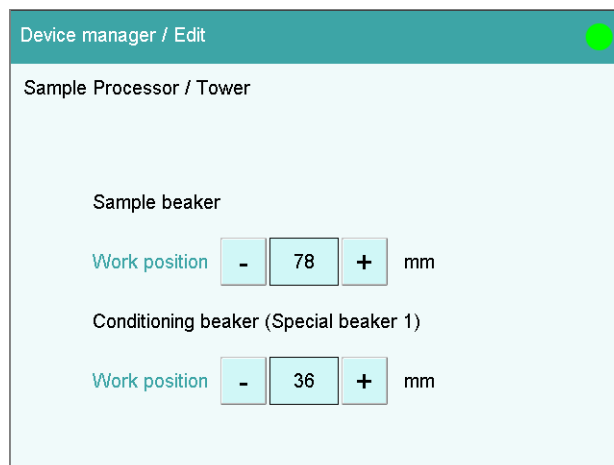

Sample beaker

Work position

At this lift position, the sample beaker is optimally placed in the oven and the needle has completely pierced the septum.

Input range	0 to 96 mm (Increment: 6)
Default value	78 mm

A lift height of 0 mm corresponds to the "home position". The lift is located at the upper stop position.

Conditioning beaker (special beaker 1)

Work position

At this lift position, the needle has completely pierced the septum of the conditioning beaker at the conditioning position. In the method run, the **[MOVE]** command is used to move the conditioning bucket with the **Objective** Special beaker 1.


Input range	0 to 96 mm (Increment: 6)
Default value	36 mm

A lift height of 0 mm corresponds to the "home position". The lift is located at the upper stop position.



10.5.3 Properties - Oven module

Instrument list: **885 Compact Oven SC** ▶ **Edit** ▶ **Oven module** ▶ **Properties**

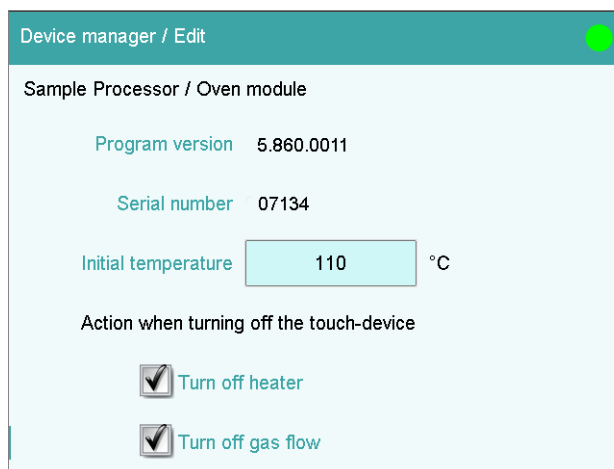

Program version

Program version of the oven module.

Serial number

Display of the serial number of the oven module.

Initial temperature

Entering an initial temperature means that the oven will heat up to the set temperature when the 917 Coulometer is switched on.

Input range	50 to 250 °C (Increment: 1)
Default value	110 °C
Selection	off

off

The oven is not switched on.

Action when switching off the touch device.

Switch off heater

Setting for the behavior of the heater of the oven module when the Coulometer is switched off.

Selection	on off
Default value	on

Switch off the gas flow

Setting for the behavior of the gas flow of the oven module when the Coulometer is switched off.

Selection	on off
Default value	on

10.6 Printer

Instrument list: **Printer ▶ Edit**

The screenshot shows a dialog box titled "Device manager / Edit" with a green close button. Below the title bar, it says "Device type: Printer". There are four input fields: "Device name" with the value "Printer", "Comment" (empty), "Printer" with a dropdown menu showing "PCL Inkjet", and "Connector" with a dropdown menu showing "USB". At the bottom, there are three buttons: "PDF settings", "Network printer", and "More parameters".

The list of devices always includes a printer, even if no corresponding device is connected. In this dialog, you also define when you would like to have a report generated as a PDF file.

Device name

Here you can enter a device name of your choice.

Entry	24 characters maximum
-------	------------------------------

Comment

Entry	24 characters maximum
-------	------------------------------

Printer

Selection of the printer type.

Selection	PCL Inkjet PCL Laser HP Epson Epson new Canon CUSTOM NEOS off
-----------	--

PCL Inkjet

for HP DeskJet printers.

PCL Laser

for HP LaserJet printers.

CUSTOM NEOS

Due to the narrow paper size, only reports marked with # in the "List of all printable reports" (*see chapter 26.3, page 209*) can be printed.

off

The reports will not be printed out on paper.

Connector

Selection of the connection type for the printer.

Selection	USB Ethernet
Default value	USB

USB

Local printer at a USB interface.

Ethernet

Network printer.

**NOTICE**

The **Coulometer** prints the reports with a fixed resolution of 300 dpi. If you are using a printer with a resolution of 360 dpi (or a multiple thereof, e.g. an Epson), then the text will be printed out somewhat smaller than with printers with a resolution of 300 dpi (or a multiple thereof, e.g. a Canon or HP).

10.6.1 PDF settings

Printer: **Edit ► PDF settings**

The settings for saving a report as a PDF file are defined in this dialog.

Memory

Memory location where the PDF file will be saved. The report will be saved in the directory *PDF_Report*. This directory is created when a PDF file is generated. The file name is appended with the date and time when the determination was started and the date and time when the PDF was generated (e.g.: **Identification 1**-20130320-141044-20130320-141220). The first date and time set indicates the start of the determination, the second date and time set indicates the generation of the PDF file. This prevents previously created PDF files from being overwritten each time a new file is created, e.g. when a result is recalculated and the report is generated again.

All three memory locations are listed as possible selections, even if they cannot be accessed at the moment.

Selection	off External memory 1 External memory 2 Shared memory
Default value	off

off

The report will not be saved as a PDF file.

Shared memory

The report will be saved in a shared directory on the network. The shared directory is selected in the **Edit device / Shared memory** dialog (see chapter 10.3.3, page 69).

Copy or extract content allowed

on | off (Default value: **on**)

If this option is activated, then content can be copied or deleted from the PDF file.

Printing allowed

on | off (Default value: **on**)

If this option is activated, then the PDF file can be printed.

Change the document allowed

on | off (Default value: **off**)

If this option is activated, then the PDF file can be edited.

Add or change comments allowed

on | off (Default value: **on**)

If this option is activated, then comments can be added to the PDF file.

10.6.2 Network printer

Printer: **Edit ► Network printer**

If you have connected your Coulometer to your network, then you can specify a network printer for your reports in this dialog.

**NOTICE**

If you have shared a memory location on a PC within your network and if you have configured the Coulometer accordingly (see chapter 10.3.3, page 69), then the settings for the parameters **Domain**, **User** and **Password** will be applied and can then no longer be modified in this dialog.



CAUTION

If the computer on which the network printer is configured does not have the same subnet as the Coulometer, then a WINS server must be present. The computer must be entered in this server. In order for the communication to work, the SMB protocol version 1 (SMBv1) must be activated. Since this function is not standard anymore in Windows 10, it has to be activated manually. From firmware version 5.917.0024, the SMB protocol version 2 and 3 is also supported.

Print server

Host name of the print server or the computer on which the network printer is configured. If you are working with Windows, then you will find the host name as follows: In the Windows **command prompt** window, enter the command **ipconfig -all**. The host name is listed together with other parameters of the print server or the computer.

Do not under any circumstances enter an IP address in this input field.

Entry	max. 32 characters
Default value	empty

Share name

Share name of the network printer on the above-specified computer. Please note that the Share name of a network printer often does not match the name of the printer. The share name can be found on the Release tab in the Properties dialog of the released network printer.

Entry	max. 32 characters
Default value	empty

Domain

Network domain in which the above-specified computer is located. If you are working with Windows, then you will find the domain name as follows: In the Windows **command window**, enter the command **ipconfig -all**. The domain is listed together with other parameters of the computer. Leave the field empty if the computer is not located within a domain.

Entry	max. 32 characters
Default value	empty

User

User name of the user authorized to access the shared network printer.

Entry	max. 32 characters
Default value	empty

Password

Password of the user configured on the computer.

Entry	max. 32 characters
Default value	empty

10.6.3 More options

Paper size

Selection of the paper size.



NOTICE

The paper size selected will be ignored when the report is printed on a **CUSTOM NEOS** printer.

Selection	A4 (210 mm x 297 mm) Letter (216 mm x 279 mm)
Default value	A4 (210 mm x 297 mm)

Color

on | off (Default value: **off**)

If this option is activated, then the report will be printed out in color.


Spooler

on | off (Default value: **on**)

If this option is activated, then you can continue to work while the printer is printing. If this option is deactivated, then the Coulometer will be blocked while the printout is being produced.

10.7 Balance

List of devices: **Balance ► Edit**

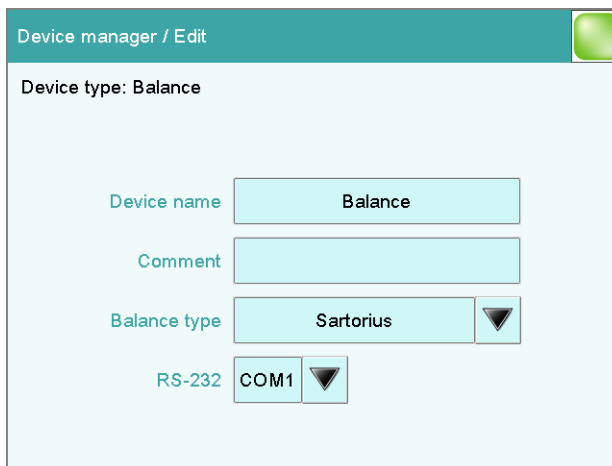

You define the type of balance and its connector in the dialog **Device manager / Edit**.

Device name

Here you can enter a device name of your choice.

Entry	24 characters maximum
-------	------------------------------

Comment

Entry	24 characters maximum
-------	------------------------------

Balance type

If you have connected a balance, then you have to specify the balance type here.

Selection	AND Mettler Mettler AT Mettler AX Ohaus Precisa Sartorius Shimadzu
Default value	Sartorius

RS-232

RS-232 interface to which the balance is connected. The interface parameters are adjusted in the **Edit device / Port parameters** dialog (see chapter 10.8, page 85).

Selection	COM 1 COM 2
Default value	COM 1

COM 2

The interface is disabled.

The following table indicates the balance type that needs to be selected for the balance model:

Balance	Balance type
AND	AND
Mettler AB, AE, AG, AM, AJ, PE, PM, PJ, PR, XP, XS	Mettler
Mettler AT	Mettler AT
Mettler AX, MX, UMX, PG, AB-S, PB-S	Mettler AX
Ohaus Voyager, Explorer, Analytical Plus	Ohaus
Precisa	Precisa
Sartorius	Sartorius
Shimadzu BX, BW	Shimadzu

10.8 USB/RS-232 adapter

Instrument list: **USB/RS-232 adapter** ► **Edit** ► **COM interface** ► **Edit**

As a rule, balances are equipped with a serial RS-232 interface. To connect a balance, you will require the 6.2148.050 cable. You can adjust the interface parameters in the dialog **Edit device / Port parameters**. These settings must match the settings on the connected device.



NOTICE

If you make changes to parameter settings in this dialog, then you must switch the Coulometer off and back on in order for the changes to take effect.

Edit device / Port parameters

Baud rate ▼

Data bits ▼

Parity ▼

Stop bits ▼

Handshake ▼

10.9 PC keyboard

Instrument list: **PC keyboard** ► **Edit**

A commercially available USB keyboard can be connected to make text and numerical input easier. It will be recognized automatically and entered in the list of devices with default settings.

The screenshot shows a dialog box titled 'Device manager / Edit'. Below the title bar, it says 'Device type: PC keyboard'. There are three input fields: 'Device name' with the value 'PC keyboard', 'Comment' which is empty, and 'Keyboard layout' with a dropdown menu showing 'English US' and a downward arrow.

Device name

Here you can enter a device name of your choice.

Entry **24 characters maximum**

Comment

Entry **24 characters maximum**

Keyboard layout

Define the country-specific keyboard layout here.

Selection **English US | German DE | French FR | Spanish ES | German CH**

Default value **English US**

To enter texts and numbers with the PC keyboard, the appropriate text and number input dialog must be opened on the Coulometer. Only the following keys on the PC keyboard have a function:

Table 6 Key assignment

Function in the editing dialog	PC keyboard key
[Cancel]	[Esc]


Function in the editing dialog	PC keyboard key
Enter respective character	Numbers, letters and special characters + Shift
[OK]	[Enter] / [↵]
[□]	[←] (backspace)
Line break (for multi-line text input)	[Enter] / [↵]
The cursor is moved up or down by one character.	[↑] [↓]
The cursor is moved to the left or to the right by one character.	[←] [→]
	[F1]
	[F2]
	[F3]
	[F4]
	[F5]
	[F6]
	[F7]

**NOTICE**

The lettering of the USB keyboard may differ from above lettering, depending on the country-specific keyboard used.

10.10 Barcode reader

List of devices: **Barcode reader** ► **Edit**

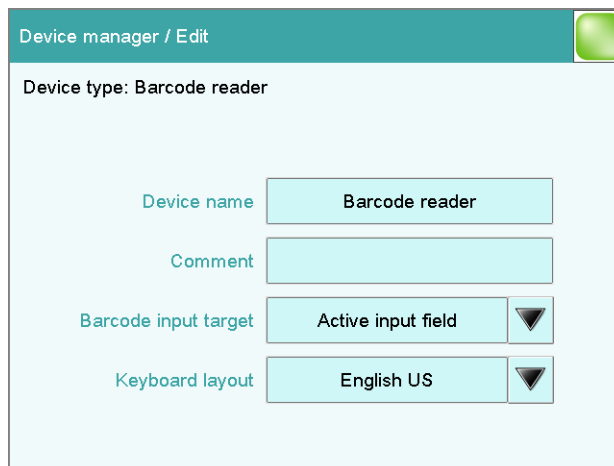

A barcode reader can be connected to read in sample data or other texts. A connected barcode reader will be recognized automatically and entered in the list of devices with default settings.

You will hear an acoustic signal as confirmation that a character string has been transmitted by the barcode reader and accepted.



NOTICE

The data will be read only if the system is in normal status, i.e. when no determination is running.

Device name

Here you can enter a device name of your choice.

Entry **24 characters maximum**

Comment

Entry **24 characters maximum**

Barcode input target

Selection of the input field for the character string read in by the barcode reader.

Selection **Active input field | Method | Identification 1 | Identification 2 | Sample size**

Active input field

The character string is entered in the input field of the opened text- or number-input dialog.

Method

The character string is entered in the input field **Method**.

**Identification 1**

The character string is entered in the input field **Identification 1**.

Identification 2

The character string is entered in the input field **Identification 2**.

Sample size

The character string is entered in the input field **Sample size**. Character strings containing characters other than numbers and decimal separators will be ignored.

Keyboard layout

Specify the country-specific keyboard layout for the emulation of the PC keyboard. This setting must match the setting on the barcode reader (see documentation for the barcode reader).

Selection	English US German DE French FR Spanish ES German CH
Default value	English US

11 File manager

Main dialog: **System ► File manager**

The saved methods, determinations, sample tables, etc. are managed in the file manager. You can also create a backup of your system (all data and settings). Also, an existing backup can be reloaded.

The file memory is organized as follows:

- **Internal memory**

The following files can be stored in the internal memory:

- Methods

- **External memory**

You can use a USB flash drive as an auxiliary storage medium, for instance. The following files can be stored on an external memory:

- Backup
- Methods
- Determinations
- Sample tables
- Result tables
- Reports as *.pdf
- PC/LIMS report

11.1 Managing files

File manager: **Internal memory / External memory 1 / External memory 2**

The saved files can be organized into groups. These groups are comparable to file directories on your PC, although unlike your PC, only one level is possible.




NOTICE

The file names must be unique for each memory location, i.e. you cannot save two files with the same name, even if they are in different groups.



NOTICE

If you use an external storage medium with the FAT or FAT32 file system, then you can save a maximum of 999 files per group. If you find it necessary to store more than 999 files in a single group, then you must reformat the storage medium with the file system **ExFAT** (see chapter 31.7.2, page 319).

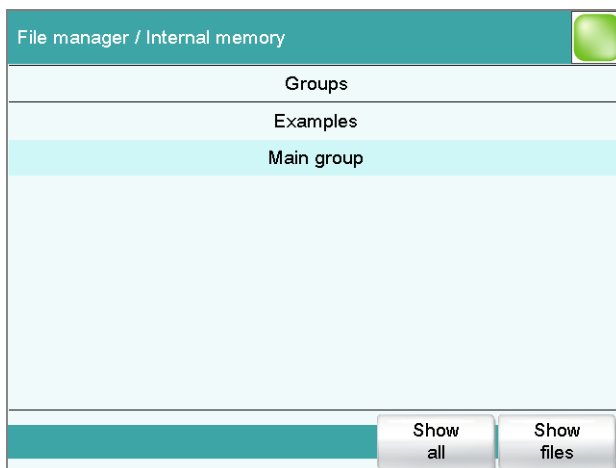


[Show all]

Display the list of all the files in the selected memory location.

[Show files]

Display the list of all the files in the selected group.

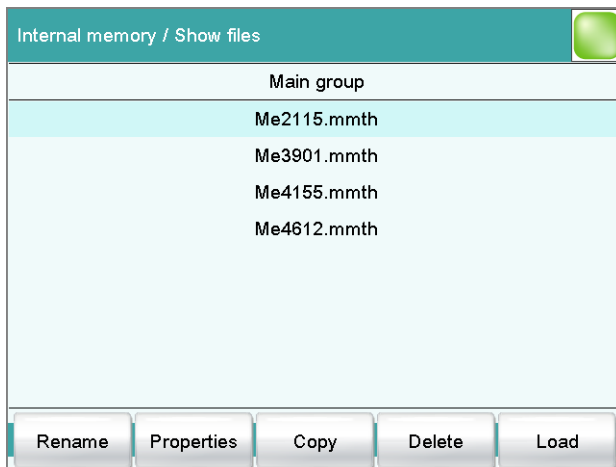

[Rename]

Rename the selected file (see chapter 11.1.2, page 93).

[Properties]

Display the file properties (see chapter 11.1.3, page 94).

[Copy]

Copy the selected file to a different memory location (*see chapter 11.1.1, page 93*).

[Delete]

Delete the selected file.

[Load]

Load the selected file.

11.1.1 Copying a file

Proceed as follows to copy a file:

1 Selecting the file

- Select the desired file.
- Tap on **[Copy]**.

The selection of memory locations is displayed.

2 Copy the file

- Tap on the desired memory location.
Only memory locations currently being accessible are active.

The file is copied. The file group is retained, i.e. the group will be created again in the new memory location if it does not already exist there.

11.1.2 Renaming a file

Proceed as follows to rename a file:

1 Selecting the file

- Select the desired file.
- Tap on **[Rename]**.


The **Show files / Rename** dialog is displayed.

2 Change the file name

- Tap on the input field **File name**.
The text editor opens.
- Enter a new file name (max. 32 characters) and confirm with **[OK]**.
- Tap on **[OK]**.

The file is saved under the new name.

11.2 External storage medium

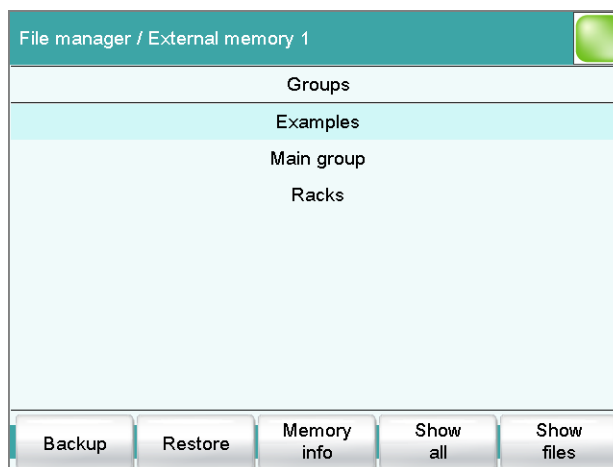

[Backup]

Creating a backup of all data and settings on this storage medium (see chapter 11.3, page 97).

[Restore]

Load the backup. This function is active only if a backup is available (see chapter 11.3, page 97).

[Memory info]

Display information on the storage medium, see following chapter.

Storage medium – Information

Detailed information about the storage medium is displayed in the dialog **External memory 1/2 / Memory info**:

- **Name**
Designation of the storage medium.
- **Write protection**
If this option is activated, then the file cannot be saved, deleted, moved or renamed. This is only an internal write protection feature and is independent of the write protection of the operating system on your computer.
- **Storage capacity**
Total capacity of the storage medium.
- **Used memory**
Storage capacity used on the storage medium.
- **Free memory**
Free storage capacity on the storage medium.

Directory structure

The directory structure appears as follows on the external storage medium:

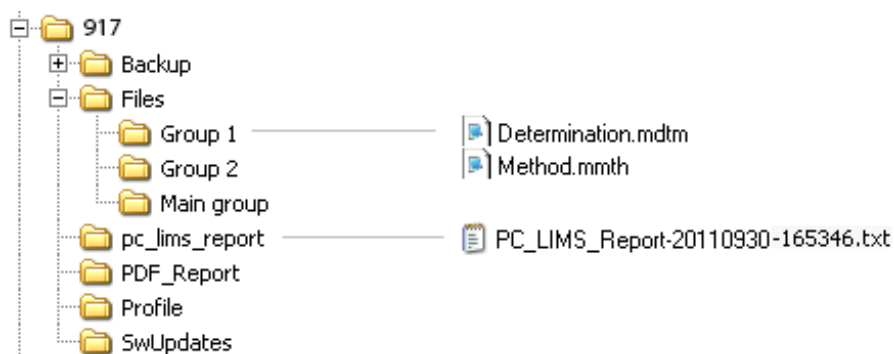


Figure 16 Directory structure on the external storage medium

- | | |
|-----------------------|--|
| Backup | All of the files of the backup are stored in this directory. The directory will be created the first time a backup is created. |
| Files | This directory contains all of the groups as well as the files contained in them (methods, determinations, etc.). |
| pc_lims_report | PC/LIMS reports are stored in this directory as TXT files. The directory will be created the first time a PC/LIMS report is printed out. |
| PDF_Report | PDF reports are stored in this directory. The directory will be created the first time a report is saved as a PDF file. |
| Profiles | The identification profile is stored in this directory. The directory will be created the first time an identification profile is generated. |
| SwUpdates | Files for software updates must be saved in this directory. |

11.3 Creating backups / Restoring data

File manager: **External memory 1 / External memory 2**

You can use the **[Backup]** function to easily create a backup containing all the data and settings of your system. You should make a backup at regular intervals in order to avoid data loss.



NOTICE

Only **one** backup can be created on a storage medium.

If a backup has already been stored on the medium, then it will be overwritten when this function is carried out again.

11.3.1 Restoring data

You can use the function **[Restore]** to restore either a complete backup or only certain data.



CAUTION

All of the methods in the internal memory will be deleted.

Backups are not backward compatible, i.e. backups of the latest version cannot be restored using earlier versions.

Proceed as follows to restore backed-up data:

1 Select the storage medium

- Connect the storage medium on which the backup is located.
- Select the storage medium under **System ► File manager**.

2 Select data

- Tap on **[Backup]**.
The selection of the data that can be restored separately is displayed (see the following).
- Deselect the data sets which are not to be restored.
- Tap on **[Load data]**.

The selected data blocks are restored.

3 Complete the restoration

- Switch the Coulometer off and on again.



Restorable data blocks

The following data blocks can be loaded individually:

- **Methods**
All of the methods stored in the internal memory.
- **Sample table**
Current sample table.
- **Result table**
Current result table.
- **Current determination data**
All the data for the current determination (including the method with which the determination was carried out).
- **Control**
Settings in the control dialog.
- **User list**
Settings for each user under **System settings / User administration**.
- **System settings / User admin.**
- **Titrants**
All titrants with all their data.
- **Sensors**
All sensors with all their data.
- **Device data**
All of the devices configured in the device manager, with all of the data.
- **GLP data**
All data from the GLP manager.
Deactivate if the backup is to be loaded onto a different system.
- **Common variables**
All common variables.
- **Templates**
- **Routine dialog settings**
Current routine dialog settings (see **Dialog options / Routine dialog**).
- **Sample racks**
The sample racks present in the properties of the USB sample processor.
- **Own subsequences**
All subsequences created and stored by the user.

12 GLP manager

Main dialog: **System ► GLP manager**

In the GLP manager you can document data for various GLP tests. The results of the automatic system tests carried out after switching on are also documented.

The following tests can be documented:

- GLP test "Measurement" (*see chapter 12.3, page 101*)
- GLP test "Titration" (*see chapter 12.3, page 101*)
- System validation (*see chapter 12.4, page 105*)

In addition, you can also:

- Create a list with your test tools (*see chapter 12.2, page 100*).
- Define a service interval for having routine maintenance carried out by a Metrohm service technician (*see chapter 12.5, page 109*).
- Define an interval for the regular performance of backups (*see chapter 12.5, page 109*).

Further information about quality management and validation is also given in the documentation series **Quality Management with Metrohm** and **Application Bulletin AB 255** (Validation of Metrohm Karl Fischer titrators in accordance with GLP/ISO 9001).

System / GLP manager		
Test	Last test	Next test
Automatic system test	2010-08-30	Power on
System validation	2010-08-02	2011-01-04
GLP test "Measurement"	2010-07-01	2011-01-03
GLP test "Titration"	2010-07-02	2011-01-04

Test tools GLP tests HW/SW System validation Monitoring View test data

The table shows the last time that each test was carried out and when the next test is to be carried out. A test is entered in the list the first time that it is documented.

[Test tools]

Configuring the test tools for GLP tests (*see chapter 12.2, page 100*).

[Delete]

Delete the selected test tool from the list.

[Edit]

Change the designation of the selected test tool.

12.3 GLP tests for measurement and titration

Main dialog: **System ► GLP manager ► GLP tests HW/SW ► GLP test "Measurement" / GLP test "Titration"**

You can document the GLP tests for measurements and titrations in the dialog **GLP manager / GLP tests Hardware/Software** and its subdialogs. The following procedure describes the GLP test "Measurement," but also applies for the GLP test "Titration".

Proceed as follows:

1 Open the properties dialog

- In the **System / GLP manager** dialog, tap on the **[GLP tests HW/SW]** button and then on the **[GLP test "Measurement"]** button.

2 Edit data

- Define the test method, the test results, etc. (see "Dialogs "GLP tests HW/SW / GLP test "Measurement" and "GLP tests HW/SW / GLP test "Titration"", page 102).

3 Define the hardware used

- Tap on **[Hardware]**.



- Select the hardware used for the test (see "Dialogs "GLP test "Measurement" / Hardware" and "GLP test "Titration" / Hardware"", page 104).
- Tap on the fixed key [].

The dialog **GLP tests HW/SW / GLP test "Measurement"** appears again.

4 Define the test interval

- Tap on [**GLP test interval**].
- Enter the time interval or the date for the next GLP test (see "Dialogs "GLP test "Measurement" / Test interval" and "GLP test "Titration" / Test interval"", page 103).
- Tap on the fixed key [].

12.3.1 Parameter description

Dialogs "GLP tests HW/SW / GLP test "Measurement"" and "GLP tests HW/SW / GLP test "Titration""

Method

Method the GLP test has been carried out with.

Entry	32 characters maximum
Selection	Selection of methods stored in the internal memory

User

User who carried out the GLP test.

Entry	24 characters maximum
Selection	Selection of configured users

Test date

Date on which the GLP was carried out.

Format: YYYY:MM:DD

Test result

Result of the test.

Selection	Test OK Test not OK
Default value	Test OK

Comment

Entry	24 characters maximum
-------	------------------------------

[GLP test interval]

Defining the time interval for the GLP test (see "Dialogs "GLP test "Measurement" / Test interval" and "GLP test "Titration" / Test interval"", page 103).

[Hardware]

Documenting the hardware with which the GLP test was carried out (see "Dialogs "GLP test "Measurement" / Hardware" and "GLP test "Titration" / Hardware"", page 104).

Dialogs "GLP test "Measurement" / Test interval" and "GLP test "Titration" / Test interval"

Test date

Date on which the GLP was carried out.

Format: YYYY:MM:DD

Monitoring

on | off (Default value: **off**)

If this parameter is activated, then the time interval after which a GLP test has to be carried out again will be monitored.

GLP test interval

If you define a time interval for the GLP test, then the date in **Next GLP test** will be tracked automatically.

Input range	1 to 999 days
Default value	999 days

Next GLP test

If you define a date for the next GLP test, then the **GLP test interval** will be tracked automatically.

Format: YYYY:MM:DD

Action

Selection of the action which is carried out when the time interval has expired.

Selection	Display message Document message Cancel determination
Default value	Display message

For all three options it is documented in the determination data (see dialog **More determination data / Messages**), that the time interval has been expired.

12.4 System validation

Main dialog: **System ► GLP manager ► System validation**

In the dialog **GLP manager / System validation** and its subdialogs, you can document the system validation results and define the time interval after which validation must be carried out again.

Proceed as follows:

1 Open the properties dialog

- In the **System / GLP manager** dialog, tap on the **[System validation]** button.

The screenshot shows the 'GLP manager / System validation' dialog. It features a teal header bar with the title 'GLP manager / System validation' and a green circle icon on the right. Below the header, there are five input fields: 'Method' (containing 'SOP_SYSVAL_02'), 'User' (containing 'chk'), 'Test date' (containing '2011-07-01'), 'Test result' (containing 'Test OK'), and 'Comment' (empty). At the bottom of the dialog, there are three buttons: 'Note (SOP)', 'Validation interval', and 'Test statistics'.

2 Edit data

- Define the method, the result, etc. (see *"GLP manager / System validation" dialog*, page 106).

3 Define the validation interval

- Tap on **[Validation interval]**.
- Enter the time interval or the date for the next system validation (see *"Dialog "System validation / Validation interval"*, page 107).
- Tap on the fixed key [**↩**].

The dialog **GLP manager / System validation** appears again.

4 Enter statistical data

- Tap on **[Test statistics]**.
- Enter the statistical data for the last system validation (see *"System validation / Test statistics" dialog*, page 108).
- Tap on the fixed key [**↩**].

The dialog **GLP manager / System validation** appears again.

5 Insert a note

- Tap on **[Note (SOP)]**.
- Enter a brief description, e.g. a summary of the SOP (standard operating procedure) according to which the system validation was carried out (see "*System validation / Note (SOP) dialog*", page 107).
- Tap on the fixed key [**↩**].

12.4.1 Parameter description

"GLP manager / System validation" dialog

Method

Method with which the system validation has been carried out.

Entry	32 characters maximum
Selection	Selection of methods stored in the internal memory

User

User who carried out the system validation.

Entry	24 characters maximum
Selection	Selection of configured users

Test date

Date on which the system validation was carried out.

Format: YYYY:MM:DD

Test result

Result of the test.

Selection	Test OK Test not OK
Default value	Test OK

Comment

Entry	24 characters maximum
-------	------------------------------

[Note (SOP)]

Enter a brief description, e.g. a summary of the SOP (standard operating procedure) according to which the system validation has been carried out.

[Validation interval]

Defining the time interval for the system validation (see "*Dialog "System validation / Validation interval"*", page 107).

[Test statistics]

Documenting the statistical data of the system validation (see *"System validation / Test statistics" dialog*, page 108).

"System validation / Note (SOP)" dialog

In this dialog, you can enter a brief text, e.g. a summary of the SOP (standard operating procedure) according to which the system validation was carried out.

Dialog "System validation / Validation interval"**Last validation**

Date on which the last system validation was carried out.

Format: YYYY:MM:DD

Monitoring

on | off (Default value: **off**)

If this parameter is activated, then the time interval after which a system validation has to be carried out again will be monitored.

Validation interval

If you define a time interval for the system validation, then the date in **Next validation** will be tracked automatically.

Input range	1 to 999 days
Default value	999 days

Next validation

If you define a time interval for the next system validation, then the **Validation interval** will be tracked automatically.

Format: YYYY:MM:DD

Action

Selection of the action which is carried out when the time interval has expired.

Selection	Display message Document message Cancel determination
Default value	Display message

For all three options it is documented in the determination data (see dialog **More determination data / Messages**), that the time interval has been expired.

**Display message**

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message

In the determination data it will be documented, that the time interval has been expired.

Cancel determination

The determination is stopped.

"System validation / Test statistics" dialog

In the dialog **System validation / Test statistics**, you can document the statistical data for the last system validation.

Number (n)

Number of determinations carried out.

Input range	1 to 999999999
Default value	empty

Mean value

Mean value out of single results.

Input range	-999999999.00000 to 999999999.00000
Default value	empty

s abs

Absolute standard deviation of the results.

Input range	0.00000 to 999999999.00000
Default value	empty

s rel

Relative standard deviation of the results.

Input range	0.00000 to 100.00000 %
Default value	empty %

d rel

Systematic deviation of the results.

Input range	0.00000 to 100.00000 %
Default value	empty %

a sys

Systematic error.

Input range	0.00000 to 999999999.00000
Default value	empty

12.5 System monitoring

12.5.1 Service interval

Main dialog: **System ► GLP manager ► Monitoring ► Service interval**

In the dialog **GLP manager / Service interval**, you can define the time interval for the maintenance of the system by the Metrohm Service department. The service interval is checked each time the system is started.

Last service

Date on which the last servicing was carried out.

Format: YYYY:MM:DD

Monitoring

on | off (Default value: **off**)

If this parameter is activated, then the time interval after which system maintenance must be carried out again will be monitored.

Service interval

If you define a time interval for the system maintenance, then the date in **Next service** will be tracked automatically.

Input range	1 to 999 days
Default value	999 days

Next service

If you specify a date for the next system maintenance, then the **Service interval** will be tracked automatically.

Format: YYYY:MM:DD

12.5.2 Backup interval

Main dialog: **System ► GLP manager ► Monitoring ► Backup interval**

In the dialog **GLP manager / Backup interval**, you can define the time interval for backups. The backup interval is checked each time the system is started.

Last backup

Date on which the last backup was created.

Format: YYYY:MM:DD



Monitoring

on | off (Default value: **off**)

If this parameter is activated, then the time interval after which a backup has to be created again will be monitored.

Backup interval

If you define a time interval for the creation of backups, then the date in **Next backup** will be tracked automatically.

Input range	1 to 999 days
Default value	999 days

Next backup

If you specify a date for the next backup, then the **Backup interval** will be tracked automatically.

Format: YYYY:MM:DD

13 Common variables

Main dialog: **System ► Common variables**

You can save 25 **method-independent variables**, or common variables. These variables can be used in future calculations (as variables **CV01... CV25**). Common variables are useful, e.g. for the following applications:

- Determination of a blank value which will be taken into account during the content determination of the sample.
- Determination of the content of a standard solution, which will be taken into account during the content determination of the sample.

System / Common variables		
CV	Name	Value
01	Blank value	0.0143 mL
02
03
04
05	Density	0.986 g/mL
06
07
08

Delete Edit

The designation and the value (including the unit) are displayed for every common variable in the list.

[Delete]

Delete the selected common variable from the list.

[Edit]

Edit the data of the selected common variable, see following chapter.

13.1 Editing common variables

List of common variables: **Common Variable ► Edit**

The common variables can be modified as follows:

- Edit manually, see the following.
- Automatic assignment from the determination run. A calculation result must be configured accordingly for this purpose (*see chapter 13.4, page 115*).

Name

Designation of the common variable.

Entry	24 characters maximum
Default value	empty

Value

Value of the common variable.

Input range	-999999999 to 999999999
Default value	empty

Unit of the common variable.

Entry	10 characters maximum
Default value	empty
Selection	% mol/L mmol/L g/L mg/L mg/mL mg/100 g ppm g mg µg mL µL mg/piece °C µg/min mL/min µL/min
Default value	%

Next assignment

This parameter is displayed only if a validity has been defined for the common variable.

Date on which the validity of the common variable expires.

Format: YYYY:MM:DD

[Validity]

Defining the time interval for the validity of the common variable (see chapter 13.3, page 114).

[Properties]

Display the properties of the common variable, see following chapter.

13.2 Properties of common variables

Common variable: **Edit ► Properties**

Detailed information concerning the common variable is displayed in the **Edit common variables / Properties** dialog:

- **Status**
Status of the common variable. If the time interval for the validity has expired, then **invalid** will be displayed.
- **Method**
Method with which the value has been assigned to the common variable. If the value was entered manually, then **manual** will be displayed.
- **Method status** (only for automatic assignment from the determination run)
- **Determination status** (only for automatic assignment from the determination run)
- **Last assignment**
Date and time of the last assignment.
- **User**
User who assigned the value to the common variable.

13.3 Monitoring validity

Common variable: **Edit ► Validity**

In the dialog **Edit common variables / Validity**, you can define the time interval after which a new value must be assigned to the common variable.

Last assignment

Date on which the common variable was last assigned a value.

Format: YYYY:MM:DD

Monitoring

on | off (Default value: **off**)

If this parameter is activated, then the time interval after which the common variable must be assigned a new value will be monitored.

Validity

If you define a time interval for the validity of the common variable, then the date in **Next assignment** will be tracked automatically.

Input range	1 to 999 days
Default value	999 days

Next assignment

If you specify a date for the next assignment, then the time interval for the **Validity** will be tracked automatically.

Format: YYYY:MM:DD

Action

Selection of the action which is carried out when the time interval has expired.

Selection	Display message Document message Cancel determination
Default value	Display message

For all three options it is documented in the determination data (see dialog **More determination data / Messages**), that the time interval has been expired.

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message

In the determination data it will be documented, that the time interval has been expired.

Cancel determination

The determination is stopped.

13.4 Assigning a result automatically to a common variable

**NOTICE**

This instruction is based on the assumption that the method contains a calculation command with a calculation.

Proceed as follows to assign a result to a common variable:

1 Opening the editing dialog of the result

- In the command list, select the command **CALC**.
- Tap on the **[Edit command]** button.
- Select the result whose value is to be assigned to a common variable and tap on **[Edit]**.

The editing dialog of the result is displayed.



2 Define result options

- Tap on the **[Result options]** button.

- Activate the parameter **Save as common variable**.
- Define **Variable = CV01...CV25**.



3 Saving the settings

Tap on the fixed keys [] or [].

In the future, the calculated result will be assigned to the selected common variable (result name, value and unit).

14 Templates

Main dialog: **System ► Templates**


You have the option of defining **system-specific templates**. You can use these templates when editing the respective data.

You can create the following templates:

- **Sample data** (see chapter 14.1, page 117)
Create sample identifications and sample assignments.
- **Custom result templates** (see chapter 14.2, page 121)
Define formulas for result calculations.
- **Input lines** (see chapter 14.3, page 125)
Define input signals on the remote interface.
- **Output lines** (see chapter 14.4, page 127)
Defining output signals on the remote interface.
- **Report header** (see chapter 14.5, page 130)
Create a system-specific report header that contains information concerning the laboratory, for example.

14.1 Sample data

Main dialog: **System ► Templates ► Sample data**

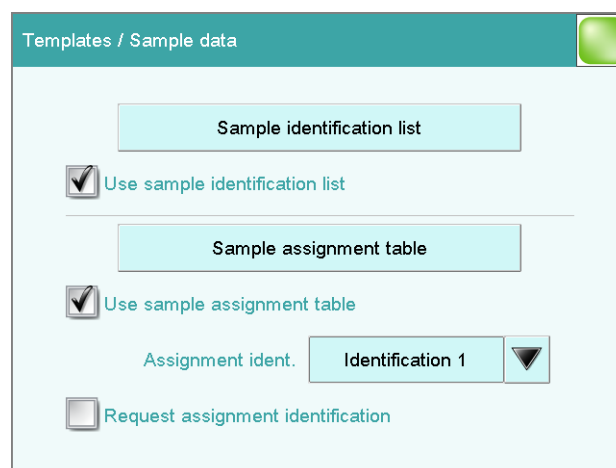

You can create the following templates in the dialog **Templates / Sample data**:

- Sample identifications (see chapter 14.1.1, page 118)
List with sample identifications. If you must enter the sample identification before starting a determination, then you can select the entries contained in this list.

- Sample assignments (*see chapter 14.1.2, page 119*)
You can assign a particular method to a sample identification.

Use sample identification list

on | off (Default value: **off**)

Activate this parameter so that the sample identifications defined in the list will be displayed as selections.

Use sample assignment table

on | off (Default value: **off**)

Activating/deactivating the use of the sample assignment table.



NOTICE

If this parameter is activated, the following will no longer be possible:

- Carrying out determinations without their sample identification being defined in the sample assignment table.
- Defining a method in the sample table. Previously defined methods will be ignored.

Assignment ident.

Identification which is used as assignment identification to load the correct method.

Selection	Identification 1 Identification 2
Default value	Identification 1


Request assignment identification

on | off (Default value: **off**)

If this parameter is activated, then the sample identification will be requested automatically at the start of a determination.

14.1.1 Sample identification list

Main dialog: **System ▶ Templates ▶ Sample data ▶ Sample identification list**

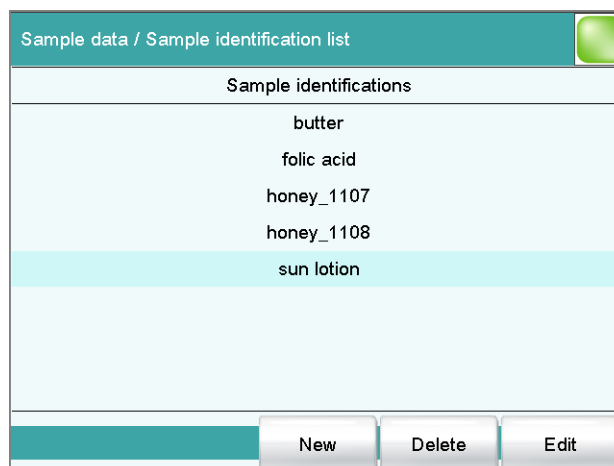

You can create a system-specific list with sample identifications in the dialog **Sample data / Sample identification list**. This list makes it easier for you to enter frequently used sample identifications, i.e. this list is available to you for entering them (main dialog, command **REQUEST**, etc.). It may be a good idea to define the unchanging part of the identification as a template and to add the variable part during sample data input.

[New]

Add a new sample identification to the list.

[Delete]


Delete the selected sample identification from the list.

[Edit]

Change the designation of the selected sample identification.

14.1.2 Sample assignment table


Main dialog: **System ► Templates ► Sample data ► Sample assignment table**

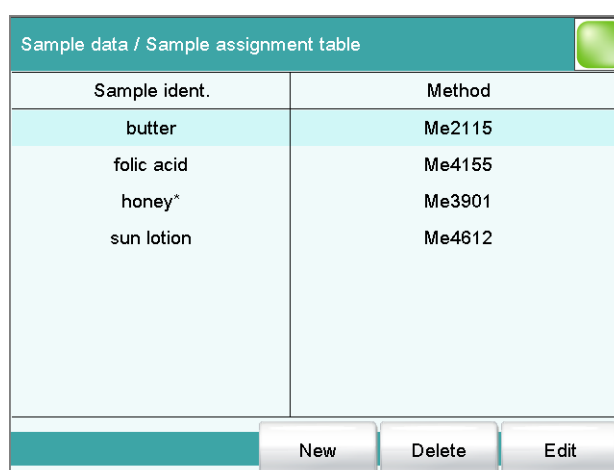

Entry	32 characters maximum
Selection	Selection of methods stored in the selected memory

14.2 Custom result templates

Main dialog: **System ► Templates ► Custom result templates**

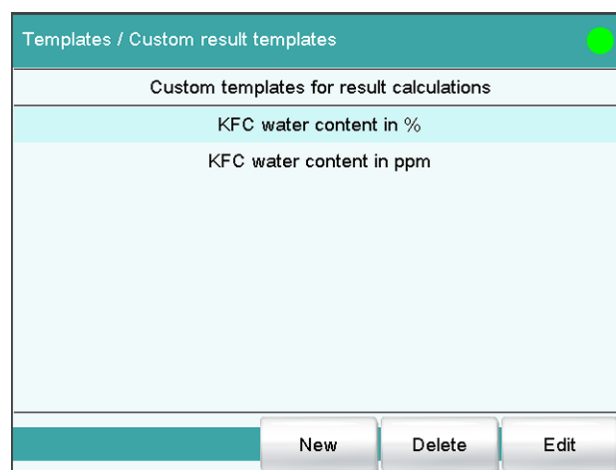

In the dialog **Templates / Custom result templates**, you can define the result calculations which can be loaded in the calculation command **CALC**. With the exception of the following points, creating a result template is identical to editing a calculation in the command **CALC** (see chapter 28.5.1.2, page 259):

- No result variable can be assigned to the template.
- The definition of limit values is not possible.



NOTICE

Up to nine **wildcards** can be inserted in the calculation formula, the **Variables F1...F9**. Use these variables, e.g. for the molar mass of your analyte. If you have loaded a result template with these wildcards in the calculation command, then you will automatically be prompted to enter the numerical values.

[New]

Add a new result template to the list, see following chapter.

[Delete]

Delete the selected result template from the list.

[Edit]

Edit the data of the selected result template, see following chapter.

Variable for mean value

If the statistic calculations have been activated (see method options), the mean value of the single results will be saved as variable SMN1 to SMN9. As default value, always the first free variable is being displayed.

Selection	SMN1 ... SMN9 off
-----------	----------------------------

off

For the result not statistic calculations are carried out.

Save as titer

The result can be saved as titer for the selected titrant.



NOTICE

The titer is assigned to the titrant defined in the last titration command before the calculation. Please note that the CALC command containing the titer assignment is inserted after the titration command with which the titer is determined.

Selection	Single value Mean value off
Default value	off

Single value

The result of the current determination is saved as titer.

Mean value

If statistic calculations are carried out for the result, the current mean value of the determination series is saved.

off

The result is not saved as titer.

Save as common variable

on | off (Default value: **off**)

The calculated result can be saved as a method-independent variable, a so-called common variable. The result is then also available in other methods for calculations. In all cases, only the single value will be saved, even if the statistics function is enabled.

Variable

Selection of the common variable to which the result is being assigned.

Selection	CV01 ... CV25
-----------	----------------------



Display result

on | off (Default value: **on**)

If you deactivate this parameter, the result is neither displayed in the result dialog nor printed in the result report. This can be advisable for intermediate results.

Save result in result table

on | off (Default value: **off**)

The calculated result can be saved in the result table. This may be advisable if e.g. the results of all determinations carried out on a particular day are to be displayed clearly. A maximum of nine results from a determination can be saved in the result table.

Precision

Setting, with which accuracy the result is used in additional calculations.

Selection	Round Truncate Full precision
Default value	Round

Round

The result is rounded to the defined number of decimal places (commercial rounding, in accordance with the US Pharmacopeia USP). If the digit at the first dropped decimal place is **1, 2, 3 or 4**, then it will be rounded off; if this digit is **5, 6, 7, 8 or 9**, then it will be rounded up. Negative digits will be rounded in accordance with their amount, i.e. away from zero.

Truncate


The result is cut to the number of decimal places defined.

Full precision

The result is used with full accuracy (floating point number either in "single precision" (32 bit) or in "double precision" (64 bit), according to the standard IEEE 754).

14.3 Input lines

Main dialog: **System ► Templates ► Input lines**

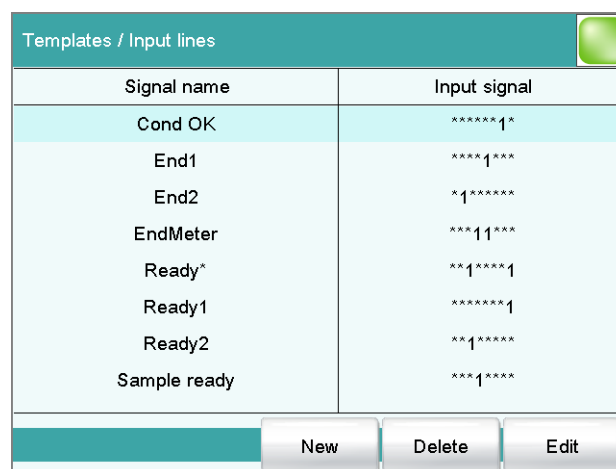

Signal name	Input signal
Cond OK	*****1*
End1	****1***
End2	*1*****
EndMeter	***11***
Ready*	**1****1
Ready1	*****1
Ready2	**1*****
Sample ready	***1****

In the dialog **Templates / Input lines**, you can define the input signals at the remote interface as a template. You can select these templates in the command **SCAN**. The list can contain a maximum of 20 templates.

Table 7 List of predefined input signals

Signal name	Input signal	Function
Cond OK	*****1*	queries "Cond OK" condition of device.
End1	****1***	waits for the EOD pulse from Device 1 (Titrimo, Titrimo).
End2	*1*****	waits for the EOD pulse from Device 2.
EndMeter	***11***	waits for the EOD pulse from 780/781 pH/Ion meter (stirrer 1 will be switched on during the waiting period).
Ready*	**1****1	queries the "Ready" condition of Devices 1 and 2 (Titrimo, Titrimo). The status of devices working in parallel can be scanned with this signal. In this case the "Ready" line of both devices must be statically set (i.e. constantly) at the end of a determination. Devices which only transmit a brief impulse cannot be controlled in parallel.
Ready1	*****1	queries the "Ready" condition of Device 1.
Ready2	**1*****	queries the "Ready" condition of Device 2.

Signal name	Input signal	Function
Sample ready	***1***	waits for a stepping pulse, of, for instance, a connected Sample Processor, as soon as it is ready.

[New]

Add a new input signal to the list, see following chapter.

[Delete]

Delete the selected input signal from the list.

[Edit]

Edit the data of the selected input signal, see following chapter.

14.3.1 Editing the input signal

List of input signals: **Input signal ► New / Edit**

Signal name

Designation of the template.

Entry	24 characters maximum
-------	------------------------------

Input signal

Entering the desired bit pattern:

Entering the bit pattern:

- 0 = line inactive
- 1 = line active
- * = retain line status

The input lines are always numbered from right to left, i.e. with the signal *******1** line 0 is expected to be active.




NOTICE

We recommend masking lines that are of no interest or for which no defined condition can be predicted with an asterisk (*).

Entry	Bit pattern of exactly 8 characters
Default value	*****

14.4 Output lines

Main dialog: **System ► Templates ► Output lines**

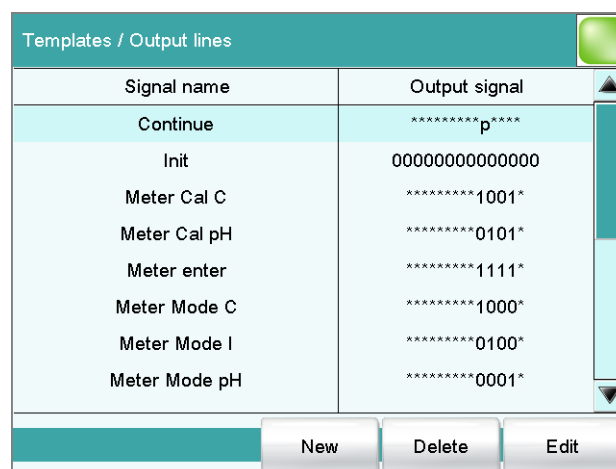

In the dialog **Templates / Output lines**, you can define the output signals at the remote interface as a template. You can select these templates in the command **CTRL**. The list can contain a maximum of 20 templates.

Table 8 List of predefined output signals

Signal name	Output signal	Function
Continue	*****p****	sends a stepping pulse to the connected Sample Processor.
Init	00000000000000	initializes the remote interface.
Meter Cal C	*****1001*	switches the 781 pH/Ion meter to concentration calibration.
Meter Cal pH	*****0101*	switches the 780/781 pH/Ion meter to pH calibration and starts the calibration.
Meter enter	*****1111*	simulates the [Enter] key of the 780/781 pH/Ion meter; mandatory for pH calibration in order to start the measurement of the second buffer.
Meter Mode C	*****1000*	switches the 781 pH/Ion meter to concentration measurement.
Meter Mode I	*****0100*	switches the 780/781 pH/Ion meter to voltametric measurement with polarization current and starts the measurement.
Meter Mode pH	*****0001*	switches the 780/781 pH/Ion meter to pH measurement and starts the measurement.

14.4.1 Editing the output signal

List of input signals: **Output signal** ► **New / Edit**

Signal name

Designation of the template.

Entry	24 characters maximum
-------	------------------------------

Output signal

Entering the desired bit pattern:

Entering the bit pattern:

- 0 = line inactive
- 1 = line active
- * = retain line status
- p = set pulse

The output lines are always numbered from right to left, i.e. with the signal *******1** line 0 is set.



NOTICE

We recommend masking lines that are of no interest or for which no defined condition can be predicted with an asterisk (*).

Entry	Bit pattern of exactly 14 characters
Default value	*****

Pulse length

Duration of the sent pulses.

Input range	100 to 1000 ms
Default value	200 ms

14.5 Report header

Main dialog: **System ► Templates ► Report header**

In the dialog **Templates / Report header**, you can create a system-specific report header that contains information concerning the laboratory, for example. The report header can contain a maximum of four lines of 46 characters each and will always be printed out in front of the standard report header. The standard report header contains the instrument type, serial number and program version of the software, instrument name and printing date.

Print logo

on | off (Default value: **off**)

If this parameter is activated, then the logo defined below will be printed out in the right margin.

Logo

Selection	Metrohm logo Custom logo
Default value	Metrohm logo

Custom logo

Select this setting if your logo is to be printed out in the report header. How to create a logo is described below.

Applying the custom logo

You can create a graphics file of your own and use it as a logo in the report header in place of the Metrohm logo.



NOTICE

Observe the size of the graphics object. The maximum size of your graphics file is 64 kB. We recommend a maximum of 500 × 200 pixels.

Proceed as follows to create your own logo:

1 Creating the logo

- Create a graphics file with the desired logo in an image processing program.
- Save the graph as a JPG file under the file name "CustomImage.jpg".

2 Copying the file to an external storage medium

- Copy the file "CustomImage.jpg" to the directory "917".
If this structure is not maintained, the file will not be found because the software directly accesses the "917" directory.
- Plug in the external storage medium at the Coulometer.

3 Loading the graphics file

- In the **Templates / Report header** dialog, tap on the **[Load]** button.
The file is copied to the internal memory.
- Switch the instrument off and back on again.

4 Applying the custom logo

- In the **Templates / Report header** dialog, activate the **Print logo** option.
- Define **Logo = Custom logo**.

Your logo will now be printed in the report header of all reports in the future.

15.2 Saving a method

If you modify method parameters, then you can save these as your own method.

To save a method, proceed as follows:

1 Open the command sequence

- In the main dialog, tap on **[Edit parameters]**.

The command sequence of the method is displayed:

Parameters / Sequence		
Current method: New method		
01	WAIT	Wait
02	REQUEST	Data request
03	KFC	Karl Fischer coulometry
04	CALC	Calculation
05	REPORT	Report
06	...	

Buttons: Save method, Method options, Insert command, Delete command, Edit command

2 Modifying/applying the method name

- Tap on **[Save method]**.
A method name will be suggested for new methods. If the method has already been saved once, then the method name will be displayed.

Sequence / Save method	
Memory	Internal memory ▼
Group	Main group ▼
File name	Me2115 ▼

Buttons: Cancel, Save

Applying the name:

- Tap on **[Save]**.

The method will be saved and the command sequence is displayed.

Entering a new name:

- Tap on the input field **File name**.
The text editor opens.
- Enter a new file name (max. 32 characters) and confirm with **[OK]**.
- Tap on **[Save]**.

The method will be saved and the command sequence is displayed.


15.3 Loading a method

Proceed as follows to load a method:

1 Opening the method table

- In the main dialog, tap on **[Load method]**.
The selection of memory locations is displayed. This selection will be skipped if only the internal memory is accessible.
- Select the memory location where the desired method is saved.
The selection of file groups is displayed (*see chapter 11.1, page 91*). If only one group is available, then the method table will be displayed directly.
- Select the group with the desired method.
- Tap on **[Show files]**.

The method table with the stored methods opens:

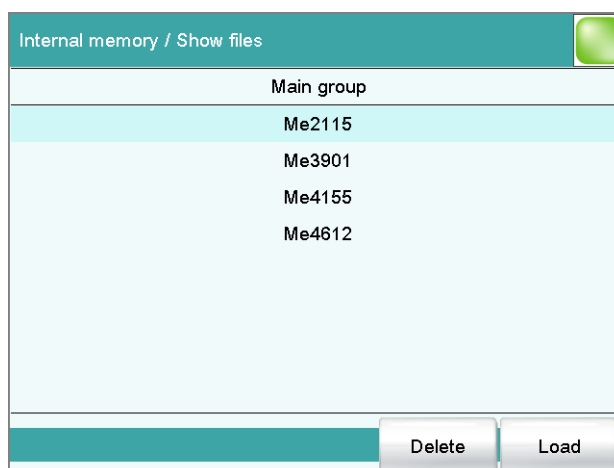

2 Selecting the method

- Select the desired method.

3 Loading the method

- Tap on **[Load]**.

The method is now loaded. The method name is displayed in the main dialog in the title line.



NOTICE

The data of the current determination will be deleted when you load a method.

15.4 Editing a method

Main dialog: **Edit parameters**

Parameters / Sequence		
Current method: New method		
01	WAIT	Wait
02	REQUEST	Data request
03	KFC	Karl Fischer coulometry
04	CALC	Calculation
05	REPORT	Report
06	...	

Save method Method options Insert command Delete command Edit command

In the dialog **Parameters / Sequence**, you can define and edit the method commands which are executed successively during a determination.

The following data is displayed in the command list for each command:

- Line number
- Name
- Comment

You can adjust the comment in accordance with your requirements.

[Save method]

Save the current method (*see chapter 15.2, page 133*).

[Method options]

Define various settings which refer to the entire method, e.g. statistics, automatic saving of determinations, settings for sample data, etc. (*see chapter 15.5, page 137*).

[Insert command]

Insert a new method command. It is inserted before the selected command.

**NOTICE**

Not all commands are available for subsequences. Commands which cannot be inserted into subsequences are disabled.

[Delete command]

Delete the selected method command.

[Edit command]

Edit the selected method command.

15.4.1 Inserting a command

A method is comprised of individual commands. When you start a determination, the commands are executed one after the other.

The following constraints apply:


- Maximum of 99 commands

Inserting a command

To insert a command in a method, proceed as follows:

1 Open the command overview

- In the main dialog, tap on **[Edit parameters]**.
- Tap on the **[Insert command]** button.

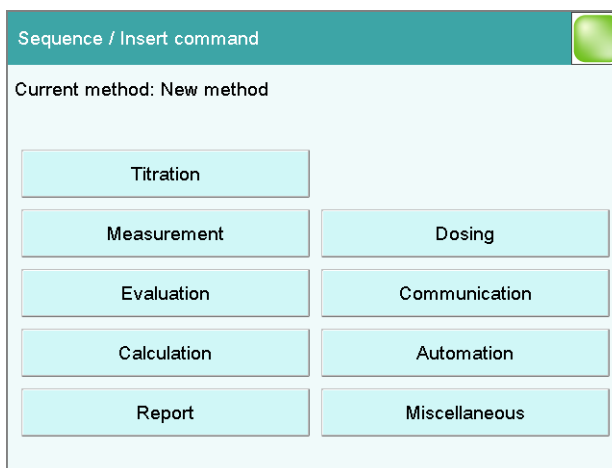

This dialog contains all of the commands which can be inserted into a method run, organized into thematic groups.

2 Select a command group

- Tap on the desired command group.

In the case of **[Report]**, the command is inserted directly into the command list. For all other command groups, the available commands will be displayed (titration / measuring commands, etc.).


3 Inserting a command

Tap on the desired command.

The dialog is closed and the command is displayed in the command list.

15.5 Method options

Main dialog: **Edit parameters ► Method options**

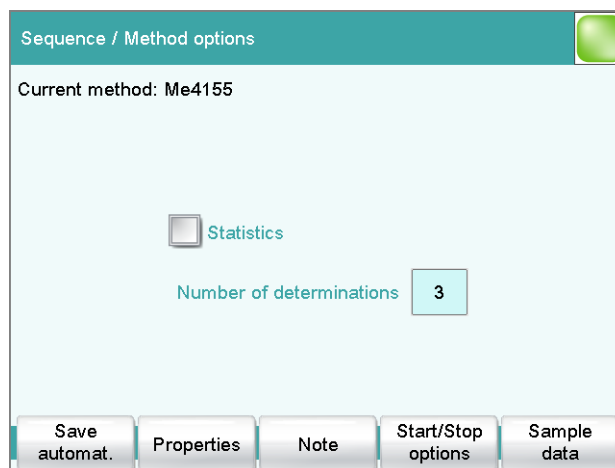

Method options are settings which apply to the method as a whole and not to an individual command.

The following settings are available:

- Activate/deactivate statistics calculations
- Save determination automatically
- Sign a method electronically
- Check the method when the determination starts
- Define a fixed sample size
- etc.

Statistics

on | off (Default value: **off**)

If this parameter is activated, then statistics calculations will be carried out for all of the defined results. The prerequisite here, however, is that a vari-

able for the mean value is defined in the result options (see "Dialog "Edit calculation / Result options"", page 261).

Number of determinations

The number of determinations that are carried out for the statistics calculations.

If an additional determination has to be added to the determination series, because one determination has been incorrect, for example, then this can be accomplished in the statistical overview (see chapter 24, page 194).

Input range	2 to 20
Default value	3

[Save automat.]

Define the settings for automatically saving the determination (see chapter 15.5.6, page 146).

[Properties]

Display the properties of the current method and sign electronically (see chapter 15.5.4, page 144).

[Note]

Enter a note on the method (see chapter 15.5.5, page 146).

[Start/Stop options]

Edit functions which are carried out immediately after the method is started or stopped, see chapter Start options, page 138 and chapter Stop options, page 140.

[Sample data]

Specify method-specific settings for the sample data, e.g. designation of the sample identifications, limits for the sample size (see chapter 15.5.3, page 141).

15.5.1 Start options

Main dialog: **Edit parameters ► Method options ► Start/Stop options ► Start options**

In the dialog **Method options / Start options**, you can configure the functions which are carried out when a determination is started.

Method check at start

on | off (Default value: **on**)

If this option is activated, then the following tests are carried out at the method start.

- Are the required control and peripheral devices connected?

- Are the titrant and the sensor on hand?
- Have the monitoring intervals expired?
- etc.



NOTICE

We recommend that the method check only be deactivated if the titration must begin immediately after the start of the determination (e.g. with very rapid reactions such as enzyme-catalyzed reactions). If the method check is deactivated, it could happen that the determination run will be interrupted, e.g. because devices, titrants or sensors are not on hand.

The following settings can be implemented for a connected S885 Compact Oven SC:

Rack reset

on | off (Default value: **off**)

If this option is activated, then the sample rack is initialized at the time of the method start. The following actions are carried out:

- The lift is moved upwards.
- The sample rack is rotated to the starting position.
- The sample variable is reset to the value 1.

Check rack

Definition of the sample rack which must be attached at the time of the method start. This ensures that the method will be carried out only with this sample rack.



NOTICE

The sample rack can, however, only be checked if the **Rack reset** option is activated.

For methods for the 885 Compact Oven SC, this parameter must be set to **no** as it has a fixed rack.

Selection	no Selection of the configured sample racks
Default value	no

Increase sample variable automatically

on | off (Default value: **on**)

The sample variable describes the current position of the sample on the rack of the Sample Processor. If this option is activated, then the sample variable is automatically increased by 1 at the end of each determination.

This option must be deactivated if the sample variable is to be altered specifically with a **SAMPLE** command.

Request rack position at start

on | off (Default value: **off**)

If this option is activated, then the rack position of the first sample to be processed is requested.

15.5.2 Stop options

Main dialog: **Edit parameters ► Method options ► Start/Stop options ► Stop options**

In the dialog **Method options / Stop options**, you can define the actions which are carried out when a method is canceled.

The method can be canceled as follows:

- Manual stop with the fixed key []
- Stop because of an error
- Stop by remote signal via the Control Remote Box

Switch off stirrers

on | off (Default value: **on**)

If this option is activated, then all of the connected stirrers will be switched off.

Set remote lines

Selection of the signal out of the templates or entering the required bit pattern. Templates are defined under **System ► Templates ► Output lines**.



NOTICE

A line set active is not being reset automatically, not even at the end of the determination.

Entering the bit pattern:

- 0 = line inactive
- 1 = line active

- * = retain line status
- p = set pulse

The output lines are always numbered from right to left, i.e. with the signal *******1** line 0 is set. With a pulse, the length is set to 200 ms. If you wish to set pulses with other lengths, you have to define a corresponding template.

Entry	Bit pattern containing exactly 14 characters or a max. of 24 characters for the name of the template
Default value	*****
Selection	Selection of the templates defined

15.5.3 Sample data

Main dialog: **Edit parameters ► Method options ► Sample data**

You can define the following method-specific settings in the dialog

Method options / Sample data:

- Modify the designation for the sample identifications
- Define a fixed sample size
- Define limits for the sample size
If you define limits, these will be monitored at the start and at the end of the determination.

Below you will find an example demonstrating the effects on the main dialog of the settings in this dialog (*see "Modifying sample data", page 142*).

Identification 1, Identification 2

on | off (Default value: **on**)

If this option is activated, then the input field for the sample identification in the main dialog will be displayed.

Designation

A designation of your own choosing can be defined for the input field for each method.

Entry	16 characters maximum
-------	------------------------------

Fixed sample size

on | off (Default value: **off**)

Activate this parameter if the same sample amount is always to be used for all determinations. If this is case, then you can define the sample size here. It will be displayed in the main dialog, but it can no longer be changed there.

Sample size

This parameter can only be edited when **Fixed sample size** is activated.

Input range	-999999999 to 999999999
Default value	1.0

Unit of sample size.

Selection	g mg µg mL µL pieces
Default value	g



NOTICE

You can also use the parameter **Fixed sample size** if you would like to establish a default for the sample size.

Example: For a particular determination, you require 10 mL of sample. Now and again, however, you have too little sample available and you must perform the determination with 5 mL, for instance. In this case, proceed as follows:

1. Activate the parameter **Fixed sample size**.
2. Define the value and unit for the sample size, in the above example, for instance, **[10 mL]**.
3. Deactivate the parameter **Fixed sample size** again.
4. Save method.

⇒ Each time you load this method, 10 mL is entered as the sample size, but in the example mentioned you can enter the lesser quantity of 5 mL.

[Sample size limits]

This button is accessible only if **Fixed sample size** is deactivated.

Define the limit values for the sample size (*see "Defining limit values for the sample size", page 143*).

Modifying sample data

This instruction is for the purpose of clarifying the settings which are possible in this dialog. The following settings should be made:

- Change the designation for the input field of the first sample identification.
- Hide the input field for the second sample identification.
- Define a fixed sample size.

Proceed as follows:

1 Change the designation for the input field

- Tap on the input field **Identification 1**.
The text editor opens.
- Change the designation in **Batch** and confirm with **[OK]**.
- Deactivate the parameter **Identification 2**.


2 Define a fixed sample size

- Activate the parameter **Fixed sample size**.
- Enter the value **10 mL** as **Sample size**.

3 Displaying the main dialog

Tap on the fixed key [].

The data is saved and the adjusted main dialog appears.

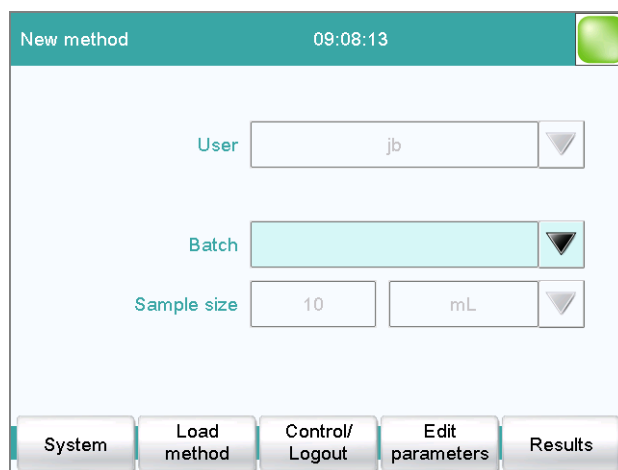

Only Identification 1 is shown with the title "Batch." The fixed sample size "10 mL" is shown, but cannot be edited.

Defining limit values for the sample size

In the dialog **Sample data / Sample size limits**, you can define the limit values for the sample size. These values are not monitored during sample data input, but rather:

- when the determination is started.
- when the determination is finished.
- upon automatic sample data query via the command **REQUEST**.
- when the determination is being recalculated.

An entry is made in the message list automatically when these limits are infringed (see chapter 23.2, page 182).

Monitoring sample size limits

on | off (Default value: **off**)

If this parameter is activated, then the sample size will be monitored.

Lower limit

Input range	-999999999 to 999999999
Default value	0

Upper limit

Input range	-999999999 to 999999999
Default value	999999999

Display message

on | off (Default value: **off**)

If this parameter is activated, a message is displayed in case the limits are infringed. You can select whether you want to continue with the determination or cancel the run.

If this parameter is deactivated, then the message will be entered only in the message list of the determination.

Timeout

The message is displayed during this time. Afterwards, the determination is continued automatically.

Input range	0 to 999999 s
Default value	30 s
Selection	off

off

The determination will not be continued until after the message has been confirmed.

15.5.4 Method properties

Main dialog: **Edit parameters ► Method options ► Properties**

Detailed information concerning the method is displayed in the **Method options / Properties** dialog and the methods can be signed electronically here:

- **Method status**

Current status of the method.

- **new**
The method has been newly created and not yet saved.
- **saved**
The current method has been saved.
- **modified**
The current method has been modified.
- **reviewed**
The current method has been signed at level 1.
- **released**
The current method has been signed at level 2.

- **Created by**

User who created the method.

- **Created on**

Date and time at which the method was created.

- **Last saved by**

User who saved the method most recently.

- **Last saved on**

Date and time at which the method was most recently saved. The version of the method is displayed as well. The version number will be increased by one each time the file is saved.

The following data is displayed only if the option **Saving modified method** under **Login options / Modification options** is activated:

- **Modified by**

User who modified and saved the method most recently.

- **Modified on**

Date and time at which the method was most recently modified and saved.

- **Reason**

Reason for the modification of the method.

- **Comment**

Comment for the modification of the method.

The following information is only shown if the method has been signed at level 1:

- **Reviewed by**

User who signed the method at level 1.

- **Reviewed on**

Date and time at which the method has been signed at level 1.

- **Reason**

Reason for the signing of the method.

- **Comment**

Comment for the signing of the method.

The following information is only shown if the method has been signed at level 2:

- **Released by**
User who signed the method at level 2.
- **Released on**
Date and time at which the method has been signed at level 2.
- **Reason**
Reason for the signing of the method.
- **Comment**
Comment for the signing of the method.

[Delete signature]

Delete all of the signatures for the method or the determination. This button is inactive if level 2 has not yet been signed (*see chapter 16.2, page 150*).

[Sign]

Sign the method or determination. This button is active only when working with activated login function and password (*see chapter 16.1, page 149*).

15.5.5 Note

Main dialog: **Edit parameters ► Method options ► Note**

In this dialog, you can enter a short text, e.g. important information about carrying out the determination.

[Display options]

Definition when the note is being displayed.

Automatically after loading the method

on | off (Default value: **off**)

If this parameter is activated, the note is displayed when loading the method. It can otherwise only be read in this dialog.

15.5.6 Saving a determination automatically

Main dialog: **Edit parameters ► Method options ► Save automat.**

In the dialog **Method options / Save automatically**, you can specify whether the determination is to be saved automatically and/or whether a PC/LIMS report is to be created.

Save determination automatically

on | off (Default value: **off**)

If this parameter is activated, then the determination will be saved automatically.

Memory

Memory location where the determination will be saved.

Selection	External memory 1 External memory 2 Shared memory
Default value	External memory 1

Shared memory

The determination will be saved in a shared directory on the network. The shared directory is selected in the **Edit device / Shared memory** dialog (see chapter 10.3.3, page 69).

Group

Directory where the determination will be saved.

Entry	32 characters maximum
Selection	Selection of available directories empty
Default value	empty

File name

File name of the determination. The file name is always extended to include the date and the time of day (YYYYMMDD-hhmmss) in order to ensure that the file names are unique for all determinations.

Entry	max. 16 characters
Selection	Identification 1 Identification 2 Method
Default value	Identification 1

Identification 1

The first 16 characters of the text which was entered in the main dialog for Identification 1 + YYYYMMDD-hhmmss.

Identification 2

The first 16 characters of the text which was entered in the main dialog for Identification 2 + YYYYMMDD-hhmmss.

Method

The first 16 characters of the text, of the method name + YYYYMMDD-hhmmss.

Write protection

on | off (Default value: **on**)

If this parameter is activated, then the file cannot be saved, deleted, or renamed. This is only an internal write protection feature and is independent of the write protection of the operating system on your computer. This write protection feature protects the saved determination data against accidental modifications or modifications by unauthorized persons.



Create PC/LIMS report

on | off (Default value: **off**)

If this parameter is activated, then a machine-readable report with all of the important data for the determination is created, a so-called PC/LIMS report.

The settings are implemented in the device manager (*see chapter 10.3.2, page 68*).

When you save the PC/LIMS report as a file, the file name will be generated automatically as follows: **PC_LIMS_Report**-first 16 characters of the value of the parameter **File name-Date-Time.txt**. If the value of the parameter **File name** contains one or more of the following characters, then these will be automatically converted in the file name into the character "_": / \ : * ? " < > |.

16 Electronic signatures

16.1 Signing methods/determinations electronically

Methods: **Main dialog ▶ Edit parameters ▶ Method options ▶ Properties ▶ Sign**

Determinations: **Results dialog ▶ More data ▶ Properties ▶ Sign**

You can sign methods or determinations in the dialog **Electronic signature**. The following information is saved along with each electronic signature:

- User (complete name)
- Date and time
- Reason
- Comment

Methods can only be signed if they have the status **saved**. The status of a method can be found in the dialog **Method options / Properties**.

Signature

Selection of the level at which the determination should be signed.

Selection	Review (signature level 1) Release (signature level 2)
	A maximum of three signatures from up to three different users can be given at each level. However, it is not possible for a user to sign at both level 1 and level 2.

Review (signature level 1)

If the method has never been signed before, then only level 1 can be selected.

Release (signature level 2)

If the method has been signed three times at level 1 or already once at level 2 then only level 2 can be selected.

User


Identification of the user who signs.

Password

Password of the user.

17 Control

Main dialog: **Control**

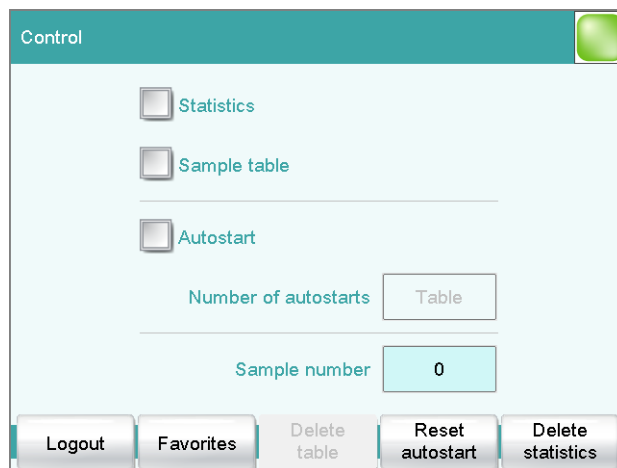

In the dialog **Control**, the settings for the execution of a single determination or of one sample series are defined.

If you work with the login function activated, you can log off the system in this dialog. The login dialog will then be shown immediately.

Statistics

on | off (Default value: **off**)

With this option, you can enable or disable the statistics calculations for individual determinations (see "Statistics", page 153).



NOTICE

The parameter **Statistics** in the method options will not be changed as a result. This setting here will be adjusted automatically to match the one in the method options.

Sample table

on | off (Default value: **off**)

If this option is activated, the sample data for a sample series can be entered in a table (see chapter 20, page 162).

Autostart

on | off (Default value: **off**)

If this option is activated, a new determination is started automatically at the end of a determination. This continues until the number specified has been reached (siehe **Number of autostarts**).



NOTICE

If you deactivate this option while a determination is running, then that determination will be ended, although the next one will however also not be started. This can be used to cancel a sample series.

Number of autostarts

Number of automatic starts.

Input range	1 to 9999
Selection	Table
Default value	Table

Table

The number of automatic starts corresponds to the number of samples in the sample table.

Sample number

The sample number is primarily used for documentation. Each time a determination is started, it is increased by one. The sample number can, however, also be entered manually. Each time the Coulometer is switched on, the value is reset to zero.

Input range	0 to 9999
Default value	0

[Logout]

User logout (visible only when work is done with activated login function).

[Favorites]

Save methods/sample tables as a favorite (*see chapter 18, page 155*). Depending on whether you are working with or without login function, user-specific or common favorites can be created. On the main dialog, a button is created for each favorite. This way, a method or a sample table can be started with one single mouse click.

[Delete table]

Delete the sample table completely.

[Reset autostart]

Reset the value of the autostart counter to zero.

[Delete statistics]

Delete all statistics data.

The statistics data should be deleted manually if a sample series has been canceled and a new sample series is to be started.

The statistics data is deleted automatically in following cases:

- when all of the determinations of the determination series have been carried out and a new determination has been started afterwards.
- when a new method is loaded.
- when a determination is loaded (the method with which the determination was carried out is loaded simultaneously with the determination).

Statistics

In the dialog **Control**, you can deactivate the statistics calculations for individual determinations. This option is required primarily in order to insert the processing of an "**urgent sample**" during the processing of a sample series when the urgent sample is not to be included in the statistics calculations.

Interrupting a determination series for an "urgent sample"

If the same method is required for the "urgent sample" as is being used for the sample series, then you need only deactivate the option **Statistics** and then reactivate it after the determination is completed. Proceed as follows if you need to process the "urgent sample" with a different method:

1 Deactivate statistics calculation

- Deactivate the option **Statistics**.

2 Saving the determination

- Save the current determination of the sample series (*see chapter 23.6, page 190*).

3 Carrying out the determination

- Load the method for the "urgent sample".
- Carry out the determination.

4 Load the last determination of the sample series

- Reload the previously saved last determination.

The determination, the method used for it and the current statistical data are loaded. The option **Statistics** is activated again and the

value of the statistics counter corresponds to that in effect before the interruption.

5 Continuing the sample series


- Carry out the remaining determinations in the sample series.

18 Favorites

Main dialog: **Control ► Favorites**

Main dialog: **System ► System settings ► User admin. ► Edit ► Favorites**

This chapter describes how you can create favorites. Favorites are comparable to the favorites/bookmarks in your Internet browser. In the main dialog, a button is created for each favorite. By pressing a single key, you can trigger an action without having to navigate through different dialogs.

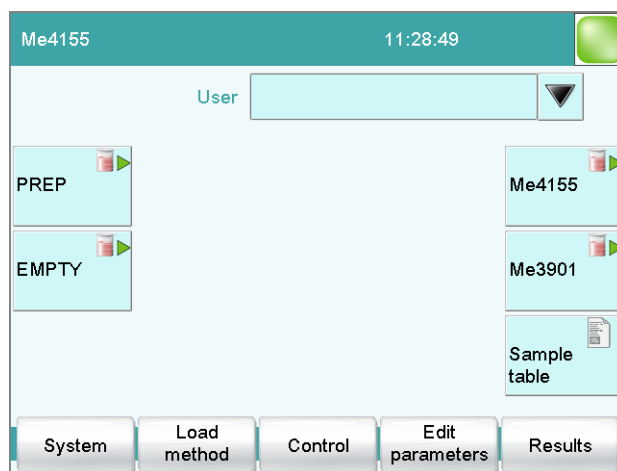

The following objects can be saved as favorites:

- Methods
- Sample tables

User-specific or common favorites can be created, depending on whether you are working with or without the login function.

- Operation with login function:
If you are working with the login function activated, **user-specific favorites** can be created. These can only be used by a certain user. User-specific favorites are created as follows:
 - in the user administration by a user with administrator rights (**System ► System settings ► User admin. ► Edit**).
 - in the dialog **Control** for the user who is logged in.
- Operation without login function:
If you are working without the login function, **common favorites** can be created. These favorites are available for all users. Common favorites are created in the **Control** dialog.

18.1 Creating favorites

Control / Favorites	
Name	Position
PREP	1
EMPTY	6
Me4155	5
Me3901	10

The list of favorites gives the designation and position of the button on the main dialog of each favorite configured.

[New]

Add a new favorite to the list, see following chapter.

[Delete]

Delete the selected favorite from the list.

[Edit]

Edit the properties of the selected favorite, see following chapter.

18.1.1 Editing favorites

Favorites / Edit	
Position	5 ▼
Name	Me4155
Type	Method ▼
Memory	Internal memory ▼
Method / Sample table	Me4155 ▼



Position

On the main dialog, a button is created for each favorite. These buttons are ordered in three rows at fixed positions. The position 1 can be found top left.

Input range	1 to 14
Default value	1

Name

The designation of the favorite is used for unambiguous identification.

Entry	24 characters maximum
-------	------------------------------

Type

Definition as to whether the favorite represents a single method or a complete sample table.

Selection	Method Sample table
Default value	Method

Memory

Memory location the method or the sample table is loaded from. Only the currently accessible memory locations are selectable.

Selection	Internal memory External memory 1 External memory 2 Shared memory
Default value	Internal memory

Shared memory

The favorite will be saved in a shared directory on the network. The shared directory is selected in the **Edit device / Shared memory** dialog (see chapter 10.3.3, page 69).

Method / Sample table

Method or sample table of the favorite.

Selection	Selection of stored methods/sample tables
-----------	--

[More options]

Parameterize the autostart function, see following chapter.

More options

The autostart function is parameterized under **[More options]**.



Autostart

on | off (Default value: **off**)

If this option is activated, a new determination is started automatically at the end of a determination. This continues until the number specified has been reached (siehe **Number of autostarts**).

Number of autostarts

This option can only be edited when **Autostart** is activated.

Number of automatic starts.

Input range	1 to 9999
Default value	1
Selection	Table

Table

The number of automatic starts corresponds to the number of samples in the sample table.

19 Sample data

You can enter the sample data (identification, sample size, etc.) in a variety of ways:

- Directly in the main dialog.
- Using the sample table. This is particularly useful with sample series. The sample table is a table in which the sample data for up to 999 samples can be entered (*see chapter 20, page 162*).
- Automatic request immediately after the start of the determination (*see chapter 19.2, page 160*).

You can also send the sample size and the unit from a connected balance in any case. With some balances, the sample identification and method can be also sent (*see chapter 31.4, page 312*).

19.1 Entering sample data in the main dialog

For a sample, you can enter the sample data directly in the main dialog, even while the determination is running (*see chapter 22, page 173*).

Identification 1

Sample identification. The sample identification can be used in calculations as the variable **CI1**. If you have defined a sample identification list (*see chapter 14.1.1, page 118*) then the entries can be selected here.

Entry	24 characters maximum
Default value	empty



Identification 2

Sample identification. The sample identification can be used in calculations as the variable **CI2**. If you have defined a sample identification list (see *chapter 14.1.1, page 118*) then the entries can be selected here.

Entry	24 characters maximum
Default value	empty

Sample size

Sample size. The value of the sample size can be used in calculations as the variable **C00**.

Input range	-99999999 to 999999999
Default value	1.0

Unit of sample size.

Entry	6 characters maximum
Selection	g mg µg mL µL pieces
Default value	g

Unit defined in method:

The unit defined in the method is ignored. The unit entered in the sample table is used.

Balance connected and configured:

Value from the connected balance.

Fixed sample size defined in method:

The value will be entered in the corresponding line when the method is loaded. The existing entry will be overwritten.

19.2 Requesting sample data at the start of the determination

The sample data can be automatically requested immediately after the start of the determination in order to ensure that the sample data entry is not forgotten. This automatic inquiry is requisite for reweighing.

Me2115

01 REQUEST Data request

Identification 1

Sample size mL

Continue

To accomplish this, you must insert and configure the command **REQUEST** in the method (see chapter 28.10.3, page 299).

If the **Hold sequence** parameter is activated, then the run will be paused and must be continued with **[Continue]** after the sample data has been entered. If the **Hold sequence** parameter is deactivated, then the determination will continue in the background until the measured data is required for further processing. This dialog will be displayed until the entering of the sample data is confirmed with **[Continue]**, even if the titration/measurement is already completed. This ensures that the sample data is available for calculations.

If a fixed sample size has been defined in the method (see chapter 15.5.3, page 141), then it will be displayed at the time of the entry but it will not be editable.

20 Sample table

20.1 General

Main dialog: **Sample table**

The sample table is a table in which the sample data for up to 999 samples can be entered. If you are processing larger-sized sample series, it makes sense to enter the sample data in this table.


Activating the sample table

Proceed as follows to enable the sample table:

1 Opening the control dialog

- In the main dialog, tap on **[Control]**.


The control dialog is displayed:

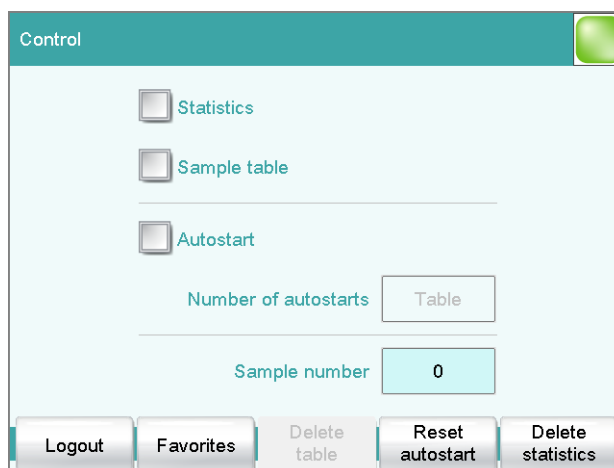

2 Activating the sample table

- Activate the **Sample table** option.
- Tap on the fixed key **[🏠]**.

The main dialog is displayed:

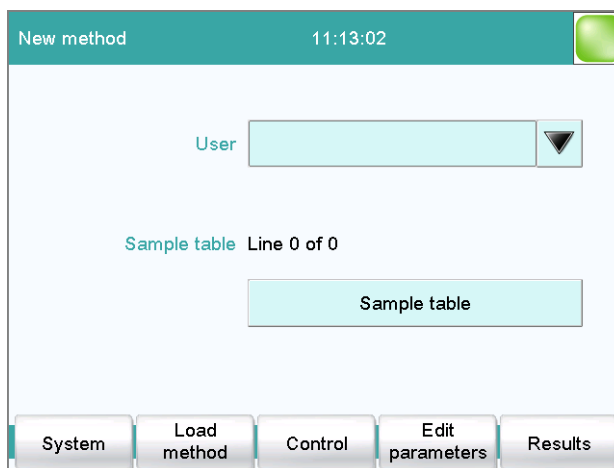

The display will show the number of determinations already carried out and the total number of sample lines containing data. The sample table is still empty in this example.

The sample table contains numbered lines; each line represents one determination. In the standard settings, the first sample identification and the sample size of each sample are displayed. In the properties, you define the columns to be displayed (*see chapter 20.3, page 167*).

Sample table		
No.	Identification 1	Sample size
1	#2370015	1.0 g
2	#2370015	1.0 g
3	#2370016	1.0 g
4	#2370016	1.0 g
5	...	

Buttons below the table: Load/Save, Properties, Insert line, Delete, Edit

[Load/ Save]

Load a saved sample table (*see "Loading a sample table", page 165*) or save the current sample table (*see "Saving a sample table", page 164*).

[Properties]

Edit the properties of the sample table (*see chapter 20.3, page 167*).

[Insert line]

Insert a new line above the line selected.

[Delete]

Delete the selected line from the sample table.

[Edit]

Edit the data of the selected line (see chapter 20.2, page 165).

Saving a sample table**NOTICE**

Sample tables can only be saved on external storage media or on a shared file system.

Proceed as follows to save a sample table:

1 Opening the dialog

- In the **Sample table** dialog, tap on the **[Load/ Save]** button.

The **Sample table / Load/Save** dialog is displayed.

If no external storage medium is plugged in or if no file system is shared, then **[Save]** is disabled.

2 Modify/apply a name

- Tap on **[Save]**.
A name is suggested for new sample tables. If the sample table has already been saved, the previous name will be displayed.

Applying the name:

- Tap on **[Save]**.

The sample table is saved.

Entering a new name:

- Tap on the input field **File name**.
The text editor opens.
- Enter a new file name (max. 32 characters) and confirm with **[OK]**.
- Tap on **[Save]**.

The sample table is saved.

Loading a sample table

Proceed as follows to load a sample table:

1 Display the list of saved sample tables

- In the **Sample table** dialog, tap on the **[Load/ Save]** button.
The **Sample table / Load/Save** dialog is displayed.
If no external storage medium is plugged in or if no file system is shared, then **[Load]** is disabled.
- Tap on **[Load]**.
The selection of memory locations is displayed. This selection will be skipped if only one memory can be accessed.
- Select the memory location where the desired sample table is stored.
The selection of file groups is displayed (*see chapter 11.1, page 91*). If only one group is available, then the list of the saved sample tables will be displayed directly.
- Select the group with the desired sample table.
- Tap on **[Show files]**.

The list with the saved sample table is opened.

2 Select the sample table

- Select the desired sample table.

3 Loading a sample table

- Tap on **[Load]**.

The sample table is now loaded.

20.2 Editing sample data

Sample table / Edit

Line number - 1 +

Method Me2115 ▼

Identification 1 #2370015

Identification 2

Sample size 1.0 g ▼



At the very top you will see the line number of the selected line. In this example, the sample data of the first line is displayed. You can scroll between individual data sets with the keys **[-]** and **[+]**.

Method

Method used for processing the sample.

If you use the sample assignment table (*see chapter 14.1.2, page 119*), then the method can no longer be defined here.

Selection	Selection of stored methods empty
Default value	empty

empty

The currently loaded method is used.

Identification 1

Sample identification. The sample identification can be used in calculations as the variable **CI1**. If you have defined a sample identification list (*see chapter 14.1.1, page 118*) then the entries can be selected here.

Entry	24 characters maximum
Default value	empty

Identification 2

Sample identification. The sample identification can be used in calculations as the variable **CI2**. If you have defined a sample identification list (*see chapter 14.1.1, page 118*) then the entries can be selected here.

Entry	24 characters maximum
Default value	empty

Sample size

Sample size. The value of the sample size can be used in calculations as the variable **C00**.

Input range	-99999999 to 9999999999
Default value	1.0

Unit of sample size.

Entry	6 characters maximum
Selection	g mg µg mL µL pieces
Default value	g

Unit defined in method:

The unit defined in the method is ignored. The unit entered in the sample table is used.

Balance connected and configured:

Value from the connected balance.

Fixed sample size defined in method:


The value will be entered in the corresponding line when the method is loaded. The existing entry will be overwritten.



NOTICE

The sample size limits defined in the method are not monitored when the sample data is being entered in the sample table.

20.3 Properties

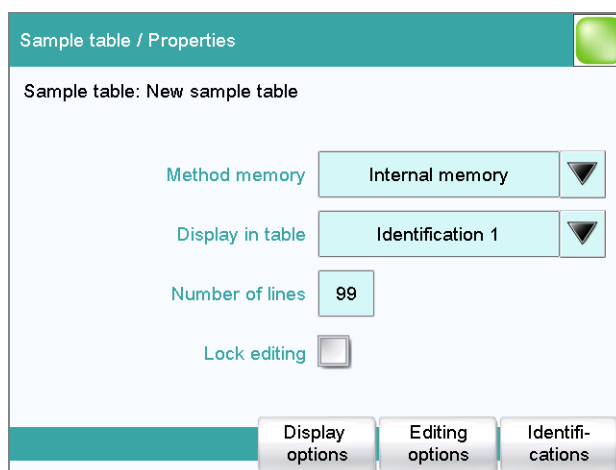

Method memory

Memory location the method is loaded from. All memory locations are selectable, even if they are currently not accessible.

Selection	Internal memory External memory 1 External memory 2 Shared memory
Default value	Internal memory

Shared memory

Shared directory in the network.

Display in table

Selection of the columns to be displayed in the sample table.

Selection	Method Identification 1 Identification 2
Default value	Identification 1



End line with

Selection as to with which input the editing dialog for the next sample is automatically displayed.

Selection	Sample size Identification 1 Identification 2 manual
Default value	Sample size

manual

The editing dialog for the next sample can be displayed with the **[New sample]** button in the editing dialog.

If, for example, you have to enter similar data for each sample, then you can automatically copy the data which is identical for each sample into the next line:

- **Method**
- **Identification 1**
- **Identification 2**
- **Sample size**

Dialog "Properties / Identifications"

Sample table: **Properties ► Identifications**

In the dialog **Properties / Identifications**, you can alter the designations for the sample identifications in accordance with specific methods.

After the confirmation of the sample size, the live curve of the titration is displayed.

Sequence of the determination

- After a determination is started, a check is made as to
 - whether all of the devices required are connected and ready for operation.
 - whether the sensor is present.
 - whether the monitoring intervals for the sensor and common variables are still valid.
 - if a sample assignment table is activated (*see chapter 14.1.2, page 119*):
 - whether the entered sample identification is defined in the sample assignment table. If this is the case, then the assigned method will be loaded automatically and the actual determination will be started.

21.2 Performing a sample series

You can use the following functions with sample series:

- **Statistics for multiple determinations**
The statistics calculations are enabled in the method options (*see chapter 15.5, page 137*).
- **Sample table**
You can enter the sample data for an entire sample series in the sample table (*see chapter 20, page 162*).

Sequence of the sample series

- After a determination is started, a check is made as to
 - whether all of the devices required are connected and ready for operation.
 -
 - whether the sensor is present.
 -
 - whether the monitoring intervals for the sensor and common variables are still valid.
 - if a sample assignment table is activated (*see chapter 14.1.2, page 119*):
 - whether the entered sample identification appears in the sample assignment table. If this is the case, then the assigned method will be loaded automatically and the actual determination will be started. If a different method has been entered in the sample table, then it will be ignored.


22 Live modifications

22.1 Editing the sample data of the running determination

The sample data can be entered or modified in the main dialog while a determination is running. The sample data entered at the end of the determination is always used in calculations.

Proceed as follows to edit the sample data:

1 Display the main dialog

- Tap on the fixed key [].

The main dialog is displayed. The determination continues to run in the background.

2 Edit the sample data

- Edit the sample data and apply with **[OK]**.

3 Display the live dialog

- Tap on **[Live display]**.

The live dialog is displayed once again.



NOTICE

Make sure that the input dialogs are closed before the determination is ended. Otherwise the determination will have to be recalculated.

22.2 Editing the sample table while a determination is running

You can insert new lines or delete existing ones or edit sample data while a determination is running.



NOTICE

To ensure that no problems arise during the run and that the current data is always available for the calculation, we recommend that the input dialogs always be closed.

Editing the sample table

Proceed as follows to edit the sample table:

1 Display the main dialog

- Tap on the fixed key [].

The main dialog is displayed. The determination continues to run in the background.

2 Select the sample data

- Tap on **Sample table**.
The sample table is displayed. The first line always contains the sample data of the running determination.
- Select the desired line.
- Tap on the button **[Edit]**.

3 Edit the sample data

- Edit the sample data and apply with **[OK]**.



NOTICE

In addition to the sample data, the method can also be modified, except in cases where the determination is running.

4 Display the live dialog

- Tap on the fixed key [].
The main dialog is displayed.
- Tap on **[Live display]**.

The live dialog is displayed once again.

22.3 Live display

Main dialog: **Live display**

The current method and the system status are displayed in the title line. The currently running command is displayed directly underneath the title line.

The following functions are available to you while a determination is running:

- **[Hold]**
Pause the determination. The label changes to **[Continue]**. This can be used to continue the run.
- **[Cancel command]**
Cancel the current method command. The next command is started.
- **[View]** (only for titrations/measurements)
To modify the settings for the curve and measured value display, see below.
- **[Sample data]** (visible only if a titration or a measurement is not currently running)
Switch to the main dialog in order to edit the sample data.
- **[Edit parameters]**
Edit the method options and live parameters (*see chapter 22.4, page 176*).
- **[Results]**
Display the results dialog (*see chapter 23, page 178*).

Dialog "View"

Main dialog: **Live display ► View**

View

Type of curve display.

Selection	Curve Curve & meas. values Measured values Curve & sample data
Default value	Curve & meas. values

Curve

Only the curve is displayed.

Curve & meas. values

The measured values are displayed to the right next to the curve.

Measured values

Only the measured values are displayed.


Curve & sample data

The sample data is displayed to the right next to the curve.

2 Edit the method parameters

- Change the desired parameters accordingly.

3 Display the live dialog


- Tap on the fixed key [].
The main dialog is displayed.
- Tap on **[Live display]**.

The live dialog is displayed once again.

23 Results and more determination data

Main dialog: **Results**

Detailed specifications concerning the current determination are displayed in the **Results** dialog:

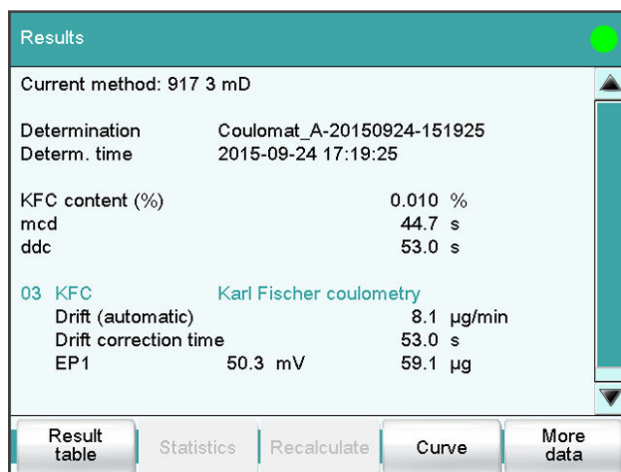

- **Current method**
Method with which the determination was carried out.
- **Determination**
Name of the determination. Displayed only if the determination has been saved.
- **Determination time**
Date and time at which the determination was carried out.
- **Results**
Calculated results. If monitoring of the result limits is activated, then the result will be highlighted in color:
 - Green lettering: The result is within the limits
 - Red lettering: The result is outside the limits
- **Commands**
List of the data-generating commands (titrations, measurements, etc.). The most important data is displayed for each command.

[Result table]

Open result table (*see chapter 25, page 199*). The result table is a table in which the results for several determinations can be saved.

[Statistics]

Display the statistical overview of a determination series (*see chapter 24, page 194*).

[Recalculate]

Recalculate the determination. The procedure will be executed immediately. Recalculation cannot be undone (*see chapter 23.8, page 193*).

[Sign]

Sign the current determination.

If the method contains a command **SIGN**, then this button will be displayed instead of **[Recalculate]**. The determination can be signed directly on level 1 (with a second command also on level 2).

[Curve]


Display the curve of the current determination (*see chapter 23.7, page 191*).

[More data]

Display detailed specifications for the determination (determination properties, measuring point list, endpoint details, etc.), save determination (*see chapter 23.1, page 179*).

23.1 More determination data

Results dialog: **More data**

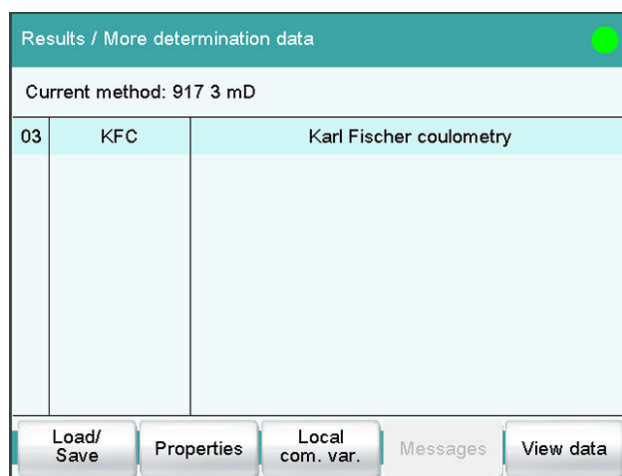

All of the commands for which you can display other data (titrations, measurements, evaluations, etc.) are listed in the **Results / More determination data** dialog.

[Load/ Save]

Load saved determinations (*see chapter 23.5, page 185*) and save current determination (*see chapter 23.6, page 190*).

[Properties]

Display the properties of the current determination (*see chapter 23.4, page 183*).

Results dialog: **Curve ► Display curve ► Measuring point list**

View data / Measuring point list		
03 KFC Karl Fischer coulometry		1M
Time [s]	Water [µg]	Meas.value [mV]
0.1	0.0	158.5
2.1	29.3	85.9
4.1	40.5	83.7
6.1	51.3	61.6
8.1	55.9	54.8
10.1	58.4	53.2
12.1	59.9	52.5

View

The individual measured values are listed in the **View data / Measuring point list** dialog. You can configure which columns are displayed (**[View]** button). These settings are saved separately for each mode (KFC, MEAS, etc.).

Column 1 / Column 2 / Column 3

Quantity displayed in the columns 1, 2 and 3. Both the selection of the quantities and the default setting depend on the mode.

Selection	Measured value Water Bromine Time Temperature Drift Index
-----------	--

Endpoint details

Results dialog: **More data ► View data ► Endpoint details**

Results dialog: **Curve ► Display curve ► Endpoint details**

Detailed information concerning each equivalence point or endpoint determined is displayed in the **View data / Endpoint details** dialog.

23.2 Messages

Results dialog: **More data ► Messages**

If messages occurred during a determination, they will be displayed in the **More determination data / Messages** dialog. The following data for the message is indicated under **[Details]**:

- **Message number**
Unique identification number.
- **Time**
Moment that the message occurred.
- **Message**
Message text.

23.3 Local common variables

Results dialog: **More data ► Local com. var.**

The list of common variables as it existed at the time of the determination is displayed in the **More determination data / Local common variables** dialog.



NOTICE

Recalculation:

The common variables can be modified retroactively. The values displayed here are used for the recalculation of a loaded determination. The list of common variables in the **System / Common variables** dialog is not changed as a result.

23.4 Determination properties

Results dialog: **More data ► Properties**

Detailed information concerning the determination is displayed in the **More determination data / Properties** dialog:

- **User**
User who carried out the determination.
- **Sample number**
Number which was entered in the dialog **Control** at the start of the determination.
- **Start counter**
This value, just like the sample number, is increased by one each time a determination starts. In contrast to the sample number, this value cannot be modified by the user.
- **Duration determ.**
Duration of the entire determination.
- **Method status**
Current status of the method.
 - **new**
The method has been newly created and not yet saved.
 - **saved**
The current method has been saved.
 - **modified**
The current method has been modified but not saved yet (e.g. live modification).
 - **reviewed**
The current method has been signed at level 1.
 - **released**
The current method has been signed at level 2.
- **Determination**
Name of the determination. Displayed only with saved determinations.
- **End of determin.**
Way in which the determination was ended.
 - **Regular without errors**
The determination was automatically ended at the end of the method.
 - **Manual stop**
The determination has been canceled with the fixed key [].
 - **Error**
The determination has been canceled due to an error.

- **Comment**

Comment for the signing of the determination.

[Delete signature]

Delete all of the signatures for the method or the determination. This button is inactive if level 2 has not yet been signed (*see chapter 16.2, page 150*).

[Sign]

Sign the method or determination. This button is active only when working with activated login function and password (*see chapter 16.1, page 149*).

23.5 Loading a determination

Proceed as follows to load a determination:

1 Opening the determination list

- In the results dialog, tap on **[More data]**.
- Tap on **[Load/ Save]**.
- Tap on **[Load]**.


The selection of memory locations is displayed. This selection will be skipped if only one external storage medium is available.

- Select the memory location where the desired determination is saved.

The selection of file groups is displayed (*see chapter 11.1, page 91*). If only one group is available, then the determination list will be displayed directly.

- Select the group with the desired determination.
- Tap on **[Show files]**.

The determination list with the stored determination opens:

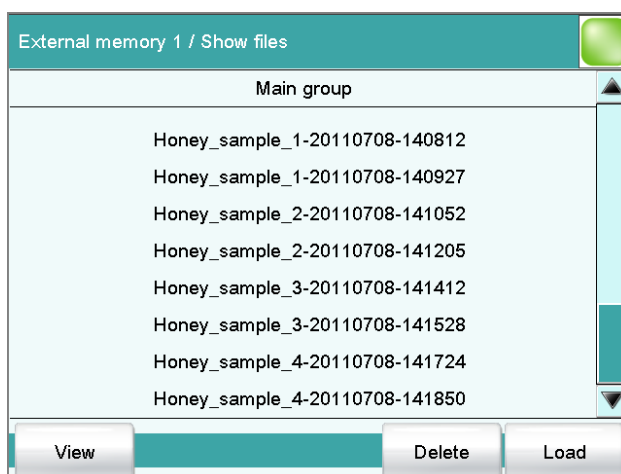

2 Selecting a determination

- Select the desired determination.

3 Loading the determination

- Tap on **[Load]**.

The determination is now loaded. The properties are displayed in the results dialog.



NOTICE

When you load a determination, the method with which the determination was carried out will be loaded automatically.

For this reason, save any modifications which may have been made to the method currently loaded before you load a determination.

23.5.1 Determination list

Results dialog: **More data ▶ Load/ Save ▶ Load ▶ Show files ▶ View**

Show files / View		
Sorted by Determ. time		27
No.	Identification 1	Result 1
20	Honey_sample_1	16.72 %
21	Honey_sample_1	16.89 %
22	Honey_sample_2	16.62 %
23	Honey_sample_2	16.70 %
24	Honey_sample_3	16.67 %
25	Honey_sample_3	16.84 %
26	Honey_sample_4	16.87 %

Previous 100 Next 100 Properties Delete Load

In the **Show files / View** dialog you can depict the individual determinations in greater detail. The sorting criterion and the number of saved determinations are displayed at the very top. The list contains numbered lines; each line represents one determination. In the default settings, the first sample identification and the first calculated result of each determination are displayed. In the properties, you define which data is displayed in the columns and the criteria according to which the determinations are sorted (see "Properties", page 187). In addition, you can define search fil-

ters to aid you in finding the desired determinations (see "Filter criteria", page 188).

[Previous 100]

Display the previous 100 determinations.

[Next 100]

Display the next 100 determinations.

[Properties]

Configure the display of the columns and define filter criteria.

[Delete]


Delete the selected file.

[Load]

Load the selected file.

Properties

Determination list: **Properties**

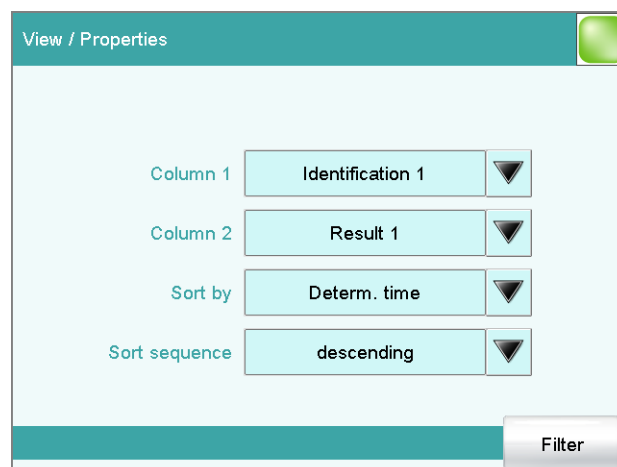

In the **View / Properties** dialog, you can define which data is displayed in the columns and the criteria according to which the determinations are to be sorted.

Column 1

Selection of what is to be displayed in the first column.

Selection	Identification 1 Identification 2 Determ. time Result name Method Sample size User
Default value	Identification 1

Column 2

Selection of what is to be displayed in the second column.

You can set filters in the **Properties / Filter** dialog. A filter is a rule defining what is to be displayed in the determination list. If a filter has been set, this is indicated with a corresponding symbol in the **Show files / View** dialog.

Filter

Define the criterion according to which the list is to be filtered. The only lines which will be displayed are those which fulfil the filter criterion.

Selection	no filter Identification 1 Identification 2 Date Result Method Sample size User
Default value	no filter

no filter

The list itself will not be filtered.

Identification 1

In the input field **is equal to**, select or enter the first sample identification according to which filtering is to be carried out.

Identification 2

In the input field **is equal to**, select or enter the second sample identification according to which filtering is to be carried out.

Date

In the input fields **Date from ... to**, specify the period of time according to which filtering is to be carried out.

Result

In the input fields **Value from ... to**, specify the value range according to which filtering is to be carried out.

Method


In the field **is equal to**, select or enter the method name according to which filtering is to be carried out.

Sample size

In the input fields **Value from ... to**, specify the value range according to which filtering is to be carried out.

23.7 Curves

Results dialog: **Curve**

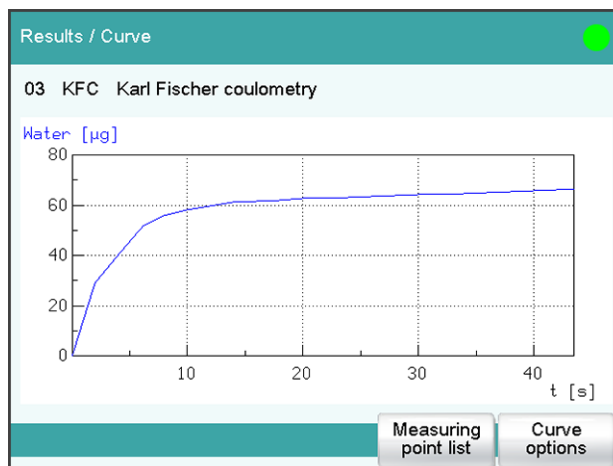

The curve of the current determination is displayed in the **Results / Curve** dialog. You can configure the curve display.

[Endpoint details]

Display details for each endpoint found (see "Endpoint details", page 181).

[Measuring point list]


Display measuring point list of titrations and measurements (see "Measuring point list", page 180).

[Curve options]

Define the settings for the curve display (see "Curve options", page 191). The settings are saved separately for each titration/measuring mode and apply not only for the live display but also for the curve display in the results dialog.

Curve options

Results dialog: **Curve** ► **Curve options**

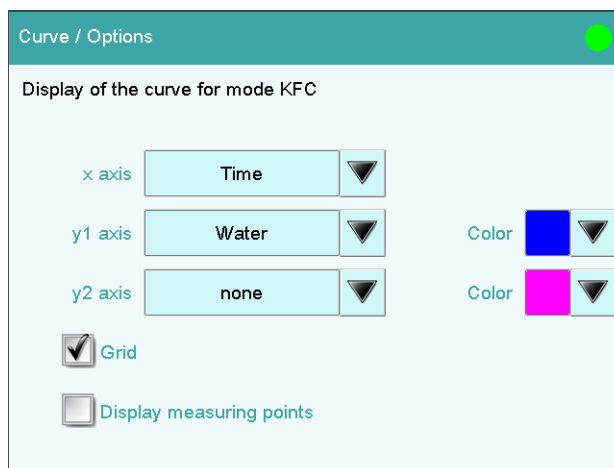

You can make the settings for the curve display in the **Curve / Options** dialog. These settings are saved separately for each mode (KFC, MEAS, etc.) and also apply for the live display. Both the selection of the quantities and the default setting depend on the mode.

x axis

Quantity which is shown on the x axis of the graph.

y1 axis

Quantity which is shown on the y1 axis of the graph.

y2 axis

Quantity which is shown on the y2 axis of the graph.

Color

Selection of a color for the curves.

Grid

on | off (Default value: **on**)

If this parameter is activated, grid lines are shown.

Display measuring points

on | off (Default value: **off**)

If this parameter is activated, the curve is shown with the single measured values.

23.8 Recalculation and reevaluation

Results dialog: **Recalculate**

You can use the **[Recalculate]** function to recalculate and reevaluate determinations. This is necessary in the following cases:

- If you have modified the evaluation parameters in a way that changes the recognized endpoints or equivalence points.
- If you have modified the calculations.
- If you have modified the variables used in calculations, e.g. sample size, titer or common variables.



NOTICE

Neither the recalculation nor the reevaluation can be undone.

We therefore recommend that the original determination be saved beforehand (*see chapter 23.6, page 190*). The best way of doing this is to use the **Save determination automatically** function (*see chapter 15.5.6, page 146*).

The following method commands are executed once again with the **[Recalculate]** function:

- EVAL commands:
All evaluations are carried out again. You can also insert new evaluations retroactively.

If a determination has been recalculated, this will be documented in the determination properties (*see chapter 23.4, page 183*).

Recalculation/reevaluation of loaded determinations

With loaded determinations, it is assumed that they were carried out a long time ago and that the **common variables** of the system no longer correspond to those of the determination. This is why the list of **common variables** is stored with each determination. This data will be used if you recalculate a loaded determination. If you wish to recalculate the result with a corrected **common variable**, you must therefore modify the respective values in the **More determination data / Local common variables** or **More determination data / View data** dialog.

If a result is assigned to a **common variable**, then the **common variable** which is currently valid in the system will not be overwritten until after a query has been confirmed.

The recalculated results of loaded determinations are not saved in the result table.

24 Statistics

Main dialog: **Results ► Statistics**


In the dialog **Results**, you can use **[Statistics]** to display the statistical overview of a determination overview. You can statistically evaluate a maximum of nine results calculated in a determination. A statistical series can contain a maximum of 20 determinations.

Activating the statistics function

Proceed as follows to activate the statistics function:

1 Editing the method options

- In the main dialog, tap on **[Edit parameters]**.
- Tap on **[Method options]**.

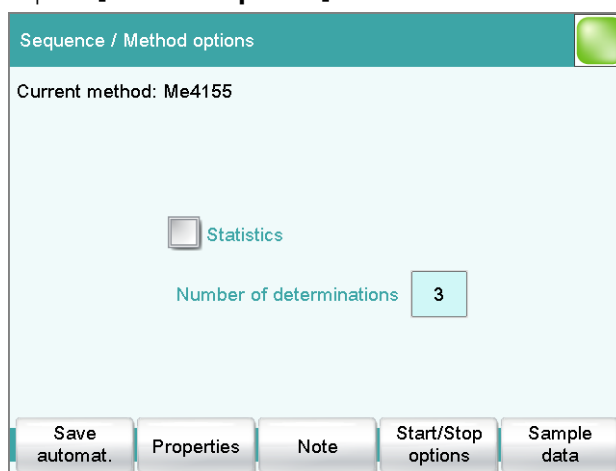

- Activate the **Statistics** option.
- Under **Number of determinations**, enter the desired number of determinations.

The results for which statistical calculations are performed are defined in the calculation command **CALC** (see "Variable for mean value", page 123).

Resultate / Statistik	
Aktuelle Methode: Me4155 3 von 3	
Resultatname	Mittelwert
KFT Titer	5.0080 mg/mL
<div style="display: flex; justify-content: space-between;"> Bestimmung hinzufügen Details </div>	

In this dialog, the method with which the determination was carried out is displayed in the first line. The second line shows how many determinations have already been carried out and the total number of determinations to be carried out. The name of the result and the mean value are shown for each result.

[Add determ.]

Add a further determination to the determination series (*see chapter 24.3, page 197*).

[Details]

Display the statistics results for the selected result and the single results from all determinations, see next chapter.



NOTICE

If you recalculate a determination, the statistics for this determination are also recalculated.

24.1 Displaying details for a result

Results dialog: **Statistics ▶ Details**

Statistics / Details		
Result name: KFC content (ppm)		SMN1
Mean value	1000 ppm	n=03
s abs	2.5 ppm	
s rel	0.25 %	
No.	Sample size	Result
1	0.9496 g	1003 ppm
2	1.7160 g	998 ppm
3	1.5813 g	1000 ppm

Sample data
 Determ. on/off
 Result on/off

In the dialog **Statistics / Details**, the following details are displayed:

- Result name
- Mean value
On the right-hand side of the dialog the number of single results from which the mean value was calculated is displayed. In this example, it is 3.
- Absolute standard deviation (**s abs**)
- Relative standard deviation (**s rel**)
- Result and sample size of each single determination
Results that could not be calculated are indicated with **invalid** and ignored.

[Sample data]

Display the sample data of the selected determination.

[Determ. on/off]

Remove all of the results of the selected determination from the statistics. All of the entries of this line are marked with an asterisk (*), the statistics will be recalculated automatically.

[Result on/off]

Remove the selected result from the statistics. The result is marked with an asterisk (*), the statistics will be recalculated automatically.

**NOTICE**

If the mean value has been assigned to a common variable (see **CALC** command), then the determination must be recalculated manually (see chapter 23.8, page 193)

24.2 Deleting statistical data

In the dialog **Control**, you can delete all of the statistics data manually with the function **[Delete statistics]**. The statistics data is deleted automatically in the following cases:

- when all of the determinations of the determination series have been carried out and a new determination has been started afterwards.
- when a new method is loaded (even if it is the same method)
If the **sample table** or the **sample assignment table** is used, the statistics data is only deleted when another method is loaded.

24.3 Adding a determination to a determination series

You can use the function **[Add determ.]** to add an additional sample to a determination series, e.g. because a determination was faulty and had to be removed from the statistics.

Resultate / Statistik	
Aktuelle Methode: Me4155 3 von 3 (+2)	
Resultatname	Mittelwert
KFT Titer	5.0080 mg/mL
<div style="display: flex; justify-content: space-between;"> Bestimmung hinzufügen Details </div>	

The number of determinations by which the set statistics counter has been increased is shown above the list in brackets. This modification is also displayed in the main dialog and in the statistics report.



NOTICE

The number of samples defined in the method options does not, however, change as a result.

25 Result table

Main dialog: **Results ► Result table**

Results / Result table		
Sorted by Determ. time		6 of 100
No.	Identification 1	Result 1
1	Composite 5	4.7159 mg/mL
2	Composite 5	4.8095 mg/mL
3	Composite 5	4.6912 mg/mL
4	Honey1 #44	16.56 %
5	Honey1 #43	16.49 %
6	Honey1 #42	16.54 %

Buttons: Load/Save, Properties, Delete all, Delete, Details

The result table is suitable for displaying in detail the results of all of the determinations performed on a single day, for instance. You can save the results of up to 200 determinations. A maximum of 9 results from a single determination can be saved in the result table. You define which results from this table are to be saved using the **CALC** command. Because the size of the display is not sufficient to display all of the data at the same time, it is practical to print out the result table.

The sorting criterion and the number of determinations from which results are saved in the result table are displayed at the top. The list contains numbered lines; each line represents one determination. In the default settings, the first sample identification and the first calculated result of each determination are displayed. In the properties, you define which data is displayed in the columns and the criteria according to which the determinations are sorted (*see chapter 25.1, page 200*). In addition, you can define search filters to aid you in finding the desired determinations (*see "Filter criteria", page 202*).

[Load/ Save]

Load a saved result table (*see chapter 25.3, page 204*) or save the current result table (*see chapter 25.2, page 204*).

[Properties]

Configure the display of the columns and define filter criteria.

[Delete all]

Delete the result table completely.

[Delete]

Delete the selected line from the result table.

[Details]

Display details concerning the selected determination.

"Result table / Details" dialog

The following determination data is displayed in this dialog:

- **User**
User who carried out the determination.
- **Method**
Method with which the determination was carried out.
- **Determ. time**
Date and time at which the determination was carried out.
- Sample data (Identification 1 and 2 and sample size).
- All of the results of this determination saved in the result table.

25.1 Properties

Result table: **Properties**

In the **Result table / Properties** dialog, you can select the data to be displayed with each determination as well as the sorting criterion for the list.

Column 1

Selection of what is to be displayed in the first column.

Selection	Identification 1 Identification 2 Determ. time Result name Method Sample size User
Default value	Identification 1

Column 2

Selection of what is to be displayed in the second column.

Selection	Result 1...9 The numbering of the results corresponds to the order in which the results were calculated in the determination run.
-----------	---

Sort by

Selection according to which criterion the list is to be sorted.

Selection	Identification 1 Identification 2 Determ. time Result Method Sample size User
Default value	Determ. time

Identification 1

Sort alphabetically according to the first sample identification.

Identification 2

Sort alphabetically according to the second sample identification.

Determ. time

Sort chronologically according to the determination time.

Result

Sort according to the numerical value of the result.

Method

Sort alphabetically according to the method name.

Sample size

Sort according to the numerical value of the sample size.

User

Sort alphabetically according to the user who carried out the determination.

Sort sequence

Selection	ascending descending
Default value	descending

Max. number of lines

If the result table contains the quantity of lines defined here, then no further results can be saved. If this is the case, save the result table and create a new one.

Input range	10 to 200
Default value	100

[Filter]

Specify the filter criteria for the list.

[Delete automat.]

Define the settings for the automatic deletion of the result table (see "Deleting the result table automatically", page 203).

Filter criteria

Result table: **Properties** ► **Filter**

You can set filters in the **Properties / Filter** dialog. A filter is a rule defining what is to be displayed in the result table. If a filter has been set, this is indicated with a corresponding symbol in the **Results / Result table** dialog.

Filter

Define the criterion according to which the list is to be filtered. The only lines which will be displayed are those which fulfil the filter criterion.

Selection	no filter Identification 1 Identification 2 Date Result Method Sample size User
Default value	no filter

no filter

The list itself will not be filtered.

Identification 1

In the input field **is equal to**, select or enter the first sample identification according to which filtering is to be carried out.

Identification 2

In the input field **is equal to**, select or enter the second sample identification according to which filtering is to be carried out.

Date

In the input fields **Date from** ... **to**, specify the period of time according to which filtering is to be carried out.

Result

In the input fields **Value from ... to**, specify the value range according to which filtering is to be carried out.

Method

In the field **is equal to**, select or enter the method name according to which filtering is to be carried out.

Sample size

In the input fields **Value from ... to**, specify the value range according to which filtering is to be carried out.

User

In the field **is equal to**, select or enter the user according to whom filtering is to be carried out.

**NOTICE**

The filter can be applied to empty fields as well, for example, if no sample identification has been entered. In this case, select only the desired filter criterion and leave the fields below empty.

Deleting the result table automatically

Result table: **Properties ► Delete automat.**

In the **Properties / Delete automatically** dialog, you define when the contents of the result table is to be automatically deleted.

Delete oldest line only

on | off (Default value: **off**)

If this option is activated at the same time that the result table is full, the oldest entry will be deleted as soon as the results of a new determination are saved.

Time when result table is deleted:**Switch on**

on | off (Default value: **off**)

If this option is activated, then the result table will be deleted each time you switch on the Coulometer.

Start a new sample series

on | off (Default value: **off**)

If this option is activated, then the result table will be deleted when you start a new sample series (with activated autostart function!).

Save result table

on | off (Default value: **on**)

If this option is activated, the result table will be deleted as soon as you have saved it.

25.2 Saving the result table



NOTICE

When you save the result table, the complete determination will not be saved with each entry. If you wish to save the determination data for each determination, use the **Save determination automatically** function (see chapter 15.5.6, page 146).

25.3 Loading the result table

Proceed as follows to load a result table:

1 Opening the current result table

- In the results dialog, tap on **[Result table]**.
- Tap on **[Load/ Save]**.
- Tap on **[Load]**.

The selection of memory locations is displayed. This selection will be skipped if only one external storage medium is available.

- Select the memory location where the desired result table is saved.

The selection of file groups is displayed (see chapter 11.1, page 91). If only one group is available, then the list of the saved result tables will be displayed directly.

- Select the group with the desired result table.
- Tap on **[Show files]**.

The list with the saved result tables opens.

2 Selecting the result table

- Select the desired result table.

3 Loading the result table


- Tap on **[Load]**.

The result table is now loaded and will be displayed immediately.

26 Printing

Reports can be printed out as follows:

- Manually with the **[Print]** fixed key in this dialog. The fixed key is disabled while a determination is running.
- Automatically in the determination run with the command **REPORT** (see chapter 28.6, page 271).

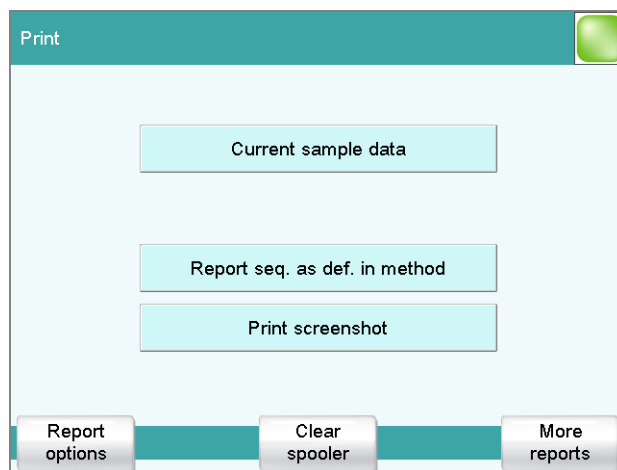

This screenshot shows the printing dialog associated with the main dialog. The sample data is entered in the main dialog, which is why you can print out the sample data from there in a context-sensitive manner with **[Current sample data]**.

The following buttons are also available in the other print dialogs:

[Report seq. as def. in method]

Print out all of the reports which are defined in the loaded method.

[Print screenshot]

Activate the "Print Screen" function. The current screen contents will be printed out the next time the fixed key **[Print]** is tapped on.

[Report options]

Defining the general settings for the report output.

[Clear spooler]

Delete all outstanding printing orders.


[More reports]

Open the list of available reports.

More reports

Fixed key **[Print]**: **More reports**

The dialog **Print / More reports** displays a list of all of the available reports.

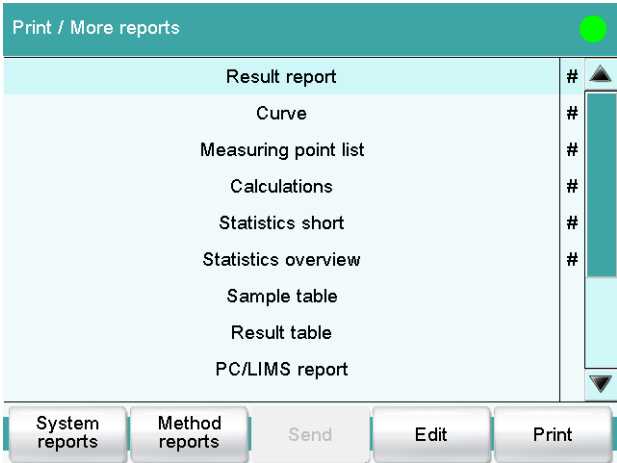

NOTICE

If you are using a **CUSTOM Neos** printer, you can print only the reports marked with a #. The designation on the reports is only visible once the **CUSTOM NEOS** printer is selected (see screenshot above).

[System reports]

Opening the list of reports for the system settings.



NOTICE

If you are using a **CUSTOM NEOS** printer, you cannot print system reports.

[Method reports]

Opening the list of reports for the current method.



NOTICE

If you are using a **CUSTOM NEOS** printer, you cannot print method reports.

[Send]

Create a PC/LIMS report.

[Edit]

Edit the settings of the selected report.

[Print]

Print the selected report.

26.1 General report options

Fixed key []: **Report options**

The settings for the report printout can be adjusted in the dialog **Print / Report options** to indicate, for example, whether a report header, a signature line or a frame is to be printed.

Report header

Defining the output of the report header. The report header contains general information such as device type, serial number, program version and the printing date.



NOTICE

You can also define your own report header, which is printed in addition to this report header (**System ► Templates ► Report header**).

Selection	off once on each page
Default value	on each page

off

The report header will not be printed.

once

The report header will only be printed on the first page.

on each page

The report header will be printed on every page.

Signature line

Output of a special line for date and signature. This line will be printed at the very bottom of every page.

Selection	off once on each page
Default value	off

off

No signature line will be printed.

once

The signature line will only be printed on the last page.

on each page

The signature line will be printed on every page.



Frame

on | off (Default value: **on**)

If this parameter is activated, a frame is printed as lateral marking.

26.2 Settings of the individual reports

Fixed key []: **More reports ► Report ► Edit**

Command REPORT: **Report ► Edit**

Settings can be edited for the following reports (for details, see online help):

- **Result report**
- **Curve**
Definition of the curve size, display of the individual measuring points and grid lines, etc.
- **Measuring point list**
Definition of the method command to which the measuring point list is to be printed.
- **Calculations**
Definition of the accuracy with which the variables used as well as the results calculated are printed.
- **Statistics short**
Definition of the time at which the report is to be printed (with each determination or only at the end of a sample series).
- **Statistics overview**
Definition of the time at which the report is to be printed (with each determination or only at the end of a sample series).
- **Sample table**
- **Result table**
Definition of the time at which the report is to be printed (with each determination, at the end of a sample series or at the end of the sample table).
- **Used devices**
- **Variables**
- **Monitoring**

26.3 List of all printable reports

The following reports can be printed out with the fixed key []:



NOTICE

If you have selected a **CUSTOM NEOS** printer, you can print only the reports that are marked with a # on the display.

Report	Contents
Result report	# Report with determination properties, sample data, calculated results, etc. If there are multiple determinations, the statistics will also be printed out.
Curve	# Curve report. The settings for the curve print-out are adopted from the curve options of the live display.
Measuring point list	# Measuring point list report.
Calculations	# Details concerning the calculations carried out (parameter settings in accordance with the CALC command and the calculated results).
Used devices	The devices used for the determination, as displayed in the More determination data / Properties dialog.
Variables	All determination variables, as displayed in the results dialog.
Monitoring	Details for the monitored quantities.
Statistics	
Statistics short	# Summary of the statistics calculations. The number of determinations, the mean value, and the absolute and relative standard deviations are printed out for each result.
Statistics overview	# Detailed statistical overview. The sample data and all individual results are printed out for each determination. The number of determinations, the mean value, and the absolute and relative standard deviations are printed out for each result.
PC/LIMS	




Report	Contents
PC/LIMS report	Machine-readable report with all of the important data for a determination. This report can be saved as a TXT or UTF-8 file or sent to a terminal program or a LIMS via an RS-232 interface. The definition is carried out in the device manager (<i>see chapter 10.3.2, page 68</i>).
Sample data	
Current sample data	Sample data of the current determination, <i>context-sensitive only</i> from the main dialog.
Sample table	List of all determinations in the sample table with the respective sample data, as entered in the sample table.
Result table	
Result table	List of all determinations in the result table with results and with the determination data, as saved in the result table.
More reports ► Method reports	
Method sequence	Method properties and the list of all method commands.
Parameters full	Method properties and options, all method commands with all parameters. All parameters which no longer have their default settings will be printed in bold . All parameters which have been modified in comparison to the stored version of the method are indicated by an *.
Titration & measuring param.	Method properties; titration and measuring commands with all parameters. All parameters which no longer have their default settings will be printed in bold . All parameters which have been modified in comparison to the stored version of the method are indicated by an *.
Modified parameters	Method properties, all method parameters which have been modified in comparison to the stored version of the method.
Non default parameters	Method properties, all method parameters which no longer have any default settings.
More reports ► System reports	
System settings	
System settings	Settings for acoustic signals and accuracy of the measured value display.

Report	Contents
Dialog options	Settings for routine dialog and expert dialog.
User list	List with all of the users configured in the system, together with their data, <i>context-sensitive only</i> from the user list.
Login options	Login options, <i>context-sensitive only</i> from the respective dialog.
Sensors	
Sensor list	List of all the sensors configured in the system.
All sensor data short	The most important sensor data for all sensors.
All sensor data full	All of the sensor data for all of the sensors (name, working life, etc.).
Sensor data short	The most important sensor data of the selected sensor, <i>context-sensitive only</i> from the editing dialog.
Sensor data full	All of the sensor data of the selected sensor, <i>context-sensitive only</i> from the editing dialog.
Device manager	
Device list	List of all devices configured in the system.
All device properties	Properties of all the devices configured in the system.
Device properties	Properties of the selected device, <i>context-sensitive only</i> from the editing dialog.
GLP manager	
GLP data	All data stored in the GLP manager.
Common variables	
Common variable list	List of all the common variables defined in the system, together with their most important data (name, value, status).
All common variable properties	Properties of all common variables (name, value, validity, status).
Common variable properties	Properties of the selected common variable (name, value, validity, status), <i>context-sensitive only</i> from the editing dialog.
Local Common variable list	List with the common variables available at the time of the determination, together with their most important data (name, value, status), <i>context-sensitive only</i> from the dialog "More determination data / Local common variables".



Report	Contents
Local common var. properties	Properties of the selected common variable (name, value, validity, status), <i>context-sensitive only</i> from the dialog "More determination data / Local common variables".
Templates	
Templates sample data	Sample identification list and sample assignment table.
Result template list	List with all of the user-generated result templates.
All result templates details	Details of the all of the user-generated result templates (calculation formula, result options, note).
Result template details	Details of the selected result template, <i>context-sensitive only</i> from the editing dialog.
Input/Output lines	List with all of the defined input and output lines at the remote interface (name, bit pattern).

27 Manual control

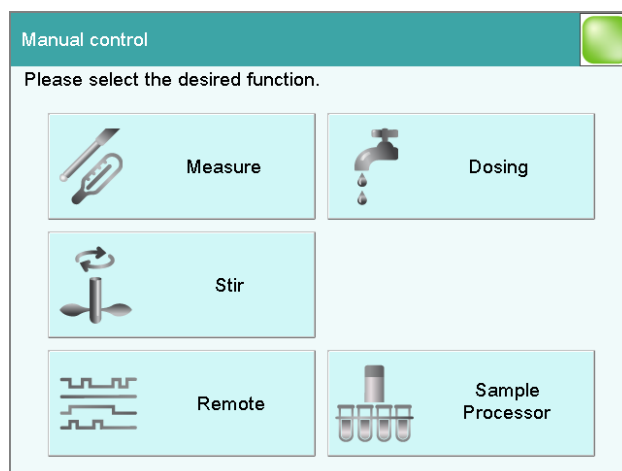

The following functions are available in the manual control:

[Measure]

Carrying out manual measurements (*see chapter 27.2, page 215*).

Measuring modes:

- **Ipol** (voltametric measurement with selectable polarization current)
- **Upol** (amperometric measurement with selectable polarization voltage)
- **T** (temperature measurement)

[Dosing]



NOTICE

Only active if a Dosino or dosing unit is connected.

Manual dosing (*see chapter 27.3, page 217*).

The following dosing functions are available:

- Preparing the exchange unit or dosing unit
- Emptying the dosing unit
- Filling the dosing cylinder of the exchange unit/dosing unit
- Dosing a specified volume
- Dosing continuously

[Stir]

Controlling the stirrer manually (*see chapter 27.4, page 223*).


[Remote]

Scanning the input lines on the remote interface and setting the output lines (see chapter 27.5, page 224).


27.1 Opening and closing the manual control

Opening the manual control

1 Opening the dialog window

- Tap on the fixed key [].

Manual control opens:

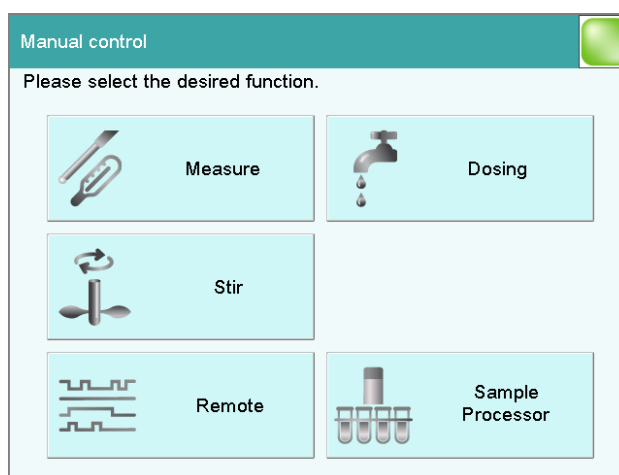

2 Select the function

- Tap on the button for the desired function.

If the function can be carried out by several devices, a dialog to select the device will appear first; otherwise the dialog for the selected function will appear immediately.

Closing the manual control

- #### 1
- Tap on the fixed key [].



NOTICE

Manual control can also be exited when a manual sequence has been started but has not yet ended. The fact that a sequence has been started in manual control can be recognized by the corresponding symbol in the title line (see chapter 6.2.3, page 34).

27.2 Measuring

Manual control ► Measure

Manual control / Measure

Sensor: Metal electrode

Measuring input: 1

Measuring mode: Ipol

Temperature: 25 °C

Pol. current: - 5.0 + μA

Info sensor Start

Manual measurements can be carried out with the function **[Measure]**.

Proceed as follows:

1 Selecting a sensor

- Select the desired sensor from the sensor list.
All of the sensors available in the sensor list are displayed. Sensors are defined under **System ► Sensors**.
- Select the measuring input at which the sensor is connected.

2 Select a measuring mode

- Select the desired measuring mode.
Only those measuring modes are displayed which are advisable for the selected sensor.

3 Enter a measuring temperature

- Enter the measuring temperature if no temperature sensor is connected. If a temperature sensor is connected, then the temperature will be measured automatically.

4 Starting the measurement

- Tap on **[Start]**.

The measured value display is opened:

The ongoing measurement is canceled with **[Stop]**. The label switches once again to **[Start]**.

27.2.1 Parameter description

"Manual control / Measure" dialog

Sensor

Selection of the sensor from the sensor list. Sensors are defined under **System ► Sensors**.

Selection	Selection of configured sensors
-----------	--

Measuring input

Selection of the measuring input the sensor is connected to. The selection depends on whether the control device has one or two measuring interfaces.

Selection	1 2
Default value	1

Measuring mode

Selection of the measuring mode. Only those measuring modes are displayed which are advisable for the selected sensor.

Selection	Ipol Upol T
-----------	------------------------

Ipol

voltametric measurement with selectable polarization current

Upol

amperometric measurement with selectable polarization voltage

T

temperature measurement

Temperature

Temperature entered manually. If a temperature sensor is connected, then the temperature will be measured continuously.

Input range	-20.0 to 150.0 °C
Default value	25.0 °C

[Info sensor]

Display information on the sensor.

27.3 Dosing

Manual control ► Dosing

The following dosing functions are available:

- Preparing the exchange or dosing unit (*see chapter 27.3.3, page 221*)
- Emptying the dosing unit (*see chapter 27.3.4, page 222*)
- Filling the dosing cylinder of the exchange unit/dosing unit (*see chapter 27.3.5, page 222*)
- Dosing a specified volume (*see chapter 27.3.2, page 220*)
- Dosing continuously (*see chapter 27.3.1, page 218*)

Selecting the dosing device

If several dosing devices are connected to a control instrument, then the dialog for selecting a dosing device will be opened.

[Dosing device X]

Selection of the dosing device. The MSB connector and the type of the connected dosing device are displayed.

Various information is shown next to the button:

- Titrant
For dosing devices of the type 8XX with an intelligent exchange unit or dosing unit the titrant name is read off the data chip.
- Cylinder volume
For dosing devices of the type 8XX with an exchange unit without data chip only the cylinder volume is displayed.
- ---
For exchange or dosing units which are not or not correctly attached.
- Manual busy
When a dosing function is being carried out.

[All dosing devices]

Carrying out the functions **[Prepare]**, **[Empty]** and **[Fill]** at the same time with several dosing devices of the control device.

Selecting the dosing function

If only one dosing device is connected, then the properties of this dosing device will be shown directly. This information varies according to the type of dosing device:

- Type 8XX with integrated data chip:
If an exchange or dosing unit with integrated data chip is attached, the data stored on the data chip is displayed.
- Type 8XX without integrated data chip:

[Empty]

Empty the cylinder and the tubings of the dosing unit (*see chapter 27.3.4, page 222*).

[Prepare]

Rinse the cylinder and the tubings of the exchange / dosing unit (*see chapter 27.3.3, page 221*).

[Fill]

Fill the cylinder of the exchange / dosing unit (*see chapter 27.3.5, page 222*).

[Dosing fixed vol.]

Dose a specified volume (*see chapter 27.3.2, page 220*).

[Dosing]

Dose continuously (*see chapter 27.3.1, page 218*).

27.3.1 Continuous dosing

Manual control: **Dosing ► Dosing**

With the function **[Dosing]**, dosing is continuous for as long as you hold the button **[Dose]** pressed down.

Proceed as follows:

1 Configure the dosing function**NOTICE**

- The dosing and filling rates should be decreased for viscous liquids.
- The maximum dosing and filling rates depend on the cylinder volume (*see chapter 31, page 310*).

- Enter the dosing rate.
- Enter the filling rate.

2 Start dosing

- Tap on **[Dose]**.

The volume dosed is displayed. When the volume of one cylinder has been dosed, the dosing cylinder will be refilled automatically.


3 Fill the cylinder

- Tap on **[Fill]**.

The dosing cylinder is filled. The displayed volume value will be reset to 0.0000 mL.



NOTICE

If you exit the dialog with the fixed key [>], the dosing cylinder is filled automatically.

Parameter description

Dosing rate

Rate at which it is dosed.

Input range	0.01 to 166.00 mL/min
Selection	maximum dynamic
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 31, page 310*).

dynamic

This setting can only be selected when the dosing cylinder is filled. The dosing is being carried out faster and faster until the maximum dosing rate is reached (starting with 1 mL/min and then doubling the dosing rate every 1.5 s).

Filling rate

Rate at which the dosing cylinder is filled.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum filling rate depends on the cylinder volume (*see chapter 31, page 310*).

27.3.2 Dosing fixed volumes

Manual control: **Dosing ► Dosing fixed vol.**

You can dose a particular volume with the **[Dosing fixed vol.]** function.

Proceed as follows:

1 Configure the dosing function



NOTICE

- The dosing and filling rates should be decreased for viscous liquids.
- The maximum dosing and filling rates depend on the cylinder volume (*see chapter 31, page 310*).

- Enter the desired volume.
- Enter the dosing rate.
- Enter the filling rate.

2 Start dosing

- Tap on **[Start]**.

The volume display is shown. When the volume of one cylinder has been dosed, the dosing cylinder will be refilled automatically.

Continuous dosing is paused with **[Hold]**. The label changes to **[Continue]**. It can be used to resume the run.

3 Fill the cylinder

With the default settings (*see "Fill automatic.", page 221*) the dosing cylinder is filled automatically.

Otherwise:

- Tap on **[Fill]**.

The dosing cylinder is filled. The displayed volume value will be reset to 0.0000 mL.

Parameter description

Volume

Volume which is dosed.

Input range	0.00000 to 99999.9 mL
Default value	0.10000 mL

Dosing rate

The rate at which dosing takes place.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 31, page 310*).

Filling rate

Rate at which the dosing cylinder is filled.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum filling rate depends on the cylinder volume (*see chapter 31, page 310*).

Fill automatic.

on | off (Default value: **on**)

If this parameter is activated, then the dosing cylinder is being automatically filled at the end of the dosing.

27.3.3 Preparing

Manual control: **Dosing ► Prepare**

The **Prepare** function is used to rinse and fill the cylinder and tubings of the exchange or dosing unit air bubble-free. You should carry out this function before the first determination or once a day.

If the titrant is selected (*see "Selecting the dosing function", page 217*), the parameters defined for the titrant for preparing/emptying and the tubing parameters will be used (dosing unit only). If the titrant is not selected, default parameters will be used (*see Chapter 31.1.2, page 310 and Chapter 31.2.2, page 311*).

The following two figures show the live display of an exchange unit and a dosing unit:

27.3.4 Emptying

Manual control: **Dosing ► Empty**



NOTICE

The function **Empty** is possible only with dosing units.

The cylinders and the tubings of the dosing unit are emptied with this function.

If the titrant is selected (see "Selecting the dosing function", page 217), the parameters defined for the titrant for preparing/emptying and the tubing parameters will be used. If the titrant is not selected, default parameters will be used (see Chapter 31.2.2, page 311 and Chapter 31.1.2, page 310).

27.3.5 Filling

Manual control: **Dosing ► Fill**

You can use the function **[Fill]** to fill the dosing cylinder of the exchange unit/dosing unit manually. The filling rate cannot be configured; the maximum filling rate will be used.

27.3.6 Replacing reagent



NOTICE

You can easily change the reagent in a dosing unit without no contact with the chemicals using the functions **Prepare** and **Empty**.


Proceed as follows:

- 1 Execute the function **[Empty]**.
- 2 Attach the dosing unit to the bottle with a suitable solvent.
- 3 Execute the function **[Prepare]**.
The last traces of the reagent in the cylinder and tubings are removed with the solvent.
- 4 Execute the function **[Empty]** again.
- 5 Attach the dosing unit to the bottle with the new reagent.

- 6** Execute the function **[Prepare]**.
The dosing unit can now be used.

27.4 Stirring

Manual control ► Stir

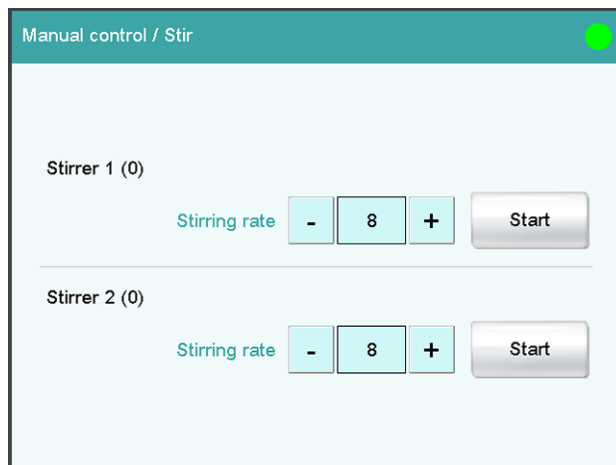

The MSB connector and type of stirrer is indicated for each connected stirrer.

You can control a connected stirrer manually with the function **[Stir]**.

Proceed as follows:

1 Setting the stirring rate

- Tap on the **[-]** or **[+]** button.
Each time one of the buttons is tapped, the stirring rate is reduced or increased by one level.

2 Switching on the stirrer

- Tap on **[Start]**.

The stirrer is started.

3 Switching off the stirrer

- Tap on **[Stop]**.

The stirrer is stopped.

Parameter description

Stirring rate

Setting the stirring rate. It can be set in steps of -15 to $+15$. The default setting **8** corresponds to 1000 rpm. The formula for calculating the rotational speed is specified in *chapter 31.3, page 311*.

The algebraic sign of the stirring rate changes the direction in which the stirring is done. When the stirrer is viewed from above, this means:

- "+": counterclockwise rotation
- "-": clockwise rotation

Input range	-15 to 15
Default value	8



NOTICE

If a stirrer is used in the determination run, it can still be manually controlled at the same time. For example, you can use the manual control to reduce the stirring rate of a stirrer which has been started in the determination run.

27.5 Remote

Manual control ► Remote

Manual control / Remote ●

	Inputs	Outputs	
Remote Box 1	00000000	00000000000000	▼
Remote Box 2	00000000	00000000000000	▼

With the function **[Remote]**, you can define manual output signals to the remote interface of a connected Remote Box. It is not possible to define input signals; the current status of the input lines is, however, displayed.

Inputs

Current status of the input lines.

Outputs

Selection of the signal out of the templates or entering the required bit pattern. Templates are defined under **System ► Templates ► Output lines**.

Entering the bit pattern:

- 0 = line inactive
- 1 = line active
- * = retain line status
- p = set pulse

The output lines are always numbered from right to left, i.e. with the signal **00000000000001** line 0 is set. With a pulse, the length is set to 200 ms. If you wish to set pulses with other lengths, you have to define a corresponding template.


Entry	Bit pattern containing exactly 14 characters or a max. of 24 characters for the name of the template
Default value	00000000000001
Selection	Selection of the templates defined

[Set]

Set a defined output signal.

27.6 885 Compact Oven SC

Manual control ► Sample Processor

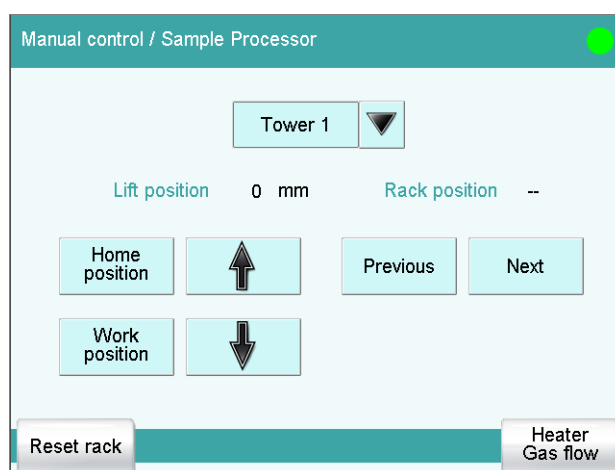


- The sample variable is reset to the value 1.

[Heater Gas flow]

Display, control and settings of the current temperature and current gas flow.

27.6.1 Controlling the heating/gas flow

Manual control: **Sample Processor ► Heater Gas flow**

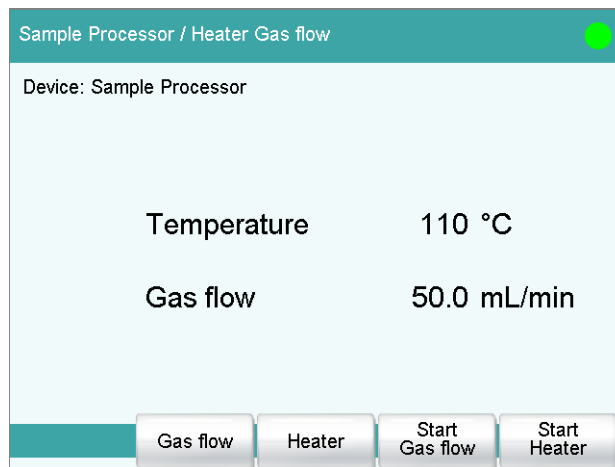

You can see, control and set the current temperature and current gas flow in the **Sample Processor / Heater Gas flow** dialog.

[Gas flow]

Edit parameters for the gas flow.

[Heater]

Edit parameters for the heater.

[Start / Stop Gas flow]

Starts or stops the gas flow.

[Start / Stop Heater]

Starts or stops the heater.

Settings for the gas flow

Flow rate

Flow rate for the gas flow.

Input range	10 to 150 mL/min (Increment: 1)
Default value	50 mL/min

Gas supply

Selection of the gas supply.



Selection	pump valve
Default value	pump

pump

The integrated air pump is used for the gas flow.

valve

The integrated inlet valve is used for the gas flow. The inlet valve in the 885 Compact Oven SC is connected to compressed air or nitrogen.

Gas type

Gas type of the carrier gas used.

Selection	air nitrogen
Default value	air

Settings for the oven temperature**Temperature**

Temperature to which the oven is heated.

Input range	50 to 250 °C (Increment: 1)
Default value	110 °C

28 Parameters



NOTICE

For most of the numerical parameters, the result of a calculation can also be utilized in place of a number. A description of this can be found in the *appendix* of the manual.

28.1 Coulometric Karl Fischer titrations (KFC)

28.1.1 Control parameters

The control parameters for the endpoint are defined under **[Control parameters]**.

Endpoint at

Measured value for the endpoint.

The values that are actually measurable may vary from the values specified here, depending on the hardware. The binding measurement value ranges can be found in the manual, in the "Technical specifications" chapter.

Input range	-1250.0 to 1250.0 mV
Default value	50.0 mV
Selection	off

Titration rate

Three predefined sets of parameters can be selected for the titration rate.

Selection	slow optimal fast user
Default value	optimal

slow

For samples with a low water content or samples which release their moisture only slowly.

optimal

For all standard titrations. The parameters have been optimized for the most frequent applications.

fast

For uncritical samples with high water content.

user

The individual titration parameters can be modified.

The settings of the individual titration rates are listed in *table 9, page 230*.

User-defined parameters

These parameters are only accessible when **Titration rate = user**.

Dynamics

This parameter defines the control range before the specified endpoint. Within the control range, the iodine is generated step by step, the generation is finely controlled. The closer the endpoint, the slower the iodine is generated until the rate defined under **Min. rate** is reached. The larger the control range, the slower the titration. Outside the control range, iodine is being continuously generated, the rate is defined under **Max. rate**.

The values that are actually measurable may vary from the values specified here, depending on the hardware. The binding measurement value ranges can be found in the manual, in the "Technical specifications" chapter.

Input range	0.1 to 1250.0 mV
Default value	70.0 mV
Selection	off

Max. rate

Rate at which iodine is generated outside the control range.

Input range	1.5 to 2241.0 µg/min
Selection	maximum
Default value	maximum

Min. rate

Rate at which iodine is generated at the very beginning of the titration and in the control range at the end of the titration. This parameter has a vital influence on the titration rate and thus also on the accuracy. The smaller the selected minimum rate, the slower the titration.

Input range	0.3 to 999.0 µg/min
Default value	15.0 µg/min

Table 9 Default values of the predefined titration rates for KFC

	Titration rate		
	slow	optimal	fast
Dynamics	120.0 mV	70.0 mV	30.0 mV
Max. rate	1000.0 µg/min	maximum	maximum
Min. rate	0.3 µg/min	15.0 µg/min	30.0 µg/min

Stop criterion

Stop criterion

The titration is stopped when the endpoint has been reached and this stop criterion has been fulfilled. If no stop criterion has been selected then the titration will not be stopped. The stop conditions (*see chapter 28.1.3, page 232*) always lead to a stop, even if the stop criterion has not been reached.

Selection	drift time rel. drift off
Default value	rel. drift

drift

The titration is stopped when the stop drift has been reached.

time

The titration is stopped if the endpoint has been exceeded during a certain time period (**Delay time**).

rel. drift

The titration is stopped when the sum of the drift at the start of the titration and the relative stop drift has been reached.

off

The titration will not be stopped until the stop conditions have been fulfilled.

Stop drift

This parameter can only be edited with **Stop criterion = drift**.

The titration is stopped when the endpoint and the stop drift have been reached.

Input range	1 to 999 µg/min
Default value	5 µg/min

Delay time

This parameter can only be edited with **Stop criterion = time**.

When the endpoint has been reached, the specified time is allowed to elapse after the last dosing and the titration is then stopped.

Input range	0 to 999 s
Default value	10 s

Relative stop drift

This parameter can only be edited with **Stop criterion = rel. drift**.

The titration is stopped when the endpoint and the sum of the drift at the start of the titration and the relative stop drift have been reached.



Input range	1 to 999 µg/min
Default value	5 µg/min

28.1.2 Titration parameters

Under [**Titration parameters**], the parameters influencing the run of the entire titration are defined.

Pause

Waiting time, for example for dissolving the sample. During this time no iodine is generated.

Input range	0 to 999999 s
Default value	0 s

Extraction time

Minimum duration of the titration. The titration will not be stopped during the extraction time, even if the endpoint has already been reached. The titration is however stopped if a stop condition is fulfilled during this time (see chapter 28.1.3, page 232). Entering an extraction time may e.g. be advisable if a Karl Fischer oven is used.

Input range	0 to 999999 s
Default value	0 s

Temperature

Temperature entered manually. The temperature is being continuously measured when a temperature sensor is connected and when **Temp. measurement = automatic** or **continuous** is defined (see sensor dialog of the command).

Input range	-20.0 to 150.0 °C
Default value	25.0 °C

Time interval MP

Time interval for entering a measuring point in the measuring point list. The measuring point list is limited to 1000 measuring points.

Input range	0.1 to 999999.0 s
Default value	2.0 s

28.1.3 Stop conditions

The conditions for canceling the titration are defined under [**Stop conditions**], if this does not occur automatically. This could be the case when the endpoint set is not reached or if the stop criterion (see "Stop criterion", page 231) is not fulfilled.

Stop time

The titration is stopped when the specified time has elapsed since the start of the titration.


Input range	1 to 999999 s
Selection	off
Default value	off

28.1.4 Conditioning

The conditions required for conditioning are defined under **[Conditioning]**.

Conditioning

on | off (Default value: **on**)

If this parameter is activated, then the first time the titration is started the working medium will be titrated to the endpoint with the specified control parameters. The status is kept stable. The actual method run does not begin until [] has been pressed once more.

Start drift

If the measured drift is smaller than this value for a certain time (so-called stabilizing time), **Conditioning OK** is displayed and the titration can be started. The stabilizing time is defined in the dialog **Conditioning / Conditioning options**.

Input range	1 to 999 µg/min
Default value	20 µg/min

Drift correction

The endpoint amount can be corrected taking the drift value into account. For this, the drift is multiplied with the drift correction time and this value is then subtracted from the endpoint amount. The drift correction time is the time interval between the end of conditioning and the end of the determination.

Selection	auto manual off
Default value	auto

auto

The value of the current drift is automatically applied at the start of the titration.

manual

If the drift is known throughout a longer period of time, this can be entered manually.

off

No drift correction takes place.

Cond. stop time

Maximum permissible time over which conditioning may take place. Conditioning is stopped when the specified time has elapsed.

Input range	1 to 999999 s
Selection	off
Default value	off

Delay reconditioning

on | off (Default value: **off**)

If this parameter is activated, the conditioning will not start before all the method commands have been processed. Otherwise, the conditioning will immediately start after the titration command.

**NOTICE**

This parameter must be activated if commands follow which must not run simultaneously with the conditioning.

Example: Emptying the titration cell followed by adding new working medium.

28.1.5 Cell

The parameters for the titration cell are defined under **[Cell]**.

Generator electrode

Type of generator electrode.

Selection	without diaphragm with diaphragm
Default value	without diaphragm

without diaphragm

Generator electrode without diaphragm.

with diaphragm

Generator electrode with diaphragm.

Generator current

Polarization current at the generator electrode.

Selection	100 mA 200 mA 400 mA auto
-----------	--

400 mA

Default value, when **Generator electrode = without diaphragm**.

- Generator electrode:
 - Is the electrode connected?
 - Is the generator system defective?
 - Is the resistance of the sample solution very high?

The electrode check is carried out continuously during the execution of this command.

Temp. measurement

Type of temperature measurement.

Selection	continuous automatic off
Default value	automatic

continuous

A temperature sensor must be connected. The temperature is measured continuously.

automatic

If a temperature sensor is connected then the temperature will be measured continuously. Otherwise, the temperature entered manually will be used (see dialog of the titration and measuring parameters).

off

The temperature will not be measured. The temperature entered manually is used (see dialog of the titration and measuring parameters).

28.1.7 Stirrer

The parameters for the stirrer are edited under **[Stirrer]**.

Stirrer

Selection of the MSB connector the stirrer is connected to. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4 off
Default value	1

off

No stirrer will be used.

Stirring rate

Setting the stirring rate. It can be set in steps of -15 to $+15$. The default setting **8** corresponds to 1000 rpm. The formula for calculating the rotational speed is specified in *chapter 31.3, page 311*. The optimum stirring rate can be tested in the manual control.

The algebraic sign of the stirring rate changes the direction in which the stirring is done. When the stirrer is viewed from above, this means:

- "+": counterclockwise rotation

User-defined parameters

These parameters are only accessible when **Titration rate = user**.

Dynamics

This parameter defines the control range before the specified endpoint. Within the control range, the bromine is generated step by step, the generation is finely controlled. The closer the endpoint, the slower the bromine is generated until the rate defined under **Min. rate** is reached. The larger the control range, the slower the titration. Outside the control range, bromine is continuously generated, the rate is defined under **Max. rate**.

The values that are actually measurable may vary from the values specified here, depending on the hardware. The binding measurement value ranges can be found in the manual, in the "Technical specifications" chapter.

Input range	0.1 to 1250.0 mV
Default value	400.0 mV
Selection	off

Max. rate

Rate at which bromine is generated outside the control range.

Input range	10.0 to 19876.0 µg/min
Default value	1000.0 µg/min
Selection	maximum

Min. rate

Rate at which bromine is generated at the very beginning of the titration and in the control range at the end of the titration. This parameter has a vital influence on the titration rate and thus also on the accuracy. The smaller the selected minimum rate, the slower the titration.

Input range	1.0 to 999.0 µg/min
Default value	25.0 µg/min

Table 10 Default values for the predefined titration rates for BRC

	Titration rate		
	slow	optimal	fast
Dynamics	400.0 mV	400.0 mV	400.0 mV
Max. rate	500.0 µg/min	1000.0 µg/min	2000.0 µg/min
Min. rate	15.0 µg/min	25.0 µg/min	50.0 µg/min

Stop criterion

Stop criterion

The titration is stopped when the endpoint has been reached and this stop criterion has been fulfilled. If no stop criterion has been selected then the titration will not be stopped. The stop conditions (*see chapter 28.2.3, page 242*) always lead to a stop, even if the stop criterion has not been reached.

Selection	drift time rel. drift off drift & time
Default value	drift & time

drift

The titration is stopped when the stop drift has been reached.

time

The titration is stopped if the endpoint has been exceeded during a certain time period (**Delay time**).

rel. drift

The titration is stopped when the sum of the drift at the start of the titration and the relative stop drift has been reached.

off

The titration will not be stopped until the stop conditions have been fulfilled.

drift & time

The titration is stopped when the stop drift has been exceeded during the time set and the endpoint has been reached.

Stop drift

This parameter can only be edited with **Stop criterion = drift** or **drift & time**.

The titration is stopped when the endpoint and the stop drift have been reached.

Input range	1 to 999 µg/min
Default value	15 µg/min

Delay time

This parameter can only be edited with **Stop criterion = time** or **drift & time**.

When the endpoint has been reached, the specified time is allowed to elapse after the last dosing and the titration is then stopped.

Input range	0 to 999 s
Default value	40 s

Relative stop drift

This parameter can only be edited with **Stop criterion = rel. drift**.

The titration is stopped when the endpoint and the sum of the drift at the start of the titration and the relative stop drift have been reached.

Input range	1 to 999 µg/min
Default value	5 µg/min

28.2.2 Titration parameters

Under **[Titration parameters]**, the parameters influencing the run of the entire titration are defined.

Pause

Waiting time, for example for dissolving the sample. During this time no bromine is generated.

Input range	0 to 999999 s
Default value	0 s

Extraction time

Minimum duration of the titration. The titration will not be stopped during the extraction time, even if the endpoint has already been reached. The titration is however stopped if a stop condition is fulfilled during this time (see chapter 28.2.3, page 242).

Input range	0 to 999999 s
Default value	0 s

Temperature

Temperature entered manually. The temperature is being continuously measured when a temperature sensor is connected and when **Temp. measurement = automatic** or **continuous** is defined (see sensor dialog of the command).

Input range	-20.0 to 150.0 °C
Default value	25.0 °C

Time interval MP

Time interval for entering a measuring point in the measuring point list. The measuring point list is limited to 1000 measuring points.

Input range	0.1 to 999999.0 s
Default value	2.0 s

auto

The value of the current drift is automatically applied at the start of the titration.

manual

If the drift is known throughout a longer period of time, this can be entered manually.

off

No drift correction takes place.

Drift value

This parameter can only be edited with **Drift correction = manual**.

Drift for manual drift correction.

Input range	0.0 to 999.9 µg/min
Default value	0.0 µg/min

Measured value display

on | off (Default value: **off**)

If this parameter is activated, the currently measured value is displayed during the conditioning.

Conditioning options [Cond. options]**Stabilizing time**

Waiting time, during which the measured drift has to be smaller than the start drift defined until **Conditioning OK** is displayed. The start drift is defined in the dialog **Edit command / Conditioning**.

Input range	0 to 999999 s
Default value	0 s

Cond. stop time

Maximum permissible time over which conditioning may take place. Conditioning is stopped when the specified time has elapsed.

Input range	1 to 999999 s
Selection	off
Default value	off

Delay reconditioning

on | off (Default value: **off**)

If this parameter is activated, the conditioning will not start before all the method commands have been processed. Otherwise, the conditioning will immediately start after the titration command.



NOTICE

This parameter must be activated if commands follow which must not run simultaneously with the conditioning.

Example: Emptying the titration cell followed by adding new working medium.

28.2.5 Cell

The parameters for the titration cell are defined under **[Cell]**.

Generator electrode

Type of generator electrode.

Selection	without diaphragm with diaphragm
Default value	with diaphragm

without diaphragm

Generator electrode without diaphragm.

with diaphragm

Generator electrode with diaphragm.

Generator current

Polarization current at the generator electrode.

Selection	100 mA 200 mA 400 mA auto
Default value	auto

auto

The current is adapted to the conductivity of the reagent and automatically reduced near the endpoint.

Reagent monitoring

Selection of the reagent from the list of reagents. The selection depends on the titration mode. Reagents are defined under **System ► Reagents**.

Selection	Selection of configured reagents off
Default value	off

off

The reagent monitoring is not active.

28.2.6 Sensor

The parameters for the sensor are edited under **[Sensor]**.

Sensor

Selection of the sensor from the sensor list. The selection depends on the measuring mode. Sensors are defined under **System ► Sensors**. You can also enter a sensor name which is not contained in the sensor list. When a determination is started there is a check whether the sensor is contained in the sensor list.

Selection	Selection of configured sensors
-----------	--

I(pol)

The polarization current is the current that is applied to a polarizable electrode during voltametric measurement.

The values that are actually adjustable may vary from the values specified here, depending on the hardware. The binding values can be found in the manual, in the "Technical specifications" chapter.

Input range	-125.0 to 125.0 μA (Increment: 2.5)
Selection	-1.0 μA 1.0 μA
Default value	1.0 μA

Electrode check

on | off (Default value: **off**)

For polarizable electrodes and the generator electrode, an electrode check can be carried out. The following will be tested:

- Polarizable electrodes
 - Is the electrode connected?
 - Is there a short circuit?

The electrode check is carried out when this command is started.

- Generator electrode:
 - Is the electrode connected?
 - Is the generator system defective?
 - Is the resistance of the sample solution very high?

The electrode check is carried out continuously during the execution of this command.

Temp. measurement

Type of temperature measurement.

Selection	continuous automatic off
Default value	automatic

28.3 Measurements (MEAS)

28.3.1 Measuring parameters

Under **[Measuring parameters]**, the parameters influencing the run of the entire measurement are defined.

Signal drift

Maximum permissible drift for the measured value acceptance, i.e. maximum change of the measured value per minute. For this value, a suitable maximum waiting time is automatically calculated except you already have modified this waiting time.



NOTICE

A constant measured value is often only reached after a certain time, as mixing and the reaction itself require a certain time. The response time of an electrode can also increase with time, i.e., reaching a constant measured value takes longer and longer. Drift-controlled measured value acceptance is particularly advisable in such cases, as the measured values are only accepted when equilibrium has almost been reached.

Measuring mode I_{pol}:

Input range	0.1 to 999.0 mV/min
Default value	10.0 mV/min
Selection	off

off

Measured value acceptance will take place after the maximum waiting time has elapsed.

Measuring mode U_{pol}:

Input range	0.01 to 99.90 µA/min
Default value	10.00 µA/min
Selection	off

off

Measured value acceptance will take place after the maximum waiting time has elapsed.

Measuring mode T:

Input range	0.1 to 999.0 °C/min
Default value	0.5 °C/min
Selection	off

**off**

Measured value acceptance will take place after the maximum waiting time has elapsed.

Min. waiting time

The measured value is not accepted until the minimum waiting time has elapsed, even if the signal drift has already been reached. The minimum waiting time is only important for drift-controlled measurements.

Input range	0 to 999999 s
Default value	0 s

Max. waiting time

If the signal drift has been switched off or has not yet been reached, then the measured value will be accepted when the maximum waiting time has elapsed. As default value, a suitable waiting time for the signal drift is automatically calculated.

Input range	0 to 999999 s
Selection	The default value depends on the measuring mode. off

off

The measurement is continued endlessly.

Stop meas. value

The measurement is canceled when the specified measured value has been reached since the start of the measurement.

The values that are actually measurable may vary from the values specified here, depending on the hardware. The binding measurement value ranges can be found in the .

Measuring mode I_{pol}:

Input range	-2000.0 to 2000.0 mV
Selection	off
Default value	off

Measuring mode U_{pol}:

Input range	-200.00 to 200.00 μA
Selection	off
Default value	off

Measuring mode T:

Input range	-20.0 to 150.0 °C
Selection	off
Default value	off

Temperature

This parameter is not available with the **MEAS T** command (temperature measurement).

Temperature entered manually. The temperature is being continuously measured when a temperature sensor is connected and when **Temp. measurement = automatic** or **continuous** is defined (see sensor dialog of the command).

Input range	-20.0 to 150.0 °C
Default value	25.0 °C

Time interval MP

Time interval for entering a measuring point in the measuring point list. The measuring point list is limited to 1000 measuring points.

Input range	0.1 to 999999.0 s
Default value	2.0 s

28.3.2 Sensor

The parameters for the sensor are edited under **[Sensor]**.

The parameters are valid for all measuring modes except for **Temp. measurement**. This parameter is not available with the **MEAS T** command (temperature measurement).

Measuring input

Selection of the measuring input the sensor is connected to. The selection is not dependent on whether the control device has one or two measuring interfaces.

Selection	1 2
Default value	1

Sensor

Selection of the sensor from the sensor list. The selection depends on the measuring mode. Sensors are defined under **System ► Sensors**. You can also enter a sensor name which is not contained in the sensor list. When a determination is started there is a check whether the sensor is contained in the sensor list.

Selection	Selection of configured sensors
-----------	--

I(pol)

The polarization current is the current that is applied to a polarizable electrode during voltametric measurement. This parameter is available only with I(pol) determinations.



The values that are actually adjustable may vary from the values specified here. The binding values can be found in the manual, in the *Technical specifications* chapter.

Input range	-125.0 to 125.0 μA (Increment: 2.5)
Default value	5.0 μA
Selection	-1.0 μA 1.0 μA

U(pol)

The polarization voltage is the voltage applied to the polarizable electrode during an amperometric measurement. This parameter is available only with U(pol) determinations.

The values that are actually adjustable may vary from the values specified here. The binding values can be found in the manual, in the *Technical specifications* chapter.

Input range	-1,250 to 1,250 mV (Increment: 25)
Default value	400 mV

Electrode check

on | off (Default value: **off**)

For the following electrodes, an electrode check can be carried out:

- Metal electrodes

A check is made that the electrode is properly connected and that no short-circuit is present. The electrode check is carried out when this command is started.

Temp. measurement

Type of temperature measurement.

Selection	continuous automatic off
Default value	automatic

continuous

A temperature sensor must be connected. The temperature is measured continuously.

automatic

If a temperature sensor is connected then the temperature will be measured continuously. Otherwise, the temperature entered manually will be used (see dialog of the titration and measuring parameters).

off

The temperature will not be measured. The temperature entered manually is used (see dialog of the titration and measuring parameters).

28.3.3 Stirrer

The parameters for the stirrer are edited under **[Stirrer]**.

Stirrer

Selection of the MSB connector the stirrer is connected to. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4 off
Default value	1

off

No stirrer will be used.

Stirring rate

Setting the stirring rate. It can be set in steps of -15 to $+15$. The default setting **8** corresponds to 1000 rpm. The formula for calculating the rotational speed is specified in *chapter 31.3, page 311*. The optimum stirring rate can be tested in the manual control.

The algebraic sign of the stirring rate changes the direction in which the stirring is done. When the stirrer is viewed from above, this means:

- "+": counterclockwise rotation
- "-": clockwise rotation

Input range	-15 to 15
Default value	8

Switch off automatically

on | off (Default value: **on**)

If this parameter is activated, the stirrer will be switched off automatically at the end of the titration, measurement, etc.

28.4.1 Fixed endpoint evaluation (EVAL FIX-EP)

For the fixed endpoint evaluation, the associated values are interpolated from the measuring point list for one quantity (measured value, volume, etc.). Up to nine fixed endpoints can be evaluated with one command.

Fixed quantity

Selection of the quantity to which the associated value is interpolated from the measuring point list.

Selection	Measured value Time Volume
Default value	Measured value

Fixed EP1 at

The fixed endpoint must lie between the first and the final entry in the measuring point list.

The values that are actually measurable may vary from the values specified here, depending on the hardware. The binding measurement value ranges can be found in the .

Measured value (measuring mode Ipol):

Input range	-2,000.0 to 2,000.0 mV
Selection	off
Default value	off

Measured value (measuring mode Upol):

Input range	-200.00 to 200.00 µA
Selection	off
Default value	off

Measured value (measuring mode T):

Input range	-20.0 to 150.0 °C
Selection	off
Default value	off

Time:

Input range	0.0 to 999999.9 s
Selection	off
Default value	off

Volume:

Input range	0.00000 to 9,999.99 mL
Selection	off
Default value	off



Fixed EP2 at...Fixed EP9 at

See **Fixed EP1 at**.

28.4.2 Minimum and maximum evaluation (EVAL MIN/MAX)

For the minimum or maximum measured value the associated volume, time and temperature are interpolated from the measuring point list. The evaluation begins as soon as the slope of the curve exceeds a particular threshold value.

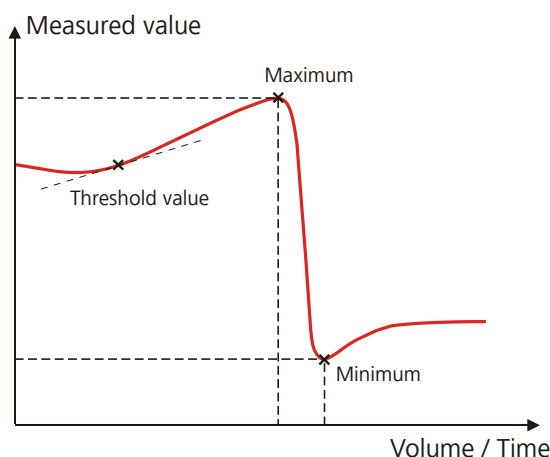


Figure 17 Evaluation of minimum and maximum

Evaluation

Selection whether the minimum or the maximum of the titration curve is being evaluated. With a command either the minimum or the maximum can be evaluated. If both values are needed, a second command has to be defined in the method.

Selection	Minimum Maximum
Default value	Maximum

Threshold value

The evaluation of the minimum or maximum begins as soon as the slope of the curve exceeds the set threshold value. Use a lower threshold value if the minimum or maximum is not found.

Measuring mode Ipol:

Input range	1.0 to 2,000.0
Default value	25.0
	With titrations the unit is mV/mL, with measurements mV/s.

Measuring mode Upol:

Input range	0.5 to 10.0
Default value	5.0
With titrations the unit is $\mu\text{A/mL}$, with measurements $\mu\text{A/s}$.	

Measuring mode T:

Input range	0.1 to 20.0 °C/s
Default value	1.0 °C/s

28.4.3 Rate evaluation (EVAL RATE)

The rate evaluation provides the median dosing rate in a defined range (so-called window). The median dosing rate is determined by means of linear regression through a minimum of three measuring points. If the dosing is carried out with only one dosing device and if more than one cylinder volume is dosed, then the median dosing rate will be reduced as a result of the filling time.

When the command is opened for the first time, a window over the entire time range is already defined. In contrary to the potentiometric evaluation of a DET/MET titration the individual windows may overlap. A maximum of 9 windows can be defined.

[New]

Define a new window.

[Delete]

Delete the selected window.

[Edit]

Edit the settings of the selected window.

Lower limit

Time for the lower limit.

Input range	0 to 999999 s
Default value	0 s

Upper limit

Time for the upper limit.

Input range	0 to 999999 s
Default value	999999 s

28.5 Calculations

The following calculation commands are available:

- | | |
|------------------|--|
| CALC | Defining calculations of a determination. |
| CALC LIVE | Defining the calculation whose result will be displayed in the live display (so-called live result) during a titration or measurement. |

28.5.1 Calculations (CALC)

Calculations are defined with the command **CALC**. A method may contain a maximum of nine calculation commands. A maximum of nine calculations can be defined per command. A series of variables (raw data from the determination, previously calculated results, etc.) is available for the calculations.

The command contains the following options, among others:

- Save result as common variable
- Save result in result table
- Define result limits

Sequence / Edit command	
02 CALC	Calculation
Result	Result name
R1	Content
<input type="button" value="New"/> <input type="button" value="Delete"/> <input type="button" value="Edit"/>	

The result name is specified in the list for each calculation.

[New]

Define a new calculation or load an existing template, see following chapter.

[Delete]

Delete the selected calculation from the list.

[Edit]

Edit the data of the selected calculation (*see chapter 28.5.1.2, page 259*).

28.5.1.1 Loading a result template


The most common calculations are already available, the so-called result templates, in order to facilitate the definition of a new calculation. You also have the option, however, of defining templates of your own (see dialog **System / Templates**).

Proceed as follows to load an existing template:

1 Select a result template


- In the **Sequence / Edit command** dialog, tap on the **[New]** button.

The list of Metrohm result templates is displayed:

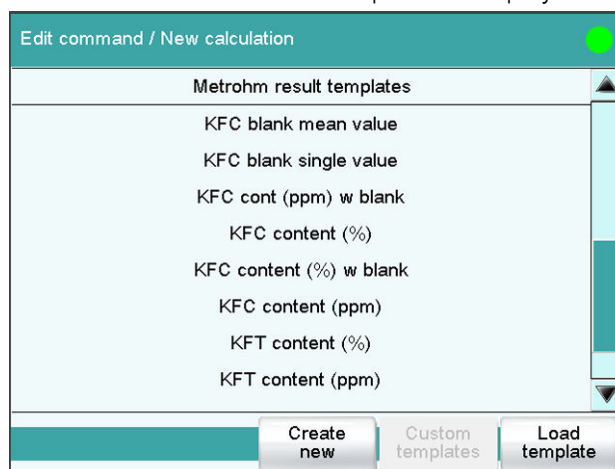

- Select the desired Metrohm result template and tap on **[Load template]**.
or
Tap on **[Custom templates]** and select a template you generated yourself.


If a statistic variable SMN# (# = 1..9) is defined in the template which is already used in another calculation, then the following message will be displayed:

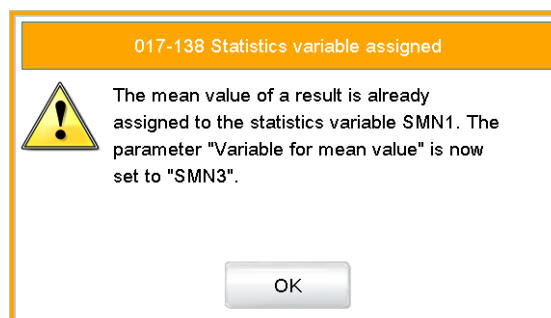

2 Confirm the message

- Tap on **[OK]**.

The Note dialog is displayed. This note is defined in the dialog **Edit result template / Note for wizard**.

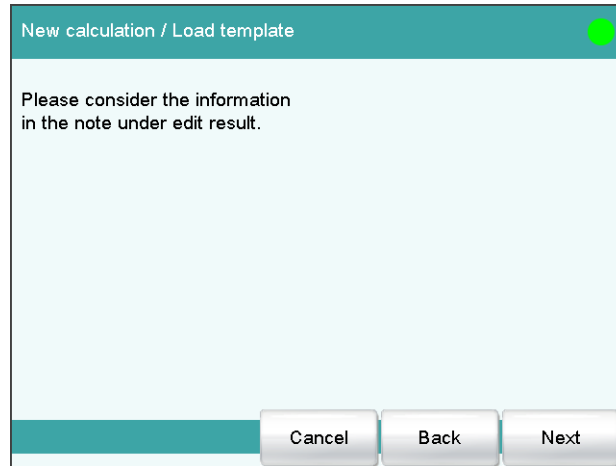

3 Enter the numerical values of the wildcards

- Tap on **[Next]**.

The list of wildcards is displayed:

- Enter the numerical values of the wildcards.
- Tap on **[Next]**.

The editing dialog of the new calculation is displayed:

The calculation can now be edited further, e.g. define result limits, see following chapter.

28.5.1.2 Editing a calculation

Result name

The result name is the text which will be shown in the result display and in the report. The default name corresponds to the result variable.

Entry	max. 24 characters
Default value	R1...R9

Calc. formula R1...R9

Shows the calculation formula. A special editor is opened for the definition (see chapter 28.5.3, page 264). The designation **R1...R9** corresponds to the result variable. Under **[Result variable]** this can later be modified.

Entry	max. 100 characters
Default value	empty

Decimal places

Number of decimal places used to display the result.

Input range	0 to 5
Default value	2

Result unit

The result unit is displayed and saved together with the result.

Entry	max. 10 characters
Selection	% mol/L mmol/L g/L mg/L mg/mL mg/100 g ppm g mg µg mL µL mg/piece °C µg/min mL/min µL/min
Default value	%

Upper limit

When this value is exceeded, the action defined in the following is triggered.

Input range	0 to 9999999999
Default value	9999999999

Action

Selection of the action, which is carried out when the result limits are exceeded.

Selection	Display message Document message Cancel determination
Default value	Display message

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message

In the determination data it will be documented, that the result limits have been exceeded.

Cancel determination

The determination is stopped.

Dialog "Edit calculation / Result options"

In the dialog **Edit calculation / Result options**, settings for how to process the calculated result are defined.

Variable for mean value

If the statistic calculations have been activated (see method options), the mean value of the single results will be saved as variable SMN1 to SMN9. As default value, always the first free variable is being displayed.

Selection	SMN1 ... SMN9 off
-----------	----------------------------

off

For the result not statistic calculations are carried out.

Save as common variable

on | off (Default value: **off**)

The calculated result can be saved as a method-independent variable, a so-called common variable. The result is then also available in other methods for calculations. In all cases, only the single value will be saved, even if the statistics function is enabled.

Variable

Selection of the common variable to which the result is being assigned.

Selection	CV01 ... CV25
-----------	----------------------

Display result

on | off (Default value: **on**)

If you deactivate this parameter, the result is neither displayed in the result dialog nor printed in the result report. This can be advisable for intermediate results.

Save result in result table

on | off (Default value: **off**)

The calculated result can be saved in the result table. This may be advisable if e.g. the results of all determinations carried out on a particular day are to be displayed clearly. A maximum of nine results from a determination can be saved in the result table.

Precision

Setting, with which accuracy the result is used in additional calculations.

Selection	Round Truncate Full precision
Default value	Round

Round

The result is rounded to the defined number of decimal places (commercial rounding, in accordance with the US Pharmacopeia USP). If the digit at the first dropped decimal place is **1, 2, 3 or 4**, then it will be rounded off; if this digit is **5, 6, 7, 8 or 9**, then it will be rounded up. Negative digits will be rounded in accordance with their amount, i.e. away from zero.

Truncate

The result is cut to the number of decimal places defined.

Full precision

The result is used with full accuracy (floating point number either in "single precision" (32 bit) or in "double precision" (64 bit), according to the standard IEEE 754).

28.5.2 Calculations (CALC LIVE)

The command **CALC LIVE** can be used to define a calculation, the result of which will be shown in the live display during a titration or measurement. The current volume or measured value, respectively, is used for this purpose. This is helpful, e.g. for Karl Fischer titrations, in order to be able to track the water content directly during the titration.

**CAUTION**

The command must be inserted directly before the corresponding titration or measuring command.

This command is identical with the **CALC** command except for the following differences:

- The number of variables is limited.
- There are no result options.
- The result cannot be monitored.
- The result variable cannot be modified.
- The result will only be displayed in the live display. It appears neither in the results dialog nor in a report.

Result name

With live calculations no custom result name can be defined, the designation **LR** cannot be modified.

Calculation formula

Shows the calculation formula. A special editor is opened for the definition (see chapter 28.5.3, page 264).

Entry	max. 100 characters
Default value	empty

Decimal places

Number of decimal places used to display the result.

Input range	0 to 5
Default value	2

Result unit

The result unit is displayed and saved together with the result.

Entry	max. 10 characters
Selection	% mol/L mmol/L g/L mg/L mg/mL mg/100 g ppm g mg µg mL µL mg/piece °C µg/min mL/min µL/min
Default value	%

[Note]

Entering a note on the calculation.



Dialog "Edit calculation / Note"

In this dialog you can enter a short text, e.g. to describe the variables used.

Displaying the live result

Proceed as follows to display the live result in the live display:

1 Starting a determination

- Tap on [▶].

The determination starts and the live display is shown:

2 Define the display options

- Tap on the **[View]** button.
- Tap on the **[Meas. value options]** button.
A maximum of three measured values can be shown in the live display.
- Select the option **Live result** at one of the three parameters.

3 Display the live display

- Tap twice on [↔].

The current result is displayed,

28.5.3 Formula editor

The formulas for the calculations are entered with the formula editor. The maximum formula length is 100 characters.

The formula editor contains buttons for entering numbers, mathematical operators and variables. The variables can be selected from lists. This is advisable in order to avoid spelling mistakes. The formula editor is equipped with an automatic syntax check, which is activated when applying the formula. The standard rules of priority apply for the calculation operations.

Edit calculation / Calculation formula

R1= (EP1-CV25)

7	8	9	/	C00	CI	EP	X
4	5	6	*	CV	R	M.	Delete entry
1	2	3	-	TITER	CONC	C.	E.
0	±	.	+	()	^	← →

Cancel Variables Math. functions OK

Key	Description
C00	Sample size
CI	Sample identification CI# (# = 1 - 2)
EP	Volume of endpoint EP# (# = 1...9)
CV	Common Variable CV# (# = 1...25)
R	Result R# (# = 1 - 9)
TITER	Titer of the titrant selected in the previous command
CONC	Concentration of the titrant selected in the previous command
□	The character in front of the cursor is deleted.
Delete entry	The content of the input field is deleted completely.
←	The cursor within the input field is shifted to the left by one character.
→	The cursor within the input field is shifted to the right by one character.

If the method contains more than one data-generating command (titration, measurement etc.) before the calculation command, then you will have to indicate the command identification in front of the variable:

Variable	Description
#M.	Titration or measuring command
#C.	Calculation command
#E.	Evaluation command

"#" stands for a sequential number that you must enter manually. The number for each command identification will be counted separately and does not necessarily correspond to the number of the command line.

Example: The variable **1M.TITER** corresponds to the titer of the titrant selected in the first titration command.



CAUTION

The command identifications are not adjusted automatically when additional commands are subsequently inserted into the method. Check the formula for correctness in such cases.

[Variables]

Opening the list of additional variables (*see chapter 28.5.3.1, page 266*).

[Math. functions]

Opening the list with mathematical functions (*see chapter 28.5.3.2, page 266*).

28.5.3.1 Variables

All of the variables which can be used for calculations are divided up according to subject in the dialog **Calculation formula / Variables**:

- **System variable**
List of the system-specific variables. These variables describe the current status of the system.
- **Result and statistics variables**
List of the variables which are required for calculating results and statistics for the current determination.
- **Common variables**
List of common variables. All of the common variables are listed, no matter whether they possess a value or not.

You will find a detailed description of all of the variables in *chapter 28.5.3.3, page 267*.

28.5.3.2 Mathematical functions

The following mathematical functions can be utilized in a calculation formula:

Function	Description
y^z	Power function Example: $4^2 = 16$

Function	Description
SQRT(X)	square root of X Example: $\sqrt{(EP1)}$
ABS(X)	absolute value of X Example: ABS(C00); in order, e.g. with reweighings, to convert the negative sample size to a positive value for later calculations
LN(X)	natural logarithm of X
LOG(X)	decimal logarithm of X
FRAC(X)	Fraction of X Example: FRAC(2.5971) = 0.5971
INT(X)	integer part of X Example: INT(2.5971) = 2
TST(X,Y)	Test function If invalid variables (e.g. missing endpoints) occur in a calculation, then these can be replaced with a valid value by using this function. In this way invalid results can be avoided. <ul style="list-style-type: none">▪ Syntax:<ul style="list-style-type: none">– X = variable to be tested– Y = substitute value

28.5.3.3 Variable list

The following table contains all of the variables which can be used for calculations. For variables having an index (e.g. 'EP1'), the index must be entered manually. In the following table, this index is characterized with the character "#".

Variable	Description
C00	Sample size
CI1, CI2	Sample identifications The sample identifications can be used in calculations only if numerical values are entered.
DD	Duration of the entire determination
Titration	
TITER	Titer of the titrant selected in the titration command
CONC	Concentration of the titrant selected in the titration command
Titration, measurements, calibrations	



Variable	Description
EP#	Volume/quantity of endpoint EP# (# = 1 - 9)
EC#	Charge at endpoint EP# (# = 1 - 9)
EM#	Measured value of endpoint EP# (# = 1...9)
EF#	ERC of endpoint EP# (# = 1 - 9)
ET#	Temperature at endpoint EP# (# = 1...9)
ED#	Time at endpoint EP# (# = 1...9)
ESI#	Recognition of endpoint EP# (# = 1 - 9) Endpoint found = 1, no endpoint = 0
FMN	Mean gas flow during MEAS T/F ON to MEAS T/F OFF
TMN	Mean temperature during MEAS T/F ON to MEAS T/F OFF
TMI	Minimum temperature during MEAS T/F ON to MEAS T/F OFF
TMA	Maximum temperature during MEAS T/F ON to MEAS T/F OFF
MIM	Initial measured value, i.e. measured value prior to the processing of the start conditions
MIT	Initial temperature, i.e. temperature prior to the processing of the start conditions
MSA	Volume for start volume
MSP	Volume for start measured value
MSS	Volume for start slope
MSV	Volume for all the start conditions
MSD	Duration of start conditions
MSM	Start measured value, i.e. measured value after the processing of the start conditions
MST	Start temperature, i.e. temperature after the processing of the start conditions
MCD	Total duration of the titration, measurement or calibration
MTM	Type of temperature measurement (Pt1000, NTC or manually) Format = Text
MDD	Duration of effective dosing, i.e. without filling times, pauses
MDC	Drift for drift correction
DDC	Time for drift correction
MCV	End volume, i.e. total dosed volume at the end of the titration
MCQ	End quantity, i.e. total amount of removed water or of generated bromine, respectively, at the end of the titration (in µg)
MCC	Final charge, i.e. total charge in mAs at the end of the titration
MCM	End measured value
MCT	End temperature
MMP	Number of measuring points in measuring point list

Variable	Description
MTS	Stop type Format = Text
LV	Currently dosed volume of the ongoing determination
LM	Current measured value of the ongoing determination
LD	Current duration of the ongoing titration or measuring command
LT	Current temperature of the ongoing determination
Sensor	
MEN	Electrode zero point pH(0) or E(0)
MSL	Electrode slope
MVA	Electrode variance (mathematically only defined for three standards or more);
MCL	Cell constant of conductivity measuring cell
Evaluations	
FP#	Volume of fixed endpoint FP# (# = 1 - 9)
FM#	Measured value of fixed endpoint FP# (# = 1 - 9)
FT#	Temperature at fixed endpoint FP# (# = 1 - 9)
FD#	Time at fixed endpoint FP# (# = 1 - 9)
HP#	Volume of pK value / half neutralization potential HP#
HM#	Measured value of pK value / half neutralization potential HP#
HT#	Temperature at pK value / half neutralization potential HP#
HD#	Time at pK value / half neutralization potential HP#
XIP	Volume/quantity of minimum measured value
XIM	Minimum measured value
XIT	Temperature at minimum measured value
XID	Time at minimum measured value
XAP	Volume/quantity of maximum measured value
XAM	Maximum measured value
XAT	Temperature at maximum measured value
XAD	Time at maximum measured value
BP#	Volume of break point BP# (# = 1 - 9)
BM#	Measured value of break point BP# (# = 1 - 9)
BT#	Temperature at break point BP# (# = 1 - 9)
BD#	Time at break point BP# (# = 1 - 9)
RD#	Mean rate in time slot # (# = 1 - 9)




Variable	Description
RDS#	Standard deviation for RD# (# = 1 - 9)
RDC#	Correlation coefficient for RD# (# = 1 - 9)
RM	Mean rate for whole range
RMS	Standard deviation for RM
RMC	Correlation coefficient for RM
Results, statistics	
R#	Result (# = 1–9)
SMN#	Mean value of result R# (# = 1–9)
SSA#	Absolute standard deviation of SMN# (# = 1 - 9)
SSR#	Relative standard deviation of SMN# in % (# = 1 - 9)
SNR#	Number of results from which the mean value SMN# was calculated (# = 1 - 9)
SSD	Number of determinations that are carried out for the statistics calculations
SNT	Statistics status (statistics activated = 1, statistics deactivated = 0)
Common variables	
CV#	Common variable (# = 1...25)
System variables	
%RN	Sample number, i.e. number that is increased by one at each start of a determination
%SC	Start counter, see dialog More determination data / Properties
%AS	Autostart status (autostart activated = 1, autostart deactivated = 0)
%AC	Autostart counter, i.e. number of autostarts already carried out
%AD	Autostart setpoint counter, i.e. number of autostarts to be carried out
%SS	Status of the sample table (sample table activated = 1, sample table deactivated = 0)
%SL	Number of the current sample table line
%SE	Last sample in the sample table processed (yes = 1, no = 0)

28.6 Reports (REPORT)

The reports to be printed out are defined with the command **REPORT**. A maximum of ten reports can be defined for each command. The following reports can be inserted, among others:

- Result report
- Calculation report
- Parameter report
- System reports (system settings, sensor list, GLP data etc.)
- Method reports (method run, method parameters etc.)

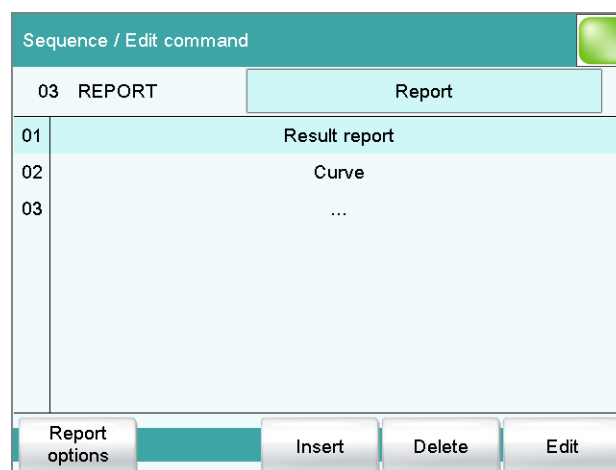

[Report options]

Defining the general settings for the report output.

[Insert]

Add a new report to the list.

[Delete]

Delete the selected report from the list.

[Edit]

Edit the settings of the selected report.

28.6.1 General report options

The settings for the report printout can be adjusted in the dialog **Edit command / Report options** to indicate, for example, whether a report header, a signature line or a frame is to be printed.

Report header

Defining the output of the report header. The report header contains general information such as device type, serial number, program version and the printing date.



NOTICE

You can also define your own report header, which is printed in addition to this report header (**System ► Templates ► Report header**).

Selection	off once on each page
Default value	on each page

off

The report header will not be printed.

once

The report header will only be printed on the first page.

on each page

The report header will be printed on every page.

Signature line

Output of a special line for date and signature. This line will be printed at the very bottom of every page.

Selection	off once on each page
Default value	off

off

No signature line will be printed.

once

The signature line will only be printed on the last page.

on each page

The signature line will be printed on every page.

Frame

on | off (Default value: **on**)

If this parameter is activated, a frame is printed as lateral marking.

28.6.2 Settings of the individual reports

Fixed key []: **More reports ► Report ► Edit**

Command REPORT: **Report ► Edit**

Settings can be edited for the following reports (for details, see online help):

- **Result report**
- **Curve**
Definition of the curve size, display of the individual measuring points and grid lines, etc.

- **Measuring point list**
Definition of the method command to which the measuring point list is to be printed.
- **Calculations**
Definition of the accuracy with which the variables used as well as the results calculated are printed.
- **Statistics short**
Definition of the time at which the report is to be printed (with each determination or only at the end of a sample series).
- **Statistics overview**
Definition of the time at which the report is to be printed (with each determination or only at the end of a sample series).
- **Sample table**
- **Result table**
Definition of the time at which the report is to be printed (with each determination, at the end of a sample series or at the end of the sample table).
- **Used devices**
- **Variables**
- **Monitoring**

28.6.3 List of reports

The following reports can be inserted in the command **REPORT**:



NOTICE

If you are using a **CUSTOM NEOS** printer, you can print only the reports that are marked with a **#** on the display.

Report	Contents
Result report	# Report with determination properties, sample data, calculated results, etc. If there are multiple determinations, the statistics will also be printed out.
Curve	# Curve report.
Measuring point list	# Measuring point list report.
Calculations	# Details concerning the calculations carried out (parameter settings in accordance with the CALC command and the calculated results).
Used devices	The devices used for the determination, as displayed in the More determination data / Properties dialog.
Variables	All determination variables, as displayed in the results dialog.
Monitoring	Details for the monitored quantities.
Statistics	

Report	Contents
Sensors	
Sensor list	List of all the sensors configured in the system.
All sensor data short	The most important sensor data for all sensors.
All sensor data full	All of the sensor data for all of the sensors (name, working life, etc.).
Device manager	
Device list	List of all devices configured in the system.
All device properties	Properties of all the devices configured in the system.
GLP manager	
GLP data	All data stored in the GLP manager.
Common variables	
Common variable list	List of all the common variables defined in the system, together with their most important data (name, value, status).
All common variable properties	Properties of all common variables (name, value, validity, status).
Templates	
Templates sample data	Sample identification list and sample assignment table.
Result template list	List with all of the user-generated result templates.
All result templates details	Details of the all of the user-generated result templates (calculation formula, result options, note).
Input/Output lines	List with all of the defined input and output lines at the remote interface (name, bit pattern).
Custom calibration buffers	Temperature tables for all defined custom calibration buffers.
Rack tables	
Sample rack list	List with all of the sample racks configured in the system, together with designation, number of positions and rack code.
Miscellaneous	
Form feed	If this entry is inserted between two reports, then these will each be printed out on a separate page.

28.7 Dosing and Liquid Handling

The following dosing commands are available:

Preparing the exchange unit or dosing unit (PREP)	Rinsing the cylinder and the tubings of the exchange unit / dosing unit.
Emptying the dosing unit (EMPTY)	Emptying the cylinder and the tubings of the dosing unit.
Dosing a fixed volume (ADD)	Dosing a specified volume.
Liquid Handling (LQH)	Carrying out complex dosing tasks with a Dosing unit.

28.7.1 Preparing an exchange or dosing unit (PREP)

The **PREP** command is used to rinse and fill air bubble-free the cylinder and tubings of the exchange or dosing unit. You should carry out this function before the first determination or once a day.

Control device

Selection	Selection of configured control devices
-----------	--

Dosing device

Selection of the MSB connector to which the dosing device is connected. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4
Default value	1

Titrant

Selection of the titrant from the titrant list. We recommend always selecting the titrant. This ensures that the parameters defined for the titrant for the preparing as well as the tubing parameters (dosing unit only) are used. Titrants are defined under **System ► Titrants**. You can also enter a name which is not contained in the titrant list. When a determination is started there is a check whether the solution is contained in the list.

For buret units with integrated data chip, a check is made in the method run to verify whether the correct titrant has been attached and whether the type of dosing drive matches. For buret units without integrated data chip, the cylinder volume and the type of dosing drive are checked. For the selected titrant, the validity of the titer, the working life of the titrant and the time interval for the GLP test of the buret unit are checked at the start of the determination.

Selection	Selection of configured titrants not defined
Default value	not defined

not defined

No check takes place. The command is carried out with default parameters (see Chapter 31.1.2, page 310 and Chapter 31.2.2, page 311).

Cylinder volume to

This parameter can only be edited with **Titrant = not defined**.

Selection of the port via which the titrant is ejected. This setting is only relevant for dosing units, in case of preparing an exchange unit, this setting is ignored.

Selection	Port 1 Port 2 Port 3 Port 4
Default value	Port 1

28.7.2 Emptying a dosing unit (EMPTY)

The **EMPTY** command is used to empty the cylinder and tubings of the dosing unit.

Control device

Selection	Selection of configured control devices
-----------	--

Dosing device

Selection of the MSB connector to which the dosing device is connected. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4
Default value	1

Titrant

Selection of the titrant from the titrant list. We recommend always selecting the titrant. This ensures that the parameters defined for the titrant for the preparing as well as the tubing parameters (dosing unit only) are used. Titrants are defined under **System ► Titrants**. You can also enter a name which is not contained in the titrant list. When a determination is started there is a check whether the solution is contained in the list.

For buret units with integrated data chip, a check is made in the method run to verify whether the correct titrant has been attached and whether the type of dosing drive matches. For buret units without integrated data chip, the cylinder volume and the type of dosing drive are checked. For the selected titrant, the validity of the titer, the working life of the titrant and the time interval for the GLP test of the buret unit are checked at the start of the determination.

Selection	Selection of configured titrants not defined
Default value	not defined

not defined

No check takes place. The command is carried out with default parameters (see Chapter 31.1.2, page 310 and Chapter 31.2.2, page 311).

Air inlet

This parameter can only be edited with **Titrant = not defined**.

Selection of the port via which air is aspirated. This setting is only relevant for dosing units, in case of preparing an exchange unit, this setting is ignored.

Selection	Port 1 Port 2 Port 3 Port 4
Default value	Port 4

28.7.3 Dosing a specified volume (ADD)

You can dose a specified volume with the command **ADD**.

28.7.3.1 Dosing parameters

The parameters for the dosing are defined under **[Dosing parameters]**.

Volume

Volume which is dosed.

Input range	0.00000 to 99999.9 mL
Default value	10.0000 mL

Dosing rate

The rate at which dosing takes place.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (see chapter 31, page 310).

Filling rate

Rate at which the dosing cylinder is filled.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum filling rate depends on the cylinder volume (*see chapter 31, page 310*).

28.7.3.2 Dosing device

The parameters for the dosing device are edited under **[Dosing device]**.

Dosing device

Selection of the MSB connector to which the dosing device is connected. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4
Default value	1

Titrant

Selection of the titrant from the titrant list. We recommend always selecting the titrant. This ensures that the correct data (titer, etc.) is always used for the calculation. Titrants are defined under **System ► Titrants**. You can also enter a name which is not contained in the titrant list. When a determination is started there is a check whether the solution is contained in the list.

For buret units with integrated data chip, a check is made in the method run to verify whether the correct titrant has been attached and whether the type of dosing drive matches. For buret units without an integrated data chip, the cylinder volume and the type of dosing drive are checked. For the selected titrant, the validity of the titer, the working life of the titrant and the time interval for the GLP test of the buret unit are checked at the start of the determination.

Selection	Selection of configured titrants not defined
Default value	not defined

not defined

No check takes place.

Tandem dosing

A second dosing device can be defined under **[Tandem dosing]** in order to enable uninterrupted dosing. Dosing is carried out with a combination of two dosing devices so that the second dosing device is dosing while the first one is being filled and vice versa.

Dosing device

Selection of the MSB connector to which the dosing device is connected. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4 off
Default value	off

Titrant

Selection of the titrant from the titrant list. We recommend always selecting the titrant. Titrants are defined under **System ► Titrants**. You can also enter a name which is not contained in the titrant list. When a determination is started there is a check whether the solution is contained in list.

For buret units with integrated data chip, a check is made in the method run to verify whether the correct titrant has been attached and whether the type of dosing drive matches. For buret units without integrated data chip, the cylinder volume and the type of dosing drive are checked. For the selected titrant, only the time interval for the GLP test of the buret unit is checked at the start of the determination.



NOTICE

The concentration, validity of the titer and the working life of the titrant are ignored.

Selection	Selection of configured titrants not defined
Default value	not defined

not defined

No check takes place.

Filling rate

Rate at which the dosing cylinder is filled.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum filling rate depends on the cylinder volume (*see chapter 31, page 310*).

However, in order to ensure uninterrupted dosing, the following points must be taken into account:

- Use the highest possible filling rate in order to keep the filling times as short as possible. The filling rate should be decreased for viscous liquids.
- When you use two buret units with different dosing cylinders, the filling rate for the larger sized cylinder must be at the minimum:

$$v_{2,Fill} \geq v_{1,Fill} \cdot \frac{V_{Cyl.2}}{V_{Cyl.1}}$$

$v_{2,Fill}$ = filling rate in mL/min for the larger sized cylinder

$v_{1,Fill}$ = filling rate in mL/min for the smaller cylinder

$V_{Cyl.2}$ = cylinder volume in mL of the buret unit of the second dosing device

$V_{Cyl.1}$ = cylinder volume in mL of the buret unit of the first dosing device

Example:

Dosing device 1: volume = 20 mL, filling rate = 50 mL/min

Dosing device 2: volume = 50 mL

Filling rate 2 \geq 50 mL/min · 50 mL / 20 mL \geq 125 mL/min

- The dosing rate must not exceed 75% of the filling rate of the smaller cylinder. These values are listed in the following table, valid at maximum filling rate:

Table 11 Maximum dosing rate for different dosing cylinders

Cylinder volume	maximum dosing rate	
	Exchange unit	Dosing unit
1 mL	2.25 mL/min	–
2 mL	–	5.00 mL/min
5 mL	11.25 mL/min	12.50 mL/min
10 mL	22.50 mL/min	25.00 mL/min
20 mL	45.00 mL/min	50.00 mL/min
50 mL	112.50 mL/min	124.50 mL/min

28.7.3.3 Stirrer

The parameters for the stirrer are edited under **[Stirrer]**.

Stirrer

Selection of the MSB connector the stirrer is connected to. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4 off
Default value	1

off

No stirrer will be used.

Stirring rate

Setting the stirring rate. It can be set in steps of -15 to $+15$. The default setting **8** corresponds to 1000 rpm. The formula for calculating the rotational speed is specified in *chapter 31.3, page 311*. The optimum stirring rate can be tested in the manual control.

The algebraic sign of the stirring rate changes the direction in which the stirring is done. When the stirrer is viewed from above, this means:

- "+": counterclockwise rotation
- "-": clockwise rotation

Input range	-15 to 15
Default value	8

Switch off automatically

on | off (Default value: **on**)

If this parameter is activated, the stirrer will be switched off automatically at the end of the titration, measurement, etc.

28.7.4 Liquid Handling (LQH)**28.7.4.1 Liquid Handling parameters**

The parameters for the Liquid Handling function are defined under [**Liquid Handling param.**].

Function

Type of Liquid Handling function.

Selection	Dose Fill Aspirate Eject Exchange position Change port Compensate End volume
-----------	---

Dose

The given volume is dosed. The dosing cylinder neither is automatically filled before nor afterwards.

Fill

The dosing cylinder is filled. The valve disc then remains at the selected port.

Aspirate

Liquid is aspirated. The dosing cylinder neither is automatically filled before nor afterwards. The aspirating volume has to be aspirated with a single piston stroke.

Eject

The whole cylinder content is ejected. The piston is, in contrast to the function **End volume**, lowered to the stop point, i.e. past the maximum volume mark. This function should only be carried out for preparing a dosing unit before the pipetting procedure.

Exchange position

The dosing cylinder is filled. This means that, for example, air can be aspirated via Port 4. The valve disc is then rotated to Port 2 and the dosing drive can be removed from the dosing unit.

Change port

The stopcock is moved to the given port, but no piston movement takes place.

Compensate

Because of the fact that the dosing units are interchangeable, the coupling of the Dosino piston rod (spindle) exhibits a low mechanical tolerance that can be noticed when the piston changes its direction of movement. This tolerance can be compensated with this function. A short piston movement is first made in the same direction as the previous movement, which is then followed by a piston movement in the reverse direction.

End volume

The whole cylinder content is ejected. The piston is lowered to the maximum volume mark i.e. until 10'000 pulses have been dosed. This command should be used for pipetting functions for emptying the cylinder.

In-/Outlet

Selection of the port via which the Liquid Handling command is carried out.

Selection	Port 1	Port 2	Port 3	Port 4
-----------	--------	--------	--------	--------

Port 1

Default value, when **Function = Dose, Aspirate** and **End volume**.

Port 2

Default value, when **Function = Fill, Exchange position** and **Change port**.

Port 4

Default value, when **Function = Eject** and **Compensate**.

Volume

This parameter can only be edited with **Function = Dose** and **Aspirate**.

Volume which is dosed or aspirated.

Selection	Selection of configured titrants not defined
Default value	not defined

not defined

No check takes place.

28.8 Communication

The following communications commands are available:

Scanning remote lines (SCAN)	Defining remote signals which are awaited before the next method command is started.
Setting remote lines (CTRL)	Defining remote signals.
Receiving RS-232 commands (SCAN RS)	Defining RS-232 commands which are awaited before the next method command is started.
Sending RS-232 commands (CTRL RS)	Defining the RS-232 commands which are sent.

28.8.1 Scanning remote lines (SCAN)

The command **SCAN** can be used to define input signals at the remote interface which are awaited before the next method command is started.

Control device

Selection	Selection of configured control devices
-----------	--

Remote Box

Selection of the MSB connector the Remote Box is connected to. The selection does not depend on the number of MSB connectors of the control instrument. The remote signal defined is requested on this Remote Box.

Selection	1 2 3 4
Default value	1

Input signal

Selection of the signal out of the templates or entering the required bit pattern. Templates are defined under **System ► Templates ► Input lines**.

Entering a bit pattern:

- 0 = line inactive
- 1 = line active

- * = retain line status

The input lines are always numbered from right to left, i.e. with the signal *******1** line 0 is expected to be active.



NOTICE

We recommend masking lines that are of no interest or for which no defined condition can be predicted with an asterisk (*).

Entry	Bit patterns containing exactly 8 characters or a max. of 24 characters for the name of the template
Default value	*****
Selection	Selection of the templates defined

Timeout

When this time interval has expired without recognizing the remote signal, the action defined in the following is triggered.

Input range	0 to 999 s
Default value	0 s

Action

Selection of the action which is carried out when the time interval has expired.

Selection	Display message Document message Cancel determination
Default value	Display message For all three options it is documented in the determination data (see dialog More determination data / Messages), that the time interval has been expired.

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message

In the determination data it will be documented, that the time interval has been expired.

Cancel determination

The determination is stopped.

28.8.2 Setting remote lines (CTRL)

The command **CTRL** can be used to define output signals at the remote interface.

Control device

Selection	Selection of configured control devices
-----------	--

Remote Box

Selection of the MSB connector the Remote Box is connected to. The selection does not depend on the number of MSB connectors of the control instrument. The remote signal defined is sent via this Remote Box.

Selection	1 2 3 4
Default value	1

Output signal

Selection of the signal out of the templates or entering the required bit pattern. Templates are defined under **System ► Templates ► Output lines**.



NOTICE

A line set active is not being reset automatically, not even at the end of the determination.

Entering a bit pattern:

- 0 = line inactive
- 1 = line active
- * = retain line status
- p = set pulse

The output lines are always numbered from right to left, i.e. with the signal *******1** line 0 is set. With a pulse, the length is set to 200 ms. If it is required to set pulses with other lengths, you have to define a corresponding template.



NOTICE

We recommend masking lines that are of no interest or for which no defined condition can be predicted with an asterisk (*).

Entry	Bit patterns containing exactly 14 characters or a max. of 24 characters for the name of the template
Default value	*****
Selection	Selection of the templates defined

28.8.3 Scanning the RS-232 interface (SCAN RS)

The command **SCAN RS** can be used to define RS-232 commands which are awaited before the next method command is started.

Serial port

Selection of the serial interface the peripheral device is connected to. The RS-232 command defined is requested on this interface.

Selection	COM1 COM2 COM3 COM4 COM5 COM6 COM7 COM8
Default value	COM1

Character string

Input of the RS-232 command as character string. All characters of the ASCII code page can be used. Control characters (Esc, FF etc.) have to be entered as a three-place, decimal ASCII code, starting with a slash. Each command is automatically terminated with the ASCII characters **CR** and **LF**. The ***** can be used for one or more characters.

Entry	ASCII string with a max. of 24 characters
Default value	!*R

Timeout

When this time interval has expired without recognizing the RS-232 command, the action defined in the following is triggered.

Input range	0 to 999 s
Default value	0 s

Action

Selection of the action which is carried out when the time interval has expired.

Selection	Display message Document message Cancel determination
Default value	Display message For all three options it is documented in the determination data (see dialog More determination data / Messages), that the time interval has been expired.

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message

In the determination data it will be documented, that the time interval has been expired.

Cancel determination

The determination is stopped.

28.8.4 Defining RS-232 commands (CONTROL RS)

The command **CONTROL RS** can be used to define the RS-232 commands which are sent.

Serial port

Selection of the serial interface the peripheral device is connected to. The RS-232 command defined is sent via this interface.

Selection	COM1 COM2 COM3 COM4 COM5 COM6 COM7 COM8
Default value	COM1

Character string

Input of the RS-232 command as character string. All characters of the ASCII code page can be used. Control characters (Esc, FF etc.) have to be entered as a three-place, decimal ASCII code, starting with a slash. Each command is automatically terminated with the ASCII characters **CR** and **LF**.

Entry	ASCII string with a max. of 24 characters
Default value	&M;\$G

28.9 Automation**28.9.1 Rotating sample rack (MOVE)**

The command **MOVE** is used to move to a rack position or an external position.

Tower

Selection of the tower, with which the command is carried out. Both Tower 1 and Tower 2 can always be selected, even if your Sample Processor has only one tower.

Selection	1 2
Default value	1

Shift direction

Direction in which the rack is moved.

Selection	auto + -
Default value	auto

auto

A shift direction, with which the shorter way has to be passed, is automatically selected.

+

Counterclockwise rotation.

-

Clockwise rotation.

Swing rate

Rate at which the robotic arm is swung.

Input range	10 to 55 °/s
Default value	55 °/s

28.9.2 Moving the lift (LIFT)

The lift of a Sample Processor is moved with the command **LIFT**. The movement can, however, only be carried out if the Sample Processor is located at a valid rack position. This will not be the case, for example, after a Rack Reset.

Tower

Selection of the tower, with which the command is carried out. Both Tower 1 and Tower 2 can always be selected, even if your Sample Processor has only one tower.

Selection	1 2
Default value	1

1

View from the front, the right tower.

2

View from the front, the left tower.

Lift position

Selecting a predefined lift position or enter any lift position. The predefined lift positions are defined in the device properties of the Sample Processor. They can, however, also be assigned in the manual control.

Input range	0 to 235 mm
Selection	Work position Shift position Rinse position Special position Home position
Default value	Work position

Home position

The lift is being located at the upper stop position, i.e. at the position 0 mm.

For the 885 Compact Oven SC, the position 0 mm is approached when **Shift position**, **Rinse position** and **Special position** is selected.

Lift rate

Speed at which the lift is moved.

Input range	5 to 25 mm/s
Default value	20 mm/s



NOTICE

This parameter is ignored with the 885 Compact Oven SC.

28.9.3 Resetting the rack (RACK)

The following actions are carried out with the command **RACK**:

- Rack, lift and robotic arm are being reset.
- The rack code of the rack attached is being read out and the corresponding rack data is transferred to the Sample Processor.
- The sample variable is reset to the value 1.

Check rack

on | off (Default value: **off**)

Activate this parameter when the rack attached is to be checked. For this you, however, have to additionally select the rack in the dialog **Method options / Start options** at the option **Check rack**.



NOTICE

This option is not available with the 885 Compact Oven SC. Therefore, this parameter must be set to **no**.

28.9.4 Defining the sample variable (SAMPLE)

The sample variable describes the current position of the sample on the rack of the Sample Processor. It is automatically increased by 1 after the end of a method run. In the following cases, the sample variable is automatically reset to the value 1:

- if the autostart counter is reset.
- if the sample rack is reset.

The sample variable can be altered specifically with the command **SAMPLE**.

Sample variable

Modifying the sample variable.

Selection	= + -
Default value	+

=

The sample variable corresponds to the number entered in the field **Value**.

+

The sample variable corresponds to the current value plus the number entered in the field **Value**.

-

The sample variable corresponds to the current value minus the number entered in the field **Value**.

Value

Value by which the current sample variable is to be modified.



NOTICE

The sample variable can only be set on rack positions which are not defined as special beakers.

Input range	1 to 999
Default value	1

28.9.5 Creating a subsequence (SUBSEQ)

It is useful, particularly with more complex tasks, e.g. rinsing sequences or Liquid Handling, to combine individual command sequences into a single subsequence. Subsequences are managed the same as method commands. A maximum of 99 commands can be combined to form a single subsequence. There is no limit to the number of subsequences within one method.

[Load/ Save]

Loading a saved subsequence or saving the current subsequence.

[Subseq. options]

Defining various settings which correspond to the entire subsequence.

[Insert command]

Insert a new method command. It is inserted before the selected command.

**NOTICE**

Not all commands are available for subsequences. Commands which cannot be inserted into subsequences are disabled.

[Delete command]


Delete the selected method command.

[Edit command]

Edit the selected method command.

28.9.5.1 Loading / saving a subsequence

[Load/ Save] is used to load stored subsequences or to save new subsequences. This means that they are available for all methods and only have to be created once.

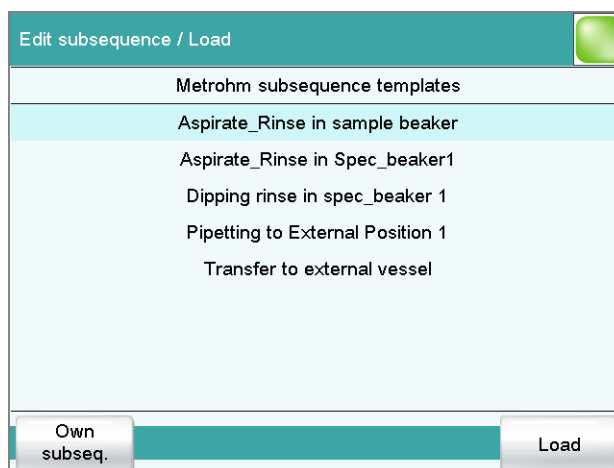Loading a subsequence**[Own subseq.]**

Opening the list of own subsequences.

[Load]

Loading the selected subsequence.

Saving a subsequence

Custom-made subsequences are saved in the internal memory.

**NOTICE**

In order to be able to use the subsequences also for other titration systems, you have to create a backup. You can only restore the subsequences from this backup on another system.

File name

File name of the subsequence.

Entry	max. 32 characters
-------	---------------------------

[Save]

Saving the subsequence to the internal memory.

28.9.5.2 Subsequence options

In this dialog, the settings can be defined which correspond to the entire subsequence.

Subsequence type

Selection	Sample sequence Start sequence End sequence Stop sequence Cond. sequence
Default value	Sample sequence

Sample sequence

The subsequence is carried out with every determination.

Start sequence

This subsequence is only carried out at the start of a sample series when the autostart counter = 1.

End sequence

This subsequence is only carried out at the last sample of a series when the autostart counter has reached the setpoint value (number of autostarts).

Stop sequence

The stop sequence is only carried out when canceling a method because of one of the following cases: Manual stop with [], stop because of an error, stop via remote signal.

Cond. sequence

This subsequence is carried out directly before the conditioning. This gives you the possibility to start e.g. a connected Polytron® high-frequency homogenizer before the conditioning as well as to set its stirring rate. Use the command **CONTROL RS** for this.

Cycles

Number of directly consecutive repetitions of the subsequence.

[Note]

Entering a note on the subsequence.

Dialog "Subsequence options / Note"

A short text can be entered in this dialog, e.g. for the description of the subsequence commands or for the application purpose.

[Display options]

Definition when the note is being displayed.

Automatically after loading the subsequence

on | off (Default value: **off**)

If this parameter is activated, the note is displayed when loading the subsequence. It can otherwise only be read in this dialog.

28.9.6 Controlling the heating (HEATER)

The heating on the 885 Compact Oven SC is controlled with the **HEATER** command.

Status

Switch the heater on or off.

Selection	on off
Default value	on

on

The heater is switched on. The command is displayed in the method run until the specified temperature is reached. The heater runs in the background until it is explicitly switched off. The current temperature can be seen in the Manual control.

off

The heater is switched off.

Temperature

Temperature to which the oven is heated.

Input range	50 to 250 °C (Increment: 1)
Default value	110 °C

28.9.7 Controlling the gas flow (FLOW)

The gas flow is controlled on the 885 Compact Oven SC with the **FLOW** command.

Status

Switch the gas flow on or off.

Selection	on off
Default value	on

on

The gas flow is switched on. The gas flow runs until it is switched off. The current gas flow can be seen in the Manual control.

off

The gas flow is switched off.

Flow rate

Flow rate for the gas flow.

Input range	10 to 150 mL/min (Increment: 1)
Default value	50 mL/min

Gas supply

Selection of the gas supply.

Selection	pump valve
Default value	pump

pump

The integrated air pump is used for the gas flow.

valve

The integrated inlet valve is used for the gas flow. The inlet valve in the 885 Compact Oven SC is connected to compressed air or nitrogen.

Gas type

Gas type of the carrier gas used.

Selection	air nitrogen
Default value	air

28.9.8 Starting the measurement for temperature and gas flow (MEAS T/F ON)

Data acquisition and evaluation of the data of the 885 Compact Oven SC is started with the **MEAS T/F ON** command.

The following values are output as raw data and are available as variables:

- Mean gas flow (evaluation starts after 10 s).
- Mean temperature
- Minimum temperature
- Maximum temperature

After a **MEAS T/F ON** command, there must also be a **MEAS T/F OFF** command present in the method run. If one of the two commands is missing or the order is not followed, the command is depicted in red. No parameters can be edited for this command.

28.9.9 Exiting the measurement for temperature and gas flow (MEAS T/F OFF)

The data acquisition evaluation of the data of the 885 Compact Oven SC is finished with the **MEAS T/F OFF** command. There must be a **MEAS T/F ON** command present before a **MEAS T/F OFF** command in the method run. The command will be depicted in red if this is not fulfilled. No parameters can be edited for this command.

28.10 Miscellaneous commands

28.10.1 Controlling a stirrer (STIR)

A stirrer is controlled with the command **STIR**.

Control device

Selection	Selection of configured control devices
-----------	---

Stirrer

Selection of the MSB connector the stirrer is connected to. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4
Default value	1

Status/Duration

Switching the stirrer on or off.

Input range	0 to 999 s (Increment: 1) During this time the stirrer is running.
Selection	on off
Default value	on

on

The stirrer is switched on. The stirrer runs until it is explicitly switched off. If the determination is canceled with the key [], the stirrer is also switched off.

off

The stirrer is switched off.

Stirring rate

Setting the stirring rate. It can be set in steps of -15 to $+15$. The default setting **8** corresponds to 1000 rpm. The formula for calculating the rotational speed is specified in *chapter 31.3, page 311*. The optimum stirring rate can be tested in the manual control.

The algebraic sign of the stirring rate changes the direction in which the stirring is done. When the stirrer is viewed from above, this means:

- "+": counterclockwise rotation
- "-": clockwise rotation

Input range	-15 to 15
Default value	8

28.10.2 Pausing the method run (WAIT)

The method run can be paused with the command **WAIT**.

Hold sequence

on | off (Default value: **off**)

If this parameter is activated, then the method run is paused until it is continued manually.

Waiting time

This parameter can only be edited when **Hold sequence** is deactivated.

The method run is automatically continued after this waiting time.

Input range	0 to 99999 s
Default value	30 s

Message

on | off (Default value: **off**)

If this parameter is activated, the text message defined in the following is displayed during the waiting time.

Message text

This parameter can only be edited when **Message** is activated.

Text which is displayed during the waiting time.

Entry	max. 28 characters
Default value	empty

28.10.3 Scan data (REQUEST)

The following data can be scanned in the method run with the command **REQUEST**:

- Sample data
 - Sample identification 1
 - Sample identification 2
 - Sample size (value and unit)
- Common variable



You can select thereby whether the sequence should be paused or resumed in the background.

Sample ident.

Selection of the sample identification that is queried in the method run.

Selection	off Identification 1 Identification 2 ID1 & ID2
Default value	off

Sample size

on | off (Default value: **on**)

If this parameter is activated, then the value for the sample size will be requested.

Sample size unit

on | off (Default value: **off**)

If this parameter is activated, then the unit for the sample size will be requested.

Common variable

Selection of the common variable which is queried in the method run.

Selection	CV01...CV25 off
Default value	off

Hold sequence

on | off (Default value: **on**)

If this parameter is activated, then the run will be paused during the request. If this parameter is deactivated, then the method continues in the background until the following measurement or titration has been completed.

28.10.4 Defining an acoustic signal (BEEP)

The **BEEP** command can be used to produce an acoustic signal.

Duration

Approximate playing time in seconds of the permanently stored melody.

Input range	1 to 9
Default value	1

28.10.5 Signing a determination (SIGN)

A determination can be signed with the **SIGN** command. The method run is automatically paused as soon as this command is reached. At the same time a switch is made to the result view. The run will not be continued until the determination has been signed or the **SIGN** command has been canceled. You can specify whether a routine user is allowed to cancel the **SIGN** command in the setting **System ► System settings ► Dialog options ► Routine dialog**. No parameters can be edited for this command. The parameters for signing determinations are described in *chapter 16.1, page 149*.



NOTICE

Methods with this command can only be started if you are working with login and password protection.

The command should always be inserted at the end directly before the **REPORT** command or at the end of the method sequence if a PC/LIMS report is to be sent automatically.

28.10.6 Canceling the method run (END)

The method run will cancel as soon as the **END** command is reached. This is a good idea if you wish to test only the first part of a method. No parameters can be edited for this command.

29.2 Maintenance

Maintenance

The electronic and mechanical functional groups of Metrohm instruments can and should be checked by specialist personnel from Metrohm as part of a regular preventive maintenance schedule. Please ask your local Metrohm representative regarding the precise terms and conditions involved in concluding a corresponding maintenance agreement.

For detailed information on this topic, please visit www.metrohm.com.

Cleaning

Cleaning the surfaces of the instrument

Prerequisites

- The instrument is disconnected from the power grid.

- 1 Clean the surfaces with a damp cloth.



NOTICE

Water or ethanol can be used as a cleaning medium.




NOTICE

The connectors at the rear of the instrument must only be cleaned with a dry cloth.


30.3 Results/Statistics

Problem	Cause	Remedy
No statistics are carried out for a result.	<i>Statistics is not activated.</i>	In the method options, activate the option Statistics (see chapter 15.5, page 137).
	<i>No statistics variable has been assigned to the calculation.</i>	In the result options (command CALC), define a variable for the mean value (see "Variable for mean value", page 123).
	<i>Statistics is not activated.</i>	In the control dialog, activate the Statistics option (see chapter 17, page 151).
A result has been removed from the statistics, the mean value has not been reassigned to the common variable or to the TITER variable.	<i>Assignment is not carried out automatically with retroactive modifications.</i>	Recalculate the determination manually (see chapter 23, page 178).
The result is not displayed in the result table.	<i>The column display in the result table is incorrectly configured.</i>	Modify the settings for the display accordingly (see chapter 25.1, page 200).
	<i>The parameter Save result in result table is not activated.</i>	In the result options (command CALC), activate the parameter (see "Save result in result table", page 124).

30.4 Printing

Problem	Cause	Remedy
Tapping on the fixed key [] has no effect.	<i>The fixed key is disabled.</i>	In the Dialog options / Fixed keys dialog, activate the Print option (see chapter 7.2, page 39).
	<i>A determination is running.</i>	Wait until the determination is finished.

30.5 Manual control

Problem	Cause	Remedy
Tapping on the fixed key [] has no effect.	<i>The fixed key is disabled.</i>	In the Dialog options / Fixed keys dialog, activate the Manual control option (see chapter 7.2, page 39).
A button in the manual control is disabled.	<i>The required hardware is either not connected or incorrectly connected.</i>	<ol style="list-style-type: none"> 1. Switch off the 917 Coulometer. 2. Connect the hardware correctly. 3. Switch the 917 Coulometer back on.
	<i>The required hardware is being used in the ongoing determination.</i>	Wait until the determination is finished.

30.6 Titration stand with pump

Problem	Cause	Remedy
Solvent is added at the same time during aspiration.	<i>The adsorber tube is blocked.</i>	<ul style="list-style-type: none"> ▪ Enlarge the hole of the drying tube cover to 2 mm at least. ▪ Refill the adsorber tube, pack not too tightly, perhaps remove some of the molecular sieve.
The pump conveys no liquid.	<i>The add/aspirate key is not pressed tightly enough.</i>	<ul style="list-style-type: none"> ▪ Usually the bottle attachments are not screwed tightly enough to the reagent and waste bottles. ▪ Check all connections for leak-tightness.
The solvent flows on into the titration vessel after the addition, without the add key being pressed.	<i>The solvent bottle is not positioned correctly.</i>	Position the solvent bottle in such a way that the liquid level in the bottle is lower than that in the titration vessel.

30.7 Coulometric Karl Fischer titration

Problem	Cause	Remedy
The drift is very high during conditioning.	<i>The titration cell is leaking.</i>	<ul style="list-style-type: none"> ▪ Check the septum and replace it if necessary. ▪ Replace the molecular sieve.

Problem	Cause	Remedy
		<ul style="list-style-type: none"> Check the ground-joint sleeves for frayed edges. If necessary, trim the edges neatly or replace the ground-joint sleeves.
	<i>Deposits containing water can be found in the titration cell.</i>	<ul style="list-style-type: none"> Shake the titration cell.
	<i>The reagent is exhausted or contaminated.</i>	<ul style="list-style-type: none"> Replace the reagent.
	<i>The catholyte is old or moist.</i>	<ul style="list-style-type: none"> The catholyte must be replaced at least once per week. Dry the catholyte with KF one-component reagent.
	<i>The diaphragm of the generator electrode is contaminated.</i>	<ul style="list-style-type: none"> Clean the diaphragm (in accordance with directions in the leaflet for the generator electrode).
	<i>A side reaction is taking place.</i>	<ul style="list-style-type: none"> Consider using a KF oven. See technical literature.
	<i>The molecular sieve on the KF oven has reached its capacity limit.</i>	<ul style="list-style-type: none"> Replace the molecular sieve.
	<i>The gas flow from the KF oven into the titration cell is too high.</i>	<ul style="list-style-type: none"> Reduce the gas flow (if you are working with the oven method with sample vials, set the gas flow between 40 and 60 mL/min).
The drift becomes greater after each titration.	<i>The sample releases water very slowly.</i>	<ul style="list-style-type: none"> Consider using a KF oven. See technical literature.
The drift is fluctuating.	<i>The sample solution is poorly stirred.</i>	<ul style="list-style-type: none"> Set the stirring rate in such a way that the sample solution will become very well mixed.
	<i>The Control parameters are set incorrectly.</i>	<ul style="list-style-type: none"> Reset the Control parameters to default values.
The titration time is too long.	<i>The drift during conditioning is not stable.</i>	<ul style="list-style-type: none"> Wait before starting the titration until the drift has become stable.
	<i>The parameter Start drift is set too high.</i>	<ul style="list-style-type: none"> Set the Start drift lower.

Problem	Cause	Remedy
	<i>The sample addition is difficult to reproduce.</i>	<ul style="list-style-type: none"> ▪ Observe the instructions in .
	<i>The drift fluctuates.</i>	<ul style="list-style-type: none"> ▪ Ensure stable drift.

31 Appendix

31.1 Dosing unit

31.1.1 Maximum dosing and filling rate

The maximum dosing rate and maximum filling rate for the dosing unit depend on the cylinder volume:

Cylinder volume	Maximum rate
2 mL	6.67 mL/min
5 mL	16.67 mL/min
10 mL	33.33 mL/min
20 mL	66.67 mL/min
50 mL	166.00 mL/min

Independent of the cylinder volume, values ranging from 0.01 to 166.00 mL/min can always be entered. When the function is carried out the rate will be, if necessary, decreased automatically to the highest possible value.

31.1.2 Default parameters for preparing (PREP) and emptying (EMPTY)

The **PREP** command is used to rinse and fill the cylinder and tubings of the dosing unit air bubble-free. You should carry out this function before the first determination or once per day.

The command **EMPTY** is used to empty the cylinders and the tubings of the dosing unit.

If no titrant is selected in the command, preparing and emptying will be carried out with the following default parameters:

- The entire cylinder volume is dosed at the maximum dosing rate through Port 1.
- The cylinder is filled at the maximum filling rate through Port 2.
- The following dimensions apply for the tubings:
 - Tubing at Port 1: length = 40.0 cm, diameter = 2 mm
 - Tubing at Port 2: length = 25.0 cm, diameter = 2 mm

31.2 Exchange unit

31.2.1 Maximum dosing and filling rate

The maximum dosing rate and maximum filling rate for the exchange unit depend on the cylinder volume:

Cylinder volume	Maximum rate
1 mL	3.00 mL/min
5 mL	15.00 mL/min
10 mL	30.00 mL/min
20 mL	60.00 mL/min
50 mL	150.00 mL/min

Independent of the cylinder volume, values ranging from 0.01 to 166.00 mL/min can always be entered. When the function is carried out the rate will be, if necessary, decreased automatically to the highest possible value.

31.2.2 Default parameters for preparing (PREP)

The **PREP** command is used to rinse and fill the cylinder and tubings of the exchange unit air bubble-free. You should carry out this function before the first determination or once per day.

If no titrant is selected in the command, preparing will be carried out with the following default parameters:

- The entire cylinder volume is dosed twice at the maximum dosing rate.
- The cylinder is filled at the maximum filling rate.

31.3 Stirring rate

The stirring rate can be adjusted in steps from -15 to +15.

The approximate rotational speed for the internal magnetic stirrer (depends on the product version) can be calculated with the following formula:

$$\text{Rotational speed/min (r/min)} = 125 \cdot \text{Stirring rate}$$

Example:

Configured stirring rate: 8

Rotational speed in revolutions per minutes = $125 \cdot 8 = 1,000$

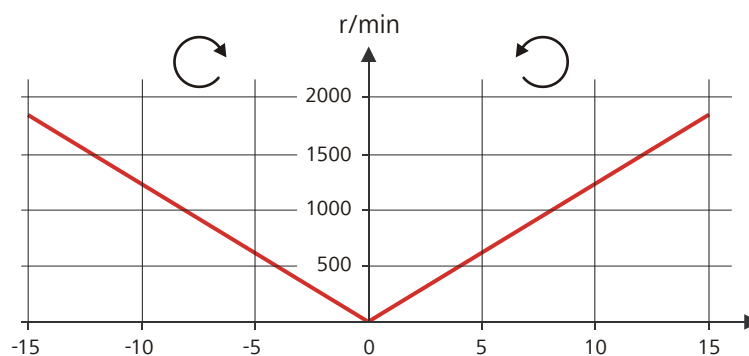


Figure 18 Rotational speed depending on the stirring rate

The information on the separately connectable 802 propeller stirrer can be found in the "802 Stirrer" manual.

31.4 Balance

The sample size and the associated unit can be sent from a connected balance. The sample size is transmitted as a number with up to ten characters (including algebraic sign and decimal point).

Sample size and unit are sent as a single character string. They are separated by a space character. The string is terminated with the ASCII characters **CR** and **LF**.

If the balance sends a negative sample size (e.g. when you are reweighing a sample), then the algebraic sign is adopted. The algebraic sign is, however, ignored for the calculations.



NOTICE

With some balances, the sample identification and the method can be sent in addition to the sample size.

Make sure that the balance does not send the sample size until the end.

Mettler AX

For the Mettler AX balance, the fields that contain the sample identification or the method must be designated as follows:

- Designation for the field with the method name: **METHOD**
- Designation for the field with sample identification 1: **ID1**
- Designation for the field with sample identification 2: **ID2**

31.5 Result variables as parameter setting

For most of the numerical parameters, the result of a calculation can also be utilized in place of a number. This result must be calculated in a preceding calculation command (command **CALC**). A typical application is the use of a relative start volume.

The following example is used to explain how this method is used to define a sample size-dependent start volume for a titration. Proceed as follows:

1 Create a new calculation

- Insert the calculation command **CALC** in front of the titration command.
- Enter the calculation formula, e.g. $R1 = C00 \cdot 3$.

The screenshot shows a software interface for editing a calculation command. The title bar reads 'Edit command / Edit calculation'. The main area displays the following settings:

- Command: 01 CALC
- Title: Calculation
- Result name: Relative start volume
- Calc. formula R1: C00*3
- Decimal places: 3
- Result unit: mL

At the bottom of the dialog, there are four buttons: 'Note', 'Result variable', 'Result limits', and 'Result options'.

2 Define a result variable as parameter setting

- In the titration command, enter the variable **R1** as the start volume.

The selection of the result variables is displayed in the number editor.



NOTICE

If the method before the command with the result variables applied contains more than one calculation commands (possibly with result variables bearing the same name), then the result variable of the preceding calculation command will always be used.

Please note that the result must be within the input range of the parameter, as otherwise the determination run will be stopped when the parameter is accessed.


31.6 Using AuditTrailViewer

When you work in accordance with the *FDA Guidance 21 CFR Part 11*, logs with detailed user activities (Audit Trail) are important. The *AuditTrailViewer* supports you in your efforts to compile, filter and export logs of this type as TXT files.

First you must install the *AuditTrailViewer* on your PC. Afterwards, you must create a backup of your Coulometer (see chapter 11.3, page 97). The Audit Trail is generated automatically when you create the backup. The file name of the Audit Trail is **log.madt**. In the following you will find detailed information about how you must proceed.

31.6.1 Installing AuditTrailViewer

You will find the file for installing the *AuditTrailViewer* on the USB flash drive provided.

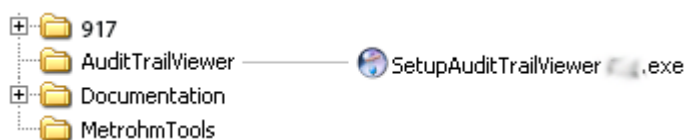

Proceed as follows:

- 1 Double-click on the file **SetupAuditTrailViewer 1.2.2.exe**.
The dialog **Welcome to the AuditTrailViewer Setup Wizard** is displayed.
- 2 Click on **[Next >]**.
The **License Agreement** dialog is displayed.
- 3 Click on **[I Agree]** to accept the contractual conditions.

The **Choose Install Location** dialog is displayed. A memory location for the *AuditTrailViewer* is suggested in this dialog.

- 4** If you accept the suggested memory location, click on **[Next >]**.

or

Select a different memory location and click on **[Next >]**.

The **Confirm Installation** dialog is displayed.

- 5** Click on **[Install]** to install the *AuditTrailViewer*.

As soon as the installation is completed, the dialog **Installation Complete** is displayed.

- 6** Click on **[Next >]**.

The dialog **Completing the AuditTrailViewer Setup Wizard** is displayed.

- 7** If the *AuditTrailViewer* should not be started immediately, deactivate the check box **Run AuditTrailViewer**.


- 8** Click on **[Finish]**.

If the preceding instruction step has been skipped, then the *AuditTrailViewer* will start.

31.6.2 Opening the Audit Trail

Create a backup of the data and of the settings of your Coulometer (see chapter 11.3, page 97).

- 1** Open the backup on your PC.

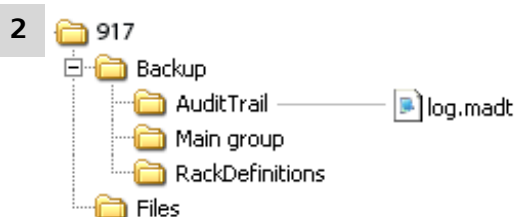

Open the file **log.madt** with the *AuditTrailViewer*.

AuditTrailViewer dialog window









File Filter Info					
No.	Date	User	Category	Action	Details
1	2011-10-31 11:30:22	Johnson	Method	New	01 Dynamic Titration pH
2	2011-10-31 11:30:34	Johnson	Method	Delete command	New method V0 03 REPORT
3	2011-10-31 11:30:40	Johnson	Method	Edit	New method V0 01 DET pH Stirrer off
4	2011-10-31 11:30:43	Johnson	Method	Edit	New method V0 01 DET pH Titrant not defined
5	2011-10-31 11:30:43	Johnson	Method	Edit	New method V0 01 DET pH Titrant
6	2011-10-31 11:30:48	Johnson	Method	Load result template	New method V0 Content (g/L)
7	2011-10-31 11:30:52	Johnson	Method	Edit	New method V0 Calc. formula EP1*CONC*TITER*3/C00
8	2011-10-31 11:31:08	Johnson	System	Message	003-908 Number of autostarts Yes/OK
9	2011-10-31 11:31:14	Johnson	Method	Start	Start key pressed
10	2011-10-31 11:31:15	Johnson	Method	Start	New method V0 1.0 g
11	2011-10-31 11:31:28	Johnson	Method	Manual stop	New method V0
12	2011-10-31 11:31:36	Johnson	System	Message	002-907 Method modified Yes/OK
13	2011-10-31 11:31:36	Johnson	Security	Logout	
14	2011-10-31 11:31:55	Meier	Security	Login message	Wrong password
15	2011-10-31 11:31:57	Johnson	System	Message	002-102 Wrong password Yes/OK
16	2011-10-31 11:31:58	Meier	Security	Log in	
17	2011-10-31 11:32:27	Meier	Method	Load	KFT 2 V1 Internal memory
18	2011-10-31 11:32:32	Meier	Security	Logout	
19	2011-10-31 11:32:45	Meier	Security	Login message	Wrong password
20	2011-10-31 11:32:47	Meier	System	Message	002-102 Wrong password Yes/OK
21	2011-10-31 11:33:01	Meier	Security	Change password	
22	2011-10-31 11:33:03	Meier	Security	Login message	Wrong password
23	2011-10-31 11:33:05	Meier	System	Message	002-102 Wrong password Yes/OK
24	2011-10-31 11:33:08	Meier	Security	Log in	
25	2011-10-31 11:33:12	Meier	Method	Start	Start key pressed
26	2011-10-31 11:33:13	Meier	Method	Start	KFT 2 V1 1.0 g
27	2011-10-31 11:33:14	Meier	System	Message	009-111 Stirrer missing Yes/OK
28	2011-10-31 11:33:15	Meier	Security	Logout	
29	2011-10-31 11:33:27	Chang	Security	Log in	
30	2011-10-31 11:33:32	Chang	Method	New	01 Dynamic Titration pH
31	2011-10-31 11:33:38	Chang	Method	Delete command	New method V0 03 REPORT
32	2011-10-31 11:33:43	Chang	Method	Edit	New method V0 01 DET pH Start volume 5 mL
33	2011-10-31 11:33:48	Chang	Method	Load result template	New method V0 Content (g/L)

31.6.3 Contents of the Audit Trail

In the following you will find a description of the dialog window and of the menu bar of the *AuditTrailViewer*.

Table 12 Dialog window

Column	Content
No.	Each entry is numbered in sequence.

Column	Content
Symbol	Classification of the entry: <ul style="list-style-type: none"> ▪  Actions that are neither relevant to safety nor alter the determination data. ▪  Actions, e.g. changes to the login options and recalculation of determinations. ▪  Errors that occur, e.g. entry of an incorrect password.
Date	Precise time of the event.
User	User who triggered the action.
Category	Category to which the entry belongs.
Action	Designation of the action.
Details	Details of the action.
Menu bar	
 Print	Print Audit Trail. An Audit Trail can be printed by means of the  symbol.
 Update	Update view. The list is updated automatically only at the time the dialog window is opened.
 Quick filter	Filter the Audit Trail.
 Show all	Display all entries once again.


31.6.4 Filtering the Audit Trail


The entries in the Audit Trail can be filtered. You can set the following filter criteria:

- **Date**
- **User**
- **Category**
- **Action**

Proceed as follows:

- 1 Click in a cell that contains the desired filter criterion (e.g. **Category** = **Method**).

2 Click on the  icon or the menu item **Filter ► Quick filter**.
The only entries which will be displayed are those which correspond to the desired filter criterion.

3 Click on the  icon or the menu item **Filter ► Show all**.
All of the entries are displayed once again.

31.6.5 Exporting the Audit Trail

You can export the Audit Trail as .txt file. This way, the Audit Trail can be archived and is available as required for inspections.

Proceed as follows:

1 Use the menu item **File ► Export** to open the dialog window for the export.

2 Enter the path and file name with the extension **.txt**.

3 Click on **[Open]** and then on **[Export]**.

The Audit Trail is exported.

31.7 Diagnosis

System ► Diagnosis

The electronic and mechanical functional groups of Metrohm instruments can and should be checked by specialist personnel from Metrohm as part of a regular maintenance schedule. Please ask your local Metrohm representative regarding the precise terms and conditions involved in concluding a corresponding maintenance agreement.

The following functions are available:

- **LCD test**
Check the display for faulty pixels (*see chapter 31.7.1, page 319*).
- **Format storage medium**
Format the external storage medium (*see chapter 31.7.2, page 319*).
- **Remove storage medium**
Remove the external storage medium safely (*see chapter 31.7.3, page 320*).
- **Touch adjustment**
Adjust the touch-sensitive screen (*see chapter 31.7.4, page 320*).

- **Touch screen test**
Test the function of the touch-sensitive screen (*see chapter 31.7.5, page 321*).
- **Software update**
Update the software of Coulometer and of the control instruments (*see chapter 31.7.6, page 322*).
- **Service**
Special functions for the service technician (*see chapter 31.7.7, page 325*).

31.7.1 LCD test

System ► Diagnosis ► LCD test

You can use the LCD test to check the display for faulty pixels. To do this, various test images are displayed one after the other.



NOTICE

The [↶] fixed key is used to display the preceding test picture; the test can be stopped at any time with the [🏠] fixed key.

Proceed as follows:

- 1 ▪ Tap on **[LCD test]**.
The entire display turns white.
- 2 ▪ Check the display for pixel errors and other irregularities.
 - Continue the test with [▶].
 Other colors and patterns will be displayed one after the other.
- 3 ▪ Repeat step 2 until the **System / Diagnosis** dialog is displayed again.
- 4 ▪ Contact your Metrohm representative in the event of faults or irregularities.

31.7.2 Formatting an external storage medium

System ► Diagnosis ► Format storage medium

With this function, you have the option of formatting an external storage medium directly on the 917 Coulometer (quick formatting).

The following file systems can be selected:

- **FAT**

- **FAT32**
- **ExFAT**



CAUTION

If you format the storage medium with the **ExFAT** file system, it is possible that it will no longer be recognized by a PC. Read the pertinent information from Microsoft Support at <http://support.microsoft.com>.

Do not use this file system unless you need to be able to save more than 999 files in a single group.

31.7.3 Removing an external storage medium

System ► Diagnosis ► Remove storage medium

As long as the stored data is not being accessed (reading/saving data), you can plug in and disconnect the storage medium without difficulty at any time. The **[Remove storage medium]** function offers additional protection. This ensures that the storage medium cannot be removed while any data is being transferred.

31.7.4 Adjusting the touch screen

System ► Diagnosis ► Touch adjustment

With time the touch-sensitive screen may not register touches at their precise position. You can readjust the screen in this dialog.

Proceed as follows:

1 In the main dialog, tap on **[System]**.

2 Tap on **[Diagnosis]**.

3 Tap on **[Touch adjustment]**.

A message in English and a cross hair will appear in the center of the screen.

4 Touch a stylus (a special pen-shaped instrument for operating devices with touch screens) to the center of the cross hair and maintain this position until the cross hair moves to a new position on the screen.

5 Repeat this process as long as the cross hair continues to be repositioned on the screen.

A message in English will appear.

6 Tap at any location on the screen to confirm the adjustment.

7 Press [] to finish the adjustment.

31.7.5 Testing the touch screen

System ► Diagnosis ► Touch screen test

You can use this test to check the function of the touch-sensitive monitor screen.

A calibration is not possible with this test. If the Touch screen does not work correctly anymore, it will need to be replaced. For that please contact your Metrohm representative.



NOTICE

You can use the [] fixed key to stop the test at any time.

Proceed as follows:


1 Tap on **[Touch screen test]**.


A red square is displayed at the upper left.

2 Tap on the square as close to the center as possible.

The actual detected point of contact will be displayed by a cross hair, and the deviation from the center will also be displayed on the lower edge of the image.

The next red square will also be displayed.

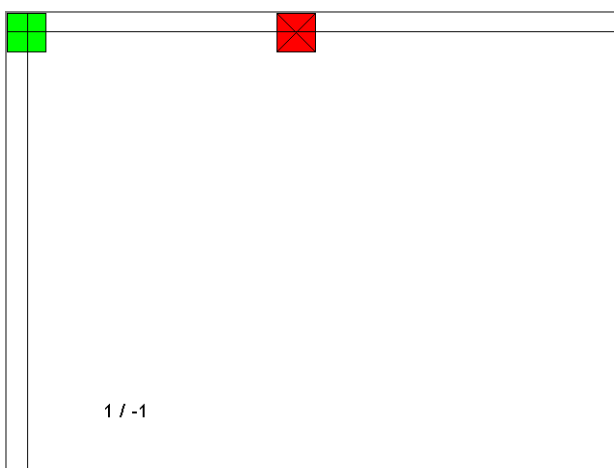

- 3 Repeat step 2 until all of the areas in the display have been tested.
A list with fixed keys will appear.
- 4 Tap on all of the fixed keys one after the other in any sequence.
The result is displayed in each case with a short information text, e.g. **[Print] key OK**.
- 5 Finish Touch screen test with [].

31.7.6 Software update (loading program versions and language files)

System ► Diagnosis ► Software update

New program versions or language files can be loaded from an external storage medium (e.g. USB flash drive). The corresponding files must be stored on the storage medium in the **917/SwUpdates** directory (see "Directory structure", page 96).



NOTICE

The memory will be reformatted when updating to software version 5.917.0023.

Before carrying out the software update, make a backup in order to save your data and the system settings (see chapter 11.3, page 97).

Use a separate USB flash drive for the software update. It must be removed from the instrument after the successful software update.

Program files

The files are instrument-specific. The file names are structured as follows:

- Control software for 917 Coulometer:
5XXXyyyy.BIN
5XXXyyyy.MBIN
 - XXX = instrument type (i.e. "917" for the 917 Coulometer)
 - yyyy = program version

Language files

Language files may have the following content:

- one or more additional dialog languages
- the online help for one or more additional dialog languages
- supplements for existing dialog languages/online help documents

They can be recognized by means of the two-digit language code in the file name. The file name has the following structure:

- **5XXXyZZZML.BIN**
5XXXyZZZML.MBIN
 - XXX = instrument type (i.e. "917" for the 917 Coulometer)
 - y = version number of the language package
 - ZZZ = program version

31.7.6.1 Carrying out a software update



CAUTION

Make sure that the power supply is guaranteed during the entire update process. Otherwise there is the chance that the instrument will no longer be able to be switched on and that it will need to be sent in for repair.



NOTICE

Make sure that no USB/RS-232 adapter is connected!

Proceed as follows:

1 Copying files to an external storage medium




NOTICE

In the case of the program versions and language files for the Coulometer, it is imperative that both the BIN file and the associated MBIN file be copied.

- Copy the files to the "SwUpdates" directory.

- Copy also the "MetrohmTools" folder with the "MUpdate-Tool.exe" file to the external storage medium. Make sure that the file is stored at the top level of the external storage medium. This also applies to the "917" folder.

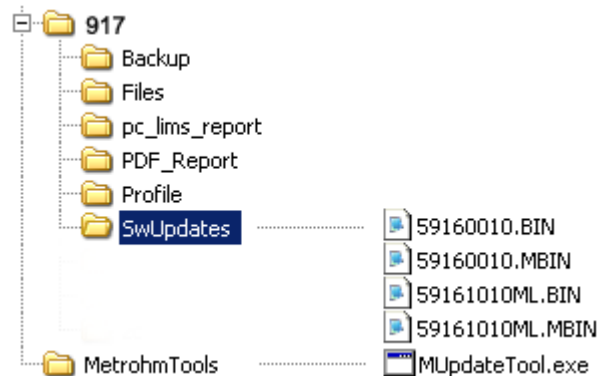

For the software update from version 5.917.0023, an additional folder that is on the same level as MetrohmTools with the name SRVPRF900 containing the file St900_SrPrf.cfg is required

If this structure is not maintained, then the files will not be found, because the directories are accessed directly during the update process.

- Plug in the external storage medium at the Coulometer.

2 Opening the update dialog

- Under **System ► Diagnosis**, tap on the button **Software update**.

3 Selecting the instrument

- Tap on the **Control device** list box and select the instrument to be updated.

4 Selecting the file

- Tap on the **Binary file** list box. The selection list with the BIN files saved in the "SwUpdates" directory is opened.
- Select the required file.
- Tap on **[Select]**.
- Update of the Coulometer: Continue with Step 5.

5 Starting the update (917 Coulometer)

- Tap on the **[Start]** button (NOT on the [▶] fixed key).

- The message **023-102 Program update** is displayed.
- Confirm the message with **[Yes]**.
The update process is started, it runs automatically. The device is automatically switched off and back on during this process, possibly several times. No user intervention is required.
- **After successfully updating the software, remove the external storage medium (USB flash drive) from the instrument.**

6 Starting the update (control instruments)

- Tap on the **[Start]** button (NOT on the [▶] fixed key).
If **Result = No errors** is displayed, then the update has been successful.
- **Remove the external storage medium (USB flash drive) from the instrument.**
- Switch the 917 Coulometer off and then back on again.

31.7.7 Service

System ► Diagnosis ► Service

Access to the service functions is password-protected and only accessible to Metrohm service technicians. We recommend that you have maintenance performed on the system regularly by the Metrohm Service Department. You can automatically monitor the time interval after which the next service visit is due (*see chapter 12.5.1, page 109*).

32.2 Measuring inputs

32.2.1 Generator electrode

One electrode connector (**Gen.**) for a generator electrode.
Iodine generation with current pulses or continuous current.

Current at the electrode **auto:** variable direct current
100, 200, 400 mA: pulsating direct current
I_{max} 400 mA
Titration rate max. 2.24 mg H₂O/min

32.2.2 Indicator electrode

One measuring input (**Ind.**) for an indicator electrode.

Measuring mode Determination with adjustable polarization current.
I_{pol} Voltametric endpoint indication with alternating current.
AC 5, 10, 20, 30 μA
DC -125 - +125 μA

32.2.3 Temperature sensor

One measuring input (**Temp.**) for a Pt1000 or NTC temperature sensor.

Measuring range
Pt1000 -150 - +250 °C
NTC -5 - +250 °C
For an NTC sensor with (R (25 °C) = 30000 Ω and B (25/50) = 4100 K)

Resolution
Pt1000 0.1 °C
NTC 0.1 °C

Measuring accuracy
Pt1000 ±0.2 °C
(Applies for measuring range -20 - +150 °C; ±1 digit; without sensor error, under reference conditions)
NTC ±0.6 °C
(Applies for measuring range +10 - +40 °C; ±1 digit; without sensor error, under reference conditions)

**32.2.4 Determination range**Recommended 10 µg to 200 mg H₂O**32.2.5 Resolution**0.1 µg H₂O**32.2.6 Reproducibility**

Sample: Standard of the reagent manufacturers

At 10 µg - 1000 µg H₂O: ±3 µgAt >1,000 µg H₂O: 0,3% or better**32.3 Internal stirrer***Rotational speed**Maximum rotational speed* ±(1700 - 1900) rpm*Rotational speed settings* ±15 steps*Change in rotational speed per step* ±(115 - 125) rpm*Power consumption* < 4 W*Fuse* Electronic overload protection**32.4 Pump***Add* > 600 mL/min (at 25 °C; standard accessories)*Aspirate* > 400 mL/min (at 25°C; standard accessories)

32.5 Interfaces and connectors

<i>Power connection</i>	For connecting the external power supply unit.
<i>MSB connector</i>	For connecting dosing devices, stirrers or a Remote Box.
<i>USB connector</i>	Type A, for connecting USB devices.
<i>Ethernet connector</i>	For connection to a data network (LAN).
<i>Sensor connectors</i>	One connector each for: <ul style="list-style-type: none"> ▪ Generator electrode ▪ Indicator electrode ▪ Temperature sensor (Pt1000 or NTC)
<i>Two connection nipples</i>	For connecting tubing for aspirating solvent and extracting the contents of the titration cell.

32.6 Power supply

<i>Instrument</i>	
<i>Input voltage</i>	24 V DC
<i>Max. power consumption</i>	2.0 A
<i>External power supply unit</i>	6.2164.010, LPS conformity according to UL60950-1
<i>Input current</i>	1.5 A
<i>Max. output current</i>	2.7 A

32.7 Declarations of Conformity

Up-to-date information on the *EU conformity* and the *requirements for electromagnetic compatibility* for your product can be found on the Internet under <https://www.metrohm.com/en-us/products-overview/>.

Using the article number, you can download the *Declaration of conformity for ...* under the documents of the product.



32.8 Ambient temperature

<i>Nominal function range</i>	+5 to +45 °C at max. 80% relative humidity, non-condensing
<i>Storage</i>	+5 to +45 °C

32.9 Reference conditions

<i>Ambient temperature</i>	+25 °C (± 3 °C)
<i>Relative humidity</i>	$\leq 60\%$
<i>Operating temperature status</i>	Instrument in operation at least 30 min
<i>Validity of the data</i>	After adjustment

32.10 Dimensions

<i>Width</i>	193 mm
<i>Height</i>	
<i>without titration vessel holder</i>	145 mm
<i>Height</i>	
<i>with titration vessel holder</i>	230 mm
<i>Depth</i>	448 mm
<i>Weight (including power supply unit)</i>	6020 g
<i>Material</i>	
<i>Housing</i>	Poly(butylene terephthalate) (PBT)
<i>Base</i>	Steel, stainless, coated

32.11 Storage capacity


<i>Random access memory</i>	128 MB
<i>Internal memory</i>	1 GB of which 900 MB available.
<i>Required memory per method</i>	3–6 KB

33 Accessories

Up-to-date information on the scope of delivery and optional accessories for your product can be found on the Internet. You can download this information using the article number as follows:

Downloading the accessories list

- 1 Enter <https://www.metrohm.com/> into your Internet browser.
- 2 Enter the article number (e.g. **917**) into the search field.
The search result is displayed.
- 3 Click on the product.
Detailed information regarding the product is shown on various tabs.
- 4 On the **Included parts** tab, click on **Download the PDF**.
The PDF file with the accessories data is created.

 **NOTICE**

Once you have received your new product, we recommend downloading the accessories list from the Internet, printing it out and keeping it together with the manual for reference purposes.

Index

- 800 Dosino 22
805 Dosimat 22
- A**
- Acoustic signals 55
ADD 278
Addition and aspiration tube
 Mount 14
Adsorber tube
 Fill 10
Audit Trail 53
 Archive 318
 Export 318
 Filter 317
 Open 315
AuditTrailViewer
 Installing 314
Automatic system test 100
Automation 289
- B**
- Backup 97
 Interval 109
Balance 25, 83, 312
Barcode reader
 Connect 26
BEEP 300
BRC 1, 30
 Parameters 238
Bromine index determination 30
- C**
- CALC 256
CALC LIVE 262
Calculation 256
 Add 256
 Delete 256
 Edit 256
 Load template 257
 Note 260
 Result options 261
Cell
 BRC 244
 KFC 235
Command
 ADD 278
 BEEP 300
 BRC 238
 CALC 256
 CALC LIVE 262
CONTROL RS 289
CTRL 287
EMPTY 277
END 301
EVAL FIX-EP 253
EVAL MIN/MAX 254
EVAL RATE 255
FLOW 296
HEATER 296
Insert 136
KFC 229
LIFT 291
LQH 282
MEAS 247
MEAS T/F OFF 298
MEAS T/F ON 297
MOVE 289
PREP 276
RACK 292
REPORT 271
REQUEST 299
SAMPLE 292
SCAN 285
SCAN RS 288
SIGN 301
STIR 298
SUBSEQ 293
WAIT 299
Command list 135
Common variable
 Delete 111
 Edit 111
 Properties 113
 Validity 114
Communication 285
Conditioning
 BRC 242
 KFC 233
Connect
 Balance 25
 Barcode reader 26
 Dosing device 22
 MSB devices 21
 PC keyboard 26
 Printer 24
 Remote Box 23
 Stirrer or titration stand 23
 USB devices 24
 USB hub 24
Control 151
Control instrument 72
Control parameters
 BRC 238
 KFC 229
Control Remote Box 66
CONTROL RS 289
Coulometer 66
Coulometric water content determination according to Karl Fischer 29
CTRL 287
Curve options 191
Curves 191
Custom logo 130
- D**
- Delete
 Signature 150
 Statistics data 153
Determination
 Cancel 172
 Carry out 170
 Data 178
 Define filters 188
 Delete signatures 150
 Endpoint details 181
 List of saved determinations 186
 Load 185
 Measuring point list 180
 Message 182
 Properties 183
 Recalculate 193
 Reevaluate 193
 Save 190
 Save automatically 146
 Send PC/LIMS report 146
 Sign 149
Determination run 170
Device
 885 Compact Oven SC 75
 Add 65
 Balance 83
 Coulometer 72
 Delete 64
 Edit 64
 New 64
 Printer 79

- USB/RS-232 adapter 85
- Device manager 64
- 885 Compact Oven SC 75
- Diagnosis 318
- Dialog language 37
- Dialog options 39
- Dialog settings
- System-specific 39
- Directory structure 96
- Display 7
- Dosing 217
- Continuous 218
- Fixed volume 220
- Dosing curve 191
- Dosing device
- Connect 22
- Parameters 279, 284
- Dosing parameters
- ADD 278
- Dosing rate
- Maximum 310, 311
- Dosing unit
- Replace reagent 222
- E**
- E-mail 67
- Electrode 60
- Connect 16
- Parameters 236, 245, 249
- Electrostatic charge 5
- EMPTY 277
- Default parameters 310
- Emptying 222
- Default parameters 310
- END 301
- Endpoint details 181
- Ethernet connector 8, 28
- EVAL FIX-EP 253
- EVAL MIN/MAX 254
- EVAL RATE 255
- Evaluation
- Fixed endpoint 253
- Maximum 254
- Minimum 254
- Rate 255
- Evaluations 252
- Expiry date
- Sensor 62
- F**
- Favorite 155
- Add 156
- Delete 156
- Edit 156
- FDA Guidance 21 CFR Part 11
..... 42, 53
- File
- Copy 93
- Properties 94
- Rename 93
- File manager 91
- Filling 222
- Filling rate
- Maximum 310, 311
- Filter
- Determination 188
- Result table 202
- Fixed endpoint
- Evaluation 253
- Fixed key
- Back 7
- Help 7
- Home 7
- Manual 7
- Print 7
- START 7
- STOP 7
- FLOW 296
- Format storage medium 319
- Formula editor 264
- G**
- Generator electrode 8, 16
- Fill 13
- GLP manager 99
- GLP test "Measurement" .. 101
- GLP test "Titration" 101
- System validation 105
- Test tool 100
- H**
- HEATER 296
- I**
- ID profile
- Create 47
- Identification profile
- Create 47
- Indicator electrode 8, 18
- Initialization 302
- Input line
- Template 125
- Instrument
- Configure 65
- Coulometer 66
- Switch off 32
- Switch on 31
- K**
- Key functions 87
- Keyboard
- Connect 26
- KFC 1, 29, 229
- L**
- LCD test 319
- LIFT 291
- Limit values
- Sample size 143
- Liquid Handling parameters 282
- Live display 175
- Live modifications 173
- Live parameters 176
- Live result
- Note 264
- Load
- Determination 185
- Sample table 165
- Load result template 257
- Login
- Options 48
- Logout automatically 49
- LQH 282
- M**
- Manual control 213
- 885 Compact Oven SC 225
- Continuous dosing 218
- Dosing 217
- Dosing fixed volume 220
- Emptying 222
- Filling 222
- Measuring 215
- Preparing 221
- Remote lines 224
- Stirring 223
- Troubleshooting 306
- Maximum evaluation 254
- MEAS 247
- MEAS T/F OFF 298
- MEAS T/F ON 297
- Measured value
- Display resolution 55
- Measurement curve 191
- Measuring 215
- Measuring parameters
- MEAS 247
- Measuring point list 180
- Message
- Determination 182
- Send e-mail 67
- Method 132
- Command list 135

- Create 132
 Delete signatures 150
 Edit 135
 Insert command 136
 Load 134
 Note 146
 Options 137
 Properties 144
 Sample data 141
 Save 133
 Save as favorite 155
 Sign 149
 Start options 138
 Stop options 140
 Troubleshooting 304
 Metrohm Serial Bus 8
 Metrohm Serial Bus MSB, see also
 "MSB" 21
 Minimum evaluation 254
 Miscellaneous commands 298
 Molecular sieve
 Replace 11
 Monitoring
 Limit values sample size 143
 MOVE 289
 MSB
 Connect devices 21
 MSB connector 8
- N**
- Network connection 28
 Note
 Method 146
 Numerical input 35
- O**
- Operation
 General 32
 Output line
 Template 127
 Oven module
 Properties 78
- P**
- Paper size 83
 Parameter
 FLOW 296
 HEATER 296
 MEAS T/F OFF 298
 MEAS T/F ON 297
 Parameters
 ADD 278
 BEEP 300
 BRC 238
 CALC 256
 CALC LIVE 262
 CONTROL RS 289
 CTRL 287
 EMPTY 277
 END 301
 EVAL FIX-EP 253
 EVAL MIN/MAX 254
 EVAL RATE 255
 KFC 229
 LIFT 291
 LQH 282
 MEAS 247
 Modify live 176
 MOVE 289
 PREP 276
 RACK 292
 REPORT 271
 REQUEST 299
 SAMPLE 292
 SCAN 285
 SCAN RS 288
 SIGN 301
 STIR 298
 SUBSEQ 293
 WAIT 299
 Password
 Forgotten 51
 Options 50
 PC keyboard
 Connect 26
 PC/LIMS report
 Save 68
 Send 146
 PDF settings 80
 Peripheral devices 64
 Power connection 8
 Power supply
 Power supply unit 15
 Power supply unit
 Connect 15
 Power switch 8
 PREP 276
 Default parameters ... 310, 311
 Preparing 221
 Default parameters ... 310, 311
 Printer 24, 79
 PDF settings 80
 Printing 205
 Troubleshooting 305
 Properties
 885 Compact Oven SC 76
 Determination 183
 File 94
 Method 144
 Oven module 78
 Tower 77
 Pump
 Troubleshooting 306
- R**
- RACK 292
 Rate evaluation 255
 Reagent 56
 Add 56
 Delete 56
 Edit 56
 Replace 222
 Reagent monitoring 57
 Reagent replacement 59
 Reason 52
 Modification option 52
 Recalculation 193
 Reevaluation 193
 Remote Box
 Connect 23
 Remote lines
 Manual control 224
 Report 205
 Print 205
 REPORT 271
 Report header 130
 REQUEST 299
 Resolution
 Measured value display 55
 Result 178
 Calculation formula 259
 Limits 260
 Save as a common variable
 115
 Template 121
 Result table 199
 Define filters 202
 Delete automatically 203
 Save 204
 Result variable 260
 As parameter setting 313
 Robotic arm
 Swing rate 291
 Routine dialog 41
 Standard configuration 41
- S**
- Safety instructions 4
 SAMPLE 292
 Sample assignment table 117
 Sample data 159
 Modify live 173
 Settings 141

Index

- Sample identification
 - Modify designation 141
 - Modify the designation 169
- Sample identification list 117
- Sample series
 - Perform 171
- Sample size
 - Fixed 141
 - Limit value 143
 - Limits 141
 - Send from a balance 312
- Sample table 162
 - Save as favorite 155
 - Troubleshooting 304
- Save
 - Determination 190
 - Result table 204
 - Sample table 164
- SCAN 285
- SCAN RS 288
- Sensor 60
 - Add 60
 - Connect 16
 - Delete 60
 - Edit 60
 - Expiry date 62
 - Parameters 236, 245, 249
 - Working life 62
- Sensor list 60
- Service 4, 325
- Set the date 38
- Set the time 38
- Sign
 - Determination 149
 - Method 149
- SIGN 301
- Signature
 - Delete 150
- Software update 322
- Spooler 83
- Start options 138
- Statistics 153, 194
 - Delete 153
 - Troubleshooting 305
- STIR 298
- Stirrer
 - Connect 23
 - Parameters 237, 246, 251, 281
- Stirring 223
- Stirring rate 311
- Stop conditions
 - BRC 242
 - KFC 232
- Stop criterion
 - BRC 240
 - KFC 231
- Stop options 140
- SUBSEQ 293
- Subsequence 293
 - Load 294
 - Note 296
 - Options 295
 - Save 294
- Supply voltage 4
- Swing rate 291
- Switch off 32
- Switch on 31
- System initialization 302
- System settings 37
- System test 100
- System validation 105
- T**
 - Tandem dosing 279
 - Temperature sensor 8, 20
 - Connect 16
 - Template
 - Input line 125
 - Output line 127
 - Report header 130
 - Result template 121
 - Sample assignment table .. 117
 - Sample identification list ... 117
- Templates 117
- Text input 35
- Time zone 38
- Titration cell
 - Equip 12
 - Fill 13
 - Insert 10
 - Set up 10
- Titration curve 191
- Titration mode
 - BRC 1
 - KFC 1
- Titration parameters
 - BRC 241
 - KFC 232
- Titration stand 7
 - Connect 23
 - Troubleshooting 306
- Touch screen 32
 - Test 321
- Tower
 - Properties 77
- Type plate 8
- U**
 - Uninterrupted dosing 279
 - Update 322
- USB
 - Connect devices 24
- USB connector 8
- USB hub
 - Connect 24
- USB/RS-232 adapter 85
- User profile
 - Create 47
- V**
 - Validation interval 107
- Validity
 - Common variable 114
- W**
 - WAIT 299
 - Working life
 - Sensor 62
 - Write protection
 - External memory 95