



# **Tin compounds in strawberry**

## **Separation of organometallic (tin) compounds**

### **Application Note**

Food Testing & Agriculture

#### **Authors**

Agilent Technologies, Inc.

#### **Introduction**

Separation of organometallic (tin) compounds is possible via gas chromatography with Agilent FactorFour VF-1701ms, using inert transfer lines. Trace analysis is done using MS detection with single ion monitoring/scanning.



**Agilent Technologies**


## Conditions

Technique : GC-capillary  
Column : Agilent FactorFour VF-1701ms, 0.25 mm x 30 m fused silica (df = 0.25 µm) (Part no. CP9151)  
Temperature : 100 °C (2 min) → 250 °C, 10 °C/min (5 min) → 280 °C, 10 °C/min (15 min)  
Carrier Gas : Helium, 1 mL/min  
Injector : Gerstel, PTV, 5 µL  
35 °C (0.5 min) → 300 °C, 12 °C/s (4 min)  
Detector : MS TIC

Courtesy : Herr DLC Godeck, Pesticide Residue Analysis, Gesellschaft für Lebensmittelforschung, Berlin

## Peak identification

1. cyhexatin
2. fentinacetate
3. fenbutatinoxide

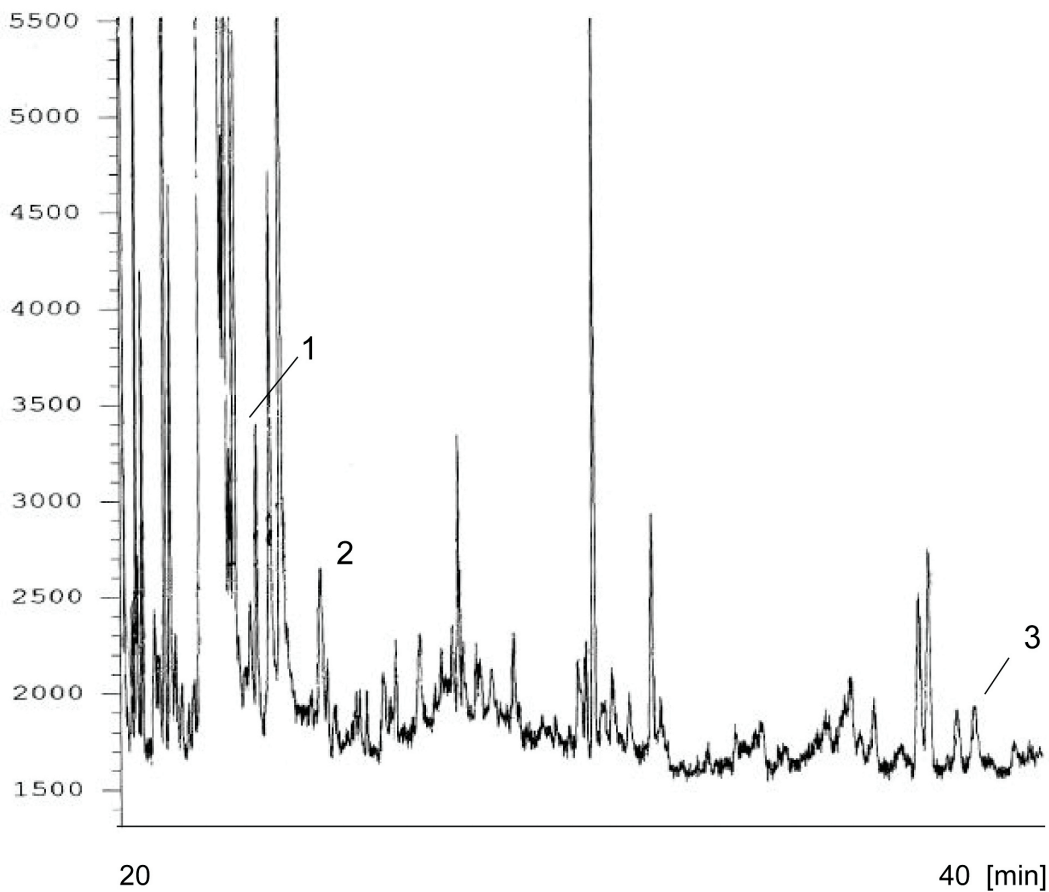

[www.agilent.com/chem](http://www.agilent.com/chem)

This information is subject to change without notice.

© Agilent Technologies, Inc. 2011

Printed in the USA

31 October, 2011

First published prior to 11 May, 2010

A02396



**Agilent Technologies**