


NOTICE: This document contains references to Varian.  
Please note that Varian, Inc. is now part of Agilent  
Technologies. For more information, go to  
[www.agilent.com/chem](http://www.agilent.com/chem).



Varian B.V.  
Herculesweg 8  
4330 EA  
Middelburg

# Micro-GC Modification Kit

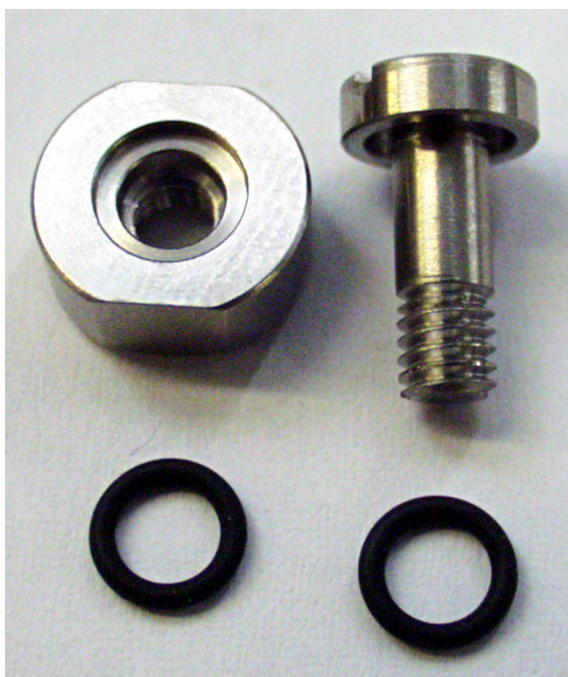

## Carriergas tube stop

North/South America  
2700 Mitchell Drive  
Walnut Creek  
94598 California, USA  
Tel: ++(1)9259392400  
Fax: ++(1)9259452360 or  
++(1)9259452344

Europe  
P.O. Box 8033  
4330 EA Middelburg  
The Netherlands  
Tel: ++(31)118671000  
Fax: ++(31)118623193

Australia/East Asia  
679 Springvale Road  
Mulgrave, Victoria 3171  
Australia  
Tel: ++(61)395607133  
Fax: ++(61)395607950

This kit CP740828 contains:

- 1) CP740029 Viton O-rings 2x
- 2) CP740209 Stop carrier tube
- 3) CP740210 Nut for stop carrier tube



- 1) Install the O-rings over the stop and nut.



- 2) Uninstall the GC-channel according the Channel exchange manual (CP501334).
- 3) Install the stop and nut on the carriergas tube.