


Maintaining Sensitivity and Reproducibility with the JetClean Self-Cleaning Ion Source for Pesticides in Food and Feed

Jessica Westland¹ and Elizabeth Almasi²

¹Agilent Technologies, Wilmington, DE 19808 USA; ²Agilent Technologies, Santa Clara, CA 95051 USA

NACRW 2016

P-064

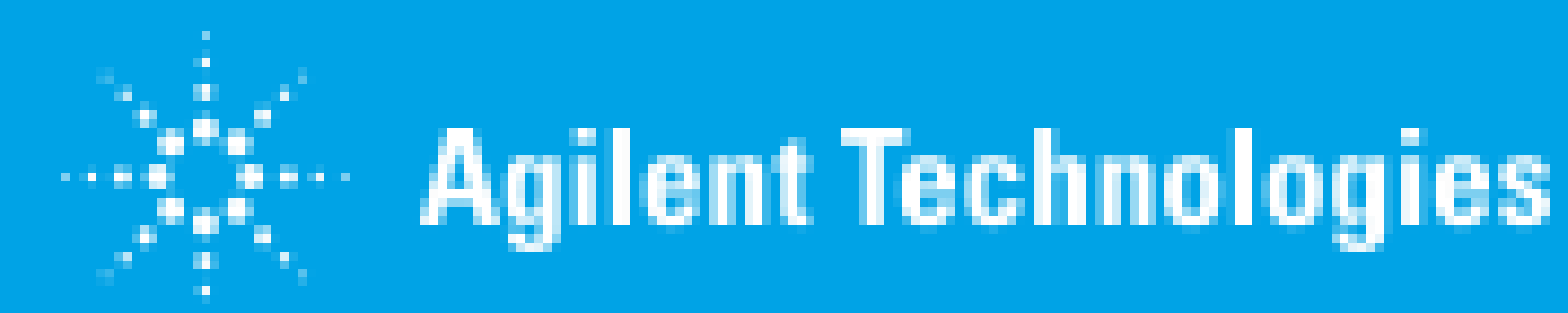

Introduction

The global agricultural industry uses over a thousand pesticides for food and foodstuffs cultivation. Producers are compelled to use pesticides to meet the growing demand for reasonably priced food, resulting in the need for pesticide residue monitoring in commodities worldwide. Concurrently, simple sample preparation practices, such as QuEChERS are routinely used for the preparation of food and feed samples, often leaving significant amount of matrix in the extracts. Analytical laboratories are challenged by these matrix residues, which with time negatively affects the response of the analyzed pesticides, and eventually requires source cleaning. Agilent's JetClean self-cleaning ion source (JetClean) reduces the need for manual source cleaning while still allowing for the analysis of complex samples without losing sensitivity and reproducibility.

JetClean

The Agilent JetClean utilizes carefully monitored hydrogen gas (H₂) introduction to the source, controlled by Agilent's MassHunter Data Acquisition Software. The appropriate H₂ flow (in the $\mu\text{L}/\text{min}$ range) generates conditions that clean the surfaces of the source, the lenses and other components. These actions aid in maintaining a stable detection environment and provide for response stability of the pesticides in difficult matrices.

JetClean has two operational modes:

1. Acquire and Clean (or on-line) mode, when H₂ is running during the analysis
2. Clean only (or off-line) mode when H₂ is introduced only post run or post sequence

Experimental

Methodology: The analysis was conducted on an Agilent 7890B GC and 7010 Series Triple Quadrupole GC/MS. See Tables 1 – 3 for method parameters. The system was configured with a Multimode Inlet, equipped with an ultra-inert liner (p/n: 5190-2293). The inlet was then connected to two HP-5ms UI columns (15 m \times 0.25 mm \times 0.25 μm ; p/n: 19091S-431 UI) coupled to each other through a purged ultimate union (PUU) for the use of backflushing (see Figure 1).

The H₂ cleaning was operated in the "Acquire & Clean" mode which allowed constant H₂ flow during the analytical runs.


Table 1. 7890B GC Method Conditions		
Injection port liner	4-mm Ultra Inert liner with wool	
Injection mode	Hot-splitless	
Injection volume	1 μL	
Inlet temperature	280 °C	
Carrier gas	He, constant flow 1.00 mL/min (column 2 = 1.20 mL/min)	
Oven program	40	60 °C 1 min
	120 °C	0 min
	5 °C/min	310 °C 0 min
MS transfer line temperature	280 °C	

Table 2. PUU Backflush Settings*	
Timing	1.5 min during post-run
Oven temperature	310 °C
Aux EPC pressure	~50 psi
Inlet pressure	~2 psi

*Backflush conditions were optimized for the application. A 1.5 min backflush duration may be too short for other methods. It can be extended up to 5 min duration.

Table 3. 7010 MS/MS Parameters	
Electron Energy	70 eV
Tune	atunes.eihs.tune.xml
EM gain	10
MS1 & MS2 resolution	Wide
Collision Cell	1.5 mL/min N ₂ & 2.25 mL/min He
Quant/Qual transitions	Matrix Optimized
Dwell times	Time Segment (TS) specific*
Source temperature	300 °C
Quad temperatures	150 °C
JetClean	0.15 mL/min H ₂


*The dwell times in each TS were the same, all with values over 10msec, resulting in a scan rate of ~5 scans/sec for the TS.

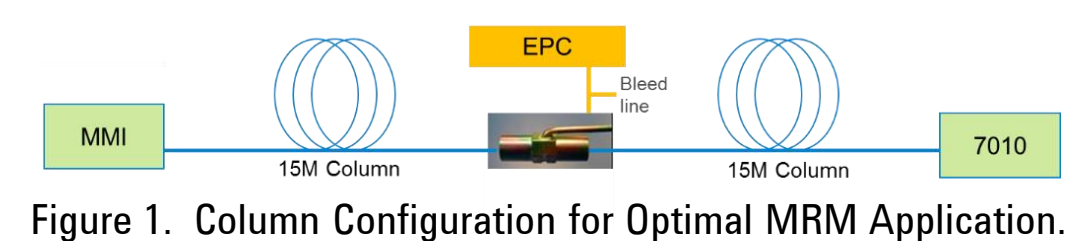

Operation

Previously the introduction of H₂ flow in to the MS source was introduced via EPC module. The next stage of Agilent's innovative technology, the JetClean SCIS, moves the control of H₂ flow to the MS. MassHunter GC/MS Data Acquisition and the MS acquisition control are shown below.

For this application JetClean was initiated in the acquire and clean mode for continuous cleaning. The MassHunter software allowed for the simple setup and operation of the process, all controlled in the MS domain.

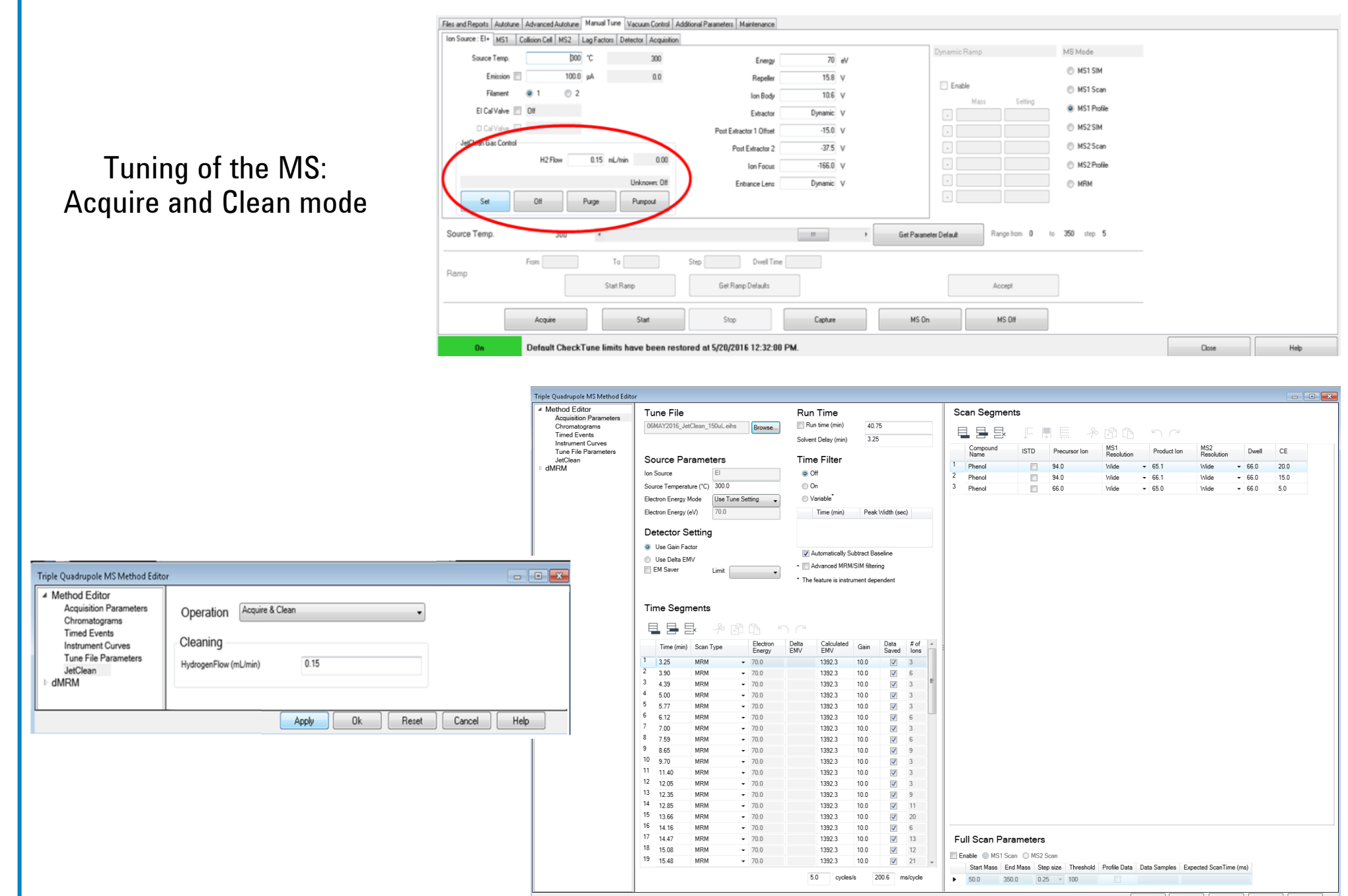


Results and Discussion

Agilent's Self-cleaning ion source has been successfully used for the extended, 64 analyte PAH analysis in environmental¹ and labs and also in food laboratories, resulting in remarkable precision, accuracy, linearity and detection levels, which were sustained for extended periods of time (several months) without manual cleaning. Considering the benefits of the Self-cleaning ion source for PAH analysis, the use of JetClean was applied for cleaning the MS source during pesticide analysis.

Chromatographic Performance: The following chromatograms show analytes at 2.5 pg concentrations, eluting at the beginning, middle and at the end of the chromatographic run. The ion plots are of target compounds and their respective matrix optimized MRM transitions in organic honey using standard source configuration and with the JetClean source. The JetClean benefits on peak shape and baseline are more obvious on the later eluting, higher MW analytes.

DBCP MW 234 Heptenophos MW 250 Hexachlorobenzene MW 282 Parathion-methyl MW 451 Endosulfan-II MW 404 Chlorobenzilate MW 324 Flucythrinate-1 MW 451

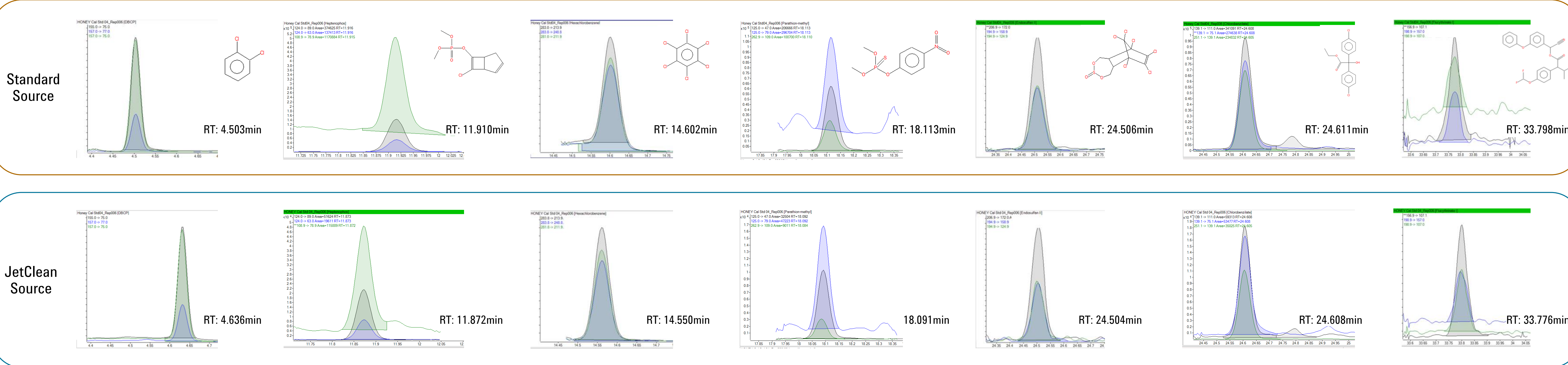


Results and Discussion


Quantitative results: Table 5 lists the R² values and the statistically derived MDLs for representative target analytes of the over 170 various pesticides tested. The calibration ranged from 0.12 pg/ μL - 50pg/ μL for the majority of the analytes, although some were not included at the lowest level. The resulting R² values were very comparable by both source type. The MDLs were calculated from 10 replicate measurement of 1.25 pg/ μL concentration spiked honey extract using 99% confidence level. Lower MDLs were obtained for the majority of the analytes using the JetClean source, with an average of 0.151 pg MDL for the standard source and 0.081pg for the JetClean source. The replicate measurements performed at 1.25pg level resulted lower %RSD using the JetClean source, although they were comparable at the 2.5 pg level.

Analyte	R ²		LOQ (pg)		MDL (pg)	
	STD	JetClean	STD	JetClean	STD	JetClean
Alachlor	0.996	0.978	0.118	0.050	0.119	0.050
Azinphos-ethyl	0.999	0.995	0.056	0.071	0.063	0.071
Bendiocarb	0.994	0.937	1.636	1.263	0.719	1.263
Benfluralin	0.991	0.994	0.173	0.042	0.143	0.042
BHC-alpha	0.995	0.995	0.414	0.098	0.221	0.098
Carboxin	0.991	0.994	0.065	0.025	0.021	0.025
Chlorfentofen	0.996	0.997	0.043	0.020	0.070	0.020
Chlorobenzilate	0.996	0.995	0.367	0.118	0.557	0.118
Chlorpyrifos-methyl	0.994	0.993	0.256	0.075	0.270	0.075
DBCP	0.995	0.986	0.364	0.208	0.155	0.208
DDT,p,p'	0.993	0.996	0.170	0.105	0.427	0.105
Dichlorobenzene, 1,2-	0.999	0.993	0.241	0.184	0.129	0.184
Dimethoate	0.999	0.992	0.143	0.028	0.150	0.028
Endosulfan I	0.995	0.996	0.045	0.032	0.027	0.032
Endosulfan II	0.991	0.998	0.034	0.040	0.037	0.040
Endosulfan sulfate	0.997	0.996	0.079	0.036	0.136	0.036
Endrin	0.992	0.994	0.045	0.025	0.045	0.025
Endrin ketone	0.993	0.996	0.022	0.016	0.041	0.016
Ethofenprox	0.996	0.992	0.293	0.036	0.161	0.036
Ethionphos	0.995	0.995	0.163	0.016	0.072	0.016
Fenitrothion	0.993	0.991	0.249	0.060	0.313	0.060
Fenitrothion sulfone	0.996	0.994	0.077	0.040	0.046	0.040
Flucythrinate I	0.997	0.992	0.066	0.023	0.040	0.023
Haloxypip-r-methyl	0.999	0.994	0.063	0.060	0.085	0.060
Heptachlor endo-epoxide	0.994	0.992	0.104	0.022	0.085	0.022
Heptachlor exo-epoxide	0.992	0.996	0.550	0.045	0.264	0.045
Hexachlorobenzene	0.999	0.997	0.606	0.116	0.227	0.116
Iprodione	0.993	0.999	0.025	0.029	0.041	0.029
Irgarol	0.993	0.997	0.100	0.037	0.117	0.037
Methoxychlor, p,p'	0.994	0.996	0.112	0.046	0.170	0.046
Nitrofen	0.990	0.995	0.118	0.030	0.097	0.030
Oxadixyl	0.994	0.994	0.076	0.030	0.245	0.030
Parathion-methyl	0.994	0.995	0.263	0.070	0.272	0.070
Pentoxazone	0.998	0.993	0.082	0.034	0.137	0.034
Permethrin, (1R)-trans-	0.997	0.968	0.061	0.105	0.033	0.105
Phosphamidon II	0.996	0.988	0.107	0.031	0.235	0.031
Pirimiphos-methyl	0.996	0.994	0.148	0.042	0.152	0.042
Profenofos	0.994	0.996	0.065	0.036	0.023	0.036
Pyrazophos	0.994	0.996	0.064	0.056	0.042	0.056
Terbufos	0.994	0.996	0.239	0.032	0.078	0.032
Tetrachlorvinphos, E-isomer	0.998	0.996	0.140	0.029	0.112	0.029
Thionazin	0.994	0.994	0.139	0.038	0.097	0.038
Trifluralin	0.997	0.994	0.189	0.033	0.088	0.033


LOQ comparison with and without JetClean obtained at the 2.5 pg level

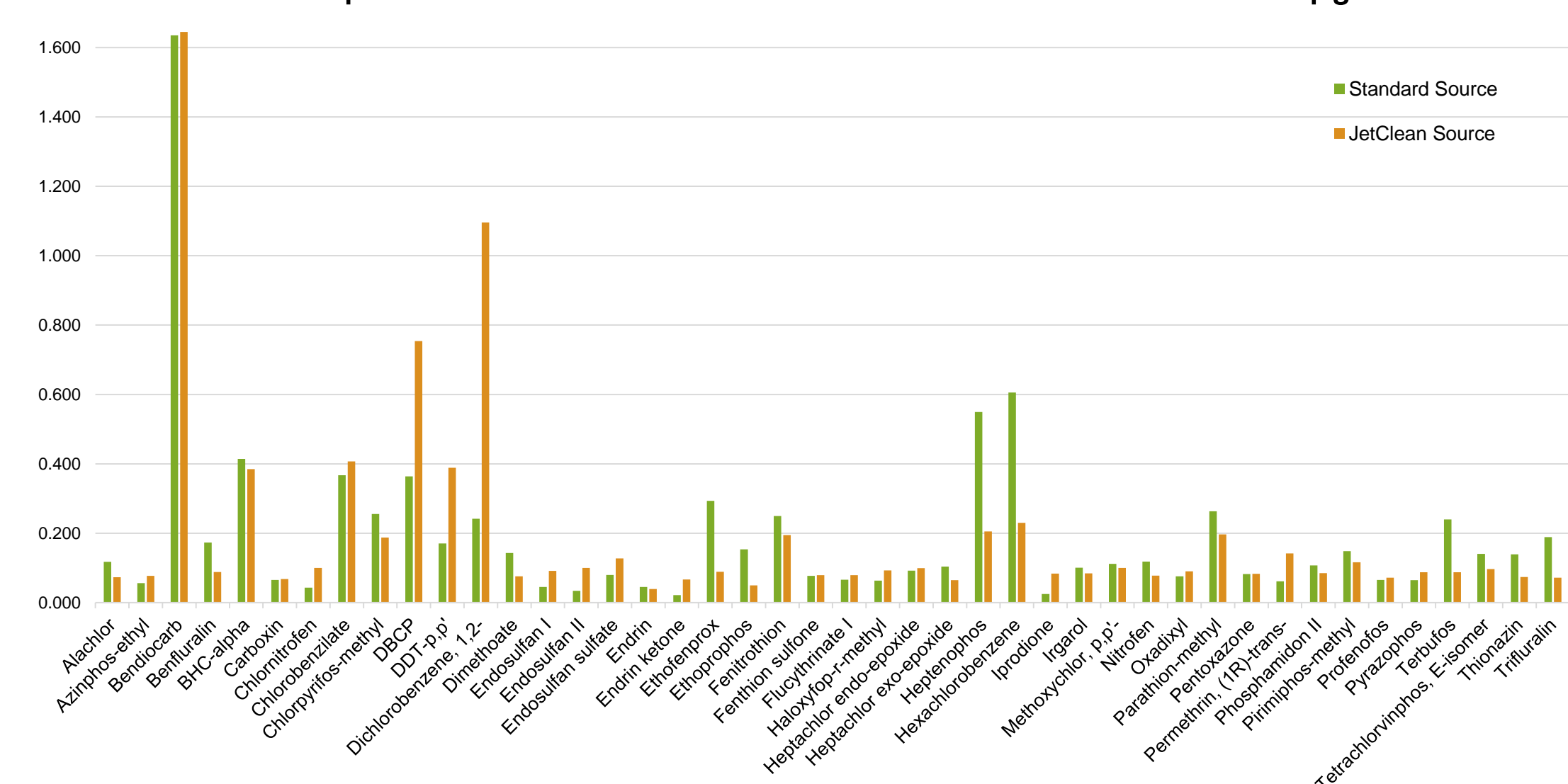

MDL comparison with and without JetClean obtained at the 2.5 pg level

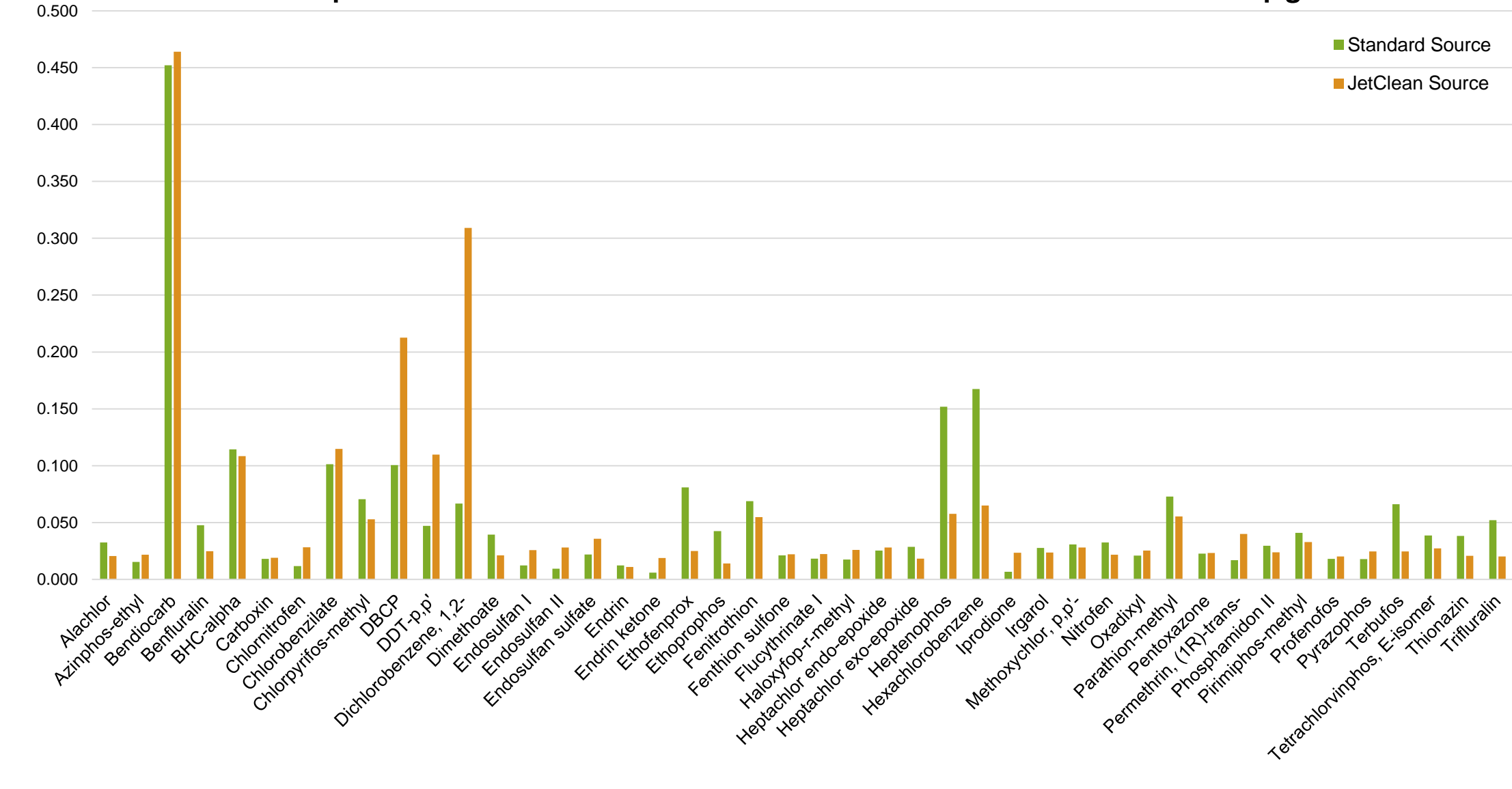

%RSD ratio of the Standard source/JetClean source results obtained at 2 lower concentration levels

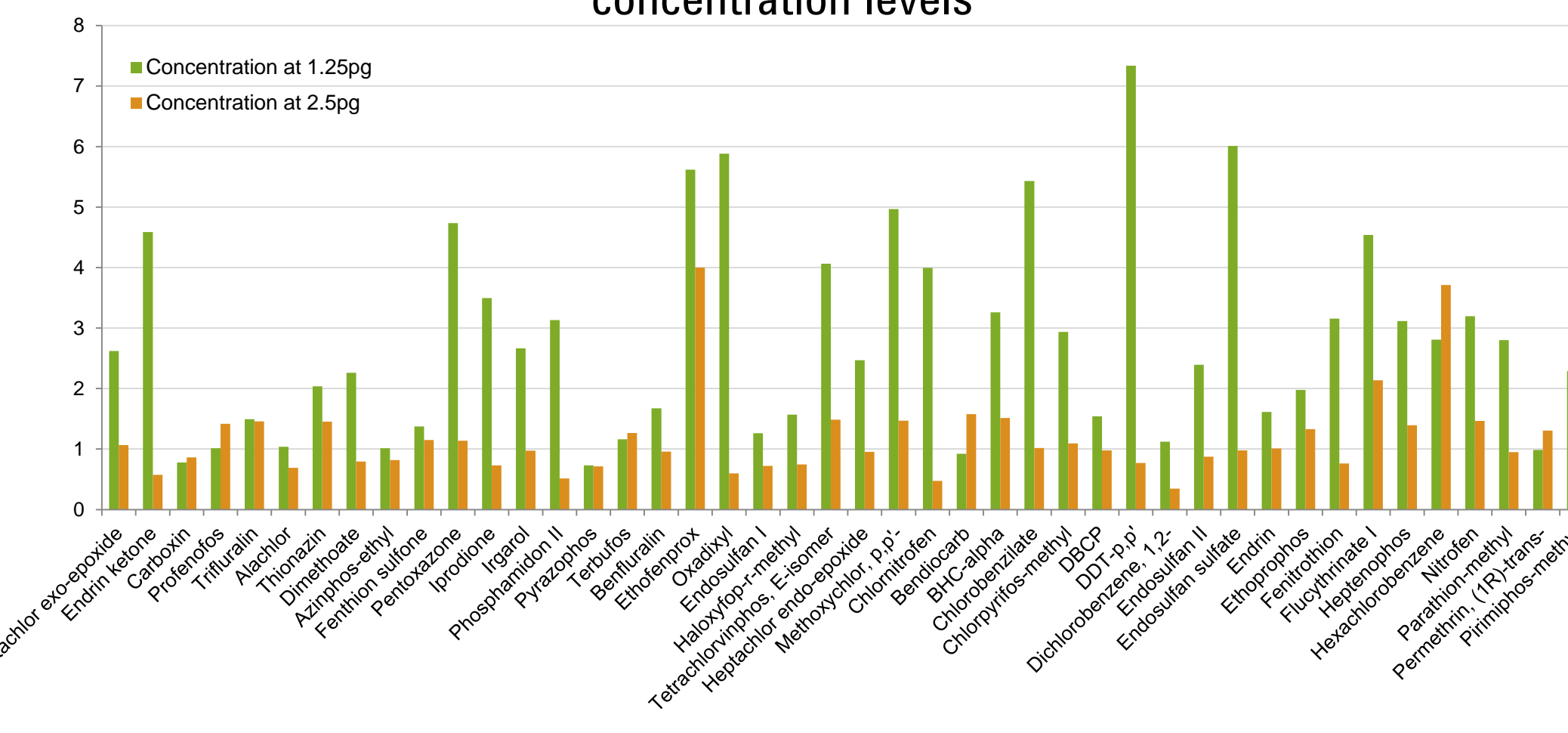

Conclusion

Approximately 170 various pesticides were analyzed in organic honey on the 7010 Series Triple Quadrupole GC/MS using standard and JetClean source in the Acquire and Clean mode, utilizing carefully introduced hydrogen flow. The JetClean control is included in the MassHunter software with easy setup and operation.

The chromatographic peak shape and baseline was improved using the JetClean source particularly for the late eluting compounds. The calibration resulted very comparable R² values by both source, while the MDLs obtained at 1.25 pg level resulted lower values using the JetClean source.

The %RSDs were comparable at higher, 2.5pg/ μL level. The results indicate that the JetClean source meets and exceed the performance delivered by the standard source. Further study is undergoing to identify how the source maintenance period is extended when JetClean is applied compared to the standard source.



Reference

¹Anderson, Kim A., et al. "Modified ion source triple quadrupole mass spectrometer gas chromatograph for polycyclic aromatic hydrocarbon analyses." Journal of Chromatography A 1419 (2015): 89-98.