

CoreFocus Report


No.498

LC-MS Reversed-phase **Shim-pack Series**

Shim-pack™ MAqC-ODS I

Analysis of Catecholamines

Keywords: Plasma, Signaling molecules

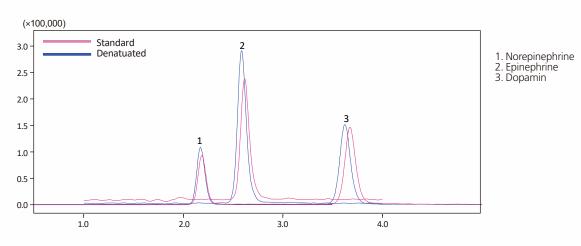

Shim-pack MAqC-ODS I Column

(150 mm L x 2.0 mm l.D., 5 μm), P/N: 228-59936-94

Mobile phase

A) 0.1 % formic acid in water
B) 0.05 % formic acid in methanol

B.Conc. 1% (0 - 0.5 min)→50% (3 min) → 99% (3.1 - 7 min) → 1% (7.1 - 12 min) Time Program

Flow rate 0.2 mL/min

Injection vol. : 5 uL

Post-column Addition : Mobile Phase B at 0.2 mL/min

LC-MS : LCMS-8600 Ionization ESI (positive mode)

Mode MRM Interface voltage Program : +0.6 kV Nebulizing gas flow : 2.2 L/min Drying gas flow 3 L/min Heating gas flow 17 L/min DL temp. 250 °C 400 °C Interface temp.

Source: Application News C114 (JP)

CoreFocus and Shim-pack are trademarks of Shimadzu Corporation or its affiliated companies in Japan and/or other countries.

Shimadzu Corporation www.shimadzu.com/an/

For Research Use Only. Not for use in diagnostic procedures.

This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of

Company names, products/service names and logos used in this publication are trademarks and trade names of Shimadzu Corporation, its subsidiaries or its affiliates, whether or not they are used with trademark symbol "TM" or "®".

Third-party trademarks and trade names may be used in this publication to refer to either the entities or their products/services, whether or not they are used with trademark symbol "TM" or "®".

Shimadzu disclaims any proprietary interest in trademarks and trade names other than its own.

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.