


# Visiprep™ Large Volume Sampler

## Product Specification

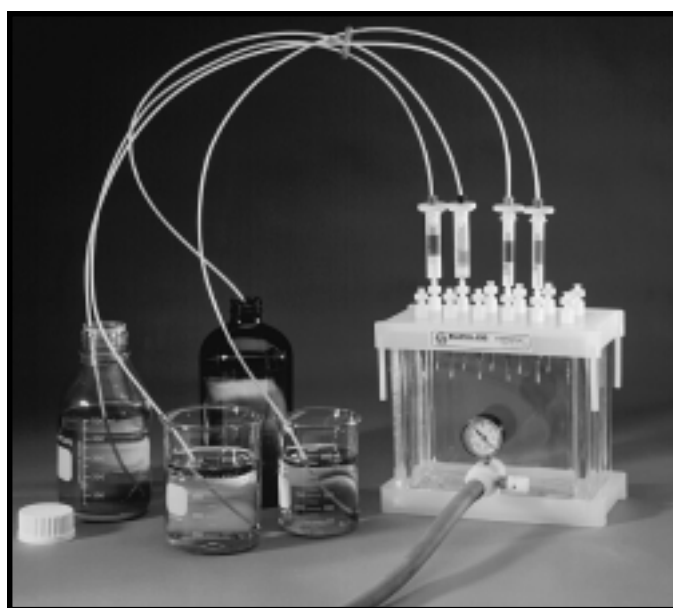

994-0086



995-0147

The Visiprep Large Volume Sampler enables you to extract multiple aqueous samples directly from any laboratory containers, using conventional solid phase extraction (SPE) tubes and a Visiprep SPE vacuum manifold.

Two versions of the large volume sampler are available:

- The Visiprep Sampler (Cat. No. 57275) fits standard 3mL and 6mL polypropylene SPE tubes. It includes four tube adaptors. In order to simultaneously extract 12 samples on a standard 12-port Visiprep manifold, three of these samplers are required.
- The Visiprep+ Sampler (Cat. No. 57272\*) is compatible with 12mL, 20mL, and 60mL SPE tubes. It contains three tube adaptors. For simultaneous extraction of 12 samples, four Visiprep+ samplers are needed.

The adaptors have replaceable screw fittings and 1/8-inch Teflon® tubing. At the free end of each piece of tubing is a reusable stainless steel weight.

To use the sampler, feed the weighted end of each Teflon tube into a sample container until it touches the bottom. Insert the adaptor ends into the conditioned SPE tubes on a Visiprep SPE vacuum manifold. When the vacuum is turned on and the manifold flow control valves opened, the samples are conveniently delivered to the SPE tubes. The flow rate of sample through each tube can be controlled by turning the flow control valves on the manifold. Both the screw fittings and the Teflon tubing are color-coded for easy sample identification.

For additional information about Visiprep SPE Vacuum Manifolds, or Supelclean™ SPE tubes, see page 2 of this publication or refer to the Supelco catalog.

### Ordering Information:

Description	Cat. No.
<b>Large Volume Sampler</b>	
Visiprep Sampler (transfers 4 samples) For 3mL and 6mL SPE tubes	57275
Visiprep+ Sampler (transfers 3 samples) For 12mL, 20mL, and 60mL SPE tubes	57272*
<b>Replacement Tubing, 1/8" Teflon</b>	
Color-coded, pk. of 4	57276
<b>Replacement Nuts and Ferrules</b>	
Color-coded, pk. of 4	57277
<b>Replacement Stainless Steel Weights</b>	
Pk. of 4	57278
<b>Visiprep SPE Vacuum Manifold</b>	
12-port model	57030-U
24-port model	57250-U
<b>Visiprep DL Disposable Liner SPE Vacuum Manifold</b>	
12-port model	57044
24-port model	57265
<b>Long Stem Flow Control Valves</b>	
Adaptable to Cat. No. 5-7030	57048

\*Restrictions for using Cat. No. 57272:

- If you are using a 12-port Visiprep Vacuum Manifold (Cat. No. 57030), and would like to use all 12 ports, you must equip alternate manifold valves with Long Stem Flow Control Valves (Cat. No. 57048).
- If you are using a Visiprep DL Vacuum Manifold or a 24-port manifold, only alternate manifold ports can be used.

## Free SPE Literature

Visiprep Solid Phase Extraction Vacuum Manifolds	494059
Visiprep DL Disposable Liner SPE Manifolds	494060
Supelclean Solid Phase Extraction Tubes	412130

## To Order Technical Literature

Supelco produces a wide variety of technical literature, all of which is free. You may request this literature several different ways:

- Use the business reply card located at the back of our catalog.
- Call our Ordering and Customer Service department (800-247-6628 or 814-359-3441).
- Use our ChromFax™ service.

### How to Use ChromFax

ChromFax is Supelco's automated digital data system, which allows you to instantly obtain technical literature any time of the day or night. Supelco technical bulletins, application notes, lab hints, and product specification sheets can be sent to your Fax machine.

1. To use ChromFax you must dial from a touch-tone phone. If you do not have a touch-tone phone, call our Ordering and Customer Service department and we will fax the publication to you.

2. Have the phone number of your Fax machine ready, and dial **800-652-8069 or 814-359-5748**. Remember you must have the *complete* Fax number (011 + country/city code + Fax number) when you are calling from outside the US or Canada, otherwise the document cannot be transmitted.
3. When ChromFax answers, follow the simple recorded instructions. Select the **ChromFax number** of the document you want. Be certain you order documents by their ChromFax numbers and NOT their literature numbers.
4. Enter your Fax number and hang up. The requested information will be sent to you within minutes. To report a problem with the ChromFax system contact our Technical Service department.

### Technical Literature Available on ChromFax

The listing of documents available on ChromFax is updated every month. We encourage you to request a list of available ChromFax documents each time you use the ChromFax system.

### Trademarks

ChromFax, Supelclean, Visiprep — Sigma-Aldrich Co.  
Teflon — E.I. du Pont de Nemours & Co., Inc.

Contact our Technical Service Department  
(phone 800-359-3041 or 814-359-3041, FAX 814-359-5468)  
for expert answers to your questions.

For more information, or current prices, contact your nearest Supelco subsidiary listed below. To obtain further contact information, visit our website ([www.sigma-aldrich.com](http://www.sigma-aldrich.com)), see the Supelco catalog, or contact Supelco, Bellefonte, PA 16823-0048 USA.

**ARGENTINA** · Sigma-Aldrich de Argentina, S.A. · Buenos Aires 1119 **AUSTRALIA** · Sigma-Aldrich Pty. Ltd. · Castle Hill NSW 2154 **AUSTRIA** · Sigma-Aldrich Handels GmbH · A-1110 Wien  
**BELGIUM** · Sigma-Aldrich N.V./S.A. · B-2880 Bornem **BRAZIL** · Sigma-Aldrich Quimica Brasil Ltda. · 01239-010 São Paulo, SP **CANADA** · Sigma-Aldrich Canada, Ltd. · 2149 Winston Park Dr., Oakville, ON L6H 6J8  
**CZECH REPUBLIC** · Sigma-Aldrich s.r.o. · 186 00 Praha 8 **DENMARK** · Sigma-Aldrich Denmark A/S · DK-2665 Vallensbaek Strand **FINLAND** · Sigma-Aldrich Finland/YA-Kemia Oy · FIN-00700 Helsinki  
**FRANCE** · Sigma-Aldrich Chimie · 38297 Saint-Quentin-Fallavier Cedex **GERMANY** · Sigma-Aldrich Chemie GmbH · D-82041 Deisenhofen **GREECE** · Sigma-Aldrich (o.m.) Ltd. · Ilioupoli 16346, Athens  
**HUNGARY** · Sigma-Aldrich Kft. · H-1067 Budapest **INDIA** · Sigma-Aldrich Co. · Bangalore 560 048 **IRELAND** · Sigma-Aldrich Ireland Ltd. · Dublin 24 **ISRAEL** · Sigma Israel Chemicals Ltd. · Rehovot 76100  
**ITALY** · Sigma-Aldrich s.r.l. · 20151 Milano **JAPAN** · Sigma-Aldrich Japan K.K. · Chuo-ku, Tokyo 103 **KOREA** · Sigma-Aldrich Korea · Seoul **MALAYSIA** · Sigma-Aldrich (M) Sdn. Bhd. · 58200 Kuala Lumpur  
**MEXICO** · Sigma-Aldrich Química S.A. de C.V. · 50200 Toluca **NETHERLANDS** · Sigma-Aldrich Chemie BV · 3330 AA Zwijndrecht **NORWAY** · Sigma-Aldrich Norway · Torshov · N-0401 Oslo  
**POLAND** · Sigma-Aldrich Sp. z o.o. · 61-663 Poznań **PORTUGAL** · Sigma-Aldrich Quimica, S.A. · Sintra 2710 **RUSSIA** · Sigma-Aldrich Russia · Moscow 103062 **SINGAPORE** · Sigma-Aldrich Pte. Ltd.  
**SOUTH AFRICA** · Sigma-Aldrich (pty) Ltd. · Jet Park 1459 **SPAIN** · Sigma-Aldrich Quimica, S.A. · 28100 Alcobendas, Madrid **SWEDEN** · Sigma-Aldrich Sweden AB · 135 70 Stockholm  
**SWITZERLAND** · Supelco · CH-9471 Buchs **UNITED KINGDOM** · Sigma-Aldrich Company Ltd. · Poole, Dorset BH12 4QH  
**UNITED STATES** · Supelco · Supelco Park · Bellefonte, PA 16823-0048 · Phone 800-247-6628 or 814-359-3441 · Fax 800-447-3044 or 814-359-3044 · email: [supelco@sial.com](mailto:supelco@sial.com)