



Mineral oil in soil and water according to DIN EN ISO 9377-2

Fast analysis of lubricating oil contamination according to DIN EN ISO 9377-2

Application Note

Environmental

Authors

Agilent Technologies, Inc.

Introduction


Fast GC analysis of lubricating oil contamination in under ten minutes using an Agilent Select Mineral Oil column according to DIN EN ISO 9377-2.



Agilent Technologies

Conditions

Technique : GC
Column : Agilent Select Mineral Oil, 0.32 mm x 15 m fused silica (optimized film thickness) (Part no. CP7491)
0.53 mm x 6 m, methyl deactivated
Temperature : 55 °C, 1.9 min → 320 °C, 80 °C/min
Carrier Gas : Nitrogen, 80 kPa
Injector : On-column
Detector : FID
Sample Size : 2 µL
Sample Size : Lubrication oil, 1.1 mg/L and 0.11 mg/L in petroleum ether
Courtesy : Thomas Karle, Chemisches Labor; Dr. Vogt, Karlsruhe, Germany

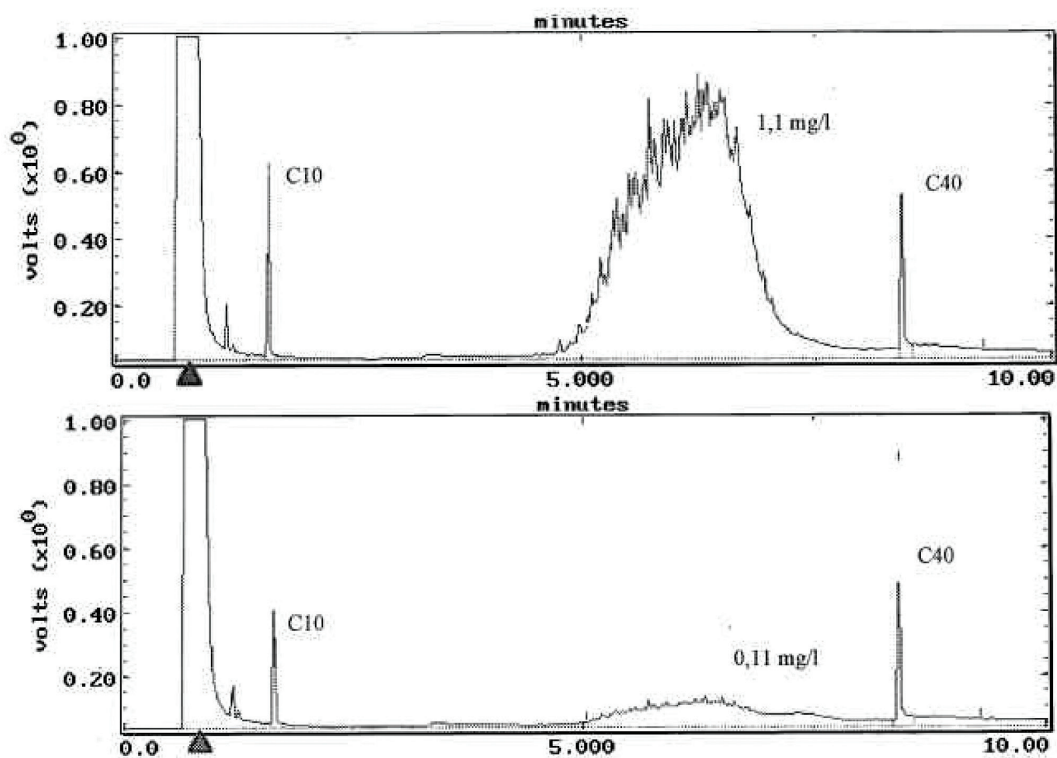

www.agilent.com/chem

This information is subject to change without notice.

© Agilent Technologies, Inc. 2011

Printed in the USA

31 October, 2011

First published prior to 11 May, 2010

A02048



Agilent Technologies