

Analysis of degradation  
products in electrolyte for  
rechargeable lithium-ion  
battery through  
high mass accuracy MS<sup>n</sup> and  
multivariate statistical  
technique

**ASMS 2012** WP26-541

Hiroki Nakajima, Satoshi Yamaki, Tsutomu Nishine,  
Masaru Furuta  
SHIMADZU CORPORATION, Kyoto, Japan

# Analysis of degradation products in electrolyte for rechargeable lithium-ion battery through high mass accuracy MS<sup>n</sup> and multivariate statistical technique

## Introduction

Rechargeable lithium-ion batteries (LiB) are one of the major power sources for portable electronic devices and electric vehicles because of their high voltage and high energy density (Fig. 1-(a)). The electrolyte of a LiB is consisting of a lithium salt in an aprotic organic solvent. The typical operational potential of a LiB is between 0 and 5 V. Therefore, solvent can be reduced or oxidized at the

negative and positive electrodes during the battery charging process. As a result, various degradation products are generated in the electrolyte and cause some problems such as a decrease in the capacitance of battery (Fig. 1-(b)). Here, we present the analysis method of degradation products generated in electrolyte using high mass accuracy MS<sup>n</sup> and multivariate statistical technique.

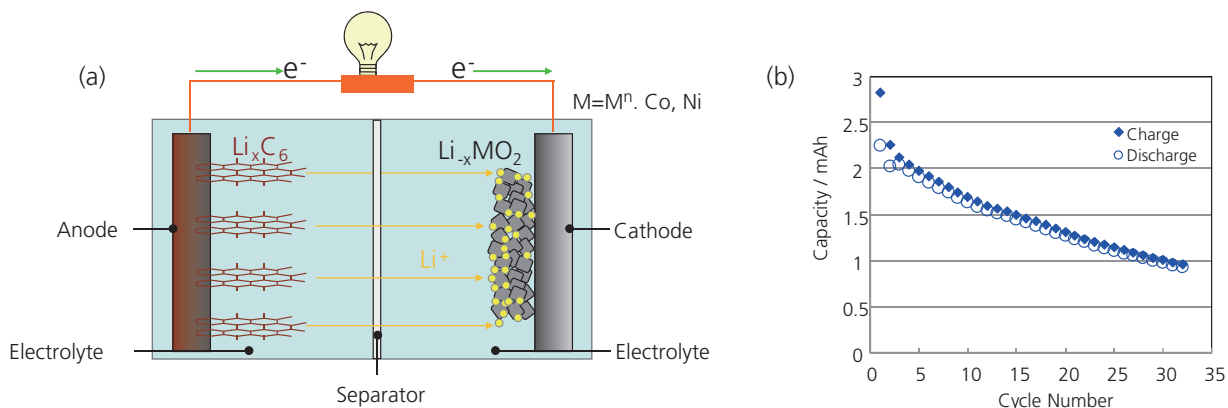


Fig. 1 Rechargeable lithium-ion battery component of lithium-ion battery (a), a decrease in the capacitance of battery (b).

## Experiment

The electrolyte was a mixture of ethylene carbonate (EC) and diethyl carbonate (DEC) (EC : DEC = 1 : 1 vol%) containing 1M lithium hexafluorophosphate (LiPF<sub>6</sub>). The electrolyte A taken from unused lithium-ion battery and the electrolyte B taken from lithium-ion battery repeated charge and discharge cycles (60°C, 30 times) were used as samples. Those samples were prepared 1/10 dilution with methanol for LCMS-IT-TOF (Shimadzu Corporation) measurement. Orthogonal Partial Least Squares


Discriminant Analysis (OPLS-DA) was performed using data acquired by LCMS-IT-TOF measurement of electrolyte A and electrolyte B (n=3) to find the compounds generated in electrolyte B. Then, those compounds were identified chemical formula using software "Formula Predictor" (Shimadzu Corporation). SIMCA-P+ (Umetrics) and Profiling Solution (Shimadzu Corporation) were used for OPLS-DA (Scheme).




1. Acquisition of the high mass accuracy MS<sup>n</sup> data (LCMS-IT-TOF)



2. Peak alignment and generation of peak list (Profiling Solution)

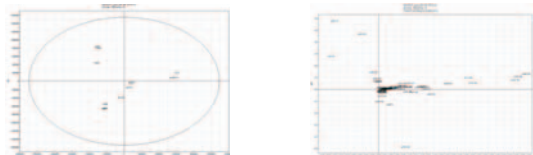

3. Searching of degradation products (SIMCA-P+)

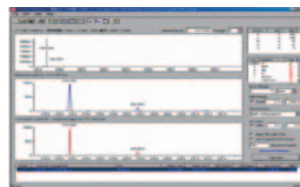

4. Prediction of chemical formula (Formula Predictor)

Structural estimation

Scheme Work flow of the analysis of degradation product in electrolyte for LiB

# Analysis of degradation products in electrolyte for rechargeable lithium-ion battery through high mass accuracy MS<sup>n</sup> and multivariate statistical technique

## Results and discussion

MS data of electrolyte A and electrolyte B were acquired using LCMS-IT-TOF under the analytical conditions shown Table 1. On the score plot of OPLS-DA, the group of electrolyte A and electrolyte B were located at left side and right side, respectively (Figure 2-(a)). 15 unique ions of electrolyte B were observed at right side on S-plot (Figure 2-(b)). And, those ions were not detected on the extracted ion chromatogram (EIC) of electrolyte A (Fig. 3). These results suggested that those ions were degradation products generated in the electrolyte of lithium-ion battery repeated charge and discharge 30 cycles.

Table 1 LCMS analytical conditions

Column	: Shim-pack FC-ODS (2.0 mmI.D.x150 mm, 3 mm)
Flow rate	: water
Column temp.	: 0.2 mL/min
Mobile phaseA	: 40°C
Mobile phaseB	: methanol
Time prog.	: 5%B (0 min) → 55%B (30 min) → 5%B (30.01 min)
Injection volume	: 1 μL
Ionization mode	: ESI(+)
Probe voltage	: 4.5 kV
CDL temperature	: 200°C
BH temperature	: 200°C
Nebulizing gas	: 1.5 L/min
Drying gas	: 0.1 MPa
Scan range	: m/z 80 - 1000

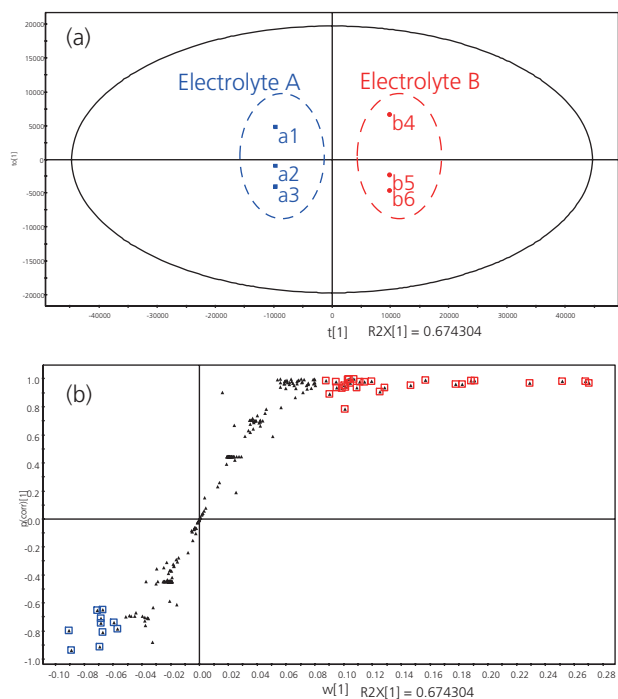


Fig. 2 The result of OPLS-DA, score plot (a), S-plot (b)

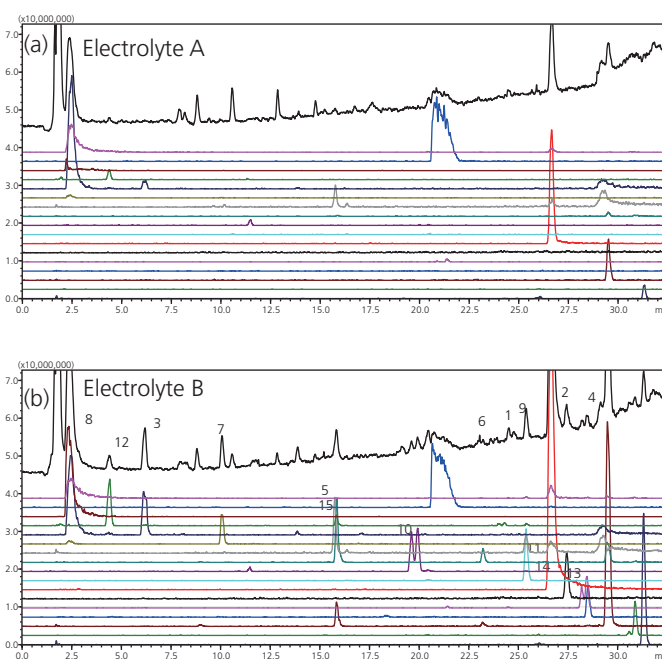


Fig. 3 EICs of ions detected in Electrolyte B (b) These ions were not detected on EIC of electrolyte A (a).

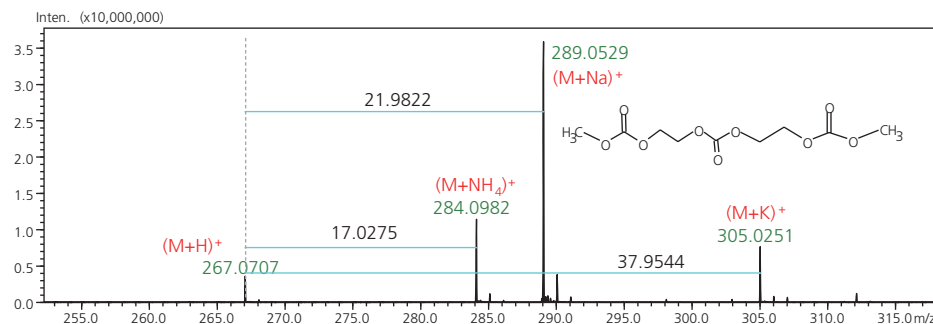


Fig. 4 MS data and predicted structure of peak number 2 (m/z 284.0982) being one of 15 unique ions of electrolyte B

# Analysis of degradation products in electrolyte for rechargeable lithium-ion battery through high mass accuracy MS<sup>n</sup> and multivariate statistical technique

The formula of peak number 2 (*m/z* 284.0982) being one of 15 unique ions of electrolyte B was predicted as C<sub>9</sub>H<sub>14</sub>O<sub>9</sub> (polycarbonate) using high mass accuracy MS data and formula predictor. Indeed, the structure of C<sub>9</sub>H<sub>14</sub>O<sub>9</sub> was predicted as H<sub>3</sub>C-(OCO<sub>2</sub>-C<sub>2</sub>H<sub>4</sub>)<sub>2</sub>-OCO-CH<sub>3</sub> referring to some articles on degradation products in electrolyte. MS<sup>n</sup>

of the ion (*m/z* 284.0982) also was measured to determine the validity of the predicted chemical structure. Each product ions and neutral loss in MS<sup>2</sup> and MS<sup>3</sup> data showed that the predicted structure of C<sub>9</sub>H<sub>14</sub>O<sub>9</sub> was correct.

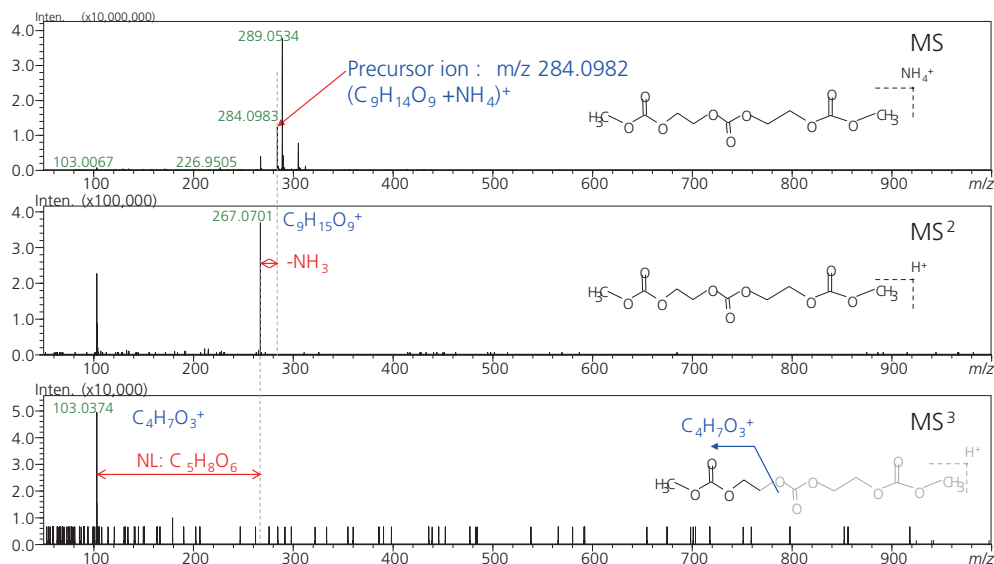


Fig. 5 MS<sup>n</sup> data of peak number 2 (*m/z* 284.0982) detected in electrolyte B

By the same method, formula of peak number 13 (*m/z* 283.0336) was identified as C<sub>3</sub>H<sub>7</sub>O<sub>4</sub>P. Structure of it was determined as phosphate shown in Fig. 6.

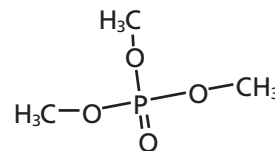


Fig. 6 of Structure peak number 13

## Conclusion

- It was clear that the chemical species of degradation products generated in the electrolyte with increasing charge and discharge cycles were carbonate and phosphate from result of this study.
- The formulae of 15 degradation products detected in Electrolyte B were identified as below.

Peak No.	<i>m/z</i>	R.T.(min)	Ion species	M.W	Predicted formula	Mass accuracy (ppm)
1	229.0678	26.645	(M+Na) <sup>+</sup>	206	C <sub>8</sub> H <sub>14</sub> O <sub>6</sub>	-2.62
2	284.0982	20.401	(M+NH <sub>4</sub> ) <sup>+</sup>	294	C <sub>9</sub> H <sub>18</sub> O <sub>9</sub>	-0.35
3	295.1032	29.482	(M+H) <sup>+</sup>	170	C <sub>11</sub> H <sub>18</sub> O <sub>9</sub>	+1.69
4	177.0512	6.178	(M+Li) <sup>+</sup>	382	C <sub>4</sub> H <sub>11</sub> O <sub>5</sub> P	+2.92
5	400.1458	31.294	(M+NH <sub>4</sub> ) <sup>+</sup>	272	C <sub>14</sub> H <sub>22</sub> O <sub>12</sub>	+2.50
6	295.056	15.833	(M+Na) <sup>+</sup>	244	C <sub>8</sub> H <sub>17</sub> O <sub>8</sub> P	-1.36
7	262.0853	25.374	(M+NH <sub>4</sub> ) <sup>+</sup>	184	C <sub>10</sub> H <sub>13</sub> O <sub>5</sub> P	+1.53
8	185.0577	10.054	(M+H) <sup>+</sup>	250	C <sub>5</sub> H <sub>13</sub> O <sub>5</sub> P	+4.32
9	251.1111	27.416	(M+H) <sup>+</sup>	268	C <sub>8</sub> H <sub>17</sub> O <sub>5</sub> F <sub>3</sub>	+2.39
10	269.0162	19.879	(M+H) <sup>+</sup>	332	C <sub>7</sub> H <sub>7</sub> O <sub>2</sub> F <sub>6</sub> P	-2.60
11	350.1003	28.426	(M+NH <sub>4</sub> ) <sup>+</sup>	138	C <sub>8</sub> H <sub>17</sub> O <sub>7</sub> F <sub>4</sub> P	+1.14
12	283.0336	4.385	(2M+Li) <sup>+</sup>	358	C <sub>3</sub> H <sub>7</sub> O <sub>4</sub> P	+4.32
13	381.0938	30.857	(M+Na) <sup>+</sup>	270	C <sub>12</sub> H <sub>23</sub> O <sub>10</sub> P	-1.57
14	293.0777	28.487	(M+Na) <sup>+</sup>	222	C <sub>9</sub> H <sub>19</sub> O <sub>7</sub> P	+1.36
15	245.0641	15.713	(M+Na) <sup>+</sup>	266	C <sub>8</sub> H <sub>14</sub> O <sub>7</sub>	-0.41



Shimadzu Corporation  
[www.shimadzu.com/an/](http://www.shimadzu.com/an/)

For Research Use Only. Not for use in diagnostic procedures.  
The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.