

Application Data Sheet

No.2

GCMS

Gas Chromatograph Mass Spectrometer

Analysis of Amino Acids Contained in Dashi (Japanese Soup Stock)

Amino acids contained in dashi were treated with EZ:faastTM (Phenomenex, Inc.), which enables easy pretreatment, and then analyzed with a GC-MS system.

Experiment

Pretreatment

Two types of dashi (katsuobushi extract (from sliced dried bonito) and kombu (from kelp of the genus Laminaria) were treated using EZ:faast. Norvaline was added as an internal standard.

Instrument

A GCMS-QP2010 Ultra (with high-power oven) was used for the measurements. The analysis conditions, shown in Table 1, were in conformity with the "Amino Acid Analysis Methods" in the "GC/MS Metabolic Components Database."

Table 1: Analysis Conditions (GC/MS Metabolic Components Database: Amino Acid Analysis Methods)

GC-MS	: GCMS-QP2010 Ultra (with high-power oven)	[MS]	
Column	: ZB-AAA (10 mL. × 0.25 mm I.D.) (Phenomenex, Inc.)	Interface temperature	: 280°C
[GC]		Ion source temperature	: 200°C
Injection quantity	: 1 μL	Solvent elution time	: 0.4 min
Vaporization chamber temperature	: 280°C	Data sampling time	: 0.5 min to 7 min
Column oven temperature	: 110°C → (30 °C /min)→ 320°C	Measurement mode	: Scan
Control mode	: Constant pressure (15 kPa)	Mass range	: <i>m/z</i> 45-450 (3,333u/sec)
Injection mode	: Split	Event time	: 0.15 sec
Split ratio	: 15		
Carrier gas	: Helium		

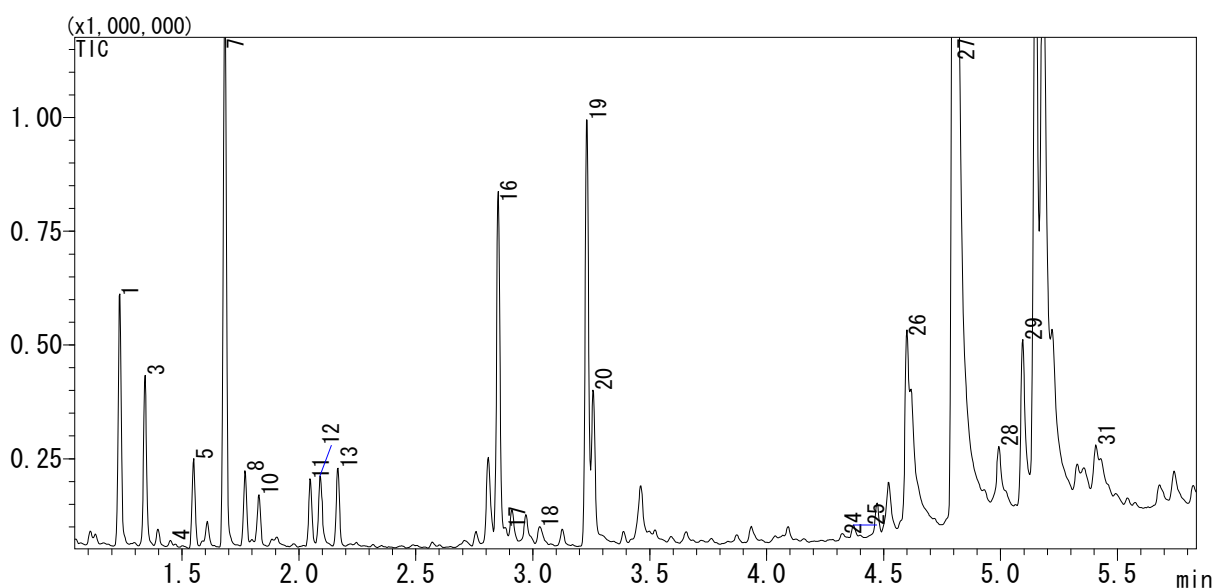


Fig. 1: Total Ion Current Chromatogram for Amino Acid Derivatives in Katsuobushi Extract

The numbers for each component follow the serial numbers in the "GC/MS Metabolic Components Database."

1 Alanine	8 Leucine	16 Asparatic acid	24 Ornithine	28 Hydroxylysine
3 Glycine	10 Isoleucine	17 Methionine	25 Glycine-proline	(2 isomers)
4 alpha-aminobutyric acid	11 Threonine	18 4-Hydroxyproline	(dipeptide)	29 Tyrosine
5 Valine	12 Serine	19 Glutamic acid	26 Lysine	31 Tryptophan
7 Norvaline (I.S.)	13 Proline	20 Phenylalanine	27 Histidine	

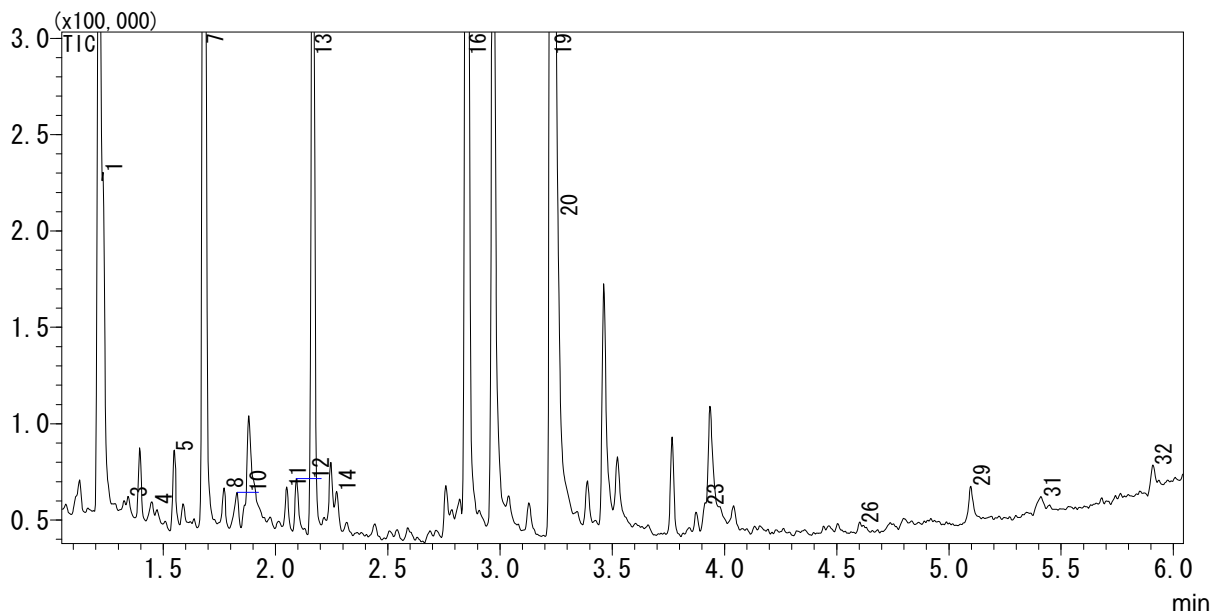


Fig. 2: Total Ion Current Chromatogram for Amino Acid Derivatives in Kombu Extract
The numbers for each component follow the serial numbers in the "GC/MS Metabolic Components Database."

1 Alanine	8 Leucine	14 Asparagine	26 Lysine
3 Glycine	10 Isoleucine	16 Aspartic acid	29 Tyrosine
4 alpha-aminobutyric acid	11 Threonine	19 Glutamic acid	31 Tryptophan
5 Valine	12 Serine	20 Phenylalanine	32 Cystathionine
7 Norvaline(I.S.)	13 Proline	23 Glutamine	

Summary

Pretreatment using the EZ:faast kit, followed by analysis using the GCMS-QP2010 Ultra, which is equipped with a high-speed scanning function, enabled rapid analysis of amino acids. With this combination, it took only 15 minutes per sample from pretreatment to analysis.

(Reference: Shimadzu Application News No. M246 Analysis of Amino Acids Using Fast-GC/MS and Metabolite Database)

For Research Use Only. Not for use in diagnostic procedures.

Shimadzu Corporation ("Shimadzu") reserves all rights including copyright in this publication. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to, or arising out of the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.

First Edition: September 2011



SHIMADZU CORPORATION. International Marketing Division

3, Kanda-Nishikicho 1-chome, Chiyoda-ku, Tokyo 101-8448, Japan

Phone: 81(3)3219-5641 Fax: 81(3)3219-5710

URL <http://www.shimadzu.com>