


Overview

- Current analytical methods for evaluating **volatile and semi volatile organic compounds (VOCs & SVOCs)** in water often require large sample volumes (1 L or more), are labor intensive, and rely upon liquid-liquid extraction into organic solvents.
- Here we present a sensitive, solvent-free method – termed **vacuum assisted sorbent extraction (VASE)** – for extraction and pre-concentration of VOCs and SVOCs in preparation for thermal desorption-GC-MS.
- In VASE, the sample (often 1 mL or less for water) is evacuated in the presence of a **headspace sorbent pen (HSP)**, which is a vacuum-tight cartridge containing sorbent. In combination with optional heat and agitation, the static reduced pressure environment promotes compounds into the headspace for adherence to an internal sorbent bed.
- Here we demonstrate the application of VASE for extraction and quantitation of several **water pollutants routinely analyzed by EPA method 8270**. Compounds with boiling points ranging from 80°C to over 550°C are examined, including 2-6-ring polycyclic aromatic hydrocarbons, phenols, pesticides, chlorinated hydrocarbons, and disinfection by-products.


Sorbent Pens & Vacuum Assisted Sorbent Extraction (VASE)

Headspace Sorbent Pen

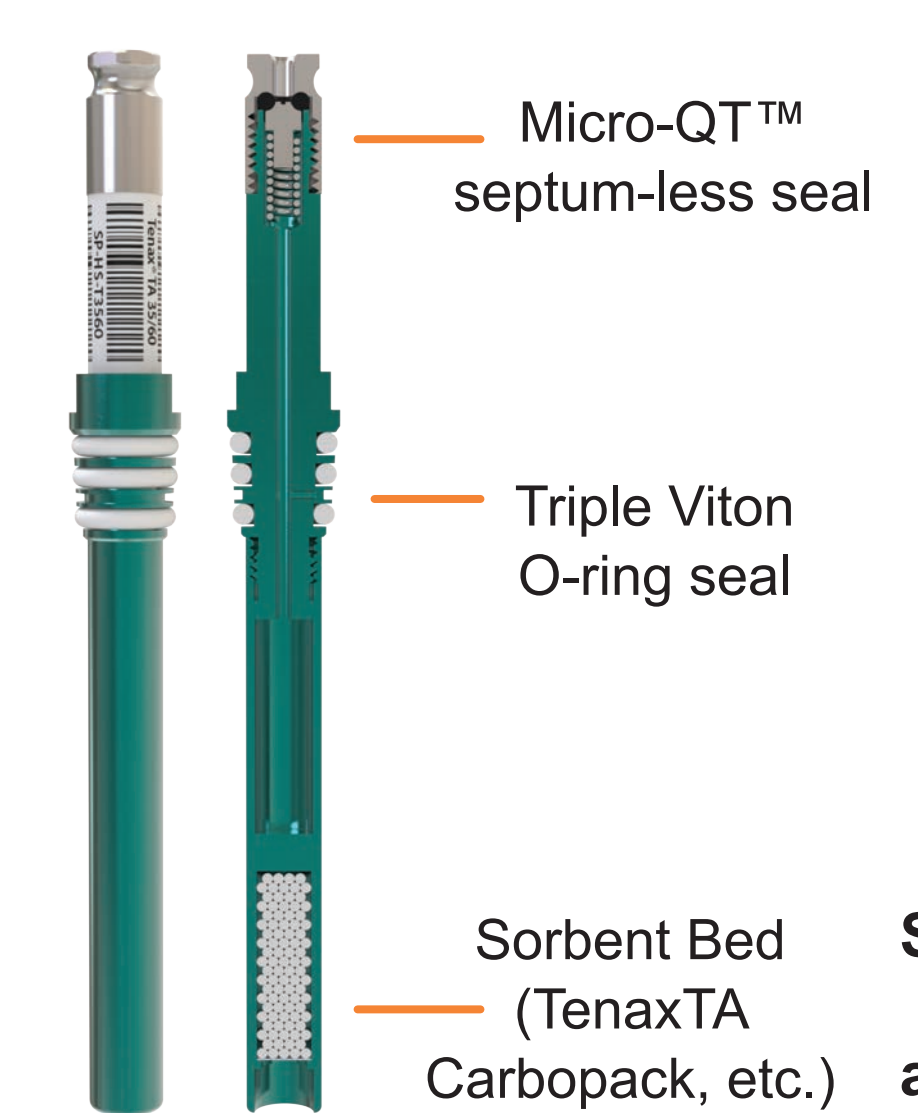


Polycyclic Aromatic Hydrocarbons (PAHs)

Composite Overlay of a 10 ppb PAH Standard Mixture (TRACE Analysis)

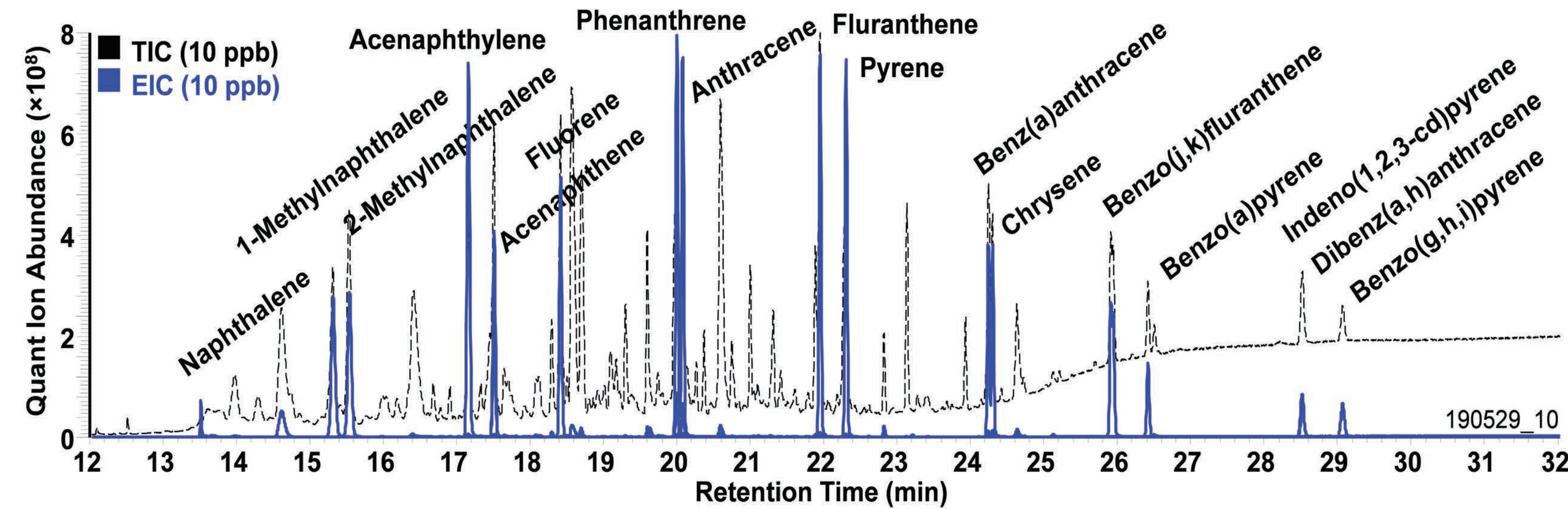


Overlay of 5 Replicate 10 ppb Desorptions (TRACE Analysis)

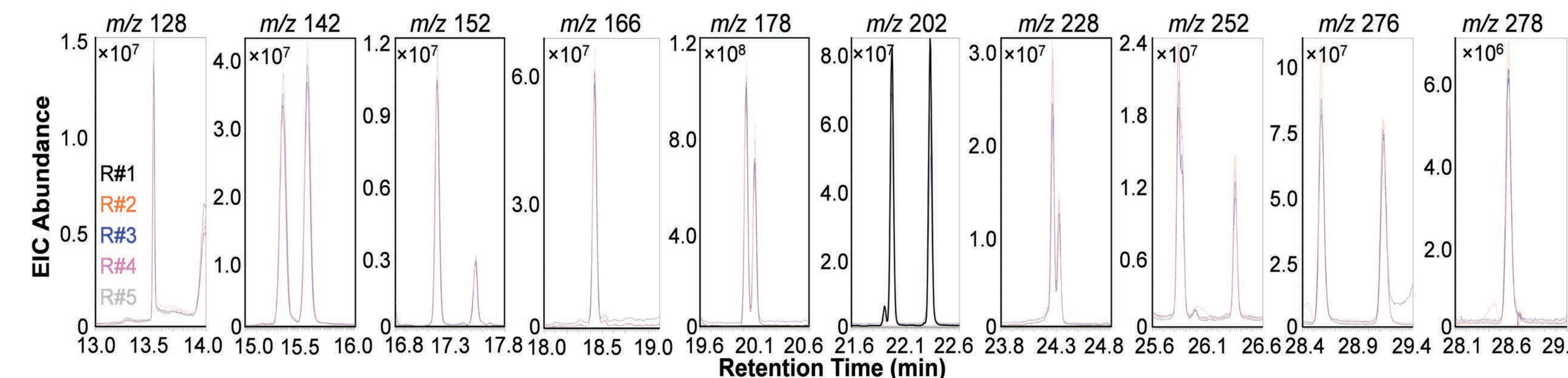

Analyte	BP (°C)	Q Ion	EPA LOQ* (ng/L)	Linear Range (ng/L) TRACE Analysis	Linear Range (ng/L) Split Analysis	R ²	RSD (Raw)	RSD (IS Normalized)
Naphthalene	218	128	10,000	10 – 10,000	500 – 100,000	0.998	4.4%	1.9%
1-Methylnaphthalene	240	142	NA	5 – 10,000	500 – 100,000	0.996	5.1%	1.8%
2-Methylnaphthalene	241	142	NA	5 – 10,000	500 – 100,000	0.997	6.8%	1.4%
Acenaphthylene	280	152	10,000	5 – 10,000	50 – 100,000	0.998	5.6%	0.8%
Acenaphthene	279	154	10,000	50 – 10,000	50 – 100,000	0.999	5.6%	2.1%
Fluorene	298	166	10,000	5 – 10,000	50 – 100,000	0.999	6.7%	2.9%
Phenanthrene	336	178	10,000	50 – 10,000	500 – 100,000	0.999	4.1%	6.1%
Anthracene	340	178	10,000	5 – 10,000	50 – 100,000	0.999	6.8%	1.1%
Fluranthene	384	202	10,000	10 – 10,000	500 – 50,000	0.985	8.9%	3.8%
Pyrene	384	202	10,000	100 – 10,000	500 – 50,000	0.999	15%	6.8%
Benzo(a)Anthracene	437	228	10,000	10 – 10,000	50 – 100,000	0.985	10%	3.0%
Chrysene	448	228	10,000	10 – 10,000	50 – 100,000	0.993	9.7%	1.3%
Benzo(k)fluoranthene	481	252	10,000	50 – 5,000	50 – 100,000	0.998	11%	5.7%
Benzo(a)Pyrene	495	252	10,000	50 – 5,000	50 – 100,000	0.997	12%	6.7%
Indeno(1,2,3-cd)Pyrene	539	276	10,000	5 – 5,000	50 – 100,000	0.991	9.9%	6.6%
Dibenz(a,h)anthracene	524	278	10,000	5 – 5,000	50 – 100,000	0.989	9.3%	8.1%
Benzo(g,h,i)Perylene	550	276	10,000	5 – 5,000	50 – 100,000	0.989	11%	7.9%

*Data from Table 2 Example lower limits of quantitation for semivolatile organics, EPA Method 8270D, Revision 4, February 2007.

Clean HSP After Desorption Enables Repeat Sampling Without Additional Cleanup Step

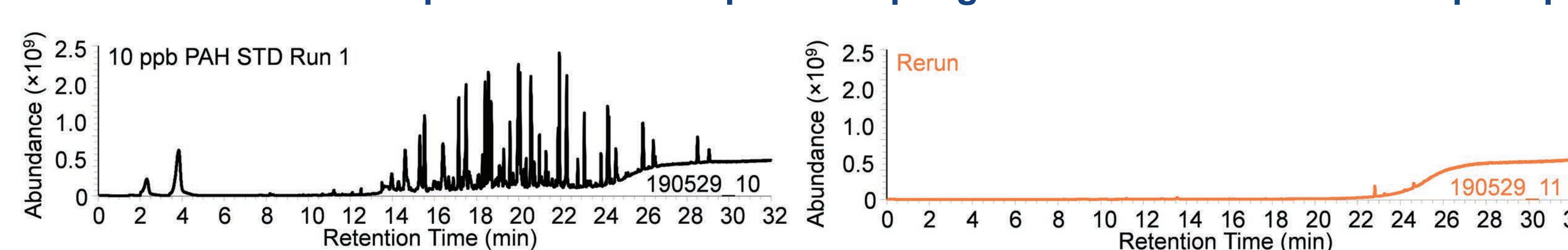

Phenols


Analyte	BP (°C)	Q Ion	EPA LOQ (µg/L)*	Linear Range (µg/L)	R ^{2**}
Phenol	182	94	10	1-50	0.997
2-Chlorophenol	175	128	10	1-50	0.998
o-Cresol	191	108	NA	1-50	0.994
m/p-Cresol	202	107	NA	1-50	1.000
2,4-Dimethylphenol	212	122	10	1-50	0.997
2,4-Dichlorophenol	210	162	10	1-50	0.995
2,6-Dichlorophenol	220	162	10	1-50	1.000
4-Chloro-3-methylphenol	235	107	20	1-50	0.995
2,4,6-Trichlorophenol	246	196	10	1-50	0.997
2,4,5-Trichlorophenol	253	196	10	1-50	1.000
2,3,4,6-Tetrachlorophenol	164	232	10	1-50	0.998
Pentachlorophenol	310	266	50	1-50	0.989
Dinoseb	332	211	20	1-50	0.994



*Data from Table 2 Example lower limits of quantitation for semivolatile organics, EPA Method 8270D, Revision 4, February 2007.

**R² refers to average R² value of n=3, 3-point calibration curves.

VASE, TD, and GC-MS Conditions

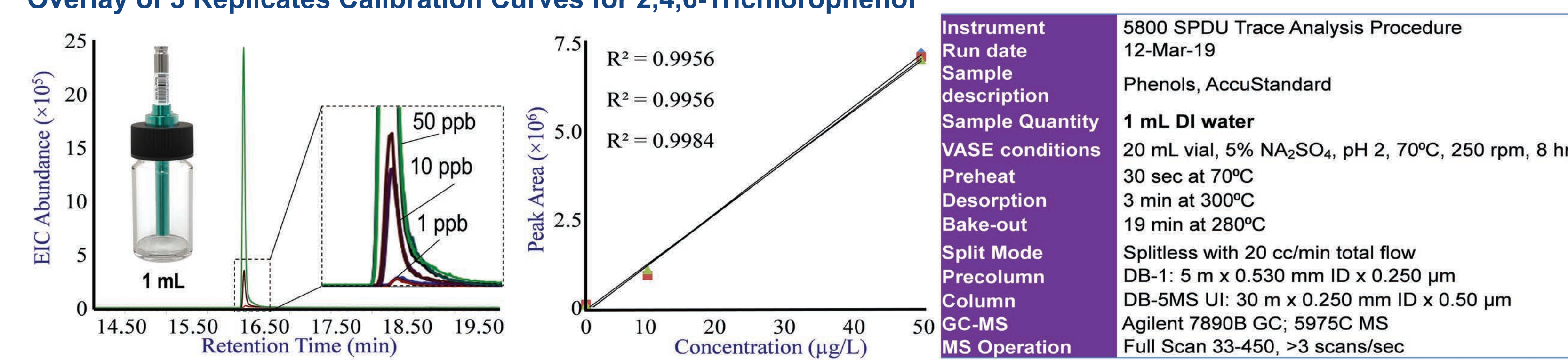


Instrument	5800 SPDU Trace Analysis Procedure
Run date	12-Mar-19
Sample description	Phenols, AccuStandard
Sample quantity	1 mL DI water
VASE conditions	20 mL vial, 5% Na ₂ SO ₄ , pH 2, 70°C, 250 rpm, 8 hr
Preheat	30 sec at 70°C
Desorption	3 min at 300°C
Bake-out	19 min at 280°C
Split Mode	Splitless with 20 cc/min total flow
Precolumn	DB-1: 5 m x 0.530 mm ID x 0.250 µm
Column	DB-5MS UI: 30 m x 0.250 mm ID x 0.50 µm
GC-MS	Agilent 7890B GC; 5975C MS
MS Operation	Full Scan 33-450, >3 scans/sec


Pesticides

Analyte	BP (°C)	Q Ion	Linear Range (µg/L)	R ^{2**}
Aldrin	145	263	0.5-10	0.998
4,4'-DDD	193	165	0.5-10	0.993
a-BHC	288	219	0.5-10	0.999
b-BHC	288	109	0.5-10	0.999
4, 4'-DDE	336	318	0.5-10	0.998
Dieldrin	385	79	0.5-10	1.000
Heptachlor	NA	272	0.5-10	0.994
Heptachlor Epoxide	NA	353	0.5-10	0.999

*R² refers to R² value of n=1, 5-point calibration curves.

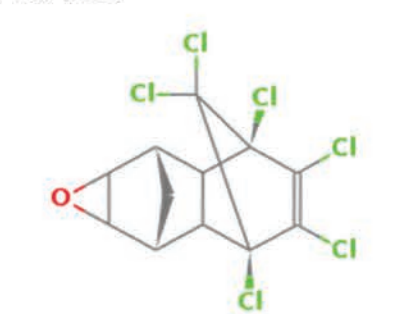


VASE, TD, and GC-MS Conditions

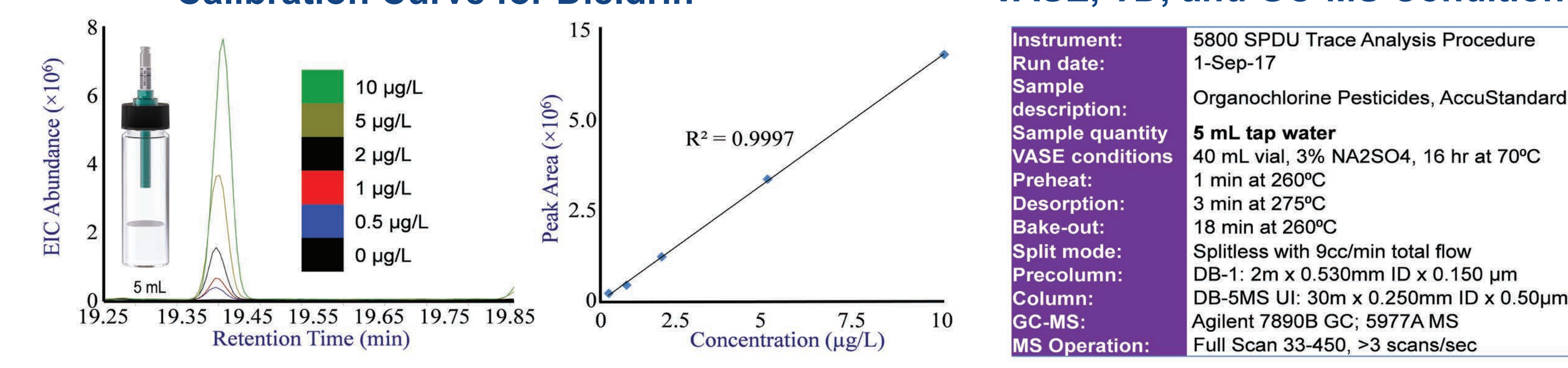

Instrument:	5800 SPDU Trace Analysis Procedure
Run date:	1-Sep-17
Sample description:	Organochlorine Pesticides, AccuStandard
Sample quantity	5 mL tap water
VASE conditions	40 mL vial, 3% Na ₂ SO ₄ , 16 hr at 70°C
Preheat:	1 min at 260°C
Desorption:	3 min at 300°C
Bake-out:	18 min at 260°C
Split mode:	Splitless with 9cc/min total flow
Precolumn:	DB-1: 2m x 0.530mm ID x 0.150 µm
Column:	DB-5MS UI: 30m x 0.250mm ID x 0.50µm
GC-MS:	Agilent 7890B GC; 5977A MS
MS Operation:	Full Scan 33-450, >3 scans/sec

Chlorinated Hydrocarbons


Analyte	BP (°C)	Q Ion	EPA LOQ (µg/L)*	Linear Range (µg/L)	R ^{2**}
Pentachloroethane	162	167	10	1-50	0.998
1,2-Dichlorobenzene	180	146	10	1-50	0.999
1,3-Dichlorobenzene	173	146	10	1-50	0.999
1,4-Dichlorobenzene	174	146	10	1-50	0.999
Hexachloroethane	187	117	10	1-50	0.999
1,2,4-Trichloroethane	214	180	10	1-50	0.999
Hexachloropropene	210	213	10	1-50	0.997
Hexachlorobutadiene	215	225	10	1-50	0.999
1,2,4,5-Tetrachlorobenzene	245	216	10	1-50	0.999
Hexachlorocyclopentadiene	239	237	10	1-50	0.990
2-Chloronaphthalene	255	127	10	1-50	0.999
Pentachlorobenzene	277	248	10	1-50	0.999
Hexachlorobenzene	322	286	10	1-50	0.999

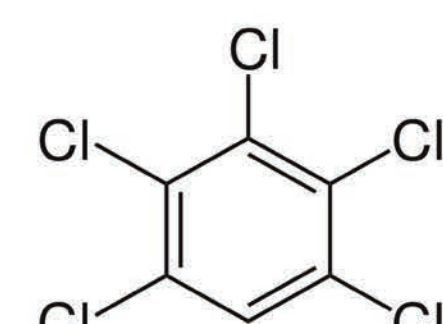

*Data from Table 2 Example lower limits of quantitation for semivolatile organics, EPA Method 8270D, Revision 4, February 2007.


**R² refers to average R² value of n=3, 3-point calibration curves.

Overlay of 3 Replicates Calibration Curves for Pentachlorobenzene

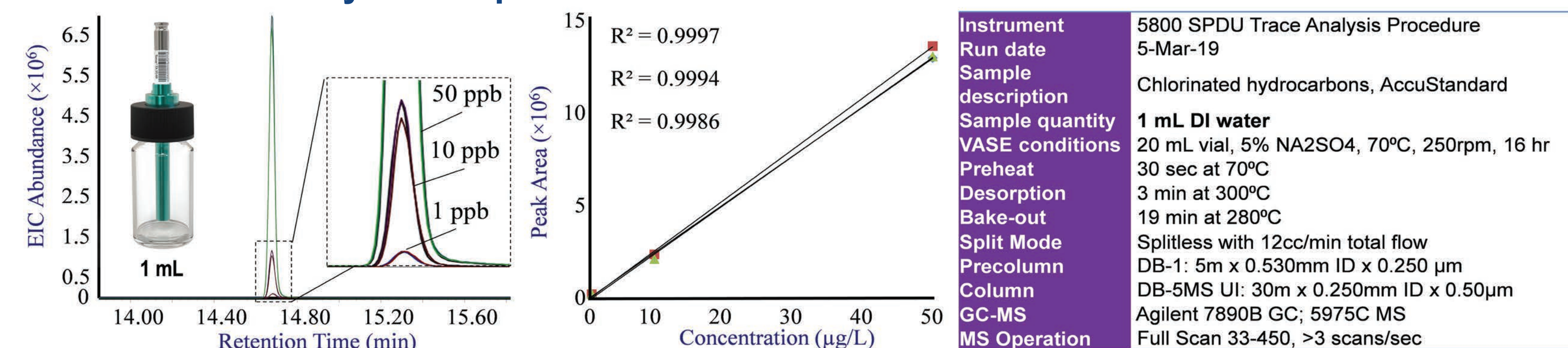


Instrument	5800 SPDU Trace Analysis Procedure
Run date	5-Mar-19
Sample description	Chlorinated hydrocarbons, AccuStandard
Sample quantity	1 mL DI water
VASE conditions	20 mL vial, 5% Na ₂ SO ₄ , 70°C, 250rpm, 16 hr
Preheat	30 sec at 70°C
Desorption	3 min at 300°C
Bake-out	19 min at 280°C
Split Mode	Splitless with 12cc/min total flow
Precolumn	DB-1: 5 m x 0.530 mm ID x 0.250 µm
Column	DB-5MS UI: 30 m x 0.250 mm ID x 0.50 µm
GC-MS	Agilent 7890B GC; 5975C MS
MS Operation	Full Scan 33-450, >3 scans/sec


Disinfectant By-products (DBPs)

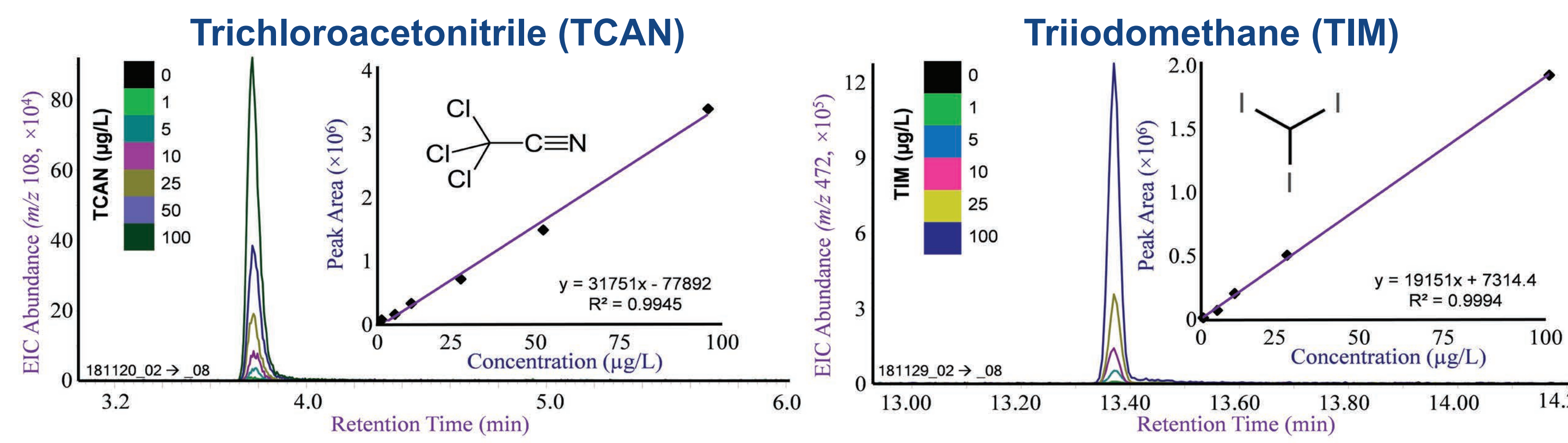

PFBHA Derivatized Dichloroacetaldehyde (DCAL)

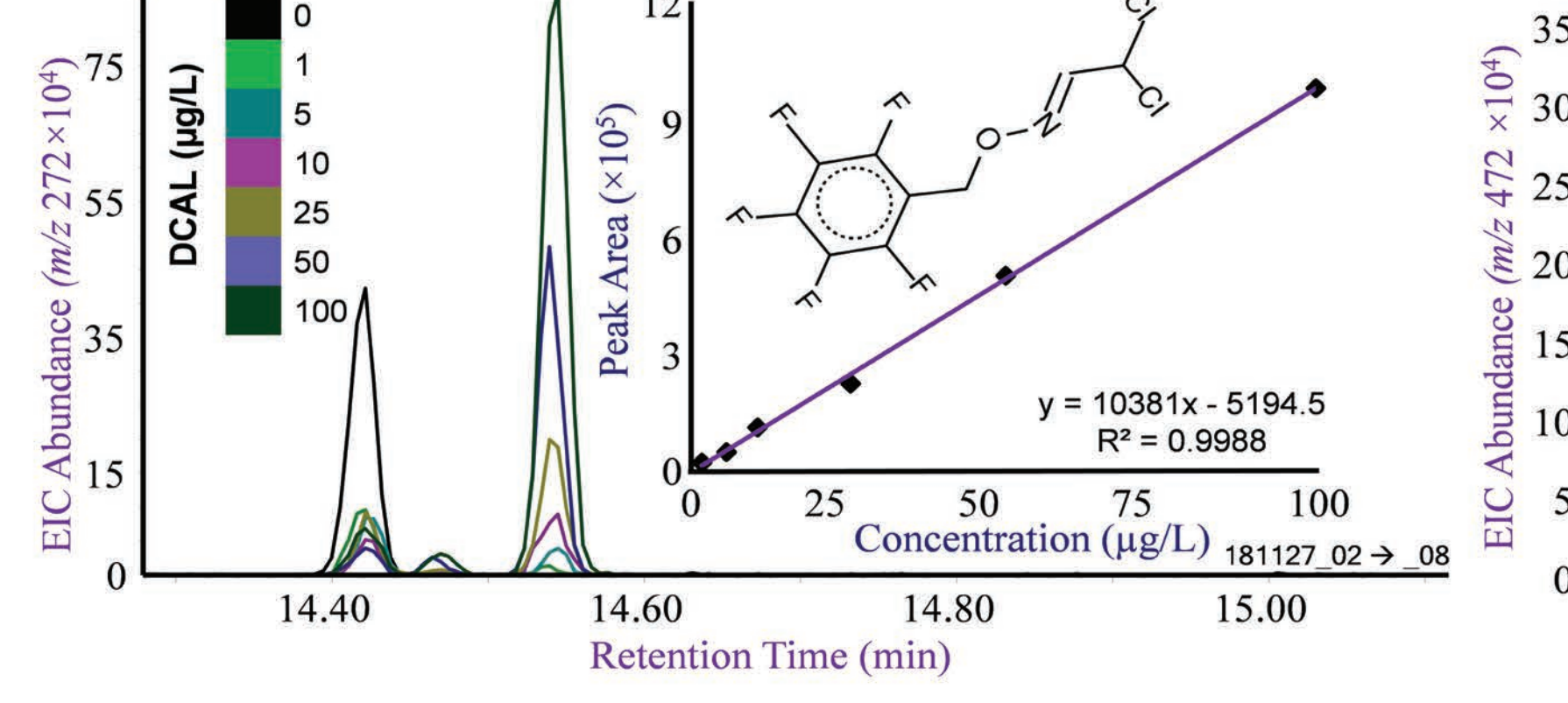

2,4,6-Triiodophenol (TIPh)

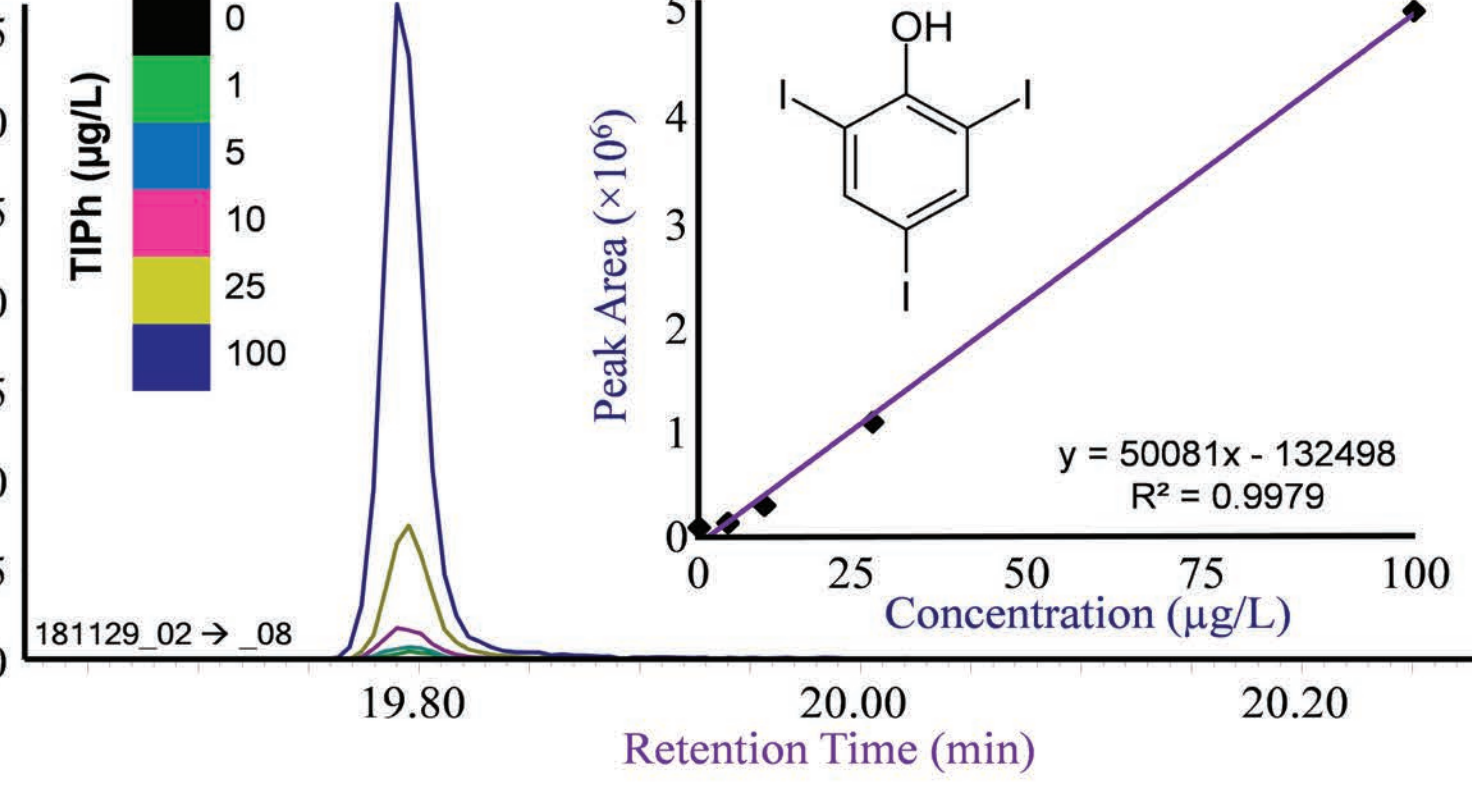

Conclusions

- The critical steps required for capturing volatile and semivolatile compounds from water via a reduced-pressure static headspace extraction technique – vacuum assisted sorbent extraction (VASE) – are shown.
- Example applications include 2 – 6-ring polycyclic aromatic hydrocarbons, phenols, pesticides, chlorinated hydrocarbons, and four classes of disinfection by-products.
- Using 1 mL of water adulterated with PAHs, LODs as low as 5 ng/L and RSDs better than 10% are obtained, demonstrating the low detection limits and exceptional repeatability of VASE.

References

U.S. EPA. 2014. "Method 8270E (SW-846): Semivolatile Organic Compounds by Gas Chromatography/ Mass Spectrometry (GC/MS)," Washington, DC.

Acknowledgments

The authors express their gratitude to the entire Entech staff, including the excellent engineers, machinists, software developers, assemblers, and graphic designers. We also thank Prof. Susan Richardson and her fantastic group of researchers for their help with the disinfectant byproducts work.