



Separation of Permanent Gases on a Liquid Phase

Separation of 5 permanent gases on a WCOT column with a liquid phase with high retention

Application Note

Authors

Rick Hamerlinck and Norbert Reuter
Agilent Technologies, Inc.

Introduction


Normally permanent gases are separated by PLOT (porous layer open tubular) columns with their high retentive phases. With WCOT (wall coated open tubular) columns sub-ambient temperatures are normally necessary. Thick films, like the 8 μm film thickness of the Agilent J&W Select CP-Sil 5CB for Formaldehyde, allow the use of high-inert liquid phases for the (pre-) separation of the standard permanent gases from carbon dioxide for possible column switching at normal ambient temperatures.



Agilent Technologies

Materials and Methods

Technique:	GC-Capillary Medium Bore	Carrier Gas:	Helium at 25 psi (170 kPa)	Detector:	Thermal Conductivity Detector at 220 °C (Filament Temp. 280 °C)
Instrument:	GC Gas Chromatograph	Temp Program:	35 °C isothermal	Sample:	All Gases 1% in Helium
Column:	CP-Sil 5 CB for Formaldehyde, 0.32 mm x 60 m, df=8 µm (part number CP7475)	Injector:	Split/Splitless-Injector (1177) at 250 °C		
		Inj Volume:	500 µL (split ratio 1:20)		

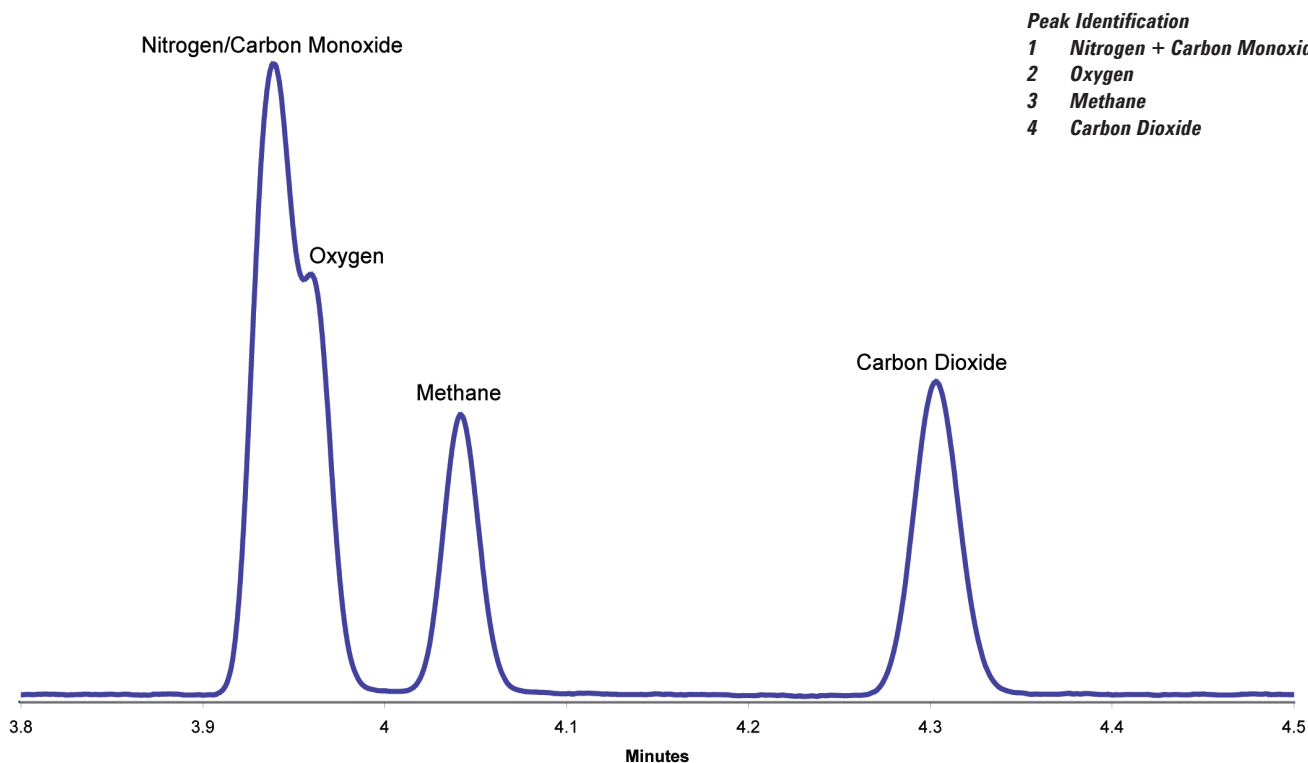

Analysis of permanent gases at 35°C

www.agilent.com/chem

This information is subject to change without notice.

© Agilent Technologies, Inc. 2010

Published in UK, October 08, 2010

SI-02166



Agilent Technologies