

Intelligent Peak and Spectrum Deconvolution Using Photodiode Array Detector

PITTCON2021

P-126

Yoshiyuki Watabe, Kosuke Nakajima, Davide Vecchiatti, Shinnosuke Horie
Shimadzu Corporation, Kyoto, Japan

Intelligent Peak and Spectrum Deconvolution Using Photodiode Array Detector

1. INTRODUCTION

A novel data analysis technique named i-PDeA II (Intelligent Peak Deconvolution Analysis II) has been developed for extracting two or more target peaks from unseparated peaks using data from three-dimensional photodiode array (PDA) detector with a multivariate curve resolution-alternating least squares (MCR-ALS). It also employs an expectation maximization (EM) algorithm with a bidirectional exponentially modified Gaussian (bemg) model function as a constraint for chromatograms and a huge number of PDA spectra aligned with the time axis. The i-PDeA II function can automatically extract respective peak profiles and absorption spectra from unseparated elution band. Through this investigation, we confirm the following issues that reveal accuracy, precision, and applicability of i-PDeA II.

- 1) Deconvolution for 2-component system
- 2) Deconvolution for 3-component system with greatly different abundance ratio
- 3) Investigation of the relationship between resolution and deconvolution accuracy
- 4) Practical applications of i-PDeA II

Intelligent Peak and Spectrum Deconvolution Using Photodiode Array Detector


2. MATHEMATICAL BACK GROUND

3D data composition


Ideal 3-compont system is expressed by vector product of 3 peak profiles and 3 spectra → Matrix expression

$$d_i^T = \alpha_i s_1^T + \beta_i s_2^T + \gamma_i s_3^T = (\alpha_i \quad \beta_i \quad \gamma_i) \begin{pmatrix} s_1^T \\ s_2^T \\ s_3^T \end{pmatrix}$$


Measured 3D data

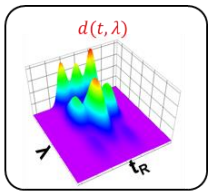


=

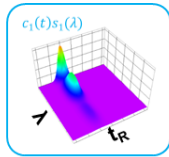

+

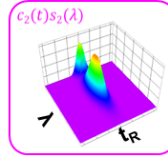


+

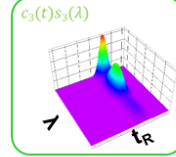

3 components system


$$D = c_1 s_1^T + c_2 s_2^T + c_3 s_3^T$$

Peak profile

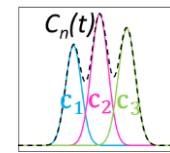

Retention time

Spectrum

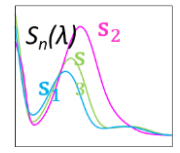

Wavelength

×



$$D = c_1 s_1^T + c_2 s_2^T + \dots + c_N s_N^T$$

N-components system

$$D = CS^T$$

Matrix expression

C: real peak profiles, **S^T**: real spectra

3D data expression by estimation

Estimation accompanies error E

$$D = C_E S_E^T + E$$

C_E: estimated peak profiles, **S_E^T**: estimated spectra, **E**: Error

$$E^2 = (D - \sum f_k S_k^T)^2$$

Minimizing square error E^2 by MCR-ALS and EM Component number is automatically optimized

S_k^T = Each spectrum

← Real spectra at the beginning and/or the ending of the elution band

f_k = $bemg(t, a, b)$

← Mathematical model for peak profile (Considering both peak tailing and leading)



$$emg(t, b) = \int_0^\infty e^{-bx} * e^{-(t-x)^2} dx$$

$$f(k) = bemg(t, a, b) = \int_{-\infty}^0 e^{ax} * emg(t-x, b) dx$$


Intelligent Peak and Spectrum Deconvolution Using Photodiode Array Detector

3. EXPERIMENTS AND RESULTS(1)

2-component system Deconvolution of Cytidine and AMP

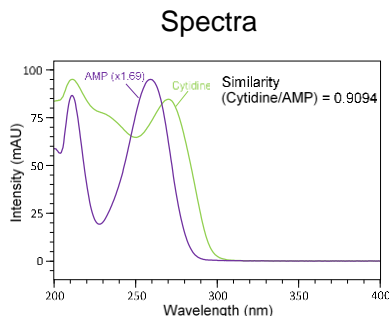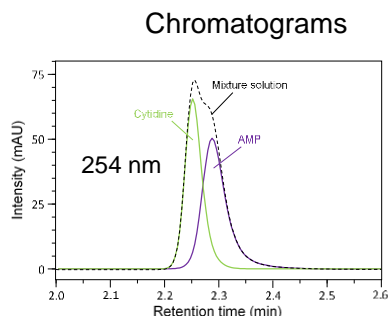


$$\text{Similarity} = \cos\theta = \frac{S_1 \cdot S_2}{|S_1||S_2|}$$

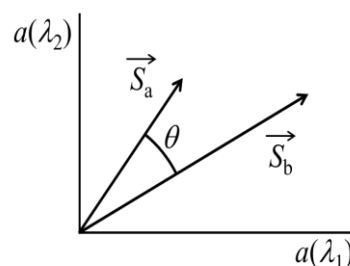

Component	Individual		Mixture (Deconvolution)		Error %	Similarity
	Area(μAU·s)	%RSD (n=6)	Area(μAU·s)	%RSD (n=6)		
Cytidine	154,514	0.14%	154,934	2.46%	0.27%	1.0000
AMP	171,643	0.22%	168,590	2.79%	-1.78%	0.9995

3-component system in different abundance ratio Deconvolution of o-, m-, p-(MAP) (Abundance ratio; 100/100/1)

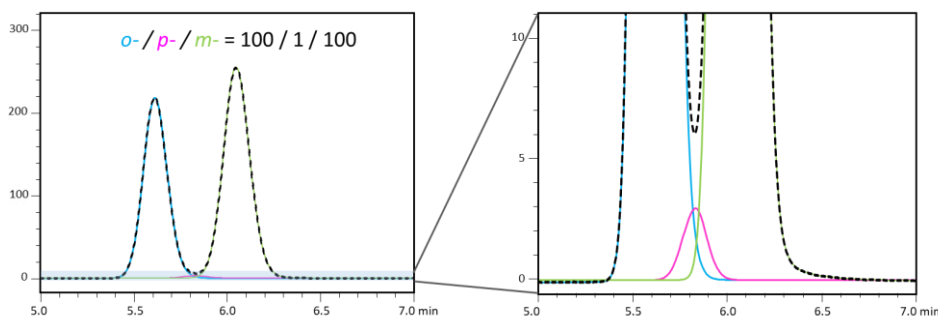


Component	Area (μAU·s)		Error %	Similarity
	Individual	Mixture		
o-MAP	2,090,806	2,080,405	-0.50%	1.0000
p-MAP	27,666	26,639	-3.71%	0.9996
m-MAP	2,658,837	2,656,836	-0.08%	1.0000

Intelligent Peak and Spectrum Deconvolution Using Photodiode Array Detector

3. EXPERIMENTS AND RESULTS(2)

Relationship between peak resolution and deconvolution accuracy Deconvolution for *o*-, *m*-, *p*-MAP

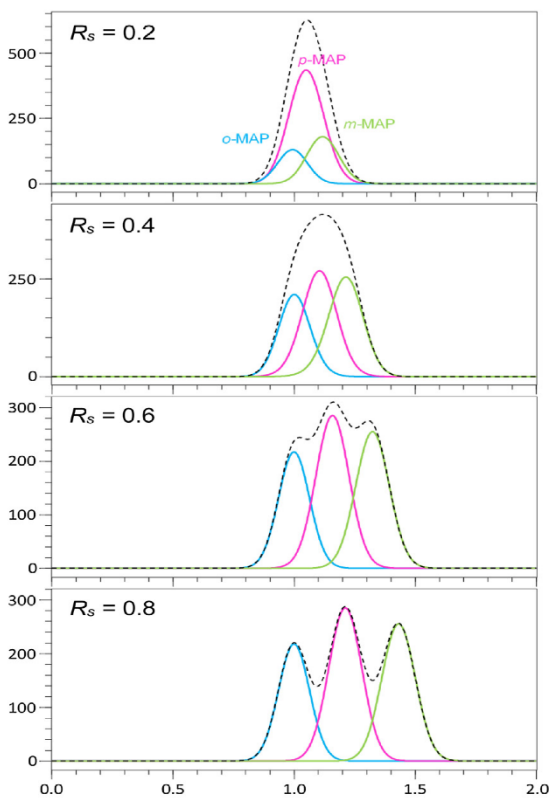

(*o*-/*m*-/*p*- : equivalent abundance ratio)



Resolution		<i>o</i> -MAP	<i>m</i> -MAP	<i>p</i> -MAP
	Individual	2.091	2.937	2.658
0.2	Area	1.204	4.698	1.784
	Error	-42.41 %	+59.98 %	-32.90 %
0.4	Area	2.046	2.900	2.740
	Error	-2.16 %	-1.23%	+3.06%
0.6	Area	2.083	2.950	2.652
	Error	-0.36 %	+0.46 %	-0.23 %
0.8	Area	2.091	2.936	2.658
	Error	0.00 %	0.00 %	0.00 %

※ Area: x 10⁶ (μV · sec)


Intelligent Peak and Spectrum Deconvolution Using Photodiode Array Detector

3. EXPERIMENTS AND RESULTS(3)

Polymer additives in SEC

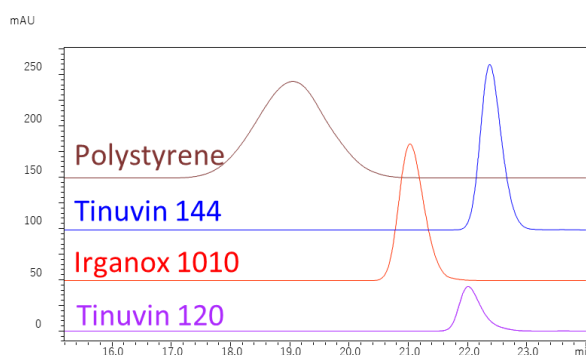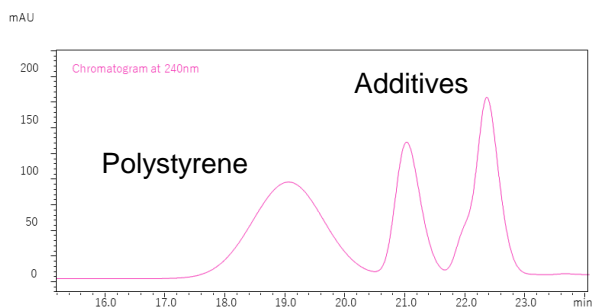

Deconvolution of three additives (PDA)


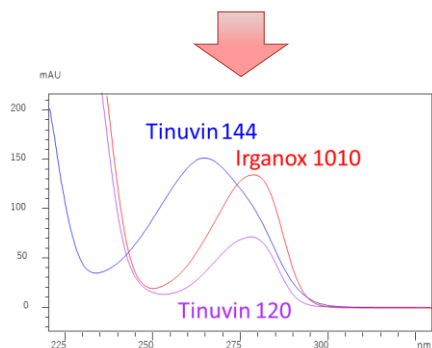


Deconvoluted Spectra

Additive	Irganox 1010	Tinuvin 144	Tinuvin 120
Linearity of calibration curve (r^2)	0.999	0.995	0.998
Determined content (mg/g)	49.2	23.1	27.4
%RSD	1.28	1.93	1.47

From 0.01 to 0.1% (w/v), n=6

Ultra high speed separation

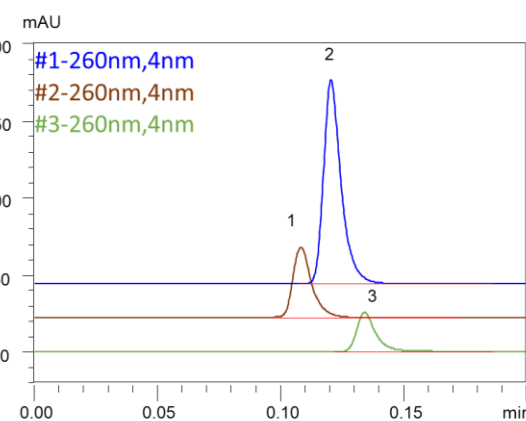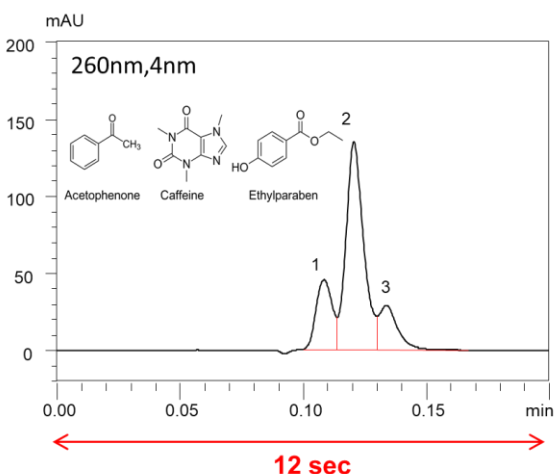

Component	t_R	Individual	Deconvolution		Vertical cut	
		Area	Area	Error	Area	Error
1.Caffeine	0.109	24680	24812	+0.5%	20486	-12.8%
2.Ethylparaben	0.121	65368	66111	+1.1%	66812	+2.9%
3.Acetophenone	0.134	13679	13912	+1.7%	15010	+13.1%

Intelligent Peak and Spectrum Deconvolution Using Photodiode Array Detector

4. CONCLUSIONS

1. i-PDeA II, an algorithm for reliable chromatographic peak deconvolution was developed by applying the MCR-ALS to PDA data.
2. Ultra fast and accurate quantitative analysis is possible at any desired wavelength even with incomplete separation.
3. i-PDeA II can be applied to analysis of isomers with identical molecular weights. It would be an alternative of MS detection for components that have identical m/z values.
4. Reliable spectral data analysis can be done even after peak deconvolution.

Reference

S. Arase et.al, *J. Chromatogr A* 1469(2016) 35

First Edition: May, 2021



For Research Use Only. Not for use in diagnostic procedure.

This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. Company names, product/service names and logos used in this publication are trademarks and trade names of Shimadzu Corporation or its affiliates, whether or not they are used with trademark symbol "TM" or "@". Third-party trademarks and trade names may be used in this publication to refer to either the entities or their products/services. Shimadzu disclaims any proprietary interest in trademarks and trade names other than its own.

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.